



# **Cedex Bio Analyzer Operator's Manual**

**Software Version 2.1** 

October 2021

For use in quality control/manufacturing process only.



# How to Use the Cedex Bio Analyzer Manual



Before reading, please review the section "Revisions" at the end of the Operator's Manual for important information.

### **Revisions**

Provides updates to the Cedex Bio Analyzer Manual, including new supplementary information and corrections to previous editions.

# **Operator's Manual**

Provides a detailed description of the Cedex Bio Analyzer, system components and all relevant software information. For installation requirements, always refer to the Operator's Manual.



# **Cedex Bio System**

*Operator's Manual Software Version 2.1* 

# **Document information**

Revision history	Manual version	Software version	Revision date	Changes
	1.0	1.0	June 2011	First publication.
	2.0	1.0	September 2011	Editorial corrections
	3.0	2.0	July 2012	Hardware improvements. Adjustments to reflect latest software version (time indication for running tests and disk space indication).
	3.1	2.1	November 2013	The <b>TRL Check</b> feature was introduced. Minor textual adjustments were made.
Edition notice	determination o optimized for sr	of chemical and bio nall throughput w	ochemical parame orkloads of appro	access analyzer intended for the ters in aqueous solutions. It is ximately 20 samples per day, ective electrodes (ISE).
	This manual is f	for users of the Ceo	<b>lex Bio</b> instrumer	nt with software version 2.1.
	manual is correc	ct at the time of pr at to make any cha	inting. However, I	ormation contained in this Roche Diagnostics GmbH thout notice as part of ongoing
	Any customer n agreement null a		instrument will r	ender the warranty or service
	Software update	es are done by serv	ice representative	S.
Intended use				access analyzer intended for the ters in aqueous solutions
	For use in qualit	ty control/manufa	cturing process or	ıly.
	It is important t system.	hat the operators r	ead this manual t	horoughly before using the
Copyrights	© 2011-2013, Ro	oche Diagnostics (	GmbH. All rights i	reserved.
Trademarks	The following tr	ademarks are acki	nowledged:	
	CEDEX, COBA	S, and ISE are trad	emarks of Roche.	
	All other traden	narks are the prope	erty of their respec	ctive owners.

# **Contact address**

Roche Diagnostics GmbH Sandhofer Strasse 116 68305 Mannheim Germany Made in Switzerland

# **Table of contents**

Document information	2
Contact address	2
Table of contents	3
Preface	5
How to use this manual	5
Online Help system	5
Conventions used in this manual	6

Part A

Part B

### **System Description**

1	Safety	
	Safety classification	A-5
	Safety information	A-5
	Data security	A-10
	Disposal recommendation	A-11
	Legal liability	A-12
	Safety labels	A-13
2	Introduction to the instrument	
	Overview	A-17
	User interface	A-20
	Wizards	A-21
	Daily operation	A-22
	Maintenance	A-33
	System status	A-34
3	Hardware	
	Covers and panels	A-39
	LEDs	A-41
	Main components	A-43
	Hardware overview	A-44
	Technical specifications	A-65
4	Software	
	Introduction	A-69
	Screen layout	A-70
	Display items	A-71
	Workflows and wizards	A-72
	Working with the user interface	A-73

working with the user interface	11 / 5
Key screens	A-82
Color interpretation for LEDs	A-117
Buttons	A-119

### Operation

5	Daily operation	
	Introduction	B-5
	Starting the shift	B-10
	Preparing the system	B-12
	Analyzing samples	B-34
	Validating sample results	B-51

	Performing calibrations	B-59
	Performing QC	B-68
	Finishing the shift	B-77
	Logging off	B-88
	Shutting down the system and switching of	f
	the instrument	B-88
	Using the barcode scanner	B-89
6	Special operations	
	Deleting sample orders	B-93
	Deleting sample results	B-94
	Calibration	B-95
	Deleting QC results	B-96
	Lot handling	B-98
	Exporting data	B-104
	Importing data	B-111
	Preparing a new disk	B-116
	Assigning tests to test tabs	B-119
	Deleting bottle sets from the <b>Inventory</b> list	B-120
	Refilling printer paper	B-122
	Removing condensation water from the	
	reagent cooler	B-124
	Replacing the probe	B-125
	Connecting and disconnecting the external	
	fluid containers	B-127
	Adjusting the touchscreen	B-131
	Cleaning the touchscreen	B-131
7	Configuration	
	Introduction	B-135
	Applications	B-137
	Configuration	B-157
Μ	aintenance	Part C
_		
8	General maintenance	<u> </u>
	Overview	C-5
	Maintenance actions	C-8
Tr	oubleshooting	Part D
9	Massages and alarms	
J	Messages and alarms	D-5
	About messages	D-5 D-5
	Message screen	
	Acoustic signals Alarm monitor	D-6
	List of alarm messages	D-6 D-10
	List of alatin messages	D-10

**Roche Diagnostics** 

10 Result flags	
About flags	D-25
Safety	D-27
List of flags	D-28
11 Troubleshooting	

D-41 D-42 D-44

D-46

roubleshooung
Introduction
Dealing with exceptional situations
Reacting to messages
Detailed procedures

ISE	Part E
12 ISE description	
Overview	E-5
Hardware	E-9
Basic operation	E-12
Technical specifications	E-13
13 ISE operation	
Daily operation	E-17
Replacing ISE fluid bottles	E-30
Replacing electrodes	E-32
Cleaning the ISE tower off the instrument	E-36
14 ISE maintenance	
Introduction	E-41
ISE maintenance actions	E-42
15 ISE troubleshooting	
Introduction	E-69
Safety	E-70
List of ISE flags	E-71
Reacting to error messages	E-80
Glossary and Index	Part F
Glossary	F-3
Index	F-11
	1 11
Revisions	Part G

# Revisions

### **18 Revisions**

## Preface

The Cedex Bio Instrument is a continuous random-access analyzer intended for the determination of chemical and biochemical parameters in aqueous solutions. It is optimized for small throughput workloads of approximately 20 samples per day, utilizing photometric analysis and a unit for ion selective electrodes (ISE).

This manual describes the **Cedex Bio** system features and general operational concepts, and it provides operating, maintenance, and emergency procedures.

<u>َ</u>ڀ

Parts B through D describe the instrument without using an ISE unit, ISE specific information is given in part E.

### How to use this manual

<u>ک</u>

- Keep this Operator's Manual in a safe place to ensure that it is not damaged and remains available for use.
- This Operator's Manual should be easily accessible at all times.

To help you find information quickly, there is a table of contents at the beginning of the manual and each chapter. In addition, a complete index can be found at the end.

### **Online Help system**

The **Cedex Bio** instrument has a context-sensitive online Help feature to aid in its operating. "Context-sensitive" means that wherever you are located within the **Cedex Bio** software, choosing Help ( ) displays Help text relating to that area of the software. The online Help offers a quick and convenient way of finding information, such as explanations of screens and dialog boxes and on how to perform particular tasks.

# **Conventions used in this manual**

Visual cues are used to help locate and interpret information in this manual quickly. This section explains the formatting conventions used in this manual.

Symbols	The following	symbols are used:
091110013	1 ne ronowing	symbols are asea.

Symbol	Used for
•	Start of procedure
•	List item
۲	Cross-reference
Ē	Call-up (software navigation path)
	Color of display item on the screen
<u>-, þ.</u>	Tip
<u>^</u>	Safety alert
X	Electrical and electronic equipment marked with this symbol are covered by the European directive WEEE.
_	The symbol denotes that the equipment must not be disposed of ir the municipal waste system.

*Screenshots* The screen representations shown in this publication are for illustrative purposes only. The screens do not necessarily show valid data.

Abbreviation	Definition
Cfas	Calibrator for automated systems
DIL	Diluent
DM	Data management
DRAM	Dynamic random access memory
e.g.	<i>Exempli gratia</i> – for example
EMC	Electromagnetic compatibility
EN	European standard
i.e.	<i>Id est</i> – that is to say
IEC	International Electrical Commission
ISE	Ion selective electrode
LED	Light-emitting diode
LIS	Laboratory information system
LLD	Liquid level detection
n/a	Not applicable
QC	Quality control
REF	Reference solution for ISE unit
ROM	Read only memory
SD	Standard deviation
SRAM	Static random access memory
TRL	Test range low (lower limit of measuring range)

Abbreviations

ons The following abbreviations are used:

Abbreviation	Description
°C	degree centigrade
μL	microliter
μm	micrometer
А	ampère
cm	centimeter
h	hour
Hz	hertz
LB	pound (weight)
in	inch
kg	kilogram
kVA	kilo volt-ampere
L	liter
m	meter
MB	megabytes
min	minute
mL	milliliter
mm	millimeter
nm	nanometer
s	second
V	volt
VA	volt-ampère
V AC	volt alternating current
V DC	volt direct current
W	watt

# A

# System Description

1	Safety	A-3
2	Introduction to the instrument	A-15
3	Hardware	A-37
4	Software	A-67

# Safety

# Protecting yourself and the environment

In this chapter, you will find information on the safe operation of the **Cedex Bio** instrument.

In this chapter	Chapter	1
Safety classification		A-!
Safety information		A-
Transport		A-
Electrical safety		A-
Optical safety		A-0
Mechanical safety		A-0
Instrument covers		A-0
Operation and maintenance		A-0
Biohazardous materials		A–0
Waste		A-'
Reagents and other working solutions		A-2
Installation		A-'
Environmental conditions		A-'
Power interruption		A-8
Electromagnetic devices		A-8
Approved parts		A-8
Third-party software		A-8
Operator qualification		A-9
Operation over an extended period of time		A-9
Cross contamination of sample		A-9
Insoluble contaminants in sample		A-9
Spillage		A-9
Data security		A-10
Disposal recommendation		A-1
Disposal label		A-1
Disposal of external components		A-1
Disposal of the instrument		
Constraint		A-12

Table of contents

Legal liability	
Safety labels	

# **Safety classification**



Caution

NOTICE

Before you attempt to use the Cedex Bio instrument, you must be fully familiar with the following symbols and their meanings:

### Warning

Indicates a hazardous situation which, if not avoided, could result in death or serious injury.

Indicates a hazardous situation which, if not avoided, may result in minor or moderate injury.

### Notice

Indicates a hazardous situation which, if not avoided, may result in property damage.

## **Safety information**

Before operating the Cedex Bio instrument, it is essential that you both read and understand the safety information listed below.

Read all Roche safety notices carefully and make sure you understand them.

### Transport



#### Injury from heavy loads

You may injure your hands, fingers, or back when putting the analyzer in place. Carry the analyzer according to the transport instructions.

### **Electrical safety**



#### Electrical shock by electronic equipment

Do not attempt to work in any electronic compartment. Installation, service, and repair must only be performed by authorized and qualified personnel.

#### **Electrical safety**

Connect the analyzer to grounded power outlets only (IEC protection class 1). All peripheral devices that are connected to the Cedex Bio instrument must comply with safety standard IEC 60950 for information technology equipment, or with IEC 61010-1, UL 61010-1 for laboratory use instruments.

### **Optical safety**



### Loss of sight

The intense light of the LEDs may severely damage you eyes. Do not stare into the LEDs. Scanning equipment using LED technology is covered by the international standard IEC 60825-1 LED Safety: Class 1.

### **Mechanical safety**



Personal injury or damage to the analyzer due to contact with instrument mechanism

Do not touch moving parts during instrument operation.

### **Instrument covers**



Personal injury or damage to the analyzer due to contact with instrument mechanism

Keep all covers closed, operate them as instructed on the screen.

### **Operation and maintenance**



Personal injury or damage to the analyzer due to contact with instrument mechanism

Do not touch any parts of the instrument other than those specified. During operation and maintenance of the instrument, proceed according to the instructions.

### **Biohazardous materials**



#### Infection by biohazardous materials

Contact with samples containing material of human origin may result in infection. All materials and mechanical components associated with samples of human origin are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

**Cedex Bio System** 

### Waste



#### Infection by waste solution

Contact with waste solution may result in infection. All materials and mechanical components associated with the waste systems are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

### **Reagents and other working solutions**



#### Injury through reagents and other working solutions

Direct contact with reagents, cleaning solutions, or other working solutions may cause personal injury.

When handling reagents, exercise the precautions required for handling laboratory reagents, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics reagents and cleaning solutions.

#### Skin inflammation caused by reagents

Direct contact with reagents may cause skin irritation, inflammation, or burns.

When handling reagents, be sure to wear protective equipment and observe the cautions given in the package insert.

### Installation



Incorrect results or damage to the analyzer due to wrong installation

Follow the specified installation instructions carefully.

### **Environmental conditions**



**Incorrect results or damage to the analyzer due to heat and humidity** Use the instrument indoor only.

 For details on the required environmental conditions, see *Environmental conditions* on page A-65. Safety information

### **Power interruption**

NOTICE

#### Data loss or damage to the system due to voltage drop

By a power failure or momentary voltage drop the operation unit or software of this system may get damaged or data loss may occur. Use only uninterruptible power supply.

### **Electromagnetic devices**



# Malfunction of instrument and incorrect results due to interfering electromagnetic fields

Devices that emit electromagnetic waves may cause the instrument to malfunction. Do not operate the following devices in the same room where the system is installed:

- Mobile phone
- Transceiver
- · Cordless phone
- Other electrical devices that generate electromagnetic waves

#### **Class B FCC rule compliance**

This equipment has been tested and found to comply with the limits for Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interferences when the equipment is operated in a residential area. However, this equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the present manual, may cause harmful interference to radio communications.

The electromagnetic environment should be evaluated prior to operation of the device.

### **Approved parts**



#### Malfunction of instrument and incorrect results due to nonapproved parts

Use of nonapproved parts or devices may result in malfunction of the instrument and may render the warranty null and void. Only use parts and devices approved by Roche Diagnostics.

### **Third-party software**



#### Malfunction of instrument and incorrect results due to third-party software

Installation of any third-party software that is not approved by Roche Diagnostics may result in incorrect behavior of the system. Do not install any nonapproved software.

### **Operator qualification**



#### Incorrect results or damage to the analyzer due to wrong operation

Operators are required to have a profound knowledge of relevant guidelines and norms as well as the information and procedures contained in the Operator's Manual.

- Do not carry out operation and maintenance unless you have been trained by Roche Diagnostics.
- Carefully follow the procedures specified in the Operator's Manual for the operation and maintenance of the system.
- Leave maintenance that is not described in the Operator's Manual to trained service representatives.
- · Follow standard laboratory practices, especially when working with biohazard material.

### Operation over an extended period of time



#### Fatigue due to long hours of operation

Looking at the monitor screen over an extended period of time may lead to fatigue of your eyes or body. Take a rest for 10 to 15 minutes every hour to relax. Avoid spending more than 6 hours per day looking at the monitor screen.

### **Cross contamination of sample**



#### Incorrect results due to carryover

Traces of analytes or reagents may be carried over one test to the next. Take adequate measures (e.g. sample aliquoting) to safeguard additional testing and to avoid potentially false results.

### Insoluble contaminants in sample



#### Incorrect results and interruption of analysis due to contaminated samples

Insoluble contaminants in samples may cause clogging or pipetting volume shortage and deterioration in measurement accuracy. When loading samples on the instrument, make sure that samples contain no insoluble contaminants such as fibrin or dust.

### Spillage

NOTICE

#### Malfunction due to spilled liquid

Any liquid spilled on the instrument may result in malfunction of the instrument. If liquid does spill on the instrument, wipe it up immediately and apply disinfectant.

# **Data security**



#### Unauthorized access and data loss due to malicious software and hacker attacks

Portable storage media can be infected with and transmit computer malware, which may be used to gain unauthorized access to data or cause unwanted changes to software.

The Cedex Bio instrument is not protected against malicious software and hacker attacks.

The customers are responsible for IT security of their IT infrastructure and for protecting it against malicious software and hacker attacks. Failure to do so may result in data loss or render the **Cedex Bio** instrument unusable.

Roche recommends the following precautions:

- Allow connection to authorized external devices only.
- Ensure that all external devices are protected by appropriate security software.
- Ensure that access to all external devices is protected by appropriate security equipment.
- Do not copy or install any software on the Cedex Bio instrument unless it is part of the system software or you are instructed to do so by a Roche service representative.
- If additional software is required, contact your Roche service representative to ensure validation of the software in question.
- Do not use the USB ports to connect other storage devices unless you are instructed to do so by official user documentation or a Roche service representative.
- Exercise utmost care when using external storage devices such as USB flash drives, CDs, or DVDs. Do not use them on public or home computers while connecting to the **Cedex Bio** instrument.
- Keep all external storage devices in a secure place and ensure that they can be accessed by authorized persons only.

• For further information, contact your Roche service representative.

# **Disposal recommendation**



Malfunction of instrument and incorrect results due to software modifications by the customer

The **Cedex Bio** instrument uses open source software. For details see document CedexBio System License Notes.

The **Cedex Bio** instrument has been designed to be operated with the unmodified software as shipped. The user assumes full responsibility for changing any part of the open source software, which excludes any liability of Roche Diagnostics Ltd.

This program is distributed without any warranty; without even the implied warranty of merchantability or fitness for a particular purpose. For details see document CedexBio System License Notes.

The portions of used open source software are part of MIKRAPs CPU-X168LCD/NET board support package and may be obtained from Mikrap AG (info@mikrap.ch).

All electrical and electronic products should be disposed of separately from the municipal waste system. Proper disposal of your old appliance prevents potential negative consequences for the environment and human health.

### **Disposal label**



Electrical and electronic equipment marked with this symbol are covered by the European directive on waste electrical and electronic equipment (WEEE) 2002/96/EC.

The symbol denotes that the equipment must not be disposed of in the municipal waste system.

### **Disposal of external components**



External components such as the scanner and the ISE power supply, which are marked with the crossed-out wheeled bin symbol, are covered by the European Directive 2002/96/EC (WEEE).

These items must be disposed of via designated collection facilities appointed by government or local authorities.

For more information about disposal of your old products, contact your city office, waste disposal service or your local service representative.

### **Disposal of the instrument**



The instrument must be treated as biologically contaminated hazardous waste. Final disposal must be organized in a way that does not endanger waste handlers. As a rule, such equipment must be sterile before it is passed on for final disposal.

For more information contact your local service representative.

Legal liability

### Constraint

It is left to the responsible laboratory organization to determine whether control unit components are contaminated or not. If contaminated, treat in the same way as the instrument.

# Legal liability

It is the users' sole responsibility to determine whether they activate the **TRL Check** feature, and if they do so to define the lower measuring range in accordance with the country specific requirements. If the users activate the **TRL Check** feature, they are also responsible for validation of the ranges they define.

# **Safety labels**

Read all safety labels on the instrument and equipment.

The following illustration shows where on the instrument labels are displayed.

A This label on the electrode block of the ISE unit indicates that there is a danger of hazardous situations arising within the vicinity of this label, which may result in death or serious injury. The relevant laboratory procedures on safe use must be observed. This label on the main cover indicates that there are potential biohazards within the vicinity of this label, which may result in death or serious injury.

The relevant laboratory procedures on safe use must be observed.



Safety labels on the Cedex Bio instrument

In addition to safety labels on the instrument, there are safety notes in the corresponding parts of the Operator's Manual.

These safety notes give more detailed information about potentially hazardous situations that may arise during daily operation or when carrying out maintenance procedures.

When working with the instrument, be sure to observe both the safety labels on the instrument and the safety notes in the Operator's Manual.

### 1 Safety

Safety labels

# Introduction to the instrument

What you need to know before you start

In this chapter, you will find basic information on the features that are relevant for working with the **Cedex Bio** instrument.

In this chapter	Chapter	2
Overview		. A–17
Principles of operation		. A–19
User interface		
Wizards		. A–21
Daily operation		. A–22
Overview		. A-22
Reagent and diluent handling		. A-24
Calibration		. A-26
Calibration type		. A-26
Calibration sequence		. A–27
Calibration status of a set		. A-28
Calibration result storage	••••••	. A–28
Validating calibration results		. A-28
Calibration procedures		. A–28
Quality control (QC)		. A-29
Sample handling		. A-30
Order handling		. A–31
Results		. A-33
Maintenance		. A-33
System status		. A–34

Table of contents

Overview

## **Overview**

The **Cedex Bio** instrument is a continuous random-access analyzer intended for metabolite analysis in mammalian cell cultures and microbial fermentation. The **Cedex Bio** instrument is a low throughput, bench top instrument used in manufacturing and quality control of pharma biotech environments. Core test parameters cover sodium, potassium, glucose, lactate, lactate dehydrogenase, ammonia, glutamine, and glutamate.

Only trained personnel working in a professional laboratory environment may operate the **Cedex Bio** instrument.



#### Incorrect results or damage to the analyzer due to wrong operation

Operators are required to have a profound knowledge of relevant guidelines and norms as well as the information and procedures contained in the Operator's Manual.

- Do not carry out operation and maintenance unless you have been trained by Roche Diagnostics.
- Carefully follow the procedures specified in the Operator's Manual for the operation and maintenance of the system.
- Leave maintenance that is not described in the Operator's Manual to trained service representatives.
- Follow standard laboratory practices, especially when working with biohazard material.

*Features* The Cedex Bio instrument offers small laboratories the following advantages:

• High analytical performance

The same bulk reagents, 12-wavelength photometer and disposable cuvettes generate results that are highly correlated to other cobas instruments.

• Efficient operation

Cooled, exchangeable reagent disks ensure economical reagent use; disposable cuvette segments allow for easy cuvette loading and removal.

High reliability, low maintenance

Innovative "low impact" instrument design and software-driven preventive maintenance improves up-time and reduces maintenance costs.

Adaptable user interface

The built-in color touchscreen, process-driven software, and reagent and sample barcode entry adapts to users of different skills and access levels.

High safety standards

Built-in safety devices, such as level detection, tube bottom detection, cuvette quality control, and ISE clot detection anticipate potential hazards during operation.

Flexible sampling


Eight on-board sample positions accommodate virtually any type of sample carrier, and enable continuous sample placing and removal during operation.

Data management

Bidirectional RS-232 and USB ports, on-board thermal printer, and drivers offer the latest in data management capabilities.

Overview

*Measuring principles* Measurements are performed by means of an absorbance photometer and an ISE (ion selective electrode) module that uses ion selective potentiometry.

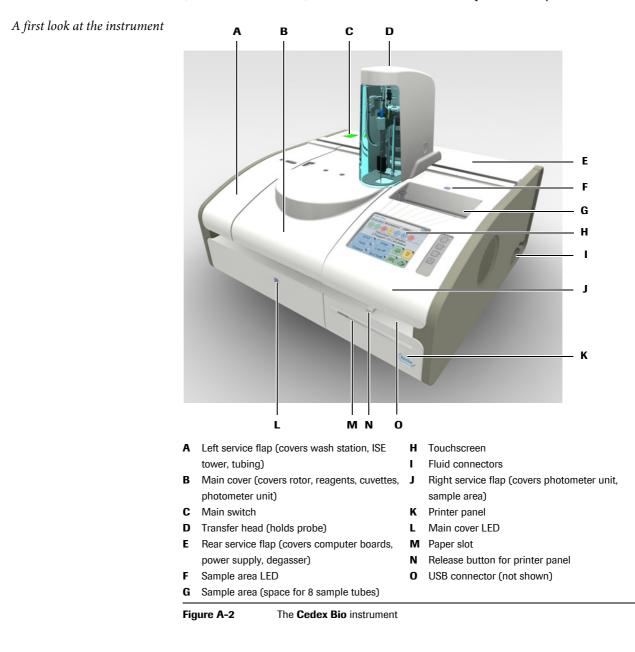

Overview

# **Principles of operation**

	The <b>Cedex Bio</b> main instrument uses absorption photometry for determining the amount of absorbance in a fluid. The absorbance is used to calculate the concentration in the solution.
Loading the sample	The operator identifies the sample, places it on the instrument, and defines the order. (If you work with a host system, the order is defined automatically.)
Measuring process	The measuring process for each test consists of forty regular cycles, each lasting 18 seconds. In each of these cycles, a measurement is taken, irrespective of what other actions take place during this cycle. The application definitions determine what is done in which cycle, and they also define which results are taken into account for the result calculation.
	With each cycle, a new test can be started.
	The basic process works as follows:
	1. Checking the cuvette.
	A measurement is taken to check the quality of the cuvette.
	2. Pipetting reagent (R1) to the cuvette.
	After each pipetting action, the system performs a wash cycle to minimize carry- over. During this cycle, the probe and tubing are flushed with water and cleaner.
	3. Wait.
	The fluid needs to reach the prescribed temperature. Such a phase can last several cycles.
	During the wait cycles, activities for other tests are performed.
	4. Pipetting the next fluid.
	Typically, this would be the sample. The details are defined in the application definitions.
	5. Wait.
	6. Pipetting the next fluid.
	7. Wait.
	8. And so on.
Calculating the results	The test result is calculated on the basis of the photometric measurement results. During this process, various checks are performed to ensure that the whole measuring process was technically correct. If values are above or below predefined limits, the test result is flagged.
	The results are stored on the system. This includes both the forty measurement results (raw data) and the calculated test result.
Sequence of processing	For a given sample, the tests are processed in the order defined by the time required for their processing (number of cycles), starting with the one that takes the longest. This order can be altered manually by defining a specific process sequence list.
Status of the measuring process	At any stage of the measuring process, the user can check its status on the screen.
Result data management	The system provides storage space for the results of one working day. For backup purposes, the results must be exported to an external storage device once a day.
<b>B</b>   <b>B</b> ' '	

### **Roche Diagnostics**

# **User interface**

The **Cedex Bio** instrument is equipped with a touchscreen, an on-screen keyboard and four global action buttons. LEDs and acoustic signals let you know when it is safe to add or remove samples, reagents and other fluids.

With buttons and other display items, "traffic light" color coding is used: Green means OK, yellow: watch out, you need to do something, and red means that your intervention is required for processing to continue.

The screens have a clear and consistent layout and are easy to use. The topics are divided in the proven work areas: **Overview** for order and fluid handling, **Workplace** for result handling and details on orders, and **Utilities** for administration tasks.

• For details on the user interface, see Chapter 4 Software.

The following is an example of a screen. It contains the full range of display items.

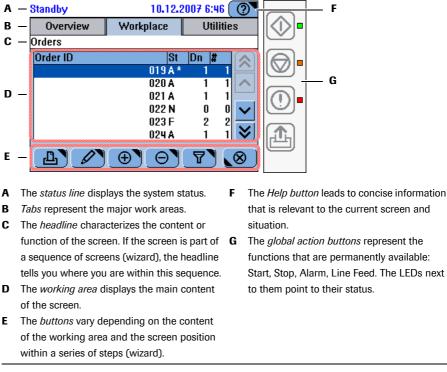


Figure A-3 Example of a screen

### Wizards

# Wizards

Screens help you perform your tasks. If not all steps of a task can be performed from one screen, the workflow is realized as a sequence of screens, a so-called wizard. **Cedex Bio** wizards do not usually *force* you to perform a task at a certain stage, they just make your work easier.

• For details on workflows, see *Workflows and wizards* on page A-72.

When intervention is required

On the screen, there are several methods of telling when your intervention is required:

• Buttons and texts are color coded.

Everything is fine.
To ensure smooth operation, you need to perform some task.
The current process or action has not started yet or stopped. You need to do something for it to start or continue.

- *Screens* can contain instructions. For example the text may ask you to place the sample on the sample area or to remove a reagent bottle from the reagent disk.
- *Messages* inform you about the status of current actions.
- A permanent *alarm monitor* alerts you to events you should know about.
- WizardsThere are three major wizards: Prepare wizard, Orders wizard, and End Shift wizard.With most tasks that involve more than one step, such as exchanging reagent or otherfluid bottles, you are supported by wizards.
- *Prepare wizard* The Prepare wizard guides you through the tasks that need to be performed at the beginning of a shift. When this wizard is done, the system is ready for processing orders.
- Orders wizard The Orders wizard guides you through the process of creating and changing orders.
- *End Shift wizard* The End Shift wizard guides you through the tasks that need to be performed at the end of the day or to prepare the instrument for handing over to another operator.

<u>ک</u>

Individual tasks can be performed outside the wizards

Most tasks that make up a workflow can be performed without using a wizard.

If you perform a task independently, you first need to navigate to the appropriate screen and then start the task from there; whereas if you use a wizard, the appropriate screen is displayed automatically.

Using the wizards also ensures that all necessary steps are performed and in the right order.

# **Daily operation**

### **Overview**

Daily operation includes the routine tasks that are required to prepare and monitor the system, and to analyze samples.

When you switch on the system, it performs several checks to make sure that all preconditions are met, for example that all covers are closed or that there are cuvettes available. It then performs self-tests to ensure that all modules function properly.

At the end of the startup phase, the screen is updated to display the current status of the system.

The following table gives an overview of the tasks you might need to perform during daily operation.

	Task	Steps		Navigation
			With wizard	As individual steps
1	Starting the system	1. Switch on the system.		
2	Logging on the system			Overview > Logon
3	Preparing the system	Start the <b>Prepare</b> wizard.	Overview >	Prepare
		1. Check the external fluid containers.		Overview > 🚵 > 🛍
		2. Perform the maintenance actions that are due.		Utilities > Maintenance
		3. Load the reagent disk.		Overview > 🕮
		4. Check the reagents.		Overview >
		5. Check the cuvettes.		Overview > 😈
		6. Perform mixing		Overview > 🔟 > test > 🖸
		7. Perform calibrations that are due.		Workplace > Calibrations > 🕀 > 💿
4	Defining orders	Start the <b>Orders</b> wizard.	Overview >	Order (or Overview > STAT)
		1. Identify the sample.		n/a
		2. Select the tests.		n/a
		3. Place the sample.		n/a
		4. Start the run.		$\bigcirc$
5	Monitoring the progress		n/a	Overview ①
		1. View results.	n/a	
6	Validating results			Workplace > Result Review
		2. Handle flagged results.	n/a	Workplace > Result Review > 🖉
				> Repeat
				> <b>Rerun</b>
		3. Accept results.	n/a	Workplace > Result Review > 🖉 >
				Accept

 Table A-1
 Overview of the daily operation tasks

Daily operation

Task	Steps		Navigation	
			With wizard	As individual steps
Performing calibrations				
Performing individual	1.	Start the wizard.	Workplace >	Calibrations > 🕀
calibrations	2.	Select the test.	_	n/a
	3.	Prepare and place the calibrators.		n/a
	4.	Start the calibration.		$\Diamond$
	5.	Validate the calibration results.		Workplace > Calibrations >
	6.	Remove the calibrators.		
Performing	1.	Start the wizard.	Workplace >	Calibrations > 🕀
all due calibrations	2.	Select all tests with due calibrations. or Select all tests with calibration due within the forecast period.	(ئ م	n/a
	3.	Prepare and place the calibrators.	_	n/a
	4.	Start the calibration.		$\diamond$
	5.	Validate the calibration results.		Workplace > Calibrations >
	6.	Remove the calibrators.		
Performing controls				
Performing	1.	Start the wizard.	Overview > 0	Order > 📲
Default QC	2.	Select a control and place the tube. Repeat until there are no controls left on the screen.	_	n/a
	3.	Start the QC measurement.		$\bigcirc$
	4.	Validate the QC results.		Workplace > QC Status >
	5.	Remove the controls.		n/a
Performing an individual	1.	Start the wizard.	Workplace >	• QC Status > 🛞
QC measurement	2.	Select a test.	_	n/a
	3.	Select a control and place the tube. Repeat until there are no controls left on the screen.		n/a
	4.	Start the QC measurement.		$\diamond$
	5.	Validate the QC results.		Workplace > QC Status >
	6.	Remove the control.		
Performing all	1.	Start the wizard.	Overview > 0	Order > 🔍 > 🕙
due QC measurements	2.	Select a control and place the tube. Repeat until there are no controls left on the screen.	-	n/a
	3.	Start the QC measurement.		$\bigcirc$
	4.	Validate the QC results.		Workplace > QC Status >
	5.	Remove the controls.		

Table A-1

Overview of the daily operation tasks (Continued)

### **2** Introduction to the instrument


Daily operation

	Task	Steps		Navigation
			With wizard	As individual steps
	Finishing the shift	1. Check for unfinished orders.		Workplace > Orders
				Choose 🍸 > Not Finished
		2. Check for non-validated results.		Workplace > Result Review
				Choose 🍸 > Not Accepted
		3. Check for non-transmitted results.		Workplace > Result Review
		(If working with a host system only.)		Choose 🍸 > Not Sent to Host
		4. Start the <b>End Shift</b> wizard.	Overview >	End Shift
		5. Perform the daily backup.		Utilities > Export > Database
		6. Export the full results.		Utilities > Export > Results
		7. Clean up the database.		Workplace > Orders > Θ
				Workplace > Result Review> Θ
				Workplace > QC Status > Θ
				Workplace > QC History > Θ
				Workplace > Calibrations > $\Theta$
		8. Perform the maintenance actions that are due.	_	Utilities > Maintenance
		9. Replace cuvettes.	_	Overview > 😈
		10. Check the external fluid containers.	_	Overview > 🔄 > 🛍
		11. Remove the reagent disk (if last shift).		Overview > 🔟 > 🖭
		12. Log off the system.		<b>Overview</b> > button with your user name
		13. Shut down the system and switch off the instrument (if last shift).	n/a	n/a

Table A-1

Overview of the daily operation tasks (Continued)

### **Reagent and diluent handling**

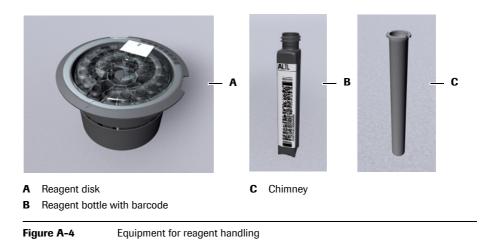

*Reagent disk* On the instrument, the reagents are stored on a reagent disk. It provides space for 27 bottles, allowing up to 14 reagent sets to be installed on the disk, assuming that most tests need two reagents. Extra diluents and cleaners are also loaded on the reagent disk.

#### **Roche Diagnostics**

You can work with up to eight different reagent disks on one Cedex Bio instrument.

You always load and remove bottles while the disk is on the instrument. (The system needs to know exactly what is loaded on the disk.)

When you finished running tests, you can remove the whole reagent disk, place it in a reagent disk container, and store it in a refrigerator.

- *Bottles* Cedex Bio reagents, diluents and extra cleaners are provided in uniform bottles. They are supplied with two dimensional barcodes and placed on the reagent disk with their cap removed.
- *Chimneys* Chimneys are bottle inserts that reduce evaporation. For reagents that are especially sensitive to concentration changes, Roche recommend using chimneys on the reagent bottles.

To generally reduce evaporation, you may use chimneys on all reagent bottles.

*Reagent set* Up to three reagents can be required to perform a certain test. These reagents are handled in reagent sets. You can define more than one reagent set for a test, but only one can be active.

A reagent set is defined as soon as its first bottle is loaded. From this moment on, whenever you remove or replace a reagent, you do so for all reagents of the set.

Each diluent or cleaner bottle is treated as a separate reagent set.

*Volume detection* For each reagent set, the number of available tests is continuously calculated.

Standby	10.12.2007 13:15 🕐
Test Status>ALTPL	
Calibration	0K
Quality control	ОК
Tests on disk 2	41
Tests ready to run	41
Test spec	cification is OK

Figure A-5 Remaining tests indication

*Periodic mixing* Reagents may have a mixing interval defined. This interval is checked by the system every 30 minutes, and mixing is performed without removing the reagent bottles from the reagent disk.

For a reagent set that contains more than one reagent for which mixing is defined, the shortest interval of all reagents of the set is used for all reagents.

Tests are blocked if any reagent they use requires mixing.

*Diluents* Both, system water and dedicated diluents are used. System water is kept in the external water container, dedicated diluents are supplied in reagent bottles and placed on the reagent disk.

*Cleaners* Both, a system cleaner and dedicated cleaners can be used. The system cleaner is kept in the external cleaner bottle, dedicated cleaners are supplied in reagent bottles and placed on the reagent disk.

**Roche Diagnostics** 

Daily operation

### Calibration

	Calibration is the process that establishes, under specified conditions, the relationship between values indicated by the analytical instrument and the corresponding known values of an analyte.
	Periodic calibration is required because the concentration of reagents can change over time.
	Reagents are typically calibrated with a two-point calibration, measuring the predefined value of a multicalibrator and of system water. Some reagents are calibrated using a set of calibrators.
	On the <b>Cedex Bio</b> instrument, reagents are handled as sets of up to three reagents. (You always load and unload all reagents of a set.) As a consequence, all reagents (bottles) of a set are calibrated when performing calibration.
	The system checks when calibration is due.
-` <b>\</b>	Each reagent set must have accepted calibration results to be available for use in tests.
	When a calibration is due depends mainly on two definitions, the calibration type and the calibration sequence.
Calibration type	
	The calibration types <i>Set</i> and <i>Lot</i> define the manner in which the system determines whether there is a valid calibration result for a particular reagent set.
Set calibration	Set calibration results are valid for the calibrated set only. They can be generated from any reagent set.
Lot calibration	Lot calibration results are valid for the reagent set they were calibrated with and for all subsequent reagent sets of the same lot. Usually, lot calibrations are generated by calibrating the first reagent set of a new lot. There can only be one accepted lot calibration result for the reagents of a given lot.
	Let us suppose that you place the first reagent set of a new lot and calibrate it straight away. Let us further assume that subsequent control measurements suggest that a new calibration is required. Within the first 24 hours of placing a set on the system, you can recalibrate it, and possibly existing lot calibration results of this set are superseded. When this period has elapsed you can no longer change the lot calibration results. (To generate new lot calibration results, you would have to delete the existing results and then calibrate a new reagent set.)
-`\$ <mark>'</mark> -	Lot calibration is relevant if you work with the calibration sequence [Each Lot and Interval].
	The following table illustrates the two calibration types in an example.
	Assumptions:
	<ul><li>Sequence: Each Lot and Interval.</li><li>Interval: 5 days.</li></ul>
	(Note that the interval (re)starts when a set is calibrated as a result of the interval expiring or a new lot being started.)

Daily operation

Day	Trigger/Event	Task	Result used	Set	Cal. type	Cal. Usage
1	NA	1. Place first reagent set of new lot.	Result 1	L1/1	Lot	Current
		2. Calibrate set L1/1.				
2	Reagent empty.	1. Remove set L1/1.	Result 1	L1/2	Set	Current
		2. Place new set L1/2.		L1/1	Lot	-
	Reagent empty.	Replace set whenever it is empty.	Result 1		Set	Current
5	Interval expired.	Calibrate current set L1/n.	Result 2	L1/n	Set	Current
				L1/n-1	Set	Obsolete
6	Reagent empty.	Place new set L1/n+1.	Result 1	L1/n+1	Set	Current
				L1/n	Set	Obsolete
8	Reagent empty.	1. Remove set L1/n+1.	Result 3	L2/1	Lot	Current
	New lot.	2. Place new set, which is the first set of a new lot		L1/1	Lot	Obsolete
		3. Calibrate set L2/1.		L1/n+1	Set	Obsolete
	Reagent empty.	Place new set L2/2.	Result 3	L2/2	Set	Current
				L2/1	Lot	-
10	Interval expired.	Calibrate current set L2/n.	Result 4	L2/n	Set	Current
				L2/n-1	Set	Obsolete
11	Reagent empty.	Place new set L2/n+1.	Result 3	L2/n+1	Set	Current
				L2/n	Set	Obsolete

Table A-2

Example for set change and calibration types

#### **Calibration sequence**

ition sequence	
	The calibration sequence is an application definition. It defines the manner in which the system determines when a calibration is due.
-`Q́-	Roche recommend not to change the calibration sequence.
	The calibration interval defines the on-board stability of a reagent.
	One of the following sequences applies to each reagent set:
No Interval	You perform calibration whenever you think fit. Use this value if you are sure that the reagent is stable until it is empty and you replace it with a new one. Calibration is due whenever a new reagent set is loaded on the instrument.
Interval Only	You perform calibration only when the interval has expired.
Each Lot and Interval	You perform calibration whenever the fist reagent of a new lot is loaded and then each time the interval has expired.
	In this case, the interval is related to the date when the lot calibration was generated, and it (re)starts whenever you calibrate a reagent set (as a result of interval expiry or starting a new lot).
	You can turn off the interval check by defining its duration as 0 (zero).
Each Set and Interval	You perform calibration whenever a new reagent is loaded and when the interval has expired.
	The interval starts again whenever you calibrate a reagent set because the interval had expired or a new lot was started.
	You can turn off the interval check by defining its duration as 0 (zero).

Daily operation

Calibration status of a set	
	Each reagent set has one of the following calibration statuses:
CU	(current) denotes that the set is on board and that its calibration results are currently used.
ОВ	(obsolete) denotes that the set's calibration results are no longer used.
	This status applies for example to the following situations:
	• The set was removed and it is empty.
	• The set was removed and it is not empty. It was removed more than 30 days ago.
SB	(standby) indicates that the set's calibration results are not currently used.
	This status applies for example to the following situations:
	• A new set was loaded and calibrated while an identical set was still in use (pre- calibration).
	• The set was removed not more than 30 days ago and it is not empty.
Calibration result storage	
	The current and up to five obsolete calibration results are stored on the system. If there are more than five obsolete calibration results, the oldest obsolete calibration results are automatically deleted as part of the daily end of shift activities.
Validating calibration results	
	Applications define checks for ranges and limits. If these are exceeded, the results are flagged.
	Each new calibration result has to be validated. If flags were generated, you must determine their cause and decide whether to accept the result, rerun the calibration or continue using the old calibration results.
	You can automatically accept unflagged results and results with flags that are contained in a specific list of flags that should be ignored.
	• See <i>Editing the acceptable flags list</i> on page B-160.
Calibration procedures	
	There are three basic procedures for performing calibration:
۲	Calibrating all reagent sets that need calibrating
*	• Calibrating all reagent sets that will need calibrating during the forecast period
<xyz></xyz>	Calibrating individual reagent sets

. .

. . 11 1

....

Forecast period	The forecast period is a configurable period of time. Calibrations that fall due within this period will be performed collectively.
	● See <i>Calibration</i> on page B-165.
	Typically you would set this period to fit your shift length, for example 8 hours. This would enable you to prepare the instrument before the work shift starts and so avoid having to interrupt sample processing for performing calibrations.
Precalibration	At any given time, there is only one accepted calibration result for each test. You can, however, install and precalibrate one reserve reagent set. This is done, for example, to ensure continuous sample processing.

### **Quality control (QC)**

QC is performed at regular intervals to check the integrity of the whole measuring system. For each test, up to three controls are defined. The results are compared against predefined ranges or values and then interpreted accordingly.

c . .

*Control* A control is a sample that has been measured using all tests it is associated with, in order to define the ranges and values that determine the correct functioning of the instrument. This is typically done both for the normal and the pathological analyte concentration.

*When QC is due* With regards to when it needs to be performed, QC is divided in the following types:

#### • QC After Cal

The QC measurement is due after calibration of the test.

Interval QC

QC is due whenever its interval has expired. QC measurements of this type are performed in a batch, typically once or twice a day.

• Default QC

QC is performed at certain times during routine operation. This is done to fit in with laboratory processes and procedures.

*Ways of performing QC* With regards to how QC is performed, the following methods are provided:

#### • Default QC

**Default QC** is an automated process for performing multiple QC measurements at the time when you define the QC orders. This is the ideal method if you want to perform QC at certain times and days.

This method only applies to tests whose controls are defined to be performed as part of **Default QC**. Therefore, if you intend to work with the **Default QC** function, you need to configure the tests accordingly.

**Default QC** follows a streamlined procedure where the necessary QC orders are automatically defined as soon as you identify a control. An order is defined for each test for which this control is defined, provided the test is currently active on the system. A wizard helps you select the controls, and a placement list supports you in preparing and loading the controls.

Daily operation

• Interval QC

This method applies to tests whose controls have an interval defined.

Interval QC is a process that is suitable both for performing a single QC measurement and for performing all QC measurements that are due. You can select all tests that require QC simply by pressing a button. (This selection also reflects QC of the type **QC After Cal**.) A wizard helps you select the controls, and a placement list supports you in preparing and loading them.

*Validating QC results* Each new QC result has to be validated. If flags were generated, you must determine their cause and decide whether to accept or ignore the result.


You can automatically accept unflagged results and results with flags that are contained in a specific list of flags that should be ignored.

• See *Editing the acceptable flags list* on page B-160.

If you ignore a QC result you exclude the result from further QC result calculations such as QC History statistics.

### Sample handling

You can place up to eight sample tubes on the sample area.

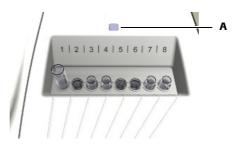

A Sample area LED. A green LED indicates that you should place a tube, a blinking yellow LED that you should keep clear of the sample area.

Figure A-6 Sample area with sample tubes

Sample types

The **Cedex Bio** instrument can process aqueous samples from mammalian and bacterial cell cultures and fermentation.



#### Incorrect results due to inadequate sample preparation

Specimen containing clots may obstruct the probe. Specimen containing bubbles or foam may cause level detection errors and air pipetting. Consequently, incorrect results may be generated.

Take adequate care when preparing the samples.

#### Incorrect results due to insufficient fluid

Insufficient fluid may lead to inaccurate pipetting and consequently to incorrect results.

Always fill the tubes with enough fluid that at least the defined dead volume of fluid is left when pipetting is complete.

See *Tubes* on page A-53.

Sample tubes	The <b>Cedex Bio</b> instrument can use both primary and secondary tubes (cups).
	You can use any type of primary tube, as long as their dimensions lay within prescribed limits. Roche recommend using approved cups only.
	• See <i>Tubes</i> on page A-53.
NOTICE	Probe damage due to not removing primary tube caps
	The probe is not designed to pierce tube caps. It can get damaged when trying to pierce tube caps.
	Always remove the caps of primary tubes before placing them on the instrument.
Sample ID	The sample ID is an identifier of up to 23 alphanumeric characters that is unique within a whole organization. It identifies the sample and is also used for host communication.
	Sample IDs are defined either by scanning a barcode or by typing them manually.
	Because there is limited space when displaying lists on screens, Roche recommend to limit the ID to 13 characters.
Sample barcode	You can use sample tubes with or without barcode.
Dilution	Pre-dilution is used when performing calibration.
	Post-dilution is used when measuring samples.
	(The dilution factor is part of the application definition and therefore does not need to be defined by the operator.)
Removing sample tubes	You can remove sample tubes as soon as pipetting is complete.

## Order handling

Order Mode	The order mode reflects the way in which you organize the tests on the test selection screen.		
	Choose <b>Easy</b> if the reagents fit on one or two reagent disks and you work with one test panel on the screen (You can fit up to 25 tests and profiles on this panel).		
Choose <b>Full</b> if you distribute the reagents across several (up to eight) reagen and if you predominantly work with specific groups of tests. You can assign tests and profiles to each panel (tab).			
Order ID	The order ID is an identifier of up to 23 alphanumeric characters that is unique within the laboratory. The order ID identifies the order and links it to the sample.		
	Order and sample IDs are often identical. Using separate IDs makes sense when working with a host system.		
	Order IDs are defined either by scanning a barcode or by typing them manually.		
	Because there is limited space when displaying lists on screens, Roche recommend to limit the ID to 13 characters.		

Host connectivity	The <b>Cedex Bio</b> instrument can be connected to an external laboratory information system (LIS) and a host computer for downloading order information and uploading results.
	If the instrument is connected to a host system, the following setups can be configured:
	Downloading order information
	When you identify a sample using the barcode scanner, the appropriate order information is automatically assigned to the order on the system. (The order information was downloaded previously.)
	Performing host queries
	When you identify a sample using the barcode scanner, a query is sent to the host, asking for the order information of the sample in question. This information is then downloaded to the <b>Cedex Bio</b> instrument and automatically assigned to the sample on the system.
	Transmitting results
	You can have results automatically transmitted to the host as soon as they are accepted.
Routine orders	Routine orders are normally defined on the <b>Overview</b> tab. The software guides you through the process of assigning the tests to the sample and placing the sample tube on the instrument.
STAT orders	STAT (short turn around time) orders are handled in the same way as routine orders, except that their tests are processed next, irrespective of the scheduling of routine order tests.
Defining orders	There must be at least one free sample tube position when defining an order. You are guided by the software when ordering the tests and placing the samples.
	There can be only one order for each test and sample.
Modifying orders	The process for changing an order is similar to that of defining it. You first identify the sample and then change the tests. You can change an order as long as its processing has not yet started.
	It is always possible to add a further test to an existing order.
Deleting orders	At the end of a shift, you should delete all orders that are defined on the system. This is to free storage space for the next shift. Deleting the orders is an integral part of the End Shift wizard. (Deleting an order also deletes the corresponding sample results.)
	You can export the data to a USB stick and store them on a computer.
Controlling the run	Controlling the execution of test runs is done via the global action buttons.
6	Press 🐼 to start the run.
	Press $\overline{\mathbf{O}}$ to stop the run.

## Results

	You can check the results on the screen as soon as they are calculated.
Units	Results are normally given in your lab units. The units can be configured.
Flags	<i>Result flags</i> are test-specific. They indicate that the limit of an internal check was exceeded or not reached.
	<i>System flags</i> point to the status of the result within the process of analysis; for example, they tell you that the result has not been accepted or that is has not been transmitted to the host successfully.
Printing results	You can print all or selected results on the built-in printer.
Validating results	All results need to be validated (result accepted, test rerun or repeated).
	Result flags help you identify critical results and point to possible actions that need to be taken.
	Each test must have accepted calibration results; tests whose associated calibration results are not accepted cannot be performed.
Ratio results	The user can manually define ratios. Ratio results are handled like any other sample result, with the exception that they cannot be accepted by the user. They are automatically accepted if all their constituent results are accepted.
Repeating and rerunning tests	If a result is flagged, you may decide to run a test again. You can either perform exactly the same test (Repeat) or perform it with a different predefined dilution (Rerun).
Storing results	The <b>Cedex Bio</b> instrument is designed to hold the sample results of one day's analyzing. Therefore, you need to back up the data regularly to an external medium. (Backing up results is an integral part of the <b>End Shift</b> wizard.)
	The QC results of the previous and the current months are stored on the system.
	Up to five calibration results are stored on the system for each test.

## Maintenance

	Completing maintenance actions correctly and on time helps to ensure smooth and uninterrupted operation of your instrument.
Maintenance scheduling	The <b>Cedex Bio</b> instrument facilitates performing the maintenance actions in bundles at the times that suit your laboratory work processes. To that purpose, you can define in the configuration settings one day of the week as your maintenance day.
	For information on scheduling maintenance actions, see <i>Scheduling maintenance actions</i> on page B-158.
	All maintenance actions can be performed any time.

System status

Interval	For most maintenance actions a fixed maintenance interval is defined. (You cannot change this interval.) This is the basis upon which the system calculates the date when the actions need to be performed.
	The interval timers and counters are reset whenever you confirm that the maintenance action has been performed.
	Maintenance actions without predefined intervals are performed whenever necessary, or they are triggered by another maintenance action.
Due date	The due date is the last possible maintenance day. This is the date you see when you check the status of maintenance actions.
Ensuring smooth operation	Performing all due maintenance actions during the daily <b>Prepare</b> or <b>End Shift</b> phase ensures that routine operation does not have to be interrupted for performing maintenance actions.

## System status

The **Cedex Bio** instrument provides several means of indicating the status of the various parts and processes:

- Color coded LEDs on the instrument inform you when and when not to open covers or place sample tubes.
  - See *Color interpretation for LEDs* on page A-117.
- The colors of buttons inform you whether you need to intervene.
  See *Color concept* on page A-72.

You can check the meaning of a button using the online Help **O**.

- Buttons on the **Overview** tab lead to detailed information on the status of selected processes and hardware items.
- Messages on the screen provide information on individual tasks and events.
- The text in the Status line provides information on the status and activities of the analyzer and photometer unit.

System status

Status	Comment	
Standby	The user and host interfaces remain active, as do the reagent cooling system, fluid system, and the cuvette heating.	
Maintenance	A maintenance action is being performed. The system is not available for performing tests.	
Diagnostics	A system diagnostics action is being performed. The system is not available for performing tests.	
Operating	Processing is in progress.	
Power Up	After switching on, the system performs initialization and functional test.	
Power Down	Regular shutdown is in progress.	
Stopped	Processing has stopped. User intervention is required to allow the system to resume regular operation.	
E-Stopped	The system has performed an emergency stop (E. Stop). This could be due to hardware failure or because any of the safety devices have requested an emergency stop.	
Table A-3	System statuses	

The following system statuses are defined:

System status

# Hardware

# The parts and how they work

In this chapter, you will find information on the main hardware components of the **Cedex Bio** instrument.

In this chapter	Chapter	3
Covers and panels		. A–39
LEDs		. A–41
Main components		. A-43
Hardware overview		. A–44
Sample area		. A–44
Fluid system		. A-45
Probe and syringe		. A–46
Syringe assembly		. A–46
Wash station		. A–47
External fluid connectors		. A–47
Transfer unit		. A–48
Fluid containers		. A–50
Waste containers		. A–50
Water container		. A–51
Cleaner bottle		. A–52
Reagent bottles		. A–53
Tubes		. A–53
Cuvettes		. A–54
Cuvette segments		. A–55
Handling cuvettes		. A–55
Reagent handling		. A–56
Rotor		. A–56
Reagent disk		. A–57
Reagent disk container		. A–58
Reagent cooler		. A–58
Cuvette ring		. A–59
Barcode scanner		. A–59
Printer		. A–61

Table of contents

Absorbance photometer	
Connectors	
Fuses	
Technical specifications	

## **Covers and panels**

	AB	C D		E F G
	A Left side panel		E	Rear service flap
	B Left service flap	)	F	Right side panel
	C Main cover		G	Right service flap
	D Transfer head o	cover	Η	Printer panel
	Figure A-7	The Cedex Bio instrument		
Transfer head cover	Provides access			
	Transfer mee	chanism		
	• Probe			
	Shut down the s	ystem and switch off the	inst	rument before you open this cover.

The following figure shows the removable panels and the lids that can be opened.

Left service flap Provides access to:

- Wash station •
- Tubing •
- Internal waste tank •
- Initialization plate •

Handle as instructed during maintenance actions, or shut down the system and switch off the instrument before handling it.

Covers and panels

Main cover Provides access to:

- Rotor
- Reagent disk
- Cuvettes
- Reagent bottles
- Photometer unit

You open this cover whenever you need to handle cuvettes, reagent bottles, or the reagent disk. A green LED indicates that you should place an item, a yellow LED that you should not open the main cover.

To open the cover, press the release button at the underside of the front of the cover.

When you should close the cover an acoustic signal is sounded and the system icon on the System Status screen turns red.

- *Rear service flap* Provides access to:
  - Computer boards
  - Power supply
  - Transfer unit
  - Degasser

This flap should be opened by service representatives only.

Shut down the system and switch off the instrument before you open this cover and remove the two side panels before opening this flap.

- *Right side panel* Provides access to:
  - Syringe assembly

Remove this panel as instructed during maintenance actions or shut down the system and switch off the instrument before you remove it.

- *Right service flap* Provides access to:
  - Photometer unit
  - Sample area
  - Touchscreen
  - Data management computer

Open this panel as instructed during maintenance actions or shut down the system and switch off the instrument before you open it.

- *Printer panel* Provides access to:
  - Printer paper.

You can open this flap any time as long as the printer is not printing.

To open the panel, press the release button.

## LEDs


	A Sample area LEI		i f Main cover LED
	Stop LED, Alarm		
	Figure A-8	The <b>Cedex Bio</b> inst	rument
Interpreting the LED colors	LED	Color	Meaning
	Main cover LED	Off	No activities in this area. You can open the main cover.
			User intervention is required, for example you are expected to place or remove a bottle.
			The system is performing some action. Do not handle the cover.
		in <b>Operating</b> stat	ll is sounded when the cover is open while the system is te. You can adjust the volume ( <b>Utilities</b> > <b>System</b> > <b>Volume</b> ).
	Sample area LED		No activities in this area. You can remove sample tubes.
		-	You are expected to place a sample tube.
		Blinking	The transfer head is approaching. Do not place
		- 22	your hand or any object in the sample area.
	Start LED	Off Off	You cannot start the measuring process.
			You can start the measuring process.
	Table A-4	LEDs and their mea	ning

LEDs are placed on key positions on the **Cedex Bio** instrument. They inform you whether you can perform certain actions.

LEDs

LED	Color	Meaning
Stop LED	Off	Pressing $\textcircled{O}$ does not have any effect.
		Pressing $\textcircled{O}$ leads to the various stop options.
Alarm LED	Off	There are no unconfirmed alarm messages.
		There is at least one unconfirmed alarm message. You need to deal with it as soon as possible.
	•	There is at least one unconfirmed alarm message. You need to deal with it immediately, processing may not be able to continue unless you do so.
	· · · ·	gnal is sounded when an alarm is generated. You can me ( <b>Utilities</b> > <b>Configuration</b> > <b>System</b> > <b>Volume</b> ).
Table A-4	LEDs and their m	eaning

## **Main components**

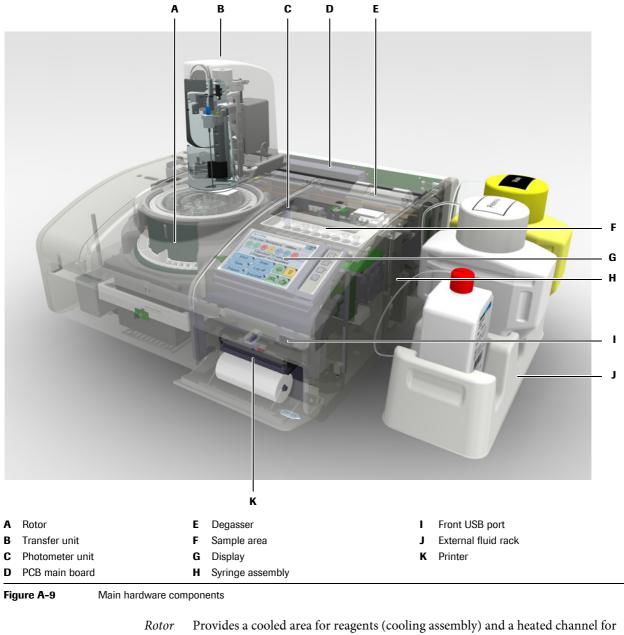

The following figure illustrates the main components of the Cedex Bio instrument.

Rotor	Provides a cooled area for reagents (cooling assembly) and a heated channel for cuvettes. It moves the containers to the correct position for loading, removal, pipetting, and measuring.
Transfer unit	Pipettes sample, reagent, and other fluids from their source to target containers such as cuvettes or the wash station.
Photometer unit	Contains the absorbance photometer used for making absorbance measurements.

*PCB main board* Controls the instrument hardware.

**Roche Diagnostics** 

Fluid system (not shown)	Transports all fluids around the instrument, including sample, reagent, calibrator, control, diluent, cleaner, system water, and waste.
Syringe assembly	Performs the aspiration and dispensing of fluids. This includes the supply of water and cleaner to clean the probes in the wash station after every pipetting action, which prevents carry-over between tests.
Degasser	Removes possible air bubbles from the system water.
Sample area	Eight positions for holding sample tubes. This area is also used for placing calibrators, controls, and auxiliary fluids.
Display	The touchscreen provides the user interface for controlling and managing the <b>Cedex Bio</b> instrument.
Front USB port	This port is used for the USB stick when backing up data or loading data on the system.

## Hardware overview

## Sample area

The sample area provides eight positions for placing sample tubes. You can place primary and secondary tubes.

See *Tubes* on page A-53.




**A** The LED indicates that the transfer tower is approaching or that you should place a sample tube.

Figure A-10	Sample area with sample tubes
inguio / io	

Place the samples when instructed by the system to do so. For calibrators, the system tells you on which position to place them, with the other fluids, you can choose any free position.

### **Fluid system**


The fluid system consists of all the valves, pumps, tubing, syringe, fluid sensors, water and waste containers, the wash station, and the probe. It transports all fluids around the instrument, including sample, reagent, calibrator, control, diluent, cleaner, system water, and waste. The fluid system also delivers the correct amounts of fluids to the cuvettes.

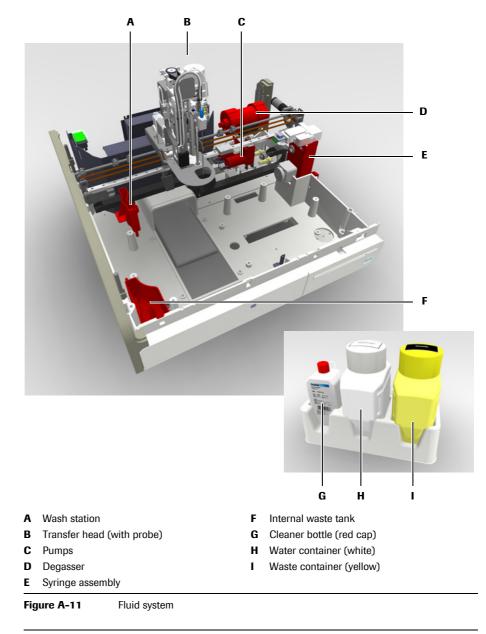

#### NOTICE

To prevent overflow of the internal waste tank when the system is in Standby status, the waste is periodically pumped to the external waste container. (Condensation can build up in the cooling assembly while the system is in **Standby** status.)

#### Probe and syringe

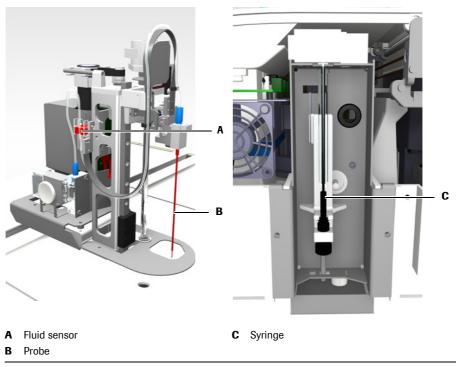




The probe is connected by tubing to:

- The syringe to ensure the pipetting of the required amounts of fluid.
- The external water container and cleaner bottle to ensure supply of fresh water and cleaner.
- See maintenance action *Clean the probe manually* on page C-12.
   See maintenance action *Deproteinize the probe* on page C-10.
   See *Replacing the probe* on page B-125.

#### Syringe assembly

The syringe assembly controls the aspiration and dispensing of fluids. It also controls the supply of water and cleaner to clean the probes in the wash station after every pipetting action, which prevents carry-over between tests.

#### Wash station

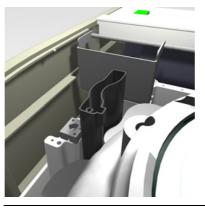




The probe is washed after each pipetting. It is lowered in the wash station and then cleaner is pumped through the probe to wash it in and outside. Next water is pumped through the probe to flush away the cleaner.

The wash station is connected by tubing to the internal waste container.

#### **External fluid connectors**

The three external fluid containers must be properly connected before you switch on the **Cedex Bio** instrument.

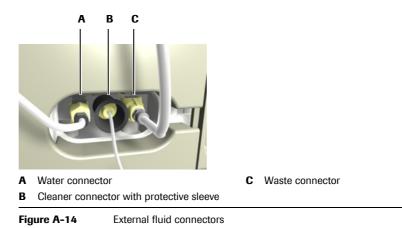

## **Transfer unit**

The transfer unit moves the probe to the correct positions for all pipetting and cleaning actions.

The following figure shows the major parts of the transfer unit.


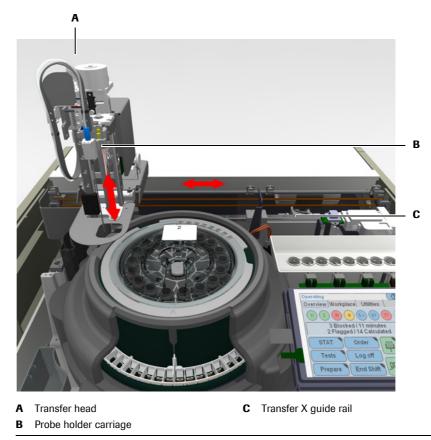


Figure A-15 Major elements of the robotic transfer unit

*Transfer head* The transfer head moves horizontally (along the X-axis); the probe moves up and down (Z-axis), and it performs a rotational movement for mixing the cuvette and reagent bottle content.

*Transfer head arrest* When the transfer head is obstructed in its horizontal movement, it immediately stops. All pipetting and processing actions stop.

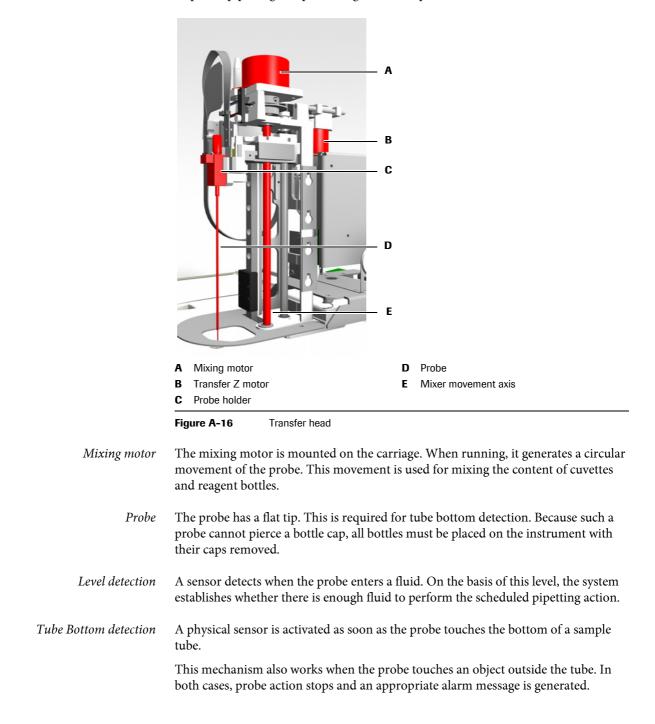

## **Fluid containers**


Fluid	Container(s)	Position
Sample	Tube	Sample area
Control	Tube	Sample area
Calibrator	Tube	Sample area
Diluent	Reagent bottle	Reagent disk
System cleaner	External cleaner bottle	External fluid rack
Cleaner	Reagent bottle	Reagent disk
Reagent	Reagent bottle	Reagent disk
Water	Bottle	External fluid rack
Waste	Bottle	External fluid rack
able A-5	Fluid containers, what they are used for a	nd where

The following table shows which container is used for which fluid:

## -`Q́-

The term tube includes all kinds of tubes, as long as their dimensions lay within prescribed limits. It also includes secondary tubes (cups). See Tubes on page A-53.

#### Waste containers

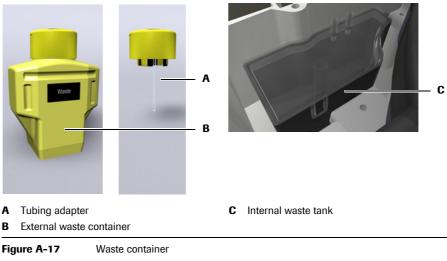

Internal waste tank

The internal waste tank collects the waste from the wash station and the ISE unit, if this is used. It also collects the condensation from the cooling assembly in the rotor.

The internal waste tank is connected by tubing to:

- External waste container •
- Wash station .
- Reagent cooler (condensation)
- ISE unit •

*External waste container* The yellow external waste container is placed on the external fluid rack. It is designed to be washed and reused.

Because the system periodically performs wash actions, an external waste container must be connected at all times. Therefore, when you empty the waste container, you immediately replace it with the spare container and then empty the original container. (The instrument is supplied with a spare waste container.)

There is no active level monitoring for the external waste container, but you are notified if the external waste container has not been emptied for more than one day.

The external waste container is connected by tubing to the internal waste tank.

See Connecting and disconnecting the external fluid containers on page B-127.
 See Checking the status of the external fluid containers on page B-15.
 See maintenance action Clean the water and waste containers on page C-16.

#### Water container

The white water container is positioned on the external fluid rack. Attached to the cap is a suction tube, which is equipped with a water filter.

There is no active level monitoring for the water container, but you are notified if the water container has not been refilled for more than one day.

The water container is designed to be washed and refilled.

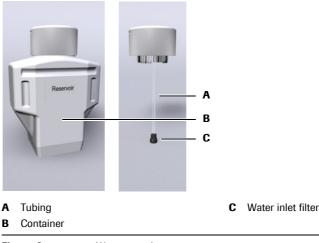


Figure A-18 Water container


The water container is connected by tubing to:

- Wash pump
- Syringe assembly
- Probe
- See Connecting and disconnecting the external fluid containers on page B-127.
   See Checking the status of the external fluid containers on page B-15.
   See maintenance action Clean the water and waste containers on page C-16.
   See maintenance action Replace water inlet filter on page C-19.

#### **Cleaner bottle**

The cleaner bottle is positioned on the external fluid rack. It is designed to be replaced when empty.

Level monitoring for the cleaner bottle is based on the number of cleaning and pipetting actions that were performed.

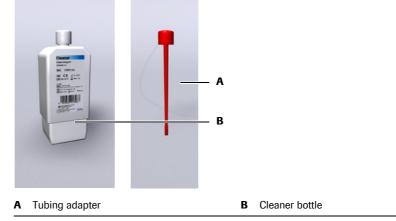



When delivered, the bottle has a white cap. During installation, this is replaced by a red cap with tubing attached to it (tubing adapter).

The cleaner bottle is connected by tubing to:

- Syringe assembly
- Probe
- See Connecting and disconnecting the external fluid containers on page B-127.
   See Checking the status of the external fluid containers on page B-15.

### **Reagent bottles**



The **Cedex Bio** instrument works exclusively with reagent bottles that are equipped with a two-dimensional barcode.

Each bottle holds up to 20 mL of fluid. The actual volume depends on the test.

Place the bottles on the reagent disk as instructed by the software.

Reagents are handled in reagent sets. A set consists of up to three reagents. You always load and replace all bottles of a set.

#### Chimneys



Chimneys are bottle inserts that reduce evaporation.

For reagents that are especially sensitive to concentration changes, Roche recommend using chimneys on the reagent bottles.

To generally reduce evaporation, you may use chimneys on all reagent bottles.



#### Incorrect results due to declining reagent quality

If the application definitions (see package insert) recommend the use of chimneys, the corresponding calibration intervals apply to conditions when working with chimneys. Roche recommend using chimneys whenever this is recommended on the test insert.

#### Tubes

The Cedex Bio instrument can use both primary and secondary tubes (cups).

You can use any type of primary tube, as long as their dimensions lay within prescribed limits.

- Maximum height (including secondary tube): 102 mm
- Minimum height: 70 mm
- Maximum outside width: 16.3 mm
- Minimum outside width: 11.8 mm



#### Incorrect results due to insufficient fluid

Insufficient fluid may lead to inaccurate pipetting and consequently to incorrect results. Always fill the tubes with sufficient fluid that at the end of the pipetting process at least the dead volume of fluid is left.

• See *Tubes* on page A-53.

WARNING

#### Incorrect results due to inappropriate tube and cup placement

Inappropriate tube and cup placement may lead to inaccurate pipetting and consequently to incorrect results.

Make sure that the primary tubes are placed centrally and perfectly vertically in the holders in the sample area and that they are inserted firmly.

Make sure that secondary tubes are placed centrally on the primary tubes and that they rest fully on them.

NOTICE

#### Probe damage due to not removing primary tube caps

The probe is not designed to pierce tube caps. It can get damaged when trying to pierce tube caps.

Always remove the caps of primary tubes before placing them on the instrument.

The following table lists a few typical tubes that are suitable, and it gives the dead volume for each of them.

Tube name	Dead volume
13 x 75 mm	500 µL
13 x 100 mm	500 µL
16 x 75 mm	700 µL
16 x 100 mm	700 µL

 Table A-6
 Typical examples of suitable tubes

Roche recommends using sample cups because they need much lower dead volume compared to primary tubes. The following table lists recommended cups.

Cup name	Dead volume	Sample volume	Placement
1.5 mL reaction tube	100 µL	1.5 mL	Using adapter cups
Roche Diagnostics Standard false bottom tube	75 μL	2.5 mL	Directly on sample area

Table A-7

Typical examples of suitable tubes

#### **Cuvettes**




All optical measurements are made using the same transparent plastic containers, called cuvettes. Samples are automatically transferred from a sample tube to the cuvettes on the cuvette ring.

Cuvettes are disposable to eliminate carry-over.

#### **Cuvette segments**

Each segment holds 10 cuvettes.

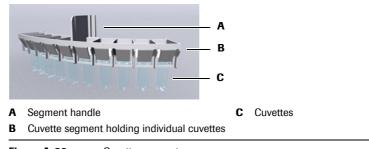



#### **Handling cuvettes**

Cuvettes are supplied in boxes containing cuvette sets. Each set contains a number of cuvette segments. This way, the cuvettes can easily be handled without touching them.

Loading and removing cuvettes is guided by the system software. When handling is required, the rotor moves the cuvette segments to the cuvette port, where you can load or remove them. You handle one segment at a time. Cuvette segments are placed in the cuvette ring of the rotor.


• See *Preparing cuvettes* on page B-28.



Incorrect results due to scratched or soiled cuvettes

Scratches and impurities on the cuvettes distort the measurements.

Do not touch the cuvettes and make sure they do not touch other items when handling them.

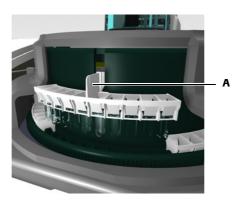


**A** Hold the segment by its handle. Make sure not to touch the cuvettes.

Figure A-21 Handling a cuvette segment

#### **Reagent handling**

Loading and removing reagents is guided by the system software. When handling is required, the rotor moves the bottles to the reagent port, where you can load or remove them. You handle one reagent bottle at a time. Reagent bottles are placed on the reagent disk.

• See *Preparing the reagents* on page B-22.

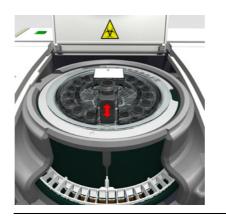




#### Rotor

The rotor provides the following features:

- Space for up to 27 reagent bottles (on the reagent disk)
- Space for up to 60 cuvettes (in the cuvette ring)
- A cooled environment for reagents (reagent cooler)
- A temperature controlled environment for samples (cuvette ring)
- A synchronized transport mechanism to move reagent bottles and cuvettes to the pipetting, loading, and measuring positions.

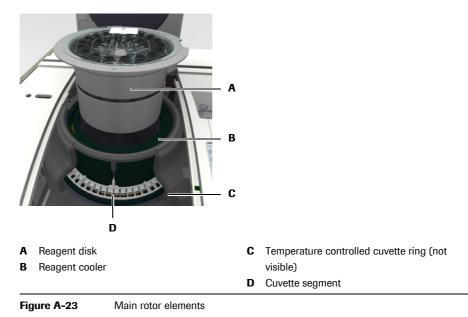

Rotational movement The bottles and cuvettes are positioned in a manner that they can be moved to the various positions by a rotational movement. There are positions for loading and removal (reagent and cuvette ports), pipetting, and measuring.

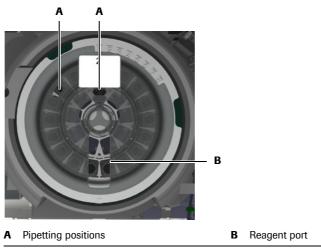


Figure A-24 Reagent port on the reagent disk

#### **Reagent disk**

The reagent disk holds up to 27 reagent bottles. It is designed to be handled as one unit, including the bottles. When not used on the instrument, the reagent disk is placed in a container and stored in a refrigerated place.



Figure A-25 Reagent disk

NOTICE


#### Damage to the reagent disk

The reagent disk is designed to handle reagents while it is loaded on the instrument. The cover is equipped with a locking mechanism.

Always remove and load reagents while the reagent disk is on the instrument and by using the software supported procedures.

• See Preparing the reagent disk on page B-21.

*Reagent disk ID* You can use up to eight different reagent disks on one **Cedex Bio** instrument. Each reagent disk is equipped with numbered tabs. For automatic disk identification by the instrument, one—and only one—of these tabs is removed. The number of this removed tab is the disk ID. When you label the disk, make sure that the number on the label corresponds to that of the removed tab.

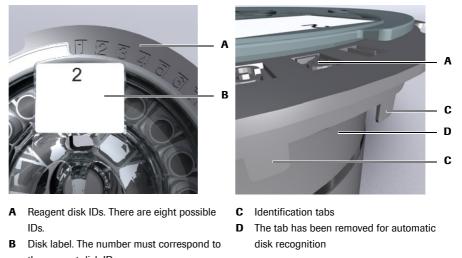


the reagent disk ID.
Figure A-26 Reagent disk ID

### **Reagent disk container**


For storage outside the instrument, the reagent disk is placed in a container. This reduces evaporation of reagents and prevents their contamination.

## **Reagent cooler**


The reagent cooler holds the reagent disk with its reagent bottles. The temperature in the cooler is kept within the range of 6 to 10°C.

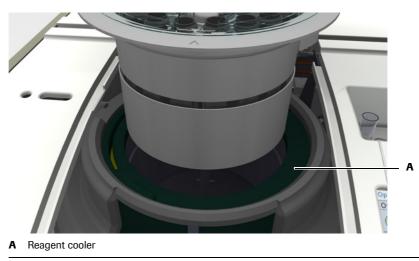



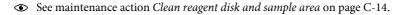

## **Cuvette ring**

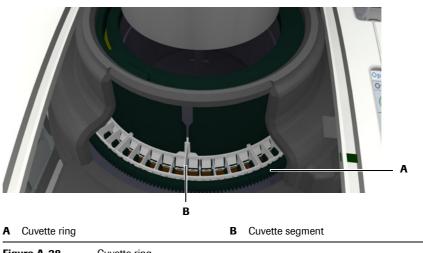

The cuvette ring holds up to 6 cuvette segments, each containing 10 cuvettes.

#### Figure A-28 Cuvette ring

Cuvettes fit neatly in the cuvette ring, without touching the walls when being moved along the ring.

See Preparing the reagents on page B-22.
 See maintenance action Clean reagent disk and sample area on page C-14.

### **Barcode scanner**



A hand-held barcode scanner is used for reading barcoded labels.

#### Loss of sight

The intense light of the LEDs may severely damage you eyes. Do not stare into the LEDs. Scanning equipment using LED technology is covered by the international standard IEC 60825-1 LED Safety: Class 1.

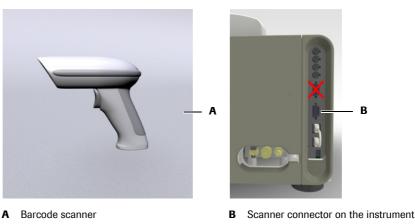


Figure A-29

Hand-held barcode scanner

-Å-	Connect to the lower COM2 port
	Always connect the barcode scanner to the lower of the two serial communication con- nectors (B).
	• For information on how to use the barcode scanner, see <i>Using the barcode scanner</i> on page B-89.
	The following containers are always supplied with barcodes:
	Reagent bottles
	Diluent bottles
	• Auxiliary fluids (diluents, cleaners etc.)
	Sample tubes can be used with or without barcoded labels.
Reagent bottle barcode	On the reagent bottles, barcode of the PDF417 format is used.
	The barcode contains the following information:
	• Part ID
	• Lot number
	Expiration date
	Reagent volume
	Serial number of bottle
	Test data
Sample barcode	The following barcode types are supported for sample tube identification:
	• Codabar
	Codabar 2 of 7
	• Code 3 of 9
	• Code 128
	• EAN
	• Interleaved 2 of 5
	• UPC (A, E)
-`\$ <u></u> '-	A sample barcode must include a checksum at the end.
E	

### **Printer**

 A Panel release button
 B

 B Paper slot

Figure A-30 Printer

The **Cedex Bio** instrument has a built-in thermal printer with a 112 mm paper roll. The printer is used for example for printing placement lists, results, maintenance action instructions, and status information on various items such as the loaded tests.

See maintenance action *Refilling printer paper* on page B-122.

See To clear the paper jam on page D-46.

### **Absorbance photometer**

The **Cedex Bio** main instrument uses the absorbance photometry measuring method.

Absorbance photometer



A Photometer unit

Figure A-31 Photometer unit

The measurements are taken without removing the cuvette from the rotor.

*Halogen lamp* The Halogen lamp is mounted on a holder for easy replacement. The system informs you when you need to replace the lamp.

• See maintenance action *Replace photometer lamp* on page C-24.

#### 3 Hardware

Hardware overview


Wavelengths for the absorbance	
photometer	

For each cuvette, the absorbance photometer measures light intensity at 12 different wavelengths:

340 nm	449 nm	520 nm	629 nm	
378 nm	480 nm	552 nm	652 nm	
409 nm	512 nm	583 nm	659 nm	

*Measurement principles* See *Principles of operation* on page A-19.

## **Connectors**

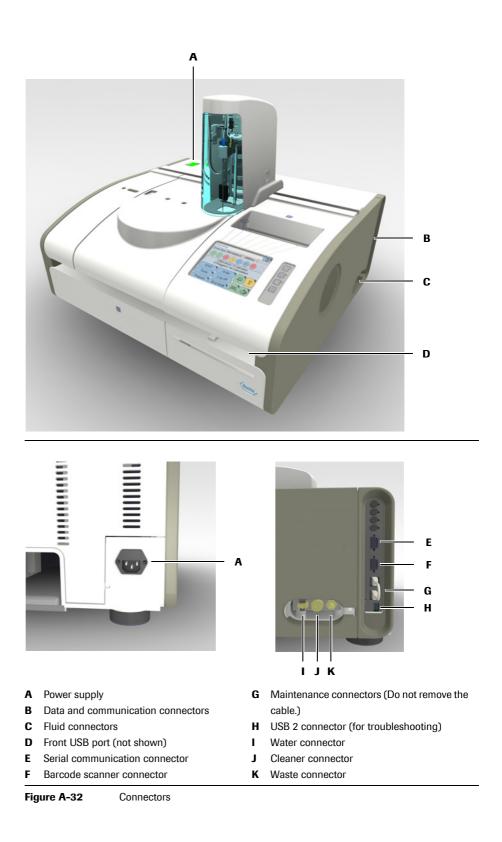

### **3 Hardware**

Hardware overview


## **Fuses**



## Electrical shock by electronic equipment

Do not attempt to work in any electronic compartment. Installation, service, and repair must only be performed by authorized and qualified personnel.

The mains fuses are situated at the rear of the instrument, above the power connector; the internal fuses are situated on the right side of the instrument, at the top of the connector panel.

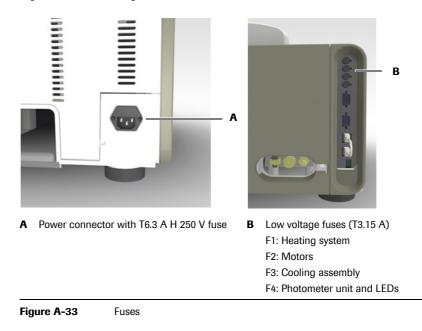

See To change the mains fuses on page D-47.
 See To change a low voltage fuse on page D-49.

# **Technical specifications**

NOTICE	Every effort has been made to ensure that all the information contained in these specifica- tions is correct at the time of printing. However, Roche Diagnostics reserves the right to make any changes necessary without notice as part of ongoing product development.					
Physical dimension	Width	720 mm				
	Depth	550 mm				
	Height	480 mm				
	Weight	Approximately 35 kg				
Power requirements	Line voltage	100-125 V and 200-240 V (-15%, +10%)				
-	Line frequency	50 Hz (±5%) and 60 Hz (±5%)				
	Power consumption	250 VA				
	Insulation coordination	Installation category II (IEC 61010-1)				
	Main fuse	T6. 3 A H 250 V				
	Low voltage fuses	T3. 15 A L 250 V				
	Battery	Lithium 3.6 V 2.3 Ah SL-360/S				
Power requirements ISE unit	Linavoltaga	100-240 V (±10%)				
	Line voltage Line frequency	50 Hz (±5%) and 60 Hz (±5%)				
	Supply voltage	19-24 V DC, Min. 2A				
	Power consumption	70 VA				
	Insulation coordination	Installation category II (IEC 61010-1)				
Measurement principles	Absorbance photometry	(enzymes, substrates, drugs of abuse)				
	Potentiometry	ISE (ion selective electrodes) Na <sup>+</sup> , K <sup>+</sup> , Cl <sup>-</sup>				
Environmental conditions	Temperature	Running conditions: 15-32°C				
		Transport and storage: -25 to +60°C				
	Humidity	Running conditions: 30-80% at 15-32°C, non condensing.				
		Transport and storage: 10-95%, non condensing				
	Pollution	Degree 2 (IEC 61010-1)				
	Altitude	Max. 2000 m above sea level				
Throughput	Photometric	85 tests/h max.; 60 tests/h consolidated (with typical test panel)				
	ISE	180 tests/h max.; 60-100 tests/h consolidated (photometric and ISE mixed)				
Samples	Sample handling	Manually by operator				
	Time to first result	5-10 min (photometric measurements)				
		2 min (ISE measurements)				

## 3 Hardware

Technical specifications

Water purity	Minimum requirements	Electrical resistivity [MΩ*cm @ 25°C]	> 1
		Electrical conductivity [µS/cm @ 25°C]	< 1
		Silicate (SiO <sub>2</sub> ) [mg/L]	< 0.1
		Particulate matter size [µm]	n/a
		Bacteria [CFU/mL]	< 1000
	Whenever the term "purified wate specified above must be used. Roche recommends using Reagen	er" is used in this document, water of at least It Grade water.	the quality
Calibrators	Roche calibrators	See the package inserts of the reagents.	
Reagent bottles	Reagent bottles	20 mL maximum	
	Identification	Barcode	
	Barcode	2-D, format PDF417	
	Number of tests	50-200 tests, depending on the test	
Cuvettes	Segments of 10 cuvettes	Manual insertion and removal of segments	
	Single use	Cuvettes are for single use only	
	Static incubation temperature in cuvette	37°C ±0.5°C	
Photometer	Absorbance photometer	20 W halogen lamp	
	12 wavelengths	340-659 nm	
	Sensor	Photosensitive diode array	
ISE unit	Ion-selective electrode	Indirect measurement	
	Sample volume	15 $\mu$ L; Dilution 1:6 (1 part sample, 5 parts v	water)
	2 Electrodes	Na, K	
	1 ISE Reference Electrode		
oftware data handling	Operating system	<ul><li>LINUX</li><li>VX Works</li></ul>	
	CPUs	Intel XScale	
	Memory system	<ul><li>Flash ROM</li><li>DRAM</li><li>SRAM</li></ul>	
Mass storage	External	USB memory stick	
-	Internal	Flash ROM	
Interfaces	USB1.1/2.0	For backing up data or loading data on the (memory stick)	system
	USB1.1/2.0	Modem	
	2 x RS232	Host, barcode scanner	
Display	Color touchscreen	5.7 inch active matrix ( $^{1}/_{4}$ VGA, 320 x 240	pixels)
Printer	Internal thermal printer	Paper width 112 mm	
	Philter	r	

# Software

# Getting the most out of the instrument

In this chapter, you will find information on how to operate the instrument by using the touchscreen. You will find out about the concept of wizards and you are introduced to the key screens.

In this chapter	Chapter	4
Introduction		A–71
Screen layout		A–72
Display items		. A-73
Color concept		. A-74
Workflows and wizards		A–74
Navigation		A–74
Working with the user interface		A–75
Adjusting the touchscreen		. A–75
Scrolling		. A–75
Expanding and collapsing lists		. A–76
Typing text		A–76
Using the filter function		. A–78
Printing information		
Using online Help		. A–81
Messages		A-82
Message screen		A-82
Alarm monitor		A-82
Key screens		. A-84
Overview tab		. A-84
Sample overview		A–86
Orders		. A-87
STAT		. A-89
Tests		A-89
Log off		A–91
Prepare		. A–91
End shift		A–91
Cuvette status		A-91

Table of contents

Disk and reagent status	
ISE status	A–93
System status	A–94
Workplace tab	A–96
Orders	A–97
Result list	A–98
QC status list	
QC history	A–101
Calibrations list	A–103
Lot data	
Lot list	
Loadlist	A–106
Worklist	
Utilities tab	
Configuration	A–108
Maintenance	
Applications related functions	
Applications	
Extra wash cycles	
Host codes	
Reagent mixing	A–113
Process sequence	A–113
System diagnostics	
Inventory	A–115
Import	A–116
Users	A–117
Export	A–117
Stopping a run	A–118
Color interpretation for LEDs	A–119
Buttons	A–121

## Introduction


To operate the **Cedex Bio** instrument, you use its touchscreen. The design and functional concept of the touchscreen support you in the way you work.

The following table lists the major items and characteristics of a **Cedex Bio** screen and describes their impact on the operation of the instrument.

Screen item	Impact on operation
Screen types	Distinct screen types make it easy for you to find your way around the user interface. For example, you immediately know whether you are on a main screen or whether you are reading a message.
Screen layout	The consistent layout allows you to quickly find the required information and to locate the various display items.
Buttons	Press a button to open and close a screen or to start a func- tion.
Colors	The color of an item on the screen points to its own status of that of the item it represents.
	The "traffic light" color scheme is used:
	<ul> <li>Green: Everything is OK.</li> <li>Yellow: The system is working, but you need to intervene for it to continue to do so.</li> <li>Red: This item does not work, you need to intervene.</li> </ul>
Wizards	A wizard denotes a predefined sequence of screens (steps) that represent a certain task, for example defining an order.
	By following the suggested sequence of steps you ensure the proper execution of the tasks and functions.
	With some critical tasks, you need to follow the wizard exactly, right through to the end. In other cases, you may skip a step and perform it later.
Table A-8	The major items and characteristics of the touchscreen

Screen layout

## **Screen layout**

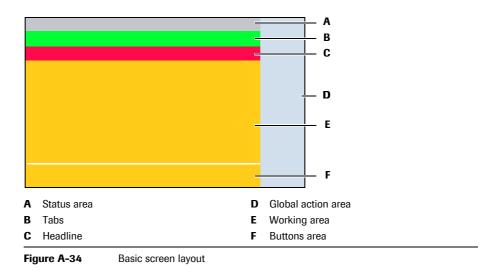


All screens are based on the following layout:



The screen representations shown in this chapter and throughout this manual are for illustrative purposes only. The screens do not necessarily show valid data.

Depending on the function of a screen, some layout items may not be displayed.

The following is an example of a screen with the full range of display items.

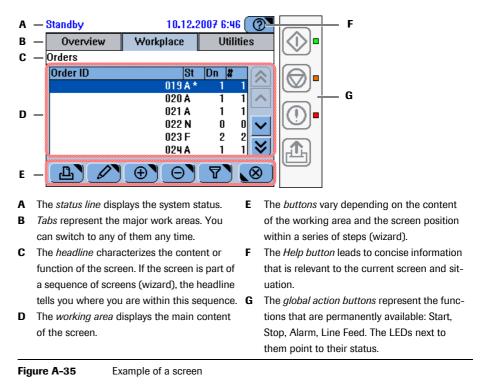

# **Display items**

The Cedex Bio screens are made up of text areas and various kinds of display items such as tabs and buttons.

The following table lists the major display items and describes their use.

Display item		Use
	Button	Press a button to start a function. In addition, many buttons also either open a new screen or close the current screen.
		A triangle in the top right corner of a button tells you that a new screen will be displayed when you press the button; a triangle in the bottom left corner that the current screen will be closed.
$\bigcirc \bigcirc \bigcirc \bigcirc \textcircled{2}$	Global action button	The global action buttons are positioned on the right side next to the screen. The LED next to each of the buttons indicates whether the button is active or not.
Standby 10.12.2007 4:22	List	Press a list item to select it. (Its color turns blue.)
Clean Works         10032           Departs However and Departs However and Departs However and Departs However and Departments However and Department However and Department However and Department However and Dep		Use the scroll buttons to display the items that are not visible.
Hace sample 12 on the instrument while the sample LED is grain. To prevent misplorement, this workflow is aborted II UK is not presed within 10 seconds.	Text	Text usually provides information or instructions. It can be color coded to indicate its importance level.
Remaining leases 0 Exploration date 01,12,2000 Lot ID 19952005 Secial number 24073		
Dender         23d(s-00) 1934 (2)           Derview         Wokalac         Utilities           Configuration         Maintenance         Applications           Applications         Utilities         Import           Users         Expert         Expert	Tab	Tabs are used to group information into units that can be displayed on one screen.
Table A-9 Ma	jor display items	

Table A-9

Workflows and wizards

## **Color concept**

The color of buttons and other display items tells you about the status of the display item or the item it represents.

The Cedex Bio instrument uses the familiar "traffic light" color scheme.

Color	Meaning for buttons				
Green	The element is OK.				
Yellow	Your intervention is required to ensure continuous operation.				
Red	Your immediate intervention is required. Operation has stopped.				
Blue	The item is selected.				

 Table A-10
 Color concept

- For details on the meaning of LED colors, see *Color interpretation for LEDs* on page A-117.
- For details on the meaning of button colors, see the explanations in the relevant operation instructions.

## Workflows and wizards


Screens and sequences of screens help you perform your tasks. If not all steps of a task can be performed from one screen, the workflow is realized as a sequence of screens, a so-called *wizard*. **Cedex Bio** wizards do not usually *force* you to perform a task at a certain stage, they just make your work easier.

## Navigation

Moving from screen to screen

Knowing where you are So

Screens on which you perform tasks provide a headline that displays the navigation path of the current screen.

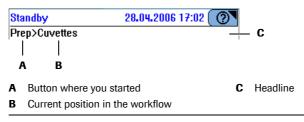

You move from screen to screen with the help of buttons.

Figure A-36Headline with navigation path

• For an overview of the navigation buttons, see *Navigation functions* on page A-121.


## Adjusting the touchscreen

With a touchscreen, it is important that the point where you press the screen corresponds exactly with its hardware equivalent. If this were not the case, pressing a screen item such as a button might not lead to the expected result.

• See Adjusting the touchscreen on page B-131.


## Scrolling

If not all text or all list elements fit on one screen or display area, use the scrolling function to display the hidden content.

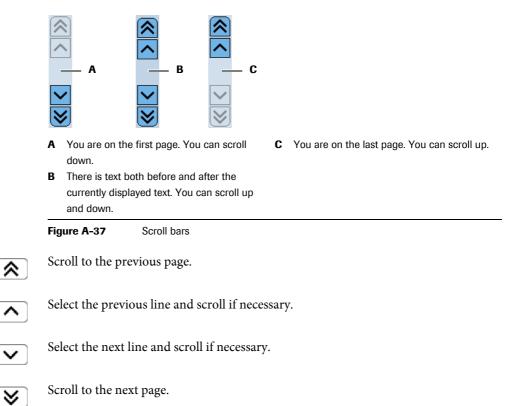

## **Expanding and collapsing lists**

In hierarchically structured lists, you initially see only the top level entries. List items that contain (but hide) lower levels of entries are marked with 🛨. List items that display lower levels of entries are marked with 🖃.

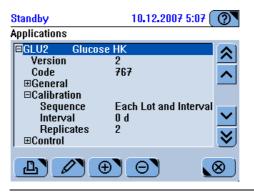



## To expand a list

- 1 Select a list item marked with 🛨.
- **2** Press **+** again or press  $\swarrow$ .
- 3 Use the scrollbar, if required, to display the items you are interested in.

## To collapse a list

- 1 Select a list item marked with  $\square$ .
- **2** Press  $\square$  again or press  $\square$ .

## **Typing text**

There are dedicated screens for typing alphanumeric and numeric characters.


You can choose from the following screens:

- Alphanumeric upper case
- Alphanumeric lower case
- Special characters
- Numeric characters

Numeric keyboard	Standby	10.12.20		
	Overview	Workplace	Utilities	
	Order>Identify S	ample		
	113			— A
	7     8     9       4     5     6       1     2     3       0     A-Z     ←       Of \$     \$		<b>»</b>	
	A Typed text			

Figure A-39 Numeric keyboard screen

- ← Delete the last character displayed in the text line.
- A-Z Switch to the uppercase alphanumeric keyboard.

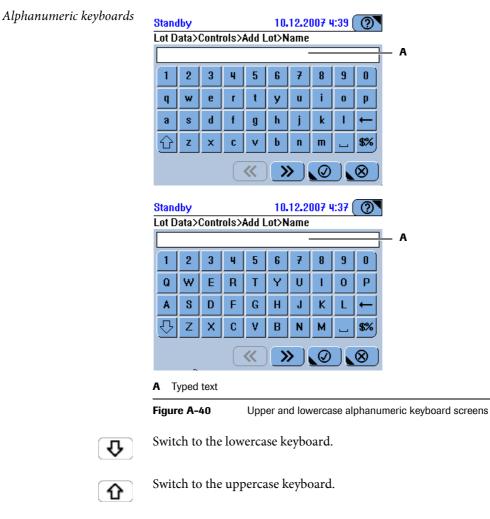

- ← Delete the last character displayed in the text line.
- **\$%** Switch to the special characters keyboard.
- Press to insert a space.

Special characters keyboard	Standby     10.12.2007 4:40       Lot Data>Controls>Add Lot>Name								
									A
	! @	# \$	%	^	&	*	ĺ	]	
	+ "	· `	<	>	]	]	{	}	
	- 1	۲ =	?	;	:	,	•	←	
	Clr	_ 1	~	Ц	ш	<b>_</b>	<b>_</b>	A-Z	
			<b>«</b>			Ø		8	
	A Typed te	ext							
	Figure A-4	1	Spec	cial c	harad	cters	keybo	oard s	screen
	Switch to	the lov	verca	ise al	lpha	num	neric	keyl	board.
A-Z	Switch to	the up	perca	ase a	lpha	nun	neric	key	board.
÷	Delete the	last c	harac	ter c	lispl	ayed	l in t	he te	ext line.
<u> </u>	Press to in	nsert a	space	е.					

## Using the filter function



In many lists you can apply a filter, that is, you can select predefined criteria for generating a selection of entries.

The way to apply a filter is the same in all screens where a filter is available. Here is an example:

## To apply a filter to a list

**1** Display the list.

Standby	10.12.2007 3:41 🕐			
Overview	Workplace	Utilitie	s	
Result Review		u		
S Test Flag				
C. HILL-021		09.12.2007		
vCH02A v	2	.38mmol/L		
Z. CASABLANC	A-023	09.12.2007		
! CA	1	.89mmol/L	$\mathbf{\sim}$	
! ->- >	1	.89mmol/L		
A LEBRUME-02	24	09.12.2007	l ≫)	
	<u>P</u> 0	<b>?</b> \	$\otimes$	

**2** Press **7**.

A screen is displayed for selecting the filter criterion.

Α	
Standby Result Review>Filter	10.12.2007 4:16 🕐
All	Flagged
STAT	Not Accepted
Not Today	Not Sent to Host
By ID	

**A** In the status line, the screen is indicated where you pressed  $\boxed{1}$ .

- **3** Choose one of the filter options.
- 4 Press 🔘.

The list is displayed again. It now contains only the entries that comply with the criterion you just applied by pressing its button.

<u>َ</u>ڳ

After you have applied a filter, the criterion name will appear as part of the **List** button, for example on screens for deleting data. If you used the **Not Accepted** filter criterion, the **List** button would be called **List [Not Accepted]**.

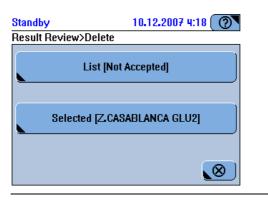


Figure A-42 List button names

## **Printing information**




On many screens, you can print the contents of the working area on the built-in printer. In many cases, a screen is first displayed for selecting the kind of data you want to print. In these cases, the print button is marked with a triangle in the top right corner **S**.

## To print information

1 Press 🖪.

If filter criteria are available, a screen is displayed for selecting what data you want to print. For example:

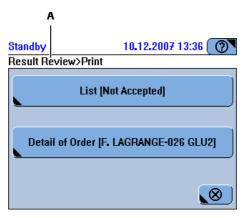

A In the status line, the screen is indicated where you pressed  $\square$ .

**2** Choose one of the print options.

The appropriate data is automatically selected and printed.



Choose  $\bigcirc$  >  $\square$  to terminate the printing task, if required.

## **Using online Help**



The Help button is a permanent feature of all screens.




**A** The Help button is always in the top right corner of the screen.

Figure A-43 Permanent Help button

## To display the Online Help

1 Press 💇.

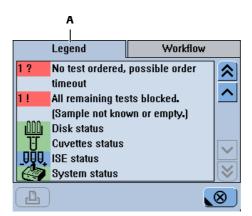

В Legend Workflow Main screen for performing the daily  $\Diamond$ routine tasks: Logging on and off, preparing the ~ system, defining orders, supervising the progress, ending the shift. Also for checking the system and module statuses. <System Status> button: If one of the underlying buttons has a status color ⊗ ப .⊗

A The Legend tab describes the buttons and B their colors.

The **Workflow** tab provides additional information on items on the screen or on actions you can perform.

Figure A-44

- **2** Use the scroll bar to display the hidden information.
  - For details on scrolling, see *Scrolling* on page A-73.

## Messages

Messages are displayed in two ways:

- Immediate feedback on user actions is displayed in a pop-up message screen.
- Information concerning a problem that occurred during operation is reported in the *alarm monitor*.

#### Message screen

Message screens are displayed automatically as soon as the message is generated.

Standby	10.12.2007 13:16 💽
Invalid Data	
	Insert all values.
	$\bigotimes$


Figure A-45 Message screen

Read the message and press  $\textcircled{\otimes}$  to close the screen.

### **Alarm monitor**

Messages concerning an irregularity that occurred during operation can be viewed in the alarm monitor. The alarm LED alerts you when such messages are generated. It is turned off when all alarm messages are confirmed.

The Alarm button 🔘 is always active, even if nobody is logged on the system.

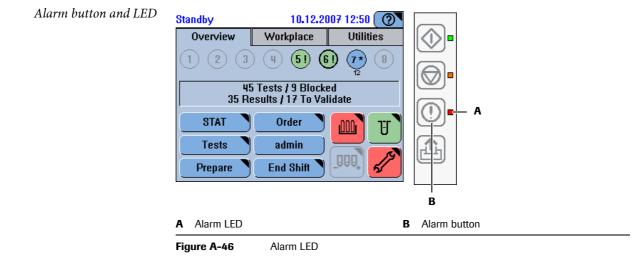

Working with the user interface Interpreting the alarm LED There are no unconfirmed alarm messages. No color, off There is at least one unconfirmed alarm message. You need to deal with it as soon as Yellow possible. There is at least one unconfirmed alarm message. You need to deal with it Red immediately, processing may not be able to continue unless you do so. An acoustic signal is sounded when an alarm is generated. You can adjust the volume Acoustic signal (Utilities > Configuration > System > Volume). To display alarm messages 1 Press ①. **2** Use T to limit the number of messages that are contained in the list. 3 Select the message you are interested in. Alarm message В A 23.06.2009 20:02 Stan łbኣ Alarr 1>20 of 20 7084: 8.100.130 [23.06.2009 18:38] С An ISE pinch valve is probably defective. Restart the system. If the error persists D

call service. Ε ዋ പ്പ  $\otimes$ Alarm ID Problem description and short remedy А D В Total number of messages in the list. suggestion Display the previous alarm message. Display the next alarm message. С Ε

Figure A-47

4 Press  $\swarrow$  to display the detailed message.

Detailed alarm message	Standby 10 Alarm>8 of 8>Detail	.12.2007 5:44 🕐	
	[8.100.130] ISE valve defec valves and tubing for prope Perform restart. If the proble call your service representa	r mounting. 🔨	
	Possible causes: - ISE valve defect, short circ connected	uit or not	
	<b>L</b>		

• For details on handling alarm messages, see *Reacting to alarm messages* on page D-7.



The screen representations shown in this chapter and throughout this manual are for illustrative purposes only. The screens do not necessarily show valid data.


The main screen is divided into tabs. These tabs represent distinct work areas.

- → The **Overview** tab is your main work area when performing the daily routine tasks.
  - Use the **Workplace** to gain information on orders and the corresponding results. You can also start the lot handling functions from this tab.
  - Use the **Utilities** tab to perform tasks that are not normally part of the routine analysis workflow. Typically, these would be administration and maintenance tasks.

The following sections describe the key screens of these tabs and point out the main tasks you can perform on them.

## **Overview tab**

The Overview tab is your main work area when performing the daily routine tasks.

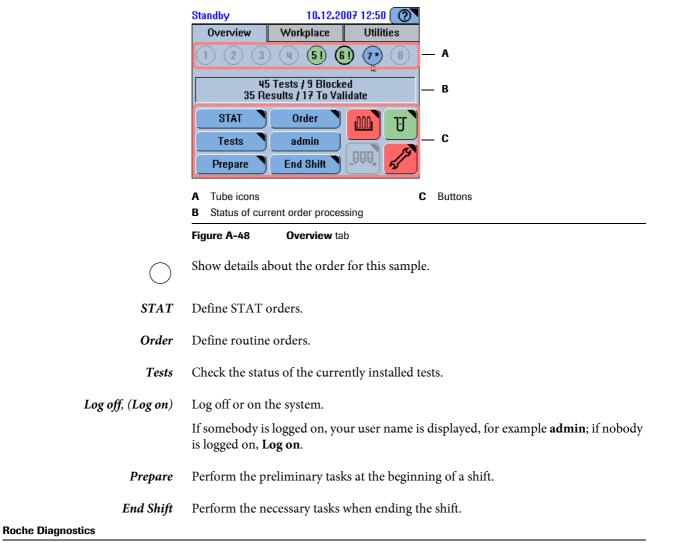

	Check the status of the reagent disk.
	The reagents are OK.
	Fewer than 10% of tests are left for a reagent set, or its expiration date has passed.
	No or unidentified disk on board.
	A reagent set is not complete or a reagent is empty.
T	Check the status of the cuvette segments currently loaded on the rotor.
	More than one segment is available.
	The last available cuvette segment is in use.
	No cuvettes are available.
System Status	The System Status button displays both the icon and the color of one of the buttons
	of the underlying system status screen (see System status on page A-92).
	of the underlying system status screen (see <i>System status</i> on page A-92). The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they are listed below.
	The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they are listed below.
	The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they are listed below. This button can show either of the following icons.
	The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they are listed below. This button can show either of the following icons.
	<ul> <li>The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they are listed below.</li> <li>This button can show either of the following icons.</li> <li>Analyzer (main cover)</li> <li>Reagent cooler and cuvette ring temperature</li> </ul>
	The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they are listed below. This button can show either of the following icons. Analyzer (main cover) Reagent cooler and cuvette ring temperature Sample area ventilation
	<ul> <li>The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they are listed below.</li> <li>This button can show either of the following icons.</li> <li>Analyzer (main cover)</li> <li>Reagent cooler and cuvette ring temperature</li> <li>Sample area ventilation</li> <li>External fluid containers</li> </ul>
	The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they are listed below. This button can show either of the following icons. Analyzer (main cover) Reagent cooler and cuvette ring temperature Sample area ventilation
	<ul> <li>The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they are listed below.</li> <li>This button can show either of the following icons.</li> <li>Analyzer (main cover)</li> <li>Reagent cooler and cuvette ring temperature</li> <li>Sample area ventilation</li> <li>External fluid containers</li> </ul>

### Sample overview

Overview	Workplace Ut	ilities 🚺	
123	) (4) (5) (6) (7		A
	Tests / 9 Blocked sults / 17 To Validate		B C
STAT	Order Drder	U	<u>ل</u>
Tests	admin 📃 💳		È
Prepare	End Shift	5	
Sample tube	buttons	C Resu	sults
Open orders			
igure A-49	Sample tube status	on the <b>Overview</b> t	r tab

1

A sample tube button with a wide edge symbolizes a STAT order.

All tests are completed and their results are accepted.

All tests are pipetted.

The number below the sample tube button indicates the estimated time to completion. You can remove the sample tube, the time to completion disappears when you do so.



All tests are completed but not yet accepted.



1

1\*

All remaining tests are blocked for one of the following reasons:

- There is not enough sample fluid. ٠
- The sample is not identified.

There is no sample on this position.



1

Tests are ordered. Processing has not yet started.



Tests are ordered, processing has started.

The number below the sample tube button indicates the estimated time to completion.



The sample is identified, but no tests were ordered yet.

If working in Order Query Mode: The order could not be obtained from the host.

## Orders


#### Overview > Order **→**

Press Order to define routine orders.

The process of defining an order, and consequently which screens are displayed, depends on how your Cedex Bio instrument is integrated in your laboratory infrastructure (barcodes, host connection).

### Identifying samples

→	Overview > Order	
---	------------------	--

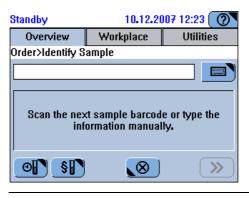

### Figure A-50



Order Interval QC.

Order Default QC.

Typing the sample ID

#### Overview > Order →

Standby	10.12.20	007 12:24 🕐
Overview	Workplace	Utilities
Order>Identify S	ample	
113		
7 8 9		
4 5 6		
1 2 3		
0 A-Z ←		
		<b>&gt;&gt;</b>

#### Figure A-51

• For information on how to use the keyboard screens, see *Typing text* on page A-74.



Order Interval QC.



Order Default QC.

#### Selecting tests, profiles, and ratios,

Standby		25.	06.2009 9	I:56( <b>⊘</b> ▼	
012>Sele	ct Tests				
ISE-I	P Dbt	P LEB	Prof1	RHbA1	
Rati1	A1W2D	A1W2I	ALBU2	ALTL	— A
ALTPL	ASTL	ASTPL	CA	CRPLX	
D-DI	GLU2				
$\mathbf{P}$			»)(«		
В					
Ī					
Standby		25.	06 <b>.</b> 2009 9	:58 🕐	
<mark>Standby</mark> 012>Sciler			1		
Standby		25. DEV DR	1	1:58 🕐	
<mark>Standby</mark> 012>Sciler			1		
Standby 012>S(let ROUT*	STAT D	DEV DR	UG 5	6	— <b>c</b>
Standby 012>Stylet ROUT* ISE-I	STAT C	DEV DR	UG 5 Prof1	6 RHbA1	— C
Standb / 012>Seler ROUT* ISE-I Rati1 ASTL	STAT C P Dbt A1W2D ASTPL	DEV DR PLEB ALBU2 CA	UG 5 Prof1 ALTL CL-1	6 RHbA1 ALTPL	— C
Standby 012>S(leo ROUT* ISE-I Rati1	STAT C P Dbt A1W2D	DEV DR PLEB ALBU2	UG 5 Prof1 ALTL	6 RHbA1 ALTPL	— C
Standb / 012>Seler ROUT* ISE-I Rati1 ASTL	STAT C P Dbt A1W2D ASTPL	DEV DR PLEB ALBU2 CA	UG 5 Prof1 ALTL CL-1 NA-1	6 RHbA1 ALTPL	— C
Standby 012>Select ROUT* ISE-I Rati1 ASTL D-DI	STAT C P Dbt A1W2D ASTPL	DEV DR PLEB ALBU2 CA K-I	UG 5 Prof1 ALTL CL-1 NA-1	6 RHbA1 ALTPL CRPLX	-

→ Overview > Order > identify sample

screen. in tabs. Tabs marked with an asterisk contain

selected tests.

В

Figure A-52 Test selection screens

Tabs are used to group information into units that can be displayed on one screen. The system administrator can define up to six test tabs, name them and assign tests, profiles, and ratios to them.

The tests, profiles, and ratios are sorted alphabetically. Profiles and ratios precede the tests, and they adopt the color of their tests.

The test is on board and ready for use.

- CA
- The test has already been pipetted.
- The test is blocked.

The expiration date of the test has passed.

There are only few tests left.

A QC is due or its result has not been accepted.

A more recent version of the application has been imported.

The test is defined but not on board.

A required diluent or cleaner is not on board.

A-86



Display a screen that contains information on the status of each test.

### STAT

## $\rightarrow$ Overview > STAT

Press STAT to define urgent (short turn-around time) orders.

The process of defining a STAT (short turn around time) order is identical to that of defining a routine order. The difference lies in the scheduling of the tasks. When a STAT order is defined, it will be the next order to be processed, irrespective of what routine orders already exist. Existing STAT orders are finished first.

#### Tests

*Displaying the test overview* 

## → Overview > Tests

Tabs

The tabs are displayed if you work with the order mode **Full**. They represent userdefined test panels. If you work with the order mode **Easy**, all tests are on one panel and there are no tabs.

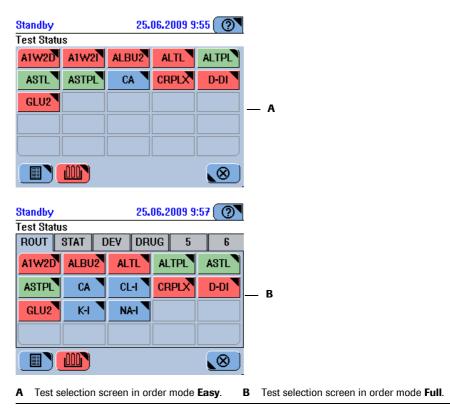


Figure A-53 Test selection screen

The color of a test button represents its status:

The test is on board and ready for use.

The test is blocked for one of the following reasons:

• The calibration is due or failed.

For the reagent set, the number of available tests is 0, or a reagent bottle is missing (incomplete reagent set).
Initial calibration is required.

The expiration date of the test has passed.

There are only few tests left.

A QC is due or its result has not been accepted.

A more recent version of the application has been imported.

The test is not on board.

A required diluent or cleaner is not on board.



Display detailed information on the status of this test.

Display a list of all defined tests, together with information on their status.



Handle the reagent disk that is currently on board.

## Displaying test details


itandby	10.12.2007 13:13 ( ②
est Status>CA	
Calibration	Due
Quality control	ОК
Fests on disk 2	94
Tests ready to run	0
	fication is OK A
	fication is OK A

*Calibration* Information on the calibration status.

*Quality control* Information on the QC status.

*Tests on disk* Total number of tests that are currently available. (There might be more than one reagent set for this test on board.)

*Tests ready to run* Number of tests that could be performed, taking into account all disks known to the system. (The reagent sets have been calibrated and are ready for use.)

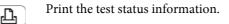

-ċ.	\$ <b>0H</b> 0	Show the fluids that were used for generating this result, together with their lot information. Profiles are user defined sets of tests. They are represented like any other test.
Log off	÷	<b>Overview</b> > button with your user name Log off the system. You can log off any time, even while the system is processing orders.
Prepare	→	<b>Overview</b> > <b>Prepare</b> Start the <b>Prepare</b> wizard to perform the preliminary tasks at the beginning of a shift.
End shift	<b>→</b>	<b>Overview</b> > <b>End Shift</b> Start the <b>End Shift</b> wizard to perform the tasks necessary for ending the shift.
Cuvette status	→	Overview > J.
		Standby     10.12.2007 11:48       Cuvettes       Select segment to replace       14 Required / 40 Available       0     0       10     10       10     10
		A Overview of required and available cuvettes
		Figure A-55Cuvette statusThe six cuvette segments are represented by buttons. The number in the button indicates how many cuvettes are free to be used.Press a segment button to exchange the corresponding segment.The segment button to exchange the defense
	_	The segment buttons are color coded: All cuvettes are used.
		Up to two cuvettes are free to be used.
		More than two cuvettes are free to be used.

Key screens

## **Disk and reagent status**

<b>→</b>	Overview > 100
	A Standby 10,12.2007,12:34 ⑦ Overview>Disk 2 ALTL ASTL BARB+ CA CHOL2 CRPLX D-DI GLUC2 My_MG THC UA2 CLEANF CLEAN NACL PYP I I III E B
	<ul> <li>A plus (+) indicates that there is already an active identical reagent set on board. (This icon is displayed as soon as the first bottle of the set is loaded.)</li> <li>B The number to the left of the button indicates the number of free positions on the disk.</li> </ul>
	Figure A-56         Reagent sets loaded on the reagent disk
	The color of a reagent set button represents the status of the set:
	The reagent set is on board, but it is blocked for one of the following reasons:
	• The number of available tests is 0.
	• The set is incomplete.
	• The test needs calibrating.
	There are fewer than 10% of tests are left for this set.
	The expiration date has passed.
	The reagent set is on board and ready for use.
	<ul><li>There is no application that uses this reagent set.</li><li>A required diluent or cleaner is not on board.</li></ul>
<xyz></xyz>	Display detailed information on the status of this test.
5	Load a reagent set. The number to the left of the button indicates the number of free positions on the disk.
	Display a list with all tests on board, together with information on their status.
	On the list, the following abbreviations are used to indicate the status of the reagent set:
	C. Calibration missing

- C: Calibration missing
- E: Empty

- I: Incomplete
- N: Not used •
- L: Low •
- X: Expired •



Handle the reagent disk.

#### **ISE status**

→ Overview > 🛄.

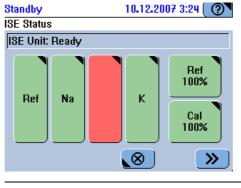


Figure A-57 ISE status

*Na*, *K*, *Ref* Display detailed information on the status of the electrode.

The expiration date of an electrode has passed.

The electrode is ready for use.

Calibration required. (Does not apply to the reference electrode.)

After installing a new electrode, the Electrode Service maintenance action is due.

*Cal*, *Ref* Display detailed information on the status of the ISE fluid bottle.

No fluid registered by a sensor. (Operation has stopped.)

Calibration required.

The fluid level in the bottle is low. (Operation will proceed until one of the sensors detects that there is no fluid.)



**》** 


There is sufficient fluid.

Display the placement list for actions that are due, for example calibration or electrode service.

#### System status


→ Overview > ( ), or ( )

The system status button on the **Overview** tab displays both the color and the icon of one of buttons of the system status screen. (The icons are first prioritized by color, first priority being red, followed by yellow and green, and then according to the sequence in which they appear on the screen.)

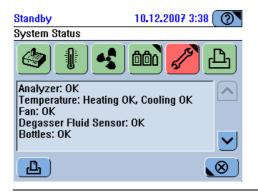



Check the texts for the status of hardware items and on IDs of installed software.

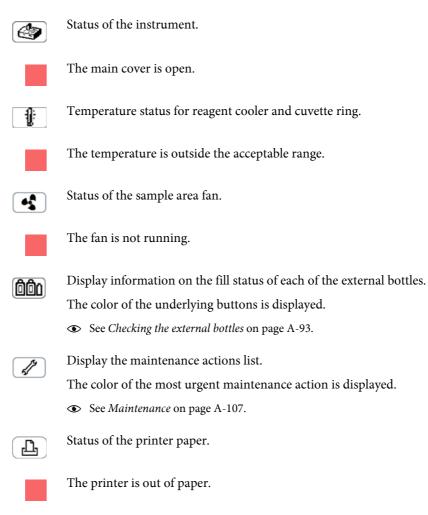

Checking the external bottles

<b>→</b>	Overview > $\bigcirc$ > $\bigcirc$ .
	Standby 10.12.2007 12:27 (?) Check Bottles
	Press fluid button when done.
	Water FullWaste EmptyCleaner 100%
	Figure A-59         Status of external containers
Water button	Press to confirm that you have refilled the external water container.
	The water container was last filled up less than one day ago.
-	The water container was last filled up more than one day ago. (The system does not monitor the filling level. It is up to the operator to fill up the water container regularly.)
	A fluid sensor detected that there is no water in the instrument tubing. No new tests can be processed, started tests may have to be restarted after refill.
Waste button	Press to confirm that you have emptied the external waste container.
	The waste container was last emptied less than one day ago.
-	The waste container was last emptied more than one day ago. (The system does not monitor the filling level. It is up to the operator to empty the waste container regularly.)
Cleaner button	Press to confirm that you have exchanged the external cleaner bottle.
	The cleaner is OK.
	The cleaner level is down to 10% or lower. See the % indication.
	A fluid sensor detected that there is no cleaner in the instrument tubing. No new tests can be processed, started tests may have to be restarted after replacing the bottle.

## Workplace tab

Standby	10.	.12.2007 3:44 🕐
Overview	Workplac	e Utilities
Orders		Result Review
QC Status		QC History
Calibration	is	Lot Data
Loadlist		Worklist

The **Workplace** tab leads to information on orders and the corresponding results.

Figure A-60	Workplace tab

Orders	View and handle orders.
Result Review	View and validate sample results.
QC Status	View and validate active QC results.
QC History	View QC results—on individual screens—of the current or the previous calendar month, or those generated before the previous month.

*Calibrations* View, validate and delete calibration results.

- *Lot Data* Define and change calibration and QC lot data by reading barcodes or typing the values.
  - *Loadlist* List of tests that are ready to be performed.
- *Worklist* Information on tubes currently placed on the sample area.

### Orders

→ Workplace > Orders


Standby	10.12.2007 6:46 🕐		
Overview	Workplace	Utilities	
Orders			
Order ID	St	Dn #	
	019 A *	1 1	
	020 A	1 1 🔨	
	021 A	1 1	
	022 N	0 0 🗸	
	023 F	2 2	
	024 A	1 1 义	
L /	<b>⊕`</b> ⊖`	7 0	

Figure A-61 Orders list

#### St column Status

Α	Accepted	В	All blocked
С	Created	D	Deleted
F	All calculated	Ν	No test order
0	+Order time-out	Р	All pipetted
R	Ready	S	Some blocked
Т	Rerun time-out	U	Unknown
11			1

blank: Done (all tests measured and results accepted)

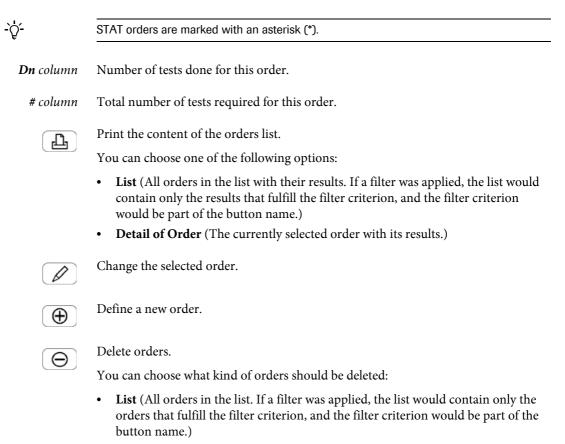

• Selected (Delete the selected order with its results.)



Apply filter criteria to the orders list.

You can choose one of the following criteria:

- All
- Incomplete
- STAT
- Not Finished
- Not Today
- By ID

## **Result list**

→ Workplace > Result Review

Reviewing sample results	Standby 10.12.2007 3:41
	Overview Workplace Utilities
	Result Review
	S Test Flag Result
	C. HILL-021 09.12.2007 VCH02A v 2.38mmol/L
	Z. CASABLANCA-023 09.12.2007
	! CA 1.89mmol/L 🗸
	! ->- > 1.89mmol/L A. LEBRUME-024 09.12.2007 ♥
	Figure A-62 Result list
	This line contains result information.
	This line contains sample information.
	This line is selected.
<b>S</b> column	Status
	!: This result has not yet been accepted.
-`Q́-	STAT orders are marked with an asterisk (*).
·	The time indication represents the time when the order was defined.
	Print results. You can choose one of the following options:
	• List (All results in the list. If a filter was applied, the list would contain only the results that fulfill the filter criterion, and the filter criterion would be part of the button name.)
	• Detail of Order (All results of the associated order)



Validate the results.

You can choose one of the following options:

- **Repeat** (Perform the same test with identical dilution.)
- Rerun (Perform the same test with different dilution.)
- Accept
- Retransmit



Display details of the selected result.



Delete results.

You can choose what kind of result should be deleted:

- List (All results in the list. If a filter was applied, the list would contain only the results that fulfill the filter criterion, and the filter criterion would be part of the button name.)
- Selected



Apply filter criteria to the results list.

You can choose one of the following criteria:

- All
- Flagged
- STAT
- Not Accepted
- Not Today
- Not Sent to Host
- By ID

### QC status list

→ Workplace > QC Status

Standby		23.06.2009 19:50 🕐			
Overview		Workplace	Utilities		
QC Status					
S Test	ID	Flag/Resul	t 🔊		
GLU2	PPU	13.14mmo	/∟		
ALTPL	PNU	52.0U/L	<b>^</b>		
! ASTL	PNU	39.4U/L			
ASTPL	PNU	37.3U/L			
I ALTPL	PPU	R1(2.5s)			
ASTL	PPU	R1(2.5s)	l ≫J		
₽₽₽₽€					

Figure A-63 QC status list

The entries are grouped first by test name, then by control.

S column Status

!: The result has not been accepted yet.

I: The result was ignored.

@: The result has not been transmitted yet.

Flag/Result column Result if no flag was generated.

Flag with highest priority, if flags were generated.

Order status if the control measurement has not been performed yet.

Previous Lot indicates that controls of more than one lot were used.



Print QC results.



Validate the QC results.

You can choose one of the following options:

- Accept
- Ignore
- Retransmit



Display details of the selected QC result.



Θ

Define a new QC order.

## Delete QC results.

You can choose what kind of result should be deleted:

- All
- All Accepted
- Selected

#### **QC** history

The QC history provides—on individual screens—information on the QC results of the current and the previous calendar month as well as on results generated before the previous month.

A graphic representation of the results provides a convenient way for comparing results over a period of time.

#### → Workplace > QC History

Standby		23.06.2009 19:46 🕐		
Overvi	iew	Workplace Utili		es
QC Histor	у			
Test	ID	Last meas	urement	
ALTPL	PNU	22.06.200	9 14:13	
ALTPL	PPU	22.06.200	9 14:13	
ASTL	PNU	22.06.200	9 14:13	
ASTL	PPU	22.06.200	9 14:13	
ASTPL	PNU	22.06.200	9 14:13	
ASTPL	PPU	22.06.200	9 14:13	I ≫∥
L		PO	7	<u>.</u>



The list contains, for each test and control, the latest QC result. The entries are grouped first by test name, then by control.

Last Measurement column

Date and time of the most recent result.

**Previous Lot** indicates that control was performed after a QC lot change.



Print QC results.



Display a graphic representation of the QC results.



Delete QC results.

You can choose what kind of result should be deleted:

- List (All results in the list. If a filter was applied, the list would contain only the results that fulfill the filter criterion, and the filter criterion would be part of the button name.)
- Older than Previous Month
- Selected


 $\left[ \begin{array}{c} \end{array} \right]$ 

Apply filter criteria to the QC results list.

You can choose one of the following criteria:

- All
- Current Lot
- Previous Lot

→ Workplace > QC History > 🔎

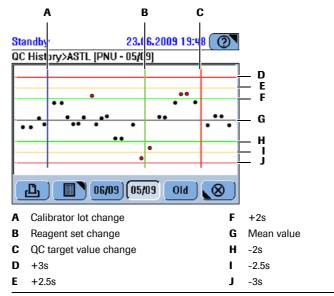





Print the graphic.



Display the results in a table.

The results are sorted chronologically, relevant events (QC lot change, reagent set change, QC target value change) precede the results.

*Month button* Display the results of the month indicated on the button.

*Old* Display the results generated before the beginning of the previous month.

#### **Calibrations list**

→ Workplace > Calibrations

Standby		10.12.2	2007 4:13 🕐
Overview		Workplace	Utilities
Calibration	is [Cui	rrent]	
Test	U	T Status	
ASTL	CU	L Blocked	
CA	CU	S 10.12.2007 1	4:36 🔨 🔨
CHO2A	CU	S 06.12.2007 1	5:24
GLU2	CU	S 03.12.2007 13	3:04 🛛 🗸 🗸
My_MG	CU	S 09.12.2007 1	D:11
UA2	CU	L Calculated	
Ŀ	Ø	<b>⊕</b> `⊖`	7 🚫



*U column* Use of calibration

CU: Current calibration

SB: Standby calibration

**OB**: Obsolete calibration

*T column* Calibration type

L: Lot calibration

S: Set calibration

is displayed.

**Status** column



Print calibration results.

You can choose one of the following options:

• List (If a filter was applied, the list would contain only the results that fulfill the filter criterion, and the filter criterion would be part of the button name.)

The date indicates when the results were accepted. If flags were generated for the result, the flag with the highest priority is displayed. In all other cases, the order status

• Detail of Calibration.



Display details of the selected calibration.



Validate the calibration.

You can choose one of the following options:

- Accept Set
- Accept Lot
- Repeat
- Use Old

Use **Use Old** to override and reset the calibration due date and to continue using the old calibration results.



Define a new calibration order.

#### 4 Software

#### Key screens

Delete the selected calibration result.



Θ


Apply filter criteria to the calibration results list.

You can choose one of the following criteria:

- All
- Current

#### Lot data

#### → Workplace > Lot Data

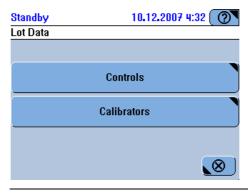



Choose the kind of material for which you want to handle lot data.


*Controls* Handle control lots.

*Calibrators* Handle calibrator lots.

Lot list

→ Workplace > Lot Data, then choose a lot type.

The content of this list depends on the selected lot type. The following screen is an example of calibrator lot data.

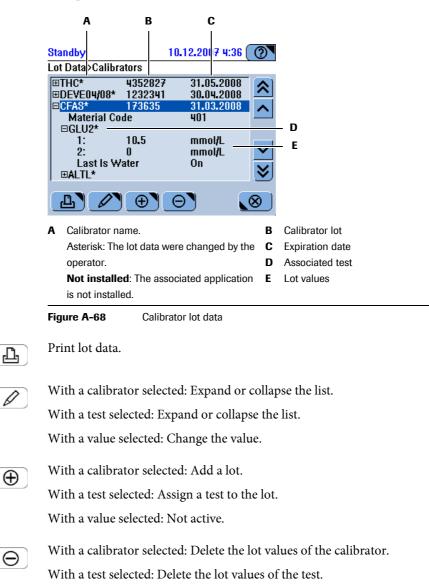

With a value selected: Not active.

#### **4 Software**

Key screens

#### Loadlist

List of tests that are ready to be performed.

#### → Workplace > Loadlist

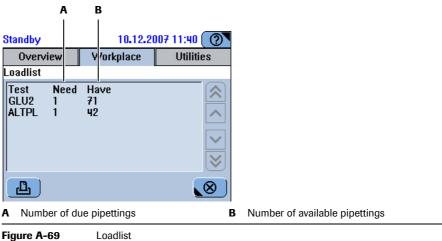


Figure A-69

Print the loadlist.

Д

#### Worklist

The worklist shows all tubes currently placed on the sample area.

ΑВ S andby 10.12.2007 11:39 🕐 Workplace Utilities (verview Y′orklist # T Material ID 5 S 111-050 6 S 222-051  $\approx$ ~  $\sim$ ப  $\otimes$ A Position on sample area E: ISE Etcher B Fluid type: P: ISE Deproteinizer A: Activator Q: Control C: Calibrator S: Sample U: Unknown D: Diluent

#### Workplace > Worklist →





Print the worklist.

#### **Utilities tab**

Use the **Utilities** tab to perform tasks that are not normally part of the routine analysis workflow. Typically, these are administration and maintenance tasks.

Standby		23.06.20	009 19:54 🕐
Overview	Workplace		Utilities
Configuration		Ma	iintenance
Applications		Diagnostics	
Inventory			Import
Users			Export



Configuration	View and change configuration values.		
Maintenance	Select and perform maintenance actions.		
Applications	<ul> <li>Handle applications and their definitions.</li> <li>Handle extra wash cycles</li> <li>Display, export and import the mapping table for test IDs of the Cedex Bio instrument and the laboratory information system.</li> </ul>		
Diagnostics	Perform diagnostics actions.		
Inventory	Display information on bottle sets that are currently defined on any disk used on the <b>Cedex Bio</b> instrument.		
Import	Import application data, software updates, the complete content of a database, certificates, extra wash cycle definitions, reagent mixing rules, or a new user interface language.		
Users	Define users and manage their user rights.		
Export	Export the complete content of the database, the full results, and log files.		

#### Configuration

→ Utilities > Configuration

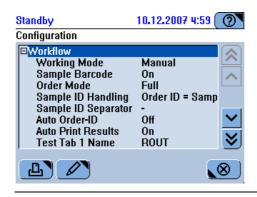


Figure A-72 Configuration table

- ☐ The list is expanded. Select the item marked with ☐ and press ☐ again to hide the items.



Print configuration data.

You can choose one of the following options:

- All
- Selected (Print the settings of the selected configuration group.)



With an item selected: configure the selected item.

With a group title selected: expand or collapse the list.

#### Maintenance

Α	В	
tandby 10.12.200	)7 4:22 🕐	
laintenance		
Clean Water and Waste	10.12. 🔊	
Replace Photometer Lamp 🛛 🗕	10.12.	- C
Clean Probe Manually	10.12. 🔨	
Deproteinize Probe	11.12.	- D
Clean Reagent Disk	07.01. 🧹	
Replace Water Inlet Filter	25.02.	
Clean Housing	02.06. Ў	
<b>L O</b>		
Maintenance action name	D	Use the scroll bar to display the
B Due dates, in ascending order		maintenance actions that are currently
Due dates, in ascending order		

 Figure A-73
 Maintenance actions list


The maintenance actions are listed according to the urgency with which they need to be performed.

No action is currently required.

**Utilities > Maintenance** 

→

This maintenance action should be performed on the next major maintenance day.

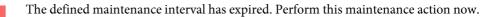



Perform the selected maintenance action.

If you interrupt the performing of a maintenance action that was due, its status will remain due, and you need to fully re-perform the action later.

#### **Applications related functions**

Standby		23.06.2009 19:55 🕜
Applications		( -
Laboratory P	Parameters	Extra Wash Cycles
Host C	odes	Reagent Mixing
Process S	equence	

Figure A-74Applications functions

Laboratory Parameters	Handle the definitions of the installed applications and import and install new applications.
Extra Wash Cycles	Handle and install extra wash cycles.
Host Codes	Display, export and import the mapping table for test IDs of the <b>Cedex Bio</b> instrument and the laboratory information system.
<b>Reagent</b> Mixing	Display, print, and delete mixing rules.
Process Sequence	Prioritize individual tests.

#### Applications

10.12.2007 5:07 0 Standby Applications **⊟**GLU2 Glucose HK 2 767 Version Code ゝ ⊞General ⊡Calibration Sequence Each Lot and Interval 0 d 2 Interval Replicates ⊞Control ⊕` Θ  $\otimes$ J

**Utilities > Applications > Laboratory Parameters** 



- ➡ There are list items of a lower level. Select the item marked with and press again to expand the list and display the items.
- ☐ The list is expanded. Select the item marked with ☐ and press ☐ again to hide the items.

Print the application list.

With an item selected that is preceded by 🛨 or 🖃: expand or collapse the list.

With any other item selected: change the selected item.

 $\bigoplus$ 

→

Import an application.

Define a new profile or a ratio. Install an application




Uninstall an application.

Delete a profile or uninstalled application.

#### Extra wash cycles

→ Utilities > Applications > Extra Wash Cycles

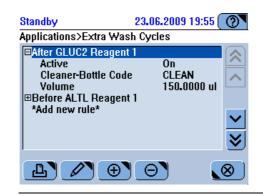





D

Θ


With an item selected that is preceded by  $\textcircled{\bullet}$  or  $\boxdot$ : expand or collapse the list.

With any other item selected: change the selected item.

Print the selected or all extra wash cycle definitions.



Import a new extra wash cycle.

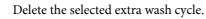

#### **Host codes**

→ Utilities > Applications > Host Codes

Standby		10.12.2007 5:3	87 ( <b>?</b> \
Host Codes	3		
;AppCode 687 767 359 686 706 685 685 684 123	HostCode 687 767 359 686 706 685 685 684 123	;ShortName ;ASTL ;GLU2 ;My_MG ;ASTPL ;CA ;ALTL ;ALTPL ;D-DI	<ul><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li><li></li></ul> <li></li>
Ъ	<b>&gt;</b>	♦ 11	$\underbrace{\otimes}$






Print the mapping table.



Import the mapping table.



Export the mapping table as a text file.

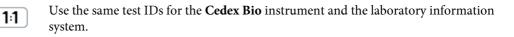

#### **Reagent mixing**



Utilities > Applications > Reagent Mixing

Figure A-78 Reagent mixing rules table



→


Print the selected or all mixing rules.



Delete the selected mixing rule.

#### **Process sequence**

Utilities > Applications > Process Sequence →

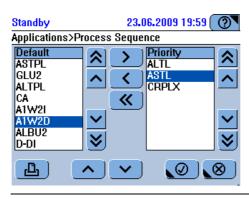






Print the process sequence list.




In the **Priority** list, move the item up.

- Š
- In the **Priority** list, move the item down.

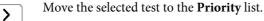



Move the selected test to the **Default** list.

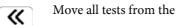

Move all tests from the **Priority** to the **Default** list.

4 Software Key screens

#### System diagnostics

#### → Utilities > Diagnostics

Diagnostics actions are designed for use by Roche representatives or when instructed by them to do so. Therefore, these functions are not described in this manual.

You need at least Lab Manager user rights to perform diagnostics functions.





L)

You can choose one of the following options:

• All

•

Selected (Print the actions of the selected action group.)



`Ų́

With a topic selected: expand or select the list.

With a diagnostic action selected: Perform the selected action.

At the end of performing a diagnostics action (system status transition from **Diagnostics** to **Ready**), the system checks that all modules are initialized, if they are not, initialization is performed.

#### Inventory

The inventory displays information on all bottle sets that are defined on any of the disks used on this **Cedex Bio** instrument, including sets that were removed from the disk, as long as they were not empty or their expiration date had not passed yet.

(Bottle sets that have been removed from the disk and whose expiration date had passed more than 30 days ago are automatically deleted from the **Inventory** list, provided there is still a valid set for the same fluid on board. If this were not the case, the set that was loaded last would remain in the inventory.)

→ Utilities > Inventory

Expiration	1 <b>#</b> C	
5702 03.07.200	9 100 3	
5702 05.06.200	978	
9701 03.07.200	9 100 3	
5751 01.12.200	6508	
5701 05.06.200	979 F	u 🗸 I
5701 08.06.200	996 F	
001 17.06.200	975 3	≫]
$\Theta$ $\nabla$		$\otimes$
	i702         03.07.200           i702         05.06.200           i701         03.07.200           i751         01.12.200           i701         05.06.200           i701         03.07.200           i701         03.07.200           i751         01.12.200           i701         05.06.200           i701         05.06.200           i701         08.06.200	5702 05.06.2009 7 8 9701 03.07.2009 100 3 5751 01.12.2006 50 8 5701 05.06.2009 79 F 5701 08.06.2009 96 F





Print the bottle set list.

You can choose one of the following options:

- List
- Detail of Bottle Set



Display details on the selected bottle set.



Delete the selected bottle set from the list.

Bottle sets must be removed from the disk before you can delete them from the list.



Apply filter criteria to the Inventory list.

You can choose one of the following criteria:

- All
- By ID

#### Import

→ Utilities > Import

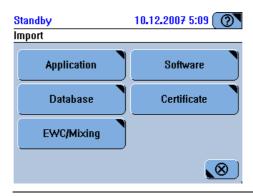



Figure A-82 Importing data

- *Application* Import application data.
  - *Software* Import software updates.
  - *Database* Import the complete content of the database.
  - *Certificate* Import certificates that authenticate reagent barcodes.
- *EWC/Mixing* Import extra wash cycle or mixing information for a particular application.

#### Users


→ Utilities > Users



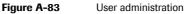


- There are list items of a lower level. Select the item marked with 
   → and press 
   →
   again to expand the list and display the items.
- ☐ The list is expanded. Select the item marked with ☐ and press ☐ again to hide the items.

Print the user data of the selected user.

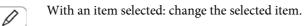

With any other item selected: expand or collapse the list.

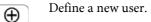


 $\Theta$  Delete the user.

(You cannot delete your own user data.)

#### Export

#### → Utilities > Export

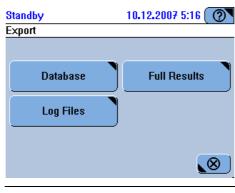



Database Export the complete content of the database.


*Full Results* Export the full result data.

*Log Files* Export the system messages, alarm logs, and possibly trace logs.

#### **Roche Diagnostics**

#### Stopping a run

 $\rightarrow$  Global action button O

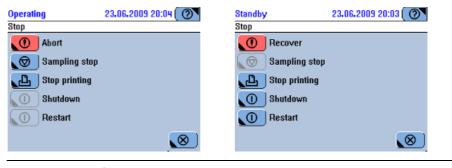



Abort When the system is in **Operating** status: Stop immediately all processing activities. Pipettings that were not finished are considered not pipetted. Measurements that have not yet yielded a valid result are considered not measured. Recover When the system is in Standby or Stopped status: Initialize all systems and modules that are currently not ready. Sampling stop Finish the current pipetting action, but do not start a new one. You can restart processing by pressing  $\mathbf{O}$ . Stop printing Stop the current printing task. (It may take a few moments before printing actually stops.) Shutdown Shut down the Cedex Bio software and the operating system. This option is active in **Standby** status only. Shut down the Cedex Bio software and automatically restart it. Restart This option is active in Standby status only. It is used in cases where a configuration

change requires restarting the software for it to become effective.

Color interpretation for LEDs

# **Color interpretation for LEDs**

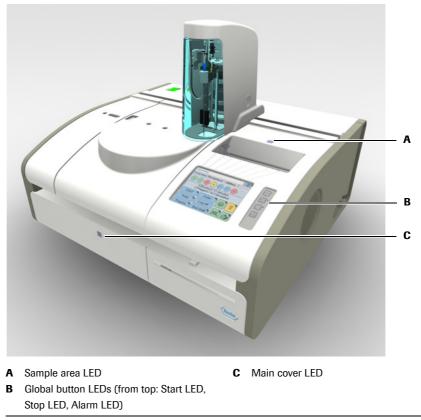


Figure A-86The Cedex Bio instrument

The following table lists the LEDs and explains what their colors mean.

LED	Color	Meaning	
Main cover LED	Off	No activities in this area. You can open the main cover.	
		User intervention is required, for example you are expected to place or remove a bottle.	
	-	The system is performing some action. Do not handle the cover.	
	in <b>Operating</b> sta	al is sounded when the cover is open while the system is te. You can adjust the volume ( <b>Utilities</b> > • <b>System</b> > <b>Volume</b> ).	
Sample area LED	Off	No activities in this area. You can remove sample tubes.	
		You are expected to place a sample tube.	
	Blinking	The transfer head is approaching. Do not place your hand or any object in the sample area.	
Start LED	Off	You cannot start the measuring process.	
		You can start the measuring process.	
Table A-11	LEDs and their meaning		

Color interpretation for LEDs

LED	Color	Meaning
Stop LED	Off	Pressing 🞯 does not have any effect.
		Pressing $\textcircled{O}$ leads to the various stop options.
Alarm LED	Off	There are no unconfirmed alarm messages.
		There is at least one unconfirmed alarm message. You need to deal with it as soon as possible.
		There is at least one unconfirmed alarm message. You need to deal with it immediately, processing may not be able to continue unless you do so.
		gnal is sounded when an alarm is generated. You can ume ( <b>Utilities</b> > <b>Configuration</b> > <b>System</b> > <b>Volume</b> ).
Table A-11	LEDs and their meaning	

# **Buttons**

The following tables list the buttons used in **Cedex Bio** screens and describe their use. The buttons are grouped according to the kind of function they represent.

General functions	lcon	Name	Use
	(⊘▼	Help	Display concise information that is relevant to the current
	0		screen and situation.
	$\bigcirc$	Start	Start processing orders.
	$\bigcirc$	Stop	Display Stop options.
	$\bigcirc$	Alarm	Look at alarms.
	£	Paper Feed	Advance the printer paper.
	Table A-	12 Global act	ion buttons
Interactive functions	lcon	Name	Use
	$( \mathbf{ \oplus } )$	Add	Add or define an item.
	Θ	Delete	Delete the selected item.
	$\bigotimes$	Cancel	Abort the operation.
	$( \blacket )$		Do not use this button while the system is performing control
			or calibration tests, this might lead to data loss
		Close	Close the screen.
	L	Print	Print the content of the current screen's working area.
		List	Display the content of the current screen's working area in a table.
	$\bigcirc$	ОК	Confirm the operation, save the data and close the screen.
	$\bigcirc$	Save	—
		Detail	Show detailed information on the selected item.
	7	Filter	Apply filter criteria to the current view.
		Edit	Modify the selected item.
	0	Perform	Perform the selected action.
	$\textcircled{\black}{\black}$	Export	Export data.
	♦	Import	Import data.
		Disk Status	Handle the reagent disk and its reagents. Show the status of the reagents.
		Cuvette Status	Show the status of the cuvette segments.
	U		Handle cuvette segments.
	Table A-	13 Interactive	buttons

con	Name	Use			
ISE Status Show the status o		Show the status of ISE electrodes and fluids.			
<u>-</u> 999 <sup>+</sup>		Exchange ISE electrodes and fluids.			
	System Status	Show the status of the main cover.			
රිරිර	Container Status	Show the status of the external fluid containers.			
		Handle the external fluid containers.			
1	Temperature Status	Show the status of the reagent cooler and cuvette ring temperature.			
Ŀ	Printer Status	Show the printer status.			
<b>4</b>	Fan Status	Show whether the sample area fan is running.			
1	Maintenance Status	Display the maintenance actions list.			
	Keyboard	Type information manually.			
₽.	Exchange Reagent Disk	Handle the reagent disk.			
100	Insert Bottle	Place a reagent bottle.			
	Remove Bottle	Remove a reagent bottle.			
0	Mixing	Perform reagent mixing.			
Replace Cuvette Confirm replacem		Confirm replacement of the cuvette segment.			
<b>I</b> ¶	Remove Cuvette	Confirm removal of the cuvette segment.			
۲	Due Calibrations	Select all tests with due calibration or QC.			
**	Forecast Calibrations	Select all calibrations that are due within the forecast period.			
\$ <b>[</b>	Default QC	Perform Default QC.			
0	Interval QC	Perform Interval QC.			
? <b>[] </b> []	Related Information	Show context information for the result. (Used fluids, together with their lot information.)			
1:1	Map Host Codes	Use identical application codes on host and <b>Cedex Bio</b> instrument.			
Ĵţ	Exchange ISE Bottle	Exchange an ISE fluid bottle.			
10	Remove ISE Bottle	Remove an ISE fluid bottle.			
<b>₽</b>	Remove Electrode	Remove an ISE electrode.			
10	Insert Electrode	Insert an ISE electrode.			
	I3 Interactive b	uttons (Continued)			

Buttons

Navigation functions	lcon	Name	Use
		Line Up	Move one line up.
		J	Select the previous line.
		Line Down	Move one line down.
	<b>_</b>	J	Select the next line.
	8	Page Up	Move one page up.
	♥	Page Down	Move one page down.
	~	То Тор	Select the first item in a list.
		To Bottom	Select the last item in a list.
	>	Forward	Open the next screen in a wizard.
		Next Step	
	<b>«</b>	Back	Open the previous screen in a wizard.
	$(\mathbf{N})$	Previous Step	
	Table A	-14 Navigatio	on buttons
ement moving functions	lcon	Name	Use
		Move Up	Move selected element one line up.
	$\mathbf{\vee}$	Move Down	Move selected element one line down.
	$\mathbf{>}$	Move Right	Move selected element to the list on the right.
		Move Left	Move selected element to the list on the left.

Table A-15

Move All Left

<

«

Element moving buttons

Move all elements in the list to the list on the left.

#### 4 Software

Buttons

# Operation B

5	Daily operation	<i>B-3</i>
6	Special operations	B-91
7	Configuration	B-133

# **Daily operation**

Performing routine tasks

In this chapter, you will find information on performing the routine tasks that are required for processing orders and keeping the system running.

In this chapter	Chapter	5
Introduction		B-5
Safety information		B-5
Overview		B-7
Working with a host system		B–9
Starting the shift		B-10
Switching on the instrument		B–10
Preconditions		B-10
Logging on to the system		B–11
Preparing the system		B-12
Short guide		B-13
Starting the Prepare wizard		B–15
Checking the status of the external fluid containers		B–15
Performing maintenance actions		
Preparing the reagent disk		
Preparing the reagents		B-22
Checking the status of reagent sets		B-23
Preparing reagents		B-25
Preparing cuvettes		B-28
Performing reagent mixing		B-29
Performing the calibrations (Prepare phase)		B-30
Analyzing samples		B-34
Safety information		B-34
Short guide		B-35
System setup and order definition workflow		B-37
Defining orders		B-38
Defining routine orders		B–39
Defining STAT orders		B-41
Making changes to an order		B-42
Starting the run		B-44

**Roche Diagnostics** 

Table of contents

	Monitoring the analysis progress	B-	-44
	Checking the sample tube status	B-	-44
	Checking the instrument status buttons	B-	-46
	Checking for alarm messages		
	Acoustic signal		
	Stopping and restarting a run		
	Removing sample tubes		
	Validating sample results		
	Repeating tests	B-	-55
	Rerunning tests	B-	-56
	Accepting results	B-	-57
	Printing sample results	B-	-58
]	Performing calibrations	B-	-59
	Safety information	B-	-59
	Short guide	B-	-60
	Process of performing calibrations	B-	-61
	Deleting ordered calibrations	B-	-63
	Validating calibration results	B-	-64
]	Performing QC	B-	-68
	Safety information		
	Short guide	B-	-69
	Performing Default QC		
	Performing interval QC measurements		
	Validating QC results		
	Interpreting the QC history		
	Performing reagent mixing		
]	Finishing the shift		
	Safety information		
	Short guide		
	Checking for unfinished tasks		
	Checking for unfinished orders		
	Checking for not accepted sample results		
	Checking the transmission of results		
	Starting the End Shift wizard		
	Performing the daily backup		
	Exporting the result data		
	Cleaning up the database		
	Performing maintenance actions		
	Replacing cuvettes		
	Dealing with the external fluid bottles		
	Emptying the waste container		
	Refilling the external water bottle		
	Removing the reagent disk		
]	Logging off		
	Shutting down the system and switching off the instrument		
	Using the barcode scanner		
	Reading reagent bottle barcodes		
	Reading barcodes from sheets		
	Reading sample tube barcodes		
	σ····		

# Introduction

Daily operation includes the routine tasks that are required to prepare and monitor the system, and to analyze samples.



The screen representations shown in this chapter and throughout this manual are for illustrative purposes only. The screens do not necessarily show valid data.

#### **Safety information**

Before you start working with the **Cedex Bio** instrument, it is essential that you both read and understand the safety information listed below.

Read carefully all safety notices given in instructions and make sure you understand them.



#### Injury through reagents and other working solutions

Direct contact with reagents, cleaning solutions, or other working solutions may cause personal injury.

When handling reagents, exercise the precautions required for handling laboratory reagents, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics reagents and cleaning solutions.

#### Infection by biohazardous materials

Contact with samples containing material of human origin may result in infection. All materials and mechanical components associated with samples of human origin are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

#### Infection by waste solution

Contact with waste solution may result in infection. All materials and mechanical components associated with the waste systems are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

# Personal injury or damage to the analyzer due to contact with instrument mechanism

Do not touch moving parts during instrument operation.

Keep all covers closed, operate them as instructed on the screen.

Introduction

#### Loss of sight

The intense light of the LEDs may severely damage you eyes. Do not stare into the LEDs. Scanning equipment using LED technology is covered by the international standard IEC 60825-1 LED Safety: Class 1.

#### Incorrect results or damage to the analyzer due to wrong operation

Operators are required to have a profound knowledge of relevant guidelines and norms as well as the information and procedures contained in the Operator's Manual.

- Do not carry out operation and maintenance unless you have been trained by Roche Diagnostics.
- Start all maintenance actions on the screen. Do not perform maintenance actions without the assistance of the user interface.
- Carefully follow the procedures specified in the Operator's Manual for the operation and maintenance of the system.
- Leave maintenance that is not described in the Operator's Manual to a trained service representative.
- Follow standard laboratory practices, especially when working with biohazard material.

#### Incorrect results or damage to the analyzer due to dust and soiling

The user can leave the main cover open while the system is in **Standby** status or while the instrument is shut down. This can cause dust and dirt being collected in the cuvette ring, which in turn might decrease the quality of the cuvettes.

Keep all covers closed. Open them only to perform operation actions.

#### **Overview**

The following table gives an overview of the tasks you might need to perform during daily operation. Roche recommend using the sequence of steps in the way given below; but you may perform the tasks differently. For details on the individual steps, see the relevant sections in this chapter.

	Task	Steps		Navigation	
			With wizard	As individual steps	
1	Starting the system	1. Switch on the system.			
2	Logging on the system			Overview > Logon	
3	Preparing the system	Start the <b>Prepare</b> wizard.	Overview > ]	Prepare	
		1. Check the external fluid containers.		Overview > 🚱 > 🛍	
		2. Perform the maintenance actions that are due.	_	Utilities > Maintenance	
		3. Load the reagent disk.		Overview >	
		4. Check the reagents.		Overview > 🕮	
		5. Check the cuvettes.		Overview > 🕑	
		6. Perform mixing		Overview > 💷 > test > 🖸	
		7. Perform calibrations that are due.		Workplace > Calibrations > 🕀 > 🕙	
4	Defining orders	Start the <b>Orders</b> wizard.	<b>Overview</b> > <b>Order</b> (or <b>Overview</b> > <b>STAT</b> )		
		1. Identify the sample.		n/a	
		2. Select the tests.		n/a	
		3. Place the sample.		n/a	
		4. Start the run.		$\Diamond$	
5	Monitoring the progress		n/a	Overview ①	
6	Validating results	1. View results.	n/a	Workplace > Result Review	
		2. Handle flagged results.	n/a	Workplace > Result Review > 🖉	
				> Repeat	
				> <b>Rerun</b>	
		3. Accept results.	n/a	Workplace > Result Review > 🖉 > Accept	

Table B-1
 Overview of the daily operation tasks

#### **Cedex Bio System**

#### 5 Daily operation

Introduction

Task	Ste	eps	Navigation		
			With wizard	As individual steps	
Performing calibrations					
Performing individual	1.	Start the wizard.	Workplace >	Calibrations > (①)	
calibrations	2.	Select the test.	_	n/a	
	3.	Prepare and place the calibrators.	_	n/a	
	4.	Start the calibration.		$\bigcirc$	
	5.	Validate the calibration results.		Workplace > Calibrations >	
	6.	Remove the calibrators.			
Performing	1.	Start the wizard.	Workplace >	- Calibrations > 🕀	
all due calibrations	2.	Select all tests with due calibrations. or Select all tests with calibration due within the forecast period.	() **	n/a	
	3.	Prepare and place the calibrators.	_	n/a	
	4.	Start the calibration.		$\bigcirc$	
	5.	Validate the calibration results.		Workplace > Calibrations >	
	6.	Remove the calibrators.			
Performing controls					
Performing		Start the wizard.	Overview > 0	Order > 📲	
Default QC	2.	Select a control and place the tube. Repeat until there are no controls left on the screen.	-	n/a	
	3.	Start the QC measurement.		$\diamond$	
	4.	Validate the QC results.		Workplace > QC Status >	
	5.	Remove the controls.		n/a	
Performing an individual	1.	Start the wizard.	Workplace >	• QC Status > 🛞	
QC measurement	2.	Select a test.	_	n/a	
	3.	Select a control and place the tube. Repeat until there are no controls left on the screen.	_	n/a	
	4.	Start the QC measurement.		$\bigcirc$	
	5.	Validate the QC results.		Workplace > QC Status >	
	6.	Remove the control.			
Performing all	1.	Start the wizard.	Overview > 0	Order > 🔍 > 🕙	
due QC measurements	2.	Select a control and place the tube. Repeat until there are no controls left on the screen.	_	n/a	
	3.	Start the QC measurement.		$\bigcirc$	
	4.	Validate the QC results.		Workplace > QC Status >	
	5.	Remove the controls.			

 Table B-1
 Overview of the daily operation tasks

	Task	Steps	Navigation	
			With wizard	As individual steps
	Finishing the shift	1. Check for unfinished orders.		Workplace > Orders
				Choose 🍸 > Not Finished
		2. Check for non-validated results.		Workplace > Result Review
				Choose Y > Not Accepted
		3. Check for non-transmitted results.		Workplace > Result Review
		(If working with a host system only.)		Choose 🍸 > Not Sent to Host
		4. Start the <b>End Shift</b> wizard.	Overview > ]	End Shift
		5. Perform the daily backup.		Utilities > Export > Database
		6. Export the full results.		Utilities > Export > Results
		7. Clean up the database.		Workplace > Orders > Θ
				Workplace > Result Review> Θ
				Workplace > QC Status > $\bigcirc$
				Workplace > QC History > Θ
			_	Workplace > Calibrations > $\bigcirc$
		8. Perform the maintenance actions that are due.		Utilities > Maintenance
		9. Replace cuvettes.	_	Overview > T
		10. Check the external fluid containers.	_	Overview > 🚱 > 🛍
		11. Remove the reagent disk (if last shift).		Overview > 🔟 > 🖭
		12. Log off the system.		<b>Overview</b> > button with your user name
		13. Shut down the system and switch off the instrument (if last shift).	n/a	n/a

Table B-1

Overview of the daily operation tasks

#### Working with a host system

Communication with the instrument is defined during installation.

• For an overview on working with a host system, see *Host connectivity* on page A-32.

Starting the shift

## Starting the shift

Starting the shift includes the tasks from switching on the instrument up to the moment when you are logged on.

The various tasks are described in the order in which they should be performed.

#### Switching on the instrument


#### Preconditions

Before you switch on the instrument, make sure the following preconditions are met:

- All covers are closed.
- The external fluid containers are connected.
- The sample area is empty.

#### To switch on the instrument

**1** Press down **I** on the power toggle switch (A).

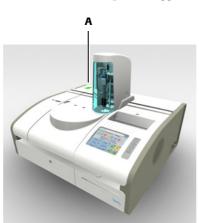

The system performs internal checks and routines.

The startup phase may take a few minutes. During this time, a splash screen is displayed.



Do not use the screen until the system is in Standby status.

When the system is ready to log on, the **Overview** tab is displayed.

**Roche Diagnostics** 

#### Logging on to the system


-χ-

#### ► To log on the system

#### 1 Press Log On.

A screen is displayed for typing your user name.

• For information on typing text, see *Typing text* on page A-74.

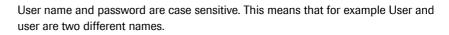

**2** Type your user name.

Use characters of the alphanumeric keyboard (upper and lower case).

As soon as you type the first characters of your user name, the system looks for a name that starts with these letters and, if it finds one, displays the complete name.

**3** Press 🔘.

A screen is displayed for typing your password.

4 Type your password.

Use characters of the alphanumeric keyboard (upper and lower case).

(For security reasons, you cannot see the actual characters when typing.)

**5** Press  $\oslash$ .

The **Overview** tab is displayed again.

Preparing the system

# Preparing the system

Before you can start analyzing samples, you need to prepare the system. Preparing the system includes both user actions and actions performed automatically by the system, such as heating up the cuvette ring or cooling down the reagent cooler. The process is designed in such a manner that user intervention is mainly required at the beginning and at the end of the **Prepare** phase. This way, the operator does not have to stand by the instrument the whole time.

The easiest way of performing the daily preparation tasks is to follow the **Prepare** wizard.



#### Preventing intervention during sample processing

Performing conscientiously all preparation steps reduces considerably the chances of you having to intervene during routine operation.

#### Follow the steps suggested by the wizard

You are guided through the preparation process by the **Prepare** wizard. This sequence of screens shows you which actions you should perform and in which order.

#### Skipping a step

You may skip a step. Before you do so, remind yourself of the consequences. In many cases, skipping a step simply means postponing the task to a more convenient time. In other cases, skipping a step could prevent the system from performing an analysis; for example if you chose not to replace a reagent that is low, a test that uses this reagent may not have enough reagent.

# Short guide

The following table provides an overview of the steps that make up the preparation process.

Ste	р	User action				
1	Start the Prepare wizard.	Standby     10.12,2007 13:46       Overview     Workplace     Utilities       1     2     3     4     5     6     7     8       0     Tests / 0     Blocked     0     Results / 0     Validate       STAT     Order     0     0     0       Tests     admin     0     0     0       Prepare     End Shift     000     0	1.	On the <b>Overview</b> tab, press <b>Prepare</b> .		
2	Check the status of the external containers.	Standby 10.12.2007 12:27 () Check Bottles Press fluid button when done. Water Waste Cleaner 100% ()	2. 3.	Refill the water container, if required, and con- firm by pressing the water button. (If you refill the water, empty the waste as well.) Empty the waste container, if required, and confirm by pressing the waste button. Replace the cleaner bottle, if required, and con- firm by pressing the cleaner button. Press D to proceed to the next stage in the <b>Prepare</b> wizard.		
3	Perform the maintenance actions that are due.	Standby       10.12.2007 4:22         Prep>Check Maintenance         Clean Water and Waste       10.12.         Replace Photometer Lamp       10.12.         Clean Probe Manually       10.12.         Deproteinize Probe       11.12.         Clean Reagent Disk       07.01.         Replace Water Inlet Filter       25.02.         Clean Housing       02.06.	2.	Check which maintenance actions are due. Perform the maintenance actions. Perform at least all red maintenance actions. Press 💓 to proceed to the next stage in the <b>Prepare</b> wizard.		
4	Prepare the reagent disk.		1. 2. 3. 4.	Open the main cover. Remove the disk from its container. Place the disk on the instrument. Close the main cover. A screen is displayed, showing the reagent sta- tus.		

Table B-2

Steps for preparing the system

# 5 Daily operation

Preparing the system

Ste	p		User action		
5	Prepare the reagents.	Standby       10.12.2007 12:34         Prep:Disk 2         ALTL       ASTL       BARBY       CA         CRPLX       D-DI       GLUC2       My_MQ       THG         UA2       CLEANY       CLEAN       NACL       PYP         UA2       CLEANY       CLEAN       NACL       PYP         Image: Comparison of the standard standar	<ol> <li>Deal with at least all red and yellow buttons.</li> <li>Press a reagent button to check the details.</li> <li>Replace the empty reagents.</li> <li>Press 10.</li> <li>Open the main cover</li> <li>Remove the bottle.</li> <li>Press 10 to confirm.</li> <li>Remove the remaining bottles of the set.</li> <li>Press 10.</li> <li>Place the new bottle.</li> <li>Press 10 to confirm.</li> <li>Load the remaining bottles of the set.</li> <li>Press 11. Press 12 to proceed to the next stage in the Prepare wizard.</li> </ol>		
6	Prepare the cuvettes.	Standby       10.12.2007 4:05         Prep>Cuvettes         Select segment to replace         0       0	<ul> <li>Replace all red cuvette segments.</li> <li>Press the cuvette button.</li> <li>Open the main cover.</li> <li>Replace the cuvette segment.</li> <li>Press ITU to confirm the replacement.</li> <li>Replace the remaining segments that need replacing.</li> <li>Press D to proceed to the next stage in the Prepare wizard.</li> </ul>		
7	Perform reagent mixing.	Standby         23.06.2009 19:44         Image: Control of the standard standar	<ol> <li>Select the test(s).</li> <li>Press  .</li> </ol>		
8	Perform calibrations.	Standby         10.12.2007 4:25 (?)           Prep>Select Cal         (*)           ROUT*         STAT*         DEV*         DRUG*         5         6           ALTL         ALTPL         ASTL         ASTPL         CA           CH02A         GLU2         UA2         (*)         (*)           (*)         **         (*)         (*)         (*)	<ol> <li>Perform the calibrations.</li> <li>Check the test selection.</li> <li>Press .</li> <li>Place the calibrators.</li> <li>On the placement screen, press .</li> <li>Press .</li> <li>Validate the results.</li> </ol>		

 Table B-2
 Steps for preparing the system

# **Starting the Prepare wizard**

- ► To start the preparation process
  - 1 Choose Overview > Prepare.


# Checking the status of the external fluid containers



Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:

- Injury through reagents and other working solutions on page B-5.
- Infection by biohazardous materials on page B-5.
- Infection by waste solution on page B-5.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page B-5.

The external fluid containers are placed on a rack.

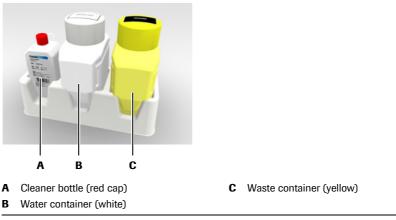



The status of all three containers is displayed on the same screen:

Standby Chack Battles	10.12.2	007 12:27 🕐		
Check Bottles				
Press fluid button when done.				
Water Full	Waste Empty	Cleaner 100%		

Figure B-2

# To check the water container

**1** Check the water button.

There is no water left. (A fluid sensor detected that there is no water coming from the water container.)
You need to refill the water now. No new tests can be processed, started tests may have to be restarted after refill.
The water container was last filled up more than one day ago. You should check the water level and refill if required.
No action is required.

`Q́

The system does not monitor the filling level. It is up to the operator to fill up the water container regularly.

**2** Refill the water, if required.





# Danger of poor measurement quality due to inadequate water quality

Inadequate water quality may lead to incorrect results. Always use purified water of the quality specified in section *Technical specifications*.

- Remove the tubing adapter from the white water container and place it on a clean surface.
- Refill the bottle with purified water.
- Insert the tubing adapter. Press down firmly.
- **3** On the screen that shows the container status, press the water button to confirm that you have refilled the water.
- **4** Roche recommend to empty the waste container whenever you refill the water container.
  - See *To check the waste container* on page B-17.



When refilling the water without using the Prepare wizard:

Choose Overview > 🎯 > 🛍.

# To check the waste container


**1** Check the waste button.

The waste container was last emptied more than one day ago.
You should check the waste level and empty if required.
No action is required.



The system does not monitor the filling level. It is up to the operator to empty the waste container regularly.

2 Empty the waste container, if required.

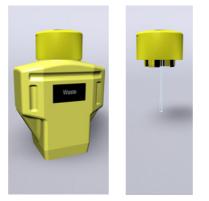

• Have the spare waste container ready.

The system periodically performs wash actions. Therefore, an external waste container must be connected at all times.

- Remove the tubing adapter from the yellow waste container and insert it in the spare container.
- Remove the bottle from the rack and place it on a firm and even surface.
- Place the spare container on the rack.
- Check that the tubing adapter is properly inserted.
- Empty the removed waste container. Treat the fluid as biohazardous waste.
- Rinse the container with water and leave it to dry.
- **3** On the screen that shows the container status, press the waste button to confirm that you have emptied the container.



When emptying the waste without using the Prepare wizard:

Choose **Overview** > 🐼 > 🛍.

# To check the cleaner bottle

**1** Check the cleaner button.

A fluid sensor detected that there is no cleaner in the instrument tubing.
You need to replace the cleaner bottle now. No new tests can be processed, started tests may have to be restarted after replacing the bottle.
The cleaner level is down to 10% or lower. See the % indication.
No action is required.

**2** Replace the cleaner bottle, if required.



- Remove the tubing adapter from the cleaner bottle and place it on a clean noncorrosive surface.
- Dispose of the bottle.
- Remove the cap from the new bottle.
- Place the new bottle on the rack.
- Insert the tubing adapter and press it down firmly.
- **3** On the screen that shows the container status, press the cleaner button to confirm that you have exchanged the cleaner bottle.

Level monitoring for the cleaner bottle is based on the number of cleaning and pipetting actions that were performed. The counter is reset when you press the cleaner button. Make sure you only press the button when you actually have replaced the bottle.

- **4** Press **>>** to proceed to the next stage in the **Prepare** wizard.



# When replacing the cleaner bottle without using the Prepare wizard:

Choose **Overview** > 🎯 > 🛍.

# **Performing maintenance actions**

Maintenance actions need to be performed periodically or after certain events.

To ensure the smooth running of the system, you should perform maintenance actions either as part of the preparation or the end of shift activities.

For information on scheduling maintenance actions, see Scheduling maintenance actions on page B-158.


Make sure that you have read and understood section Safety information on page C-8.

If you interrupt the performing of a maintenance action that was due, its status will remain due, and you need to fully re-perform the action later.

# To perform maintenance actions

Checking what maintenance actions are due

1 Check the colors of the maintenance actions.

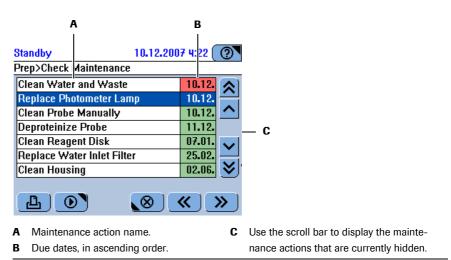

#### Figure B-3

The maintenance actions are sorted by the maintenance date by when they should be performed. Use these dates for planning the maintenance actions, for example for ordering the required materials.

Interpreting the colors	The defined maintenance interval has expired. Perform this maintenance action now. The system may not be able to operate until this action is performed.
	This maintenance action should be performed on the next major maintenance day.
	No action is currently required.
	This maintenance action is selected.

Performing the maintenance actions



**2** Select the maintenance action you want to perform.

The selected line turns blue.

# Incorrect results or processing stop due to skipping maintenance actions

Not performing maintenance actions that are due may lead to situations where the system cannot continue processing orders, or it may lead to incorrect results. If at all possible, perform the maintenance actions when they are due.

3 Press 💽.

The maintenance definition screen is displayed.

- For information on how to perform individual maintenance actions, see Chapter 8 General maintenance.
- **4** Press 🔘 to start the action.
- **5** Follow the instructions on the screen.
- 6 Perform the next maintenance action that is due.
- 7 When you have finished performing maintenance actions, press ≫ to proceed to the next stage in the wizard.



When performing maintenance actions without using the Prepare wizard: Choose Utilities > Maintenance.

# Preparing the reagent disk

The reagent disk holds the reagent and diluent bottles. During periods when you are not performing tests, for example at night or during holidays, the disk is placed in a reagent disk container and stored in a refrigerated place. Reagents should be stored at temperatures in the range of +2 to +8°C.



Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:

- Injury through reagents and other working solutions on page B-5.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page B-5.



### Spillage through tipping reagent disk

The reagent disk container can slip off or tip over if it is not placed on an even horizontal surface.

When storing the reagent disk container, make sure to place it on a firm, even, horizontal surface that is easily accessible.

When handling the reagent disk, make sure not to tilt it.

#### To load the reagent disk

- 1 Get the reagent disk from the refrigerator and remove it from its container.
- **2** Open the main cover.

A screen is displayed, asking you to insert the disk.

**3** Place the reagent disk in the reagent cooler.

Make sure the reagent port faces the front and align the cut-outs with their counterparts on the reagent cooler.

The system automatically detects that a disk was inserted.

A screen is displayed, asking you to close the main cover.

**4** Close the main cover.

At this stage, the system identifies the disk.

A screen is displayed that shows the status of the reagent sets.



When handling the reagent disk without using the Prepare wizard: Choose Overview > ()).

# Preparing the reagents

Reagents are handled in sets. A set consists of up to three reagents. If for example one reagent bottle of a set is empty, all reagents of the set need to be replaced. The system only uses reagents of complete sets.

On the screen, each set is represented by a separate button.



### Use reagents of the same lot, if possible

When replacing a new reagent set, try to use one of the same lot as the old one. This way you avoid having to perform a calibration.

### **Diluent and cleaner bottles**

An extra diluent or cleaner bottle is considered one set. Extra cleaners and diluents are handled in the same type of bottles that are used for reagents. Their handling works the same way. In the following, the process is described for reagents.



Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:

- Injury through reagents and other working solutions on page B-5.
- Infection by biohazardous materials on page B-5.
- Loss of sight on page B-6.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page B-5.



#### Incorrect results due to declining reagent quality

Due to the declining quality of reagents, the quality of measurement may decline. This may cause incorrect results.

Make sure to use reagents whose expiration date has not passed yet.

#### Incorrect results due to reagent dilution through condensation

In conditions of high temperature and humidity, condensation can build up inside the reagent bottles. This leads to dilution of the reagent.

In such conditions and when not performing tests, especially if you observe condensation on the reagent disk or in the reagent cooler, be sure to remove the reagent disk from the instrument and place it in the reagent disk container. Close the container with its lid and place it in a refrigerator.

#### Incorrect results due to condensation in the reagent cooler

In conditions of high temperature and humidity, condensation can build up in the reagent cooler. Water can spill on the analyzer unit and get into the cuvettes when removing the reagent disk.

In conditions of high temperature and humidity, be sure to periodically wipe up the condensation water in the reagent cooler.

#### Incorrect results due to inappropriate reagent handling

Removing and loading reagents while the reagent disk is outside the instrument may lead to inconsistencies between the recorded and the physically loaded reagents and consequently to incorrect results.

Always remove and load reagents while the reagent disk is on the instrument and by using the software supported procedures.

NOTICE

#### Damage to the reagent disk

The reagent disk is designed to handle reagents while it is loaded on the instrument. The cover is equipped with a locking mechanism.

Always remove and load reagents while the reagent disk is on the instrument and by using the software supported procedures.

#### Checking the status of reagent sets

There are two main ways of checking the status of reagent sets:

- To get a general overview of the status of all reagents that are defined on any disk, use **Utilities** > **Inventory**.
  - For information on deleting reagent sets from the Inventory list, see Deleting bottle sets from the Inventory list on page B-120.
- To check the status of individual reagent sets, use Overview > Tests > test icon.

### ► To gain a general overview of the currently defined reagent sets

1 Choose Utilities > Inventory.

The **Bottle sets** list is displayed. It contains all bottle sets that are defined on any of the disks used on this **Cedex Bio** instrument, including sets that were removed from the disk, as long as they were not empty or their expiration date has not passed yet.

			1	A	В	
-	Standby Bottle Se Set A1C-2 A1C-2 ALBT2 ALP2S ALTL	ts 60165702 60165702 69409701 67505751 60915701	23.06.2009	2 0:01 00 3 00 3		— C
	ALTL	60915701 60630001	08.06.2009 90	6 1	R	U
		20030001		<u> </u>	⊗	
ŀ		per of tests le			С	<b>R</b> indicates that the set was removed from
E	define	0	on which the se	et is		the disk.

### Figure B-4

**2** Select a set and press  $\square$ .

A dialog box is displayed that contains detailed set information.


3 Press M.

A dialog box is displayed that contains set, calibration, and QC information.

- **4** Press 🛞 to close the dialog boxes until the **Utilities** tab is displayed.

# ► To check the status of individual reagent sets

- **1** Choose **Overview** > **(11)**.
- **2** Check the colors of the reagent set buttons.

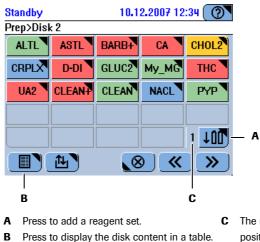

The number indicates the number of free Press to display the disk content in a table. positions on the disk.

### Figure B-5

Interpreting the colors

Color	Meaning	Possible actions
	No tests can be performed with this reagent set.	
	The number of available tests is 0.	Replace the reagent set.
	The set is incomplete.	Add the missing reagent.
	The test needs calibrating.	Perform calibration.
	There are fewer than 10% of tests left for this set.	Load a new reagent set as soon as pos- sible.
	The expiration date has passed.	
	A QC is due or its result has not been accepted.	Perform QC as soon as possible or accept the result
	A more recent version of the applica- tion has been imported.	Install the new version of the applica- tion to start working with it.
	This development channel test is run	Activate the extra wash cycle.
	without extra wash cycles.	Roche strongly recommends to always use extra wash cycles with develop- ment channel applications and also to always load extra cleaner when tests with extra wash cycles are used.
	Ready for use.	No action is currently required.
	There is no application that uses this reagent set.	Add the missing application.
	A required diluent or cleaner is not on board.	Load the missing fluid.

Displaying detailed information on a reagent set **3** Press a reagent button.

A screen is displayed that shows the details of the selected reagent.

Standby	10.12.2007 12:30 🕐	
Overview>Disk 2>BA	IRB	
Remaining tests	0	
Expiration date	07.06.2008	
Lot ID	DCCI-TS1	
Serial number	11/11	
Bottle	e set is empty	_ A
L	11 200 🚫	
A Status description	1	

Figure B-6

# **Preparing reagents**

During preparation, you typically replace reagents that are empty or whose expiration date has passed, or you add new reagents.



#### Incorrect results due to impurities and carryover

Traces of analytes or reagents may be carried over one test to the next when reusing bottle caps.

Do not remove reagent bottles that are not empty with the purpose of loading them again later.

### To prepare the reagents

1 Choose Overview > Tests.

*Preparing the reagent set* **2** Obtain the reagent sets you want to add or replace.

When replacing a reagent set, try to use one of the same lot as the old one. (This way, if you work with the calibration sequence **Each Lot and Interval**, you avoid having to perform a calibration.)

### *Removing a reagent set* **3** Press **1**.


- **4** Select the reagent you want to remove.
- **5** Press **1**.

The system moves the first bottle of the set to the reagent port.

6 Wait for the main cover LED to turn green.

A message is displayed, asking you to open the main cover and remove the bottle.

7 Open the main cover and remove the bottle.

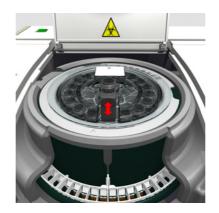

**8** Press 🖉 to confirm the removal.

The system does not check whether you have actually removed the bottle.

# **9** Do one of the following:

lf	Do this		
There is another bottle belonging to the set:	The system moves this bottle to the reagent port. A screen is displayed, asking you to remove the bottle.		
	<ol> <li>Remove the bottle.</li> <li>Press O to confirm the removal.</li> <li>Start adding the new reagent set.</li> </ol>		
All bottles of the set are removed:	1. Start adding the new reagent set.		

The screen with the reagent set buttons is displayed again. The button for the removed set is no longer present. If you did not remove all bottles of the set, the reagent set would be disabled, its button would be red.

Addin <u>reagent</u> set

Before you add a set, check the number to the left of the button, it indicates the number of free positions on the disk. If this number is smaller than the number of bottles in the set you need to add, remove a set first. For the number of bottles in a set refer to the respective package insert.

- **10** Open the main cover if required.
- 11 Press 100.

A screen is displayed, asking you to scan the bottle barcode.

**12** Scan the barcode on the bottle.



# When scanning, bear this in mind:

- Scanning its barcode is the only way of identifying a reagent, diluent, and cleaner bottle.
- If there are not enough free positions for all bottles of the set, a message will inform you.
- You cannot re-insert empty bottles that were previously removed from the system. If you scan such bottle, a message will inform you of this fact.

A screen is displayed, asking you to place the bottle on the reagent disk.



WARNING

### Time limit for placing reagent bottles

- The system assumes that you place the bottle that you just scanned.
- You need to place the reagent bottle on the reagent disk within 15 seconds of scanning the reagent barcode. Failing to confirm the placement within 15 seconds cancels the current identification process. You will be asked to scan the bottle again.

#### Incorrect results due to not placing the identified reagent

The system assumes that the operator places the reagent that was just identified. Failing to do so might lead to wrong results.

- **13** Remove the cap from the bottle and place it in the reagent bay.
- **14** Press 🞯 to confirm the insertion.

If you confirm without placing the bottle, the system assumes that the bottle is placed.

If you press 🛞 after placing the bottle, the position is deemed empty.

When the first reagent bottle is loaded, the reagent set is defined. From this moment on, the reagents are handled as part of the set. You no longer handle them as individual reagents.

# **15** Do one of the following:

lf	Do this
There is another bottle	A screen is displayed, asking you to scan the bottle.
belonging to the set:	1. Scan the bottle
	2. Insert the bottle.
	3. Press 🖉 to confirm the insertion.
All bottles of the set are	1. Close the main cover.
inserted:	On the screen with the reagent set buttons, the button
	for the new set is now present.

**16** On the screen with the reagent set buttons, press the button for the set you just inserted.

A screen is displayed that shows details of the set.

The status description points to the action you need to take. If you just inserted a reagent set, you may have to perform the initial calibration and QC.

Standby	10.12.2007 12:29 🕐	
Overview>Disk 2>CA		
Remaining tests	94	
Expiration date	02.03.2008	
Lot ID	12345678	
Serial number	24283/24293	
Missing	ı calibration: CA	_ A
<b>L</b>	11 ?00 🚫	
A Status description	1	

Figure B-7

# 17 Press 🛞.

**18** When you have finished adding and replacing reagent sets, press *≫* to proceed to the next stage in the **Prepare** wizard.



When handling the reagent sets without using the Prepare wizard: Choose **Overview** > (11).

# **Preparing cuvettes**

Cuvettes are supplied and handled in cuvette segments. Each segment contains ten cuvettes. The segments are placed on the cuvette ring of the rotor.

For information on cuvette segments, see *Cuvette segments* on page A-55.
 For information on the cuvette ring, see *Cuvette ring* on page A-59.

Each segment on the cuvette ring is represented by a button on the screen.

Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:


- Personal injury or damage to the analyzer due to contact with instrument mechanism on page B-5.
- Infection by waste solution on page B-5.

### To prepare the cuvettes

### *Checking the cuvette status*

1 Check the color of the cuvette segment buttons.

You can refresh the cuvette screen by pressing O. (If there were pending orders, this would also start their processing.)

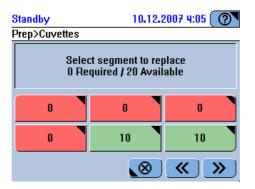

Interpreting the screen

The number in the button indicates how many cuvettes are free to be used.

Icon	Meaning	Possible actions
	All cuvettes are used.	Replace the segment.
	Up to two cuvettes are free to be used.	
	More than two cuvettes are free to be used.	No action is currently required.

#### **Roche Diagnostics**

Replacing cuvette segments	2	Press a segment button.		
Reputeting envente segments	-	e	at the system is ready for cuvette handling.	
	•			
	3	Wait for the main cover LED to turn gr	een.	
	4	Open the main cover.		
	5	Remove the segment and treat it as biohazardous waste.		
		Incorrect results due to scratched or s	soiled cuvettes	
		Scratches and impurities on the cuvettes distort the measurements.		
WARNING		Do not touch the cuvettes and make sure they do not touch other items when handling them.		
	6	6 Insert a new cuvette segment.		
	<b>7</b> Press <b>I</b> U to confirm the replacement.			
		(Press <i>IV</i> if you removed the segment	t without replacing it.)	
	8	Do one of the following:		
		lf	Do this	
		You want to exchange another segment:	Perform steps 2 through 7.	
		This was the last segment you wanted to exchange:	Close the main cover.	
	9	When you have finished replacing cuve stage in the <b>Prepare</b> wizard.	ettes, press $>>$ to proceed to the next	
-ò-	w	hen handling cuvette segments without	t using the Prepare wizard:	
E	Ch	boose <b>Overview</b> > T. The system must b	pe in Standby status.	

# Performing reagent mixing

Reagents that contain Latex granules require periodic mixing. The mixing interval is part of the application definitions and cannot be changed or deleted by the user. The system checks every 30 minutes for reagent sets that require mixing.

Within the **Prepare** phase a screen is displayed that lists all reagent sets for which a mixing interval is defined.

# To perform the mixing

1 Select the reagent sets you want to mix.

If working with the **Prepare** wizard, the reagent sets that need mixing are automatically selected.


**2** Press **O**.

Mixing starts. A screen is displayed informing you about the progress of the mixing action.

(If there is more than one bottle in the set with a mixing interval defined, it will be mixed as well, irrespective of being due or not.)

**3** When mixing is finished, press  $\bigotimes$  to close the screen.

# **Roche Diagnostics**

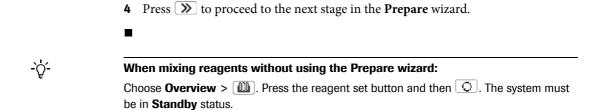

# Performing the calibrations (Prepare phase)

At this stage, the system checks for all calibrations that are due.

For an overview on calibration, see *Calibration* on page A-26.
 For performing individual calibrations, see *Performing calibrations* on page B-59.

By default, all calibrations that are due or will be due within the forecast period are taken into account during the **Prepare** phase.



Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:

- Injury through reagents and other working solutions on page B-5.
- Infection by biohazardous materials on page B-5.



### Skin inflammation caused by reagents

Direct contact with reagents may cause skin irritation, inflammation, or burns.

When handling reagents, be sure to wear protective equipment and observe the cautions given in the package insert.

#### Incorrect results due to expired calibration

Calibrations are performed to compensate for changes over time in reagents and in the measurement systems. Failing to perform calibrations when they are due may lead to incorrect results.

Make sure to perform calibrations when they are due.

### Incorrect results due to wrong tube placement

Be sure to place the calibrators in the specified positions.

#### To perform the calibrations ►


A screen is displayed that shows all active tests.

Checking which calibrations are due

Interpreting the colors

Modifying selections

1 Check the colors of the test buttons.

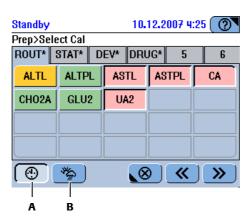

- A Press to cancel the selection of all tests that need calibrating
- Press to cancel the selection of all tests that need calibrating within the forecast period.

Color	Meaning	Possible actions
	No sample tests can be measured. Possible reasons:	
	Initial calibration is required.	Perform the calibration.
	The calibration failed.	Check the flag to find out why the cal- ibration failed. Repeat the calibration if necessary.
	The calibration was successfully per- formed, but its results have not been accepted yet.	Validate the calibration results.
	QC is due.	Perform QC as soon as possible.
	The calibration is OK. Its results have been accepted.	No action is currently required.
	The calibration is OK. It applies to a reagent set that is not active.	No action is currently required.

В

period are automatically selected. You can cancel these selections by pressing 🕲, 🤔 or a test button.


If you do not perform a calibration that is due, the affected test is blocked; you cannot perform this test.

# Performing calibrations

-Ò́-

**2** On the screen that shows the tests, press  $\gg$ .

A screen is displayed that lists all required calibrators, and it shows on which position to place them.

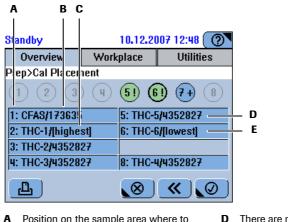

A Position on the sample area where to place the calibrator.

Calibrator with highest concentration

- There are not enough empty positions, the system suggests using occupied positions.
- **E** Calibrator with lowest concentration

### Figure B-9

Lot number

R

С

- The system first uses the free positions, if there are not enough free positions, occupied positions are suggested. For these you would have to replace the currently loaded tubes with calibrator tubes.
- With absorbance tests that require several calibrators, the calibrators are placed according to their concentration, starting with the highest concentration.

Calibrators with several concentrations are listed according to their concentration, starting with the one with the highest concentration. They are displayed as follows:

Number of calibrators	Displayed information	
1	Position, name, lot number	
2	<ol> <li>Position, name, highest</li> <li>Position, name, lot number</li> </ol>	
More than 2	<ol> <li>Position, name, highest</li> <li>Position, name, lot number</li> <li></li> <li>Position, name, lowest</li> </ol>	

- **3** Prepare the calibrators.
- 4 Place the calibrators on the sample area positions indicated on the placement list.
- **5** Press 🕖 to confirm the placement.
- **6** Press 0 to start the calibration.
- 7 Choose Workplace > Calibrations to check the status of the calibrations.
- *Validating the results* **8** Validate the results.

• See Validating calibration results on page B-64.



When performing calibrations without using the Prepare wizard: Choose Workplace > Calibrations >  $\textcircled{\oplus}$ .

End of **Prepare** phase

The instrument is now ready for analyzing samples.

You may want to perform QC before analyzing samples.

See *Performing QC* on page B-68.

# **Safety information**



Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:

- Infection by biohazardous materials on page B-5.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page B-5.
- Loss of sight on page B-6.



### Incorrect results due to inadequate sample preparation

Specimen containing clots may obstruct the probe. Specimen containing bubbles or foam may cause level detection errors and air pipetting. Consequently, incorrect results may be generated.

Take adequate care when preparing the samples.

### Incorrect results due to insufficient fluid

Insufficient fluid may lead to inaccurate pipetting and consequently to incorrect results.

Always fill the tubes with enough fluid that at least the defined dead volume of fluid is left when pipetting is complete.

• See *Tubes* on page A-53.



# Time limit for placing samples

The system assumes that you place the sample tube that you just identified.

You need to place the sample on the sample area within 10 seconds of confirming the test selection. Failing to place the sample tube within 10 seconds cancels the current identification process. You will be asked to identify the sample tube again.

# Short guide

The following table gives an overview of the tasks you typically perform when analyzing samples.

Ste	p		User action
1	Identify the sample.	Standby     10.12.2007 12:23       Overview     Workplace       Utilities       Order>Identify Sample       Scan the next sample barcode or type the information manually.       Oli Si No.	If working with sample barcodes: 1. Scan the barcode using the barcode scanner.
		Standby         10.12.2007 12:24         Image: Constraint of the state of the st	<ul> <li>If working without sample barcodes, type the sample ID by using the screen keyboards.</li> <li>1. Press .</li> <li>2. Type the ID.</li> <li>3. Press .</li> </ul>
2	Select the tests, ratios and profiles. This step is not required if the workflow setting <b>Working</b> <b>Mode</b> is <b>Host</b> . (Utilities > Con- figuration > Workflow.)	Standby     25.06,2009 9:58       D12>Select Tests       ROUT*     STAT       DEV     DRUG       5     6       ISE-I     P Dbt       P LEB     Prof1       RHbA1       Rati1     A1W2D       ALBU2     ALTL       ASTL     ASTPL       CA     CL-I       CPDI     GLU2       K-I     NA-I	<ol> <li>Press a tab to select the test panel (if required).</li> <li>Press an active test, ratio or profile to select it. Press a selected item to cancel its selection.</li> <li>Press D.</li> </ol>
3	Place the sample.	Standby 10.12.2007 12:25 ( 113>Select Tests>Place Sample Place sample 113 on the instrument while the sample LED is green. To prevent misplacement, this workflow is aborted if no tube placement is detected within 10 seconds.	<ol> <li>Place the sample on any free position on the sample area.</li> <li>You need to place the sample within 10 seconds of pressing . Failing to place the sample tube within 10 seconds cancels the current identification process. You will again be asked to identify the sample tube.</li> </ol>

Table B-3 St

Steps for analyzing samples

# 5 Daily operation

Analyzing samples

Ste	p		User action
5	Validate the results.	Standby         10.12.2007 3:41         Image: Constraint of the state of the sta	<ol> <li>Choose Workplace &gt; Results.</li> <li>Validate the result (You can accept the result or perform the test again.)</li> </ol>
6	Remove the samples.	Standby     10.12.2007 12:50       Overview     Workplace       Utilities       1     2       3     4       5     6       7     8       45     5       75     7       8     35       75     8       75     7       76     7       8     7       76     8       77     7       76     7       8     7       76     8       77     7       76     7       8     7       76     8       77     8       76     7       8     7       76     8       77     8       77     8       77     8       77     8       77     8       78     8       79     8       70     9       70     9       77     8       77     8       77     8       77     8       77     8       77     8       78     8       79     8	<ol> <li>Press the <b>Overview</b> tab.</li> <li>Check the status of the sample tube button that corresponds to the sample you want to remove.</li> <li>If the sample tube button is green, remove the tube.</li> </ol>

Table B-3

Steps for analyzing samples

# System setup and order definition workflow

	The workflow for defining orders depends on a few configuration settings:
$\rightarrow$	Utilities > Configuration > Workflow
Working Mode	The Working Mode defines the way you select tests when defining orders.
	• Use Manual if you use the Cedex Bio instrument as a stand-alone system.
	During order definition, the test selection screen is displayed, allowing you to select the tests or to make changes to the selection.
	• Use <b>Host</b> if the <b>Cedex Bio</b> instrument is connected to a host computer.
	The orders are automatically defined, the test selection screen is not displayed. After identifying the sample, you are asked to place the sample.
Sample Barcode	Use <b>On</b> if you exclusively work with barcoded sample tubes. When defining orders, the screen for typing the sample ID manually is skipped.
Order Mode	The order mode reflects the way in which you organize the tests on the test selection screen.
	Use <b>Easy</b> if the reagents fit on one or two reagent disks and you work with one test panel on the screen (You can fit up to 25 tests and profiles on this panel).
	Use <b>Full</b> if you distribute the reagents across several (up to eight) reagent disks and if you predominantly work with specific groups of tests. You can define up to 20 tests and profiles on each panel (tab).
	• See Assigning tests to test tabs on page B-119.
Sample ID Handling	The <i>sample ID</i> is an identifier of up to 23 alphanumeric characters that is unique within a whole organization. It cannot be changed once the order is saved. This ID is used for communication with a laboratory information system.
	The <i>order ID</i> is an identifier of up to 23 alphanumeric characters that is unique within the laboratory. In practice, this could for example be an autonumber followed by the tested badges description.
-, Å	Because there is limited space when displaying lists on screen, Roche recommend to limit the IDs to 13 characters.
	If you work with one ID only, the sample ID is used as the order ID as well. There exists a fixed relation between the sample and order IDs. Make sure this ID is unique within the <b>Cedex Bio</b> instrument.
	• Use <b>Order ID</b> = <b>Sample ID</b> to have the system automatically define the order ID identical to the sample ID.
	• Use <b>Independent IDs</b> to define the sample and order IDs independently of each other.
	<ul> <li>Use Grouped Sample ID to use identical order and sample IDs as a constant part of any order for a give sample, and to let the system automatically append a running number to the ID for each order.</li> </ul>
Auto Order-ID	Use <b>On</b> to automatically increment the order ID number by one whenever you define a new order. (You only need to define the number of the first order of your shift.)
	If you use <b>On</b> for this feature, use <b>Off</b> for <b>Order ID</b> = <b>Sample ID</b> .

**Roche Diagnostics** 

# **Defining orders**

Preconditions • All prepa

- All preparation tasks are complete.
  - The required tests are installed on the system and ready for use (calibration and QC performed).
  - The system status is either Standby or Operating.
  - At least one sample position on the sample area is free.
  - There can be only one order per test and sample.
  - Possible post-dilution is predefined in the application definition.



**Incorrect results due to declining sample quality** Evaporation of sample fluid may lead to incorrect results. In ambient temperatures of more than 25°C, be sure to start processing immediately after placing the sample and defining the order. When processing of the order is finished, be sure to immediately remove the sample from the sample area.



# Incorrect results due to insufficient fluid

Insufficient fluid may lead to inaccurate pipetting and consequently to incorrect results.

Always fill the tubes with enough fluid that at least the defined dead volume of fluid is left when pipetting is complete.

• See *Tubes* on page A-53.

### Incorrect results due to inappropriate tube and cup placement

Inappropriate tube and cup placement may lead to inaccurate pipetting and consequently to incorrect results.

Make sure that the primary tubes are placed centrally and perfectly vertically in the holders in the sample area and that they are inserted firmly.

Make sure that secondary tubes are placed centrally on the primary tubes and that they rest fully on them.

# *NOTICE* Probe damage and instrument malfunction due to not removing primary tube caps

The probe is not designed to pierce tube caps. It can get damaged when trying to pierce tube caps, which can lead to instrument malfunction.

Always remove the caps of primary tubes before placing them on the instrument.

# **Defining routine orders**

### ► To define an order

# 1 Choose Overview > Order.


A screen is displayed, asking you to identify the sample.

**2** Identify the sample. Do one of the following:

lf	Do this
You work with sample barcodes:	Use the barcode scanner to scan the bar- code.
You do not work with sample barcodes or if the barcode could not be read for some rea- son:	
	Note: This step is only required if the configuration setting <b>Order ID = Sam ple ID</b> is not selected.

A screen is displayed that shows all active tests, profiles, and ratios. Note that all applications of a profile or ratio must be installed for it to be active.

(If the workflow setting **Working Mode** is **Host**, this screen will not be displayed. **Utilities** > **Configuration** > **Workflow**.)

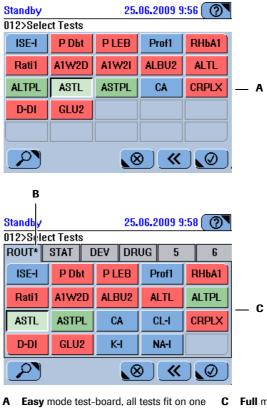

**Full** mode test-board. The tests are distributed across several panels (tabs).

Figure B-10

selected tests.

screen.

Tabs marked with an asterisk contain

В

# **5 Daily operation**

Analyzing samples

Interpreting the colors	<ul> <li>The test is blocked for one of the following reasons:</li> <li>The calibration failed.</li> <li>Initial calibration is required.</li> <li>For the reagent set, the number of available tests is 0, or a reagent bottle is missing (incomplete reagent set).</li> </ul>
	The expiration date of the test has passed. There are only few tests left. A QC is due or its result has not been accepted. A more recent version of the application has been imported.
	The test is on board and ready for use.
	The test is defined but not on board. A required diluent or cleaner is not on board.

The tests are sorted alphabetically. Profiles precede the tests. Profiles display the color of their tests.

**3** Select one or several of the tests, profiles, and ratios.

You can select items from more than one tab. (An item can be contained in more than one tab. If it is selected in one, it will automatically be selected in the others.)

Tabs with selected items are marked with an asterisk (\*).

- **4** Press 🔘 to confirm your selection.
- **5** Wait for the Sample Area LED to turn green. (Make sure to keep clear of the sample area while the LED is yellow and blinking.)

A screen is displayed, asking you to place the sample on the instrument.

6 Place the sample on any free position on the sample area.



# Time limit for placing samples

You need to place the sample on the sample area within 10 seconds of confirming the test selection. Failing to place the sample tube within 10 seconds cancels the current identification process. You will be asked to identify the sample tube again.



# Incorrect results due to not placing the identified sample

The system assumes that the operator places the sample that was just identified. Failing to do so might lead to wrong results.

The system registers where you placed the sample and associates this position with the order you just defined.

The screen for identifying samples is displayed again. You can now start defining the next order. (If there were no free sample positions, the screen for selecting the tests would be displayed instead.)


**7** Do one of the following:

If	Do this
There is another order to be defined:	Identify the sample and repeat the order definition process.
There are no more orders to define:	Press $\textcircled{\otimes}$ to close the screen.

	8 Press the $\textcircled{0}$ global action button to start processing.
	See Starting the run on page B-44.
	If you defined the order while the system was processing orders, new orders will automatically be processed, without pressing $\textcircled{O}$ .
-`\ <u>\</u> '-	Roche strongly recommends to always load extra cleaner when tests with extra wash cycles are used.
Order of processing	When you first start processing orders, the order of the sample placed on the leftmost position on the sample area is processed first. The others follow in sequence from left to right. Once processing is in progress, the orders are processed according to the sequence in which they were defined.
	Repeats and reruns of routine orders are performed before routine orders.
	<ul> <li>See Repeating tests on page B-55.</li> <li>See Rerunning tests on page B-56.</li> </ul>
Defining STAT orders	
	Short turn around time (STAT) orders are defined in the same way as routine orders.
►	To define a STAT order
	1 On the <b>Overview</b> tab, press <b>STAT</b> .
	<b>2</b> Continue as if defining a routine order.
	• See <i>Defining routine orders</i> on page B-39.
-`位҉-	If there is no space on the sample area
F	Remove any sample whose tube button is green 🦲.
Order of processing	When you have defined a STAT order and started the run, the system reacts as follows:
	• Existing STAT orders are finished first.
	• Started tests of routine orders are finished.
	• Repeats and reruns of STAT orders are treated as normal STAT orders. (The
	order that was defined first will be performed first.)

Recognizing STAT orders on the screens

On the **Overview** tab, the sample buttons for STAT orders are marked with a wide edge.

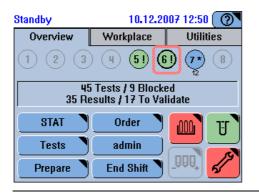



In the orders list (**Workplace** > **Orders**), STAT orders are marked with an asterisk (\*).

# Making changes to an order

You can make changes to orders that either have not yet been processed or that are fully processed. (You can add further tests to an order any time.)

The process of changing an order depends on whether the sample tube is still on the instrument or not.

Making a change to an order when the sample tube is no longer on the sample area is basically the same as defining a new order.

# ► To change an order

**1** Identify the sample.

If the sample is still on the instrument

• On the **Overview** tab, press the sample tube button of the order you want to change.

A screen is displayed, that shows details on the order and the sample.

• Press 🛞.

A screen for selecting tests is displayed.

<i>If the sample is no longer on the</i>	•	• Do one of the following:				
instrument		If the Starting point was	Perform these steps			
		Overview > Order	A screen is displayed, asking you to scan the sample or to type its ID manually.			
			1. Identify the sample either by scanning its barcode or by typing the sample ID.			
			If the same barcode was scanned before (during the current day) or you use the same sample ID as in the original order, the system recognizes the original order and displays its associated information.			
			2. Press 🔊.			
			A screen for selecting tests is displayed.			
		Workplace > Orders	A screen is displayed that lists all orders.			
			<ol> <li>Select the order you want to change.</li> <li>Press </li> </ol>			
			A screen for selecting tests is displayed.			
	2 3	Select the tests.				
	•	<ul> <li>Press an available test to select it.</li> <li>Press a selected test to cancel its selection. (Tests that are already scheduled to be performed cannot be cancelled.)</li> </ul>				
	•					
	<b>3</b> I	Press 🖉 to confirm your	selection.			
<i>If the sample is still on the instrument</i>	1	A message is displayed show	ving the position the sample is placed on.			
	•	• Press 🞯 to confirm the position				
		The screen with details on the order and the sample is displayed again.				
	•	Press $\textcircled{\begin{subarray}{c} \ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \ \$	creen.			
<i>If the sample is no longer on the instrument</i>	1	A message is displayed, aski	ing you to place the sample.			
		Place the sample on the sample area.				
		The orders list is display	-			
			•			
	4	i ne system reschedules the	order and processes it as normal.			

# Starting the run

Το	start	processing	an	order
•••	Sturt	processing	un	oraci

**1** Press the 0 global action button.

The system checks whether there are enough cuvettes and whether all required reagents are on the reagent disk.

If any of this is not the case, a screen is displayed informing you of what is missing.

*Order of processing* When you first start processing orders, the order of the sample placed on the leftmost position on the sample area is processed first. The others follow in sequence from left to right. Once processing is in progress, the orders are processed according to sequence in which they were defined.

Repeats and reruns of routine orders are performed before routine orders.

# Monitoring the analysis progress

During operation, you should regularly check the following items:

- The status of the buttons on the Overview tab
- The Alarm Monitor LED

# Checking the sample tube status

# To check the sample status

1 Press the **Overview** tab.

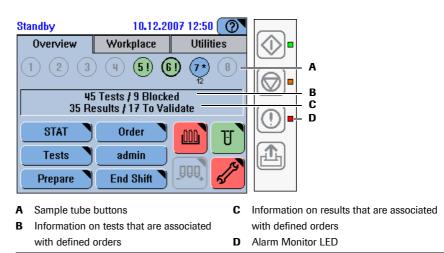


Figure B-12


Interpreting the sample tube *buttons* 

lcon	Meaning	Possible action
1	The number in the button denotes the position on the sample area.	n/a
0	A sample tube button with a wide edge symbolizes a STAT order.	n/a

lcon	Meaning	Possible action	
1	All tests are accepted.	You can remove the sample tube.	
1* 10	All tests are pipetted.	You can remove the sample tube, the time to completion disappears when you do so.	
1!	All tests are completed but not yet accepted.	Validate the results.	
1	All remaining tests are blocked because:		
	There is not enough sample fluid.	1. Remove the sample tube and add fluid.	
		Do not delete the old order!	
		<ol> <li>Scan the sample barcode or type the same sample ID as before.</li> <li>Reinsert the sample tube.</li> </ol>	
		Processing continues where it stopped.	
		(Removing a sample and then placing it again means defining a new order. The system does not remember where the sample was placed in the previous order.)	
	The sample is not identified.	<ol> <li>Remove the sample.</li> <li>Scan the sample barcode or type the same sample ID as before.</li> <li>Reinsert the sample on any posi- tion.</li> </ol>	
		(Removing a sample and then placing it again means defining a new order. The system does not remember where the sample was placed in the previous order.)	
1	There is no sample on this position.	You can place a sample tube on this position.	
1+	Tests are ordered. Processing has not yet started.	You can still cancel ordered tests and add additional tests to the order.	
<b>1</b> * 12	Tests are ordered. Processing has started.	You can no longer cancel ordered tests, but you can add additional tests	
- -	The number below the sample tube button indicates the estimated time to completion.	to the order.	
1?	The sample is identified, but no tests were ordered yet.	This should be a temporary status. No action is required.	
	If working in <b>Order Query Mode</b> : The order could not be obtained from the host.		

**2** Press the sample tube button.


A screen is displayed that shows details about the status of the sample tube.

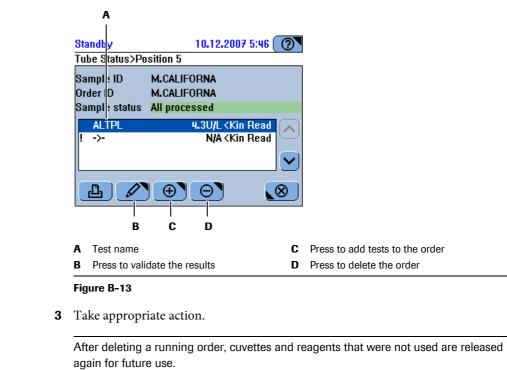

<del>`</del>گ

**Checking the instrument status buttons** 

- ► To check the instrument status
  - **1** Press the **Overview** tab.

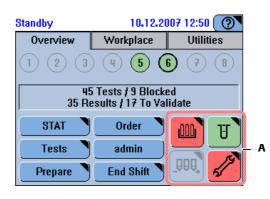

A Instrument status buttons

Figure B-14

Interpreting the instrument	Icon	Meaning	Possible action
status buttons		The reagents are OK.	No action is required.
		There is no disk on board.	Load a disk.
		Fewer than 10% of tests are left for a reagent set, or its expiration date has passed.	Load a new reagent set.
	Ī	A reagent set is not complete or a reagent is empty.	Complete the reagent set or replace it.
		A reagent set is blocked because it requires calibration or mixing.	Perform calibration or mixing.
		The disk could not be identified.	Remove the disk and make sure that one and only one ID tab is removed. Reinsert the disk.
	T	There is more than one segment available.	No action is required.
	ļ	The last available segment is in use.	Replace the used cuvette segments as soon as possible.
	ļ	There are no empty cuvettes avail- able.	Replace the cuvette segments.
	Status I	The system status button displays both the buttons of the underlying system status so by color, first priority being red, followed according to the sequence in which they a	creen. (The icons are first prioritized by yellow and green, and then
		Analyzer (main cover)	Press the button and, on the Sys- tem Status screen, check the text about the status of the main cover
	1	Reagent cooler and cuvette ring temperature	Press the button and, on the Sys- tem Status screen, check the text about the temperature.
		Sample area ventilation	Press the button and, on the Sys- tem Status screen, check the text about the ventilation status.
	ÔÔO	External fluid containers	Press the button and, on the Sys- tem Status screen, press it again to display the screen for handling the external fluid containers.
	A	Maintenance	Press the button and, on the Sys- tem Status screen, press it again to display the maintenance actions
			list.

# **Checking for alarm messages**

# ► To check for problems during processing

**1** Observe the Alarm LED.

LED	Meaning	Possible action
No color (off)	There are no unconfirmed alarm messages.	No action is required.
Yellow	There is at least one unconfirmed alarm message.	User intervention is required as soon as possible. Processing can continue for the time being.
		Check the details of the message.
Red	There is at least one unconfirmed alarm message.	Immediate user intervention is required. Processing may not be able to continue without it.
		Check the details of the message.

An acoustic signal is sounded when an alarm is created. You can adjust the volume (Utilities > Configuration > System > Volume).

**2** Take appropriate action.

● For details on dealing with alarm messages, see *Alarm monitor* on page D-6.

**Acoustic signal** 

An acoustic signal informs you of the fact that all tests are finished and the system status has changed to **Standby**.

# Stopping and restarting a run

## To stop a run

**1** Press the 🞯 global action button.

A screen is displayed that offers several kinds of stopping, each of them representing a certain level of interrupt.

Operating	23.06.2009 20:04 💽
Stop	
Abort	
🧑 Sampling stop	
C Stop printing	
Shutdown	
Restart	
	$\bigotimes$

#### Figure B-15

Interpreting the screen

Abort	When the system is in <b>Operating</b> status: Stop immediately all pro- cessing activities.			
	Pipettings that were not finished are considered not pipetted.			
	Measurements that have not yielded a valid result are considered not measured.			
Recover	When in <b>Standby</b> or <b>Stopped</b> status: Initialize all systems and modules that are currently not ready.			
Sampling stop	Finish the current pipetting action, but do not start a new one.			
	You can restart processing by pressing the $\textcircled{O}$ global action button.			
Stop Printing	Stop the current printing task. (It may take a few moments before printing actually stops.)			
Shutdown	Shut down the Cedex Bio software and the operating system.			
	This option can only be performed in <b>Standby</b> status.			
Restart	Shut down the Cedex Bio software and automatically restart it.			
	This option is available in <b>Standby</b> status only. It is used in cases where a configuration change requires restarting the software for it to become effective.			

Analyzing samples

# **Removing sample tubes**

You can remove a sample when its tube button is green —.



All tests are accepted.
All tests are pipetted.



If you remove a sample tube before the pipetting is complete, the tests that were pipetted will be performed as normal. The order remains unfinished.

You can check the resulting order details in Workplace > Order >  $\swarrow$ .

• See *Checking the sample tube status* on page B-44.

The Cedex Bio instrument provides several aids for validating results:

- In the results list, results that fall outside predefined technical ranges are flagged.
- You can display detailed result information that allows you to make a considered decision.
- Non flagged results can be accepted automatically.
- You can print the results.
- You can export the results and process them on an external computer.

The following ways of dealing with results are available:

- Accept the result.
- Re-perform the identical test (**Repeat**).
- Re-perform the test using a predefined different dilution (Rerun).

Results must be accepted before they can be transmitted to a host or before they can be automatically printed.



Results must be accepted before they can be transmitted to a host or before they can be automatically printed.

Ratio results cannot be manually accepted. They are automatically accepted if their constituent results are accepted.

The effect of flagged results depends on the configuration (**Utilities** > **Configuration** > **Result Handling**). The following table shows how.

Configuration setting	Effect
Sample Auto Accept: On	Results that do not contain a flag are auto- matically accepted. Results with flags that are marked in a predefined list of flags that should be ignored are accepted as well.
Sample Auto Accept: Off	All results need to be manually accepted.
Sample Accept Flags	Flags that are marked in this list will be ignored by the system.

 Table B-4
 Sample flag configuration and its effect

• See *Editing the acceptable flags list* on page B-160.


There are two main approaches to validating results:

- To validate sample results of an order whose sample is still on board, press the tube button on the **Overview** tab.
- In all other situations, choose Workplace > Result Review.

## ► To validate sample results of on-board samples

**1** On the **Overview** tab, press the sample button.

A screen is displayed that lists the results of the order that is associated with the sample.

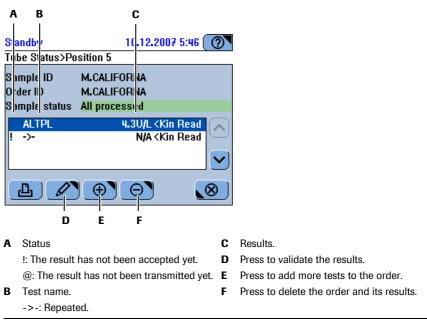

#### Figure B-16

If **TRL Check** is activated, the result might be displayed for example as <0.5, and the **<Test Rng** flag is added. This means that the measured value is below the defined TRL value.

- 2 Select the result, if there is more than one test.
- **3** Continue with *Validating the results* on page B-53.

#### ► To validate sample results from the results list

1 Choose Workplace > Result Review.

A screen is displayed that lists the results.

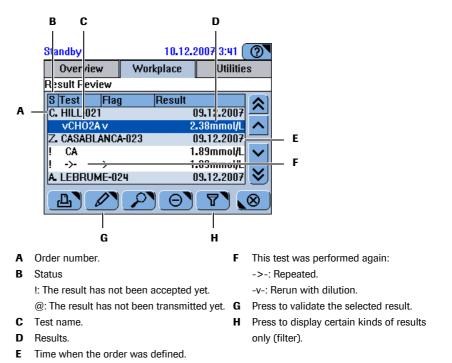


Figure B-17

If **TRL Check** is activated, the result might be displayed for example as <0.5, and the **<Test Rng** flag is added. This means that the measured value is below the defined TRL value.

<u>\</u>	
-()-	
¥	

- To check when the result was generated press  $\mathcal{P}$ .
- For results, that were not generated on the current day, only their date of generation is displayed. You can look at the results by pressing
- To display context information of the fluids used to arrive at this result, press 
   P,
   then on the detail screen press 
   Imp.

Displaying details of a result

- **2** Select the result.
- **3** Press  $\swarrow$  to display detailed information on the selected result.
- **4** Press  $\bigotimes$  to close the screen.

The screen with the result list is displayed again.

**5** Select a result.

Validating the results The system can be set up to automatically accept results that are not flagged. In addition, it can be set up to accept results with flags that are marked in an editable list of flags that should be ignored, which is particularly suitable if the flags are assessed on the host system anyway.

See Result handling on page B-165.
 See Editing the acceptable flags list on page B-160.

6 Press 🖉.

A screen is displayed for selecting your reaction.

- **7** Do one of the following:
  - Press **Repeat** to re-perform the test using the same dilution.
    See *Repeating tests* on page B-55.
  - Press Rerun to re-perform the test using a predefined (different) dilution.
    See *Rerunning tests* on page B-56.
  - Press Accept to accept the result.
     See Accepting results on page B-57.
  - Press **Retransmit** to send the result to the host again. Use this function if you suspect that the result has not been stored properly on the host system, for example because of a communication problem.

# **Repeating tests**

*Repeat* Performing a repeat means reperforming the same test with identical dilution. Typically, you perform repeats if the result is flagged and you want to confirm the result.



## Incorrect results due to insufficient fluid

Insufficient fluid may lead to inaccurate pipetting and consequently to incorrect results. Always fill the tubes with enough fluid that at least the defined dead volume of fluid is left when pipetting is complete.

• See *Tubes* on page A-53.

#### To repeat a test

- **1** Validate the result.
  - If the sample is still on board, see *To validate sample results of on-board samples* on page B-52.
  - If the sample is no longer on board, see *To validate sample results from the results list* on page B-53.

#### 2 Press Repeat.

The system automatically creates a new order and selects the test. It then performs the test.

In the results list, the result of the repeat is shown on a separate line:

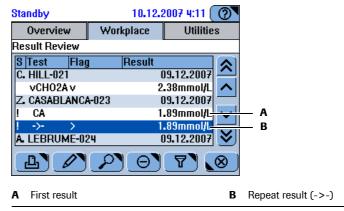


Figure B-18

# **Rerunning tests**

Rerun

## Incorrect results due to insufficient fluid

Insufficient fluid may lead to inaccurate pipetting and consequently to incorrect results.

Performing a rerun means reperforming the same test with a different predefined dilution. Typically, you perform reruns if the result is outside the test range.

Always fill the tubes with enough fluid that at least the defined dead volume of fluid is left when pipetting is complete.

• See *Tubes* on page A-53.

## To rerun a test

- 1 Validate the result.
  - If the sample is still on board, see *To validate sample results of on-board samples* on page B-52.
  - If the sample is no longer on board, see *To validate sample results from the results list* on page B-53.

#### 2 Press Rerun.

The system automatically creates a new order and selects the test. (The dilution factor is part of the test definition and therefore automatically selected. It cannot be changed.) It then performs the test.

In the results list, the result of the rerun is shown on a separate line:

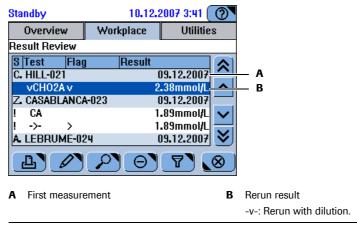


Figure B-19

# **Accepting results**

Results must be accepted before they can be printed, sent to the host, or deleted.

Automatic acceptance

The system can be set up to automatically accept results that are not flagged. In addition, it can be set up to accept results with flags that are marked in an editable list of flags that should be ignored (**Utilities** > **Configuration** > **Result Handling**).

See Result handling on page B-165.
 See Editing the acceptable flags list on page B-160.

## To accept a result

- **1** Validate the result.
  - If the sample is still on board, see *To validate sample results of on-board samples* on page B-52.
  - If the sample is no longer on board, see *To validate sample results from the results list* on page B-53.
- → Press Accept.

The result list is displayed again.

<u>`</u>Q́-

Ratio results cannot be manually accepted. They are automatically accepted if all their constituent results are accepted.

**Printing sample results** 

## ► To print results:

**1** Do one of the following:

lf	Do this
You want to print all results of an order:	Choose Workplace > Orders.
You want to print individual results:	Choose Workplace > Result Review.

## **2** Press **(**.

A screen is displayed for selecting which results should be printed.

**3** Press one of the buttons.

Press **List** to print the items currently displayed in the list. If a filter was applied, the list would contain only the items that fulfill the filter criterion, and the filter criterion would be part of the button name.

Press **Detail of Order** to print all results of the associated order.

For repeated or rerun orders, all results of this order are printed, including those that were accepted earlier.

#### 

Automatic printing

You can set up the system to automatically print results as soon as all results of an order are accepted (**Configuration** > **Workflow** > **Auto Print Results**).



Choose  $\bigcirc$  >  $\square$  to terminate the printing task, if required.

• For an overview on calibration concepts, see *Calibration* on page A-26.

For performing calibrations in the **Prepare** phase, see *Performing the calibrations (Prepare phase)* on page B-30.

For information on configuring calibration, see *Defining calibrator definitions and lots* on page B-150.

Tests with a due calibration are blocked.

• For validating calibration results see Validating calibration results on page B-64.

# **Safety information**



Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:

- Injury through reagents and other working solutions on page B-5.
- Infection by biohazardous materials on page B-5.



Skin inflammation caused by reagents

Direct contact with reagents may cause skin irritation, inflammation, or burns.

When handling reagents, be sure to wear protective equipment and observe the cautions given in the package insert.

#### Incorrect results due to expired calibration

Calibrations are performed to compensate for changes over time in reagents and in the measurement systems. Failing to perform calibrations when they are due may lead to incorrect results.

Make sure to perform calibrations when they are due.

#### Incorrect results due to wrong tube placement

Be sure to place the calibrators in the specified positions.

#### Incorrect results due to inappropriate tube and cup placement

Inappropriate tube and cup placement may lead to inaccurate pipetting and consequently to incorrect results.

Make sure that the primary tubes are placed centrally and perfectly vertically in the holders in the sample area and that they are inserted firmly.

Make sure that secondary tubes are placed centrally on the primary tubes and that they rest fully on them.

# Short guide

The following table provides an overview on the steps that make up the calibration process.

te	p		User action
	Perform the calibrations.		
	Perform all due calibrations.	Standby     10.12.2007 4:25       Calibrations [Current]>Add>Select Cal       ROUT*     STAT*       DEV*     DRUG*       ALTL     ALTPL       ASTL     ASTPL       CH02A     GLU2       UA2	<ol> <li>Choose Workplace &gt; Calibrations &gt; ⊕.</li> <li>Press ⊕ to select all tests that currently need calibrating.</li> <li>Check and, if required, print the placement list.</li> <li>Prepare and place the calibrators according to the placement list.</li> <li>Press ⊕.</li> </ol>
	Perform all calibrations that will be due within the forecast period.		<ol> <li>Choose Workplace &gt; Calibrations &gt; ⊕.</li> <li>Press ⇒ to select all tests that will require calibrating within the forecast period.</li> <li>Check and, if required, print the placement list.</li> <li>Prepare and place the calibrators according to the placement list.</li> <li>Press Φ.</li> </ol>
	Perform an individual calibra- tion.	-	<ol> <li>Choose Workplace &gt; Calibrations &gt; ⊕.</li> <li>Select the test.</li> <li>If there is more than one reagent set on board for this test, choose whether to calibrate the cur- rent set or to precalibrate a standby set.</li> <li>Check and, if required, print the placement list.</li> <li>Prepare and place the calibrators according to the placement list.</li> <li>Press <sup>(</sup>D).</li> </ol>
	Validate the results.	Standby 10.12.2007 4:27 () Calibrations [Current]>Detail>Edit UA2 CU L Calculated Accept Set Accept Lot Repeat Use Old	<ol> <li>Choose Workplace &gt; Calibrations.</li> <li>Select the calibration result.</li> <li>Press P to look at result details.</li> <li>Press I to validate the calibration.</li> <li>Remove the calibrators.</li> </ol>

 Table B-5
 Steps for performing calibrations

# **Process of performing calibrations**

The process of defining calibration orders depends on what you want to achieve:

- Calibrating all tests that need calibrating
- Calibrating all tests that will need calibrating during the forecast period
- Calibrating individual tests
  - Calibrating the current set
  - Pre-calibrating a standby set
- For information on performing calibrations in the **Prepare** phase, see *Performing the calibrations (Prepare phase)* on page B-30.

## To perform calibrations

1 Choose Workplace > Calibrations.

A screen is displayed that lists all tests and their currently valid calibration.

**2** Press 🕀.

A screen is displayed that shows all active tests.

Standby 10.12.2007 4:25 (2) Calibrations [Current]>Add>Select Cal							
ROUT*	STAT*	DEV*	DEV* DRUG* 5			6	
ALTL	ALTP	'L AS	TL	AS	TPL	CA	
CH02A	CHO2A GLU2		UA2				
( the second sec	**	[	8			<b>&gt;</b>	ĥ
				_			_
A	в						

A Press to select all tests that now need calibrating.

**B** Press to select all tests that need calibration within the forecast period.

#### Figure B-20

Interpreting the colors

Color	Meaning	Possible actions		
	Calibration is due.	Perform the calibration.		
	Calibration failed.	Check the result flag to find out why the calibration failed. Repeat the cali- bration if necessary.		
	The calibration was successfully per- formed, but its result has not been accepted yet.	Validate the calibration results.		
	Not relevant for calibration.			
	The calibration is OK. Its result has been accepted.	No action is currently required.		
	The calibration is OK. It applies to a reagent set that is not active.	No action is currently required.		

#### **Roche Diagnostics**

lf	Do this			
You want to perform all due calibrations:		Press 💮. All tests with due calibrations are selected. (The cur- rently active lots and sets are used for the calibration.)		
	2.	Press 🔊.		
You want to perform all calibrations that fall due during the forecast period:		Press S. All tests are selected whose calibration will be due within the forecast period. (The currently active lots and sets are used for the calibration.) Press S.		
You want to calibrate the current set:	1. 2.	Select the test.		

**3** To select the tests, do one of the following:

A screen is displayed that lists the calibrators that are required for the ordered calibrations, and it shows on which sample position to place them.

	A B C		
P (	Andby       Overview     Work       ep>Cal Pla cement       1     2       3     4       1: CFAS/17363       2: THC-1/[highest]	10.12.2007       12:48       (2)         place       Utilities         (5)       (6)       (7 + )       (8)         5:       THC-5/4352827       (6)       (7 + )       (6)         6:       THC-6/[lowest]       (7 + ) <th>— D — E</th>	— D — E
	3: THC-2/4352827 4: THC-3/4352827	8: THC-4/4352827	
(	品		
A	Position on the samp	le area where to <b>D</b>	There are not enough empty positions, the
	place the calibrator.		system suggests using occupied positions.
В	Lot number	E	Calibrator with lowest concentration
ſ	Calibrator with highe	st concentration	

**C** Calibrator with highest concentration

# Figure B-21

- <u>`</u>\\_\_\_
- The system first uses the free positions, if there are not enough free positions, occupied positions are suggested. For these you would have to replace the currently loaded tubes with calibrator tubes.
- With absorbance tests that require several calibrators, the calibrators are placed according to their concentration, starting with the highest concentration.
- 4 Prepare the calibrators.
- 5 Place the calibrators on the sample area positions indicated on the placement list.
- **6** Press 🔘 to confirm the placement.
- **7** Press 0 to start the calibration.

You can delete a running calibration order. After deleting the order, the calibration retains the **Due** status if it was due before



- 8 Validate the calibration results.
  - See Validating calibration results on page B-64.
- **9** Remove the calibrator tubes.

#### To replace the current lot calibration

1 Choose Workplace > Calibrations.

A screen is displayed that lists all tests and their currently valid calibration.

- 2 Select the test whose lot calibration want to replace.
- **3** Press  $\bigcirc$  to delete the result.
- **4** Press 🔘 to confirm the deletion.
- 5 Load the new reagent set. (See Adding a reagent set on page B-26.)

You can only generate a new lot calibration result when calibrating a new reagent set.

- **6** Choose Workplace > Calibrations.
- **7** Press ⊕.
- 8 Select the test.
- **9** Press **>>**.

If a screen is displayed for selecting the calibration type, choose Lot Master.

- **10** Prepare the calibrators.
- 11 Place the calibrators on the sample area positions indicated on the placement list.
- **12** Press 🔘 to confirm the placement.
- **13** Press 0 to start the calibration.

## **Deleting ordered calibrations**

You can delete a calibration order while it is being processed.

#### To delete an ordered calibration

- 1 Choose Workplace > Calibrations.
- **2** Select the order.
- **3** Press  $\Theta$ .

A confirmation dialog box is displayed.

**4** Press 🔘 to confirm the deletion.

The processing activities for this order are stopped. Measurements that have not yielded a valid result are considered not measured.



After deleting the order, the calibration retains the **Due** status if it was due before.

#### **Roche Diagnostics**

# Validating calibration results

For a calibration to become active, you need to accept its result.

Flagged calibration results

If flags are generated, you must determine their cause and decide whether to accept the result, continue working with the old results, or to rerun the calibration.

The effect of flagged calibration results depends on the configuration.

Configuration setting	Effect		
Cal Auto Accept: On	Results that do not contain a flag are auto- matically accepted. Results with flags that are marked in a predefined list of flags that should be ignored are also automatically accepted.		
Cal Auto Accept: Off	All results need to be manually accepted.		
Cal Acceptable Flags	Flags that are marked in this list will be ignored by the system.		

 Table B-6
 Calibration flag configuration and its effect

• See *Result handling* on page B-165.

See Editing the acceptable flags list on page B-160.

## ► To validate the calibration results

#### 1 Choose Workplace > Calibrations

A screen is displayed that lists all tests and their currently valid calibration.

	A	В	C	D			
St	an Iby		111-1		2007 4:1	3 🕐	
C	Overvie alit ration:			place		nues	
A () () () () ()	Test ASTL CA CHO2A GLU2 My_MG JA2		S 06.1 S 03.1	ked 2.2007 1 2.2007 1 2.2007 1 2.2007 1 2.2007 1	5:24 3:04	< 	
A	Test ID.					E	Press
В	Calibrat	ion u	se.			F	Press
С	Calibrat		•			G	Press
D	Status i	ndicat	tion.				and i
						н	Press

- Press to validate the result.
- F Press to define a new calibration order.
- G Press to delete the selected calibration and its results.
- **H** Press to display all or only the current results.

#### Figure B-22

Interpreting the screen	ltem	Value	Interpretation
	U (Use)	CU	Current.
			The calibration results of this set are currently used.
		SB	Standby.
			The calibration results of this set are not currently used. (An identical set is currently in use or the set was removed and is not empty.)
		OB	Obsolete.
			The calibration results of this set are no longer used. (The set was empty and removed or it was removed more than 30 days ago).
	Т (Туре)	L	Lot calibration (first calibration of a lot, also applies to sub- sequent sets of the lot).
		S	Set calibration (applies to calibrated set only).
	Status		The date indicates when the results were accepted. If flags were generated for the result, the flag with the highest pri- ority is displayed. In all other cases, the order status is dis- played.

• For details on calibration types, see *Calibration type* on page A-26.

**2** Select the result.

- в с A D 23.06.2009 19:49 Stai idby Calibrations (Current)>Detail 11.05.2009 16:20 F 25.05.2009 ALTL SB L Accepted F admin, 11.05.2009 16:26 G Accepted by н Flags Cal Exp R0 -0.000665665 F 5624.88 I ?**(∏**]  $\mathbf{N}$ പ്പ Κ L Test ID. н Flags. А Calibration results. В Calibration use. н С Calibration type. н Press to validate the results. Calibration status. Press to display result details. D κ Ε Calculation date and time. Press to display context information of the L Calibrator, lot, expiration date. fluids used to arrive at this result F G Accepted by: User name. \$SYS\$ means automatically accepted.
- **3** Press  $\swarrow$  to display information on the selected calibration and its results.

Figure B-23

4 Press 🖉.

A screen is displayed for selecting your decision.

- **5** Do one of the following:
  - Press Accept Set to accept the set calibration results for the selected test.
  - Press Accept Lot to accept the lot calibration results for the selected test.

This button is only active if the calibration sequence **Each Lot and Interval** is defined for the application and if no flags were generated for the result.

- For information on lot and set calibration, see *Calibration type* on page A-26.
- Press Repeat to have the calibration performed again.
- Press **Use Old** to discard the new result, reset the calibration due date, and to continue using the old calibration results.

This possibility is available if you selected an accepted result. When you press **Use Old**, a copy of the old result is made and a new entry is displayed in the calibration results list. Note that the intervals are reset as if a new calibration result were generated.



#### Incorrect results when using Use Old

Calibrations are performed to compensate for changes over time in reagents and in the measurement systems. Failing to perform calibrations when they are due may lead to incorrect results.

- Roche recommend performing a QC measurement before you continue working with the old calibration results.
- In the application definitions, choose On for QC After Cal.
   (Utilities > Applications > Laboratory Parameters > Control > QC After Cal)



## Accepting calibration results

Calibration results will be used by the system only when you have accepted them.

## There can be only one unaccepted calibration result for a test.

If there already exists a not accepted calibration for a test and you place another calibration order for this test, the order will be blocked.

QC is performed at regular intervals to check the integrity of the whole measuring system.

*Control* A control is a sample that has been measured using all tests it is associated with, in order to define the ranges and values that determine the correct functioning of the instrument. This is typically done both for the normal and the pathological analyte concentration.

For an overview on QC, see *Quality control (QC)* on page A-29.
 For information on QC configuration, see *Defining control definitions and lots* on page B-149.

There are two basic ways of performing QC on the Cedex Bio instrument:

Default QC

**Default QC** is an automated process for performing QC measurements in one request. This is the ideal method if you want to perform QC on certain days or at certain times.

This process only applies to tests whose controls are defined to be performed as part of **Default QC**. Therefore, if you intend to work with **Default QC**, you need to configure the tests accordingly.

• See Defining control definitions and lots on page B-149.

Performing **Default QC** follows a streamlined procedure whereby QC orders are automatically defined as soon as you identify a control. An order is defined for all tests for which this control is defined as the **Default QC**, provided the test is currently active on the system.

Interval QC



Interval QC is a process that is suitable both for performing a single QC measurement and for performing all QC measurements that are due. You can select all tests that require QC simply by pressing a button (I). (This selection also reflects QC of the type **QC after Cal**.) A wizard helps you select the controls, and a placement list supports you in preparing and loading them.

# **Safety information**



Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:

- Injury through reagents and other working solutions on page B-5.
- Infection by biohazardous materials on page B-5.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page B-5.



#### Incorrect results due to inappropriate tube and cup placement

Inappropriate tube and cup placement may lead to inaccurate pipetting and consequently to incorrect results.

Make sure that the primary tubes are placed centrally and perfectly vertically in the holders in the sample area and that they are inserted firmly.

Make sure that secondary tubes are placed centrally on the primary tubes and that they rest fully on them.

# Short guide

Ste	eb (1997)		User action
	Perform QC		
	Perform <b>Default QC</b> .	Standby       10.12.2007 15:25 ()         Overview       Workplace       Utilities         Order>Identify       Sample>Select Control         P_NORM	<ol> <li>Choose Overview &gt; Order.</li> <li>Press II.</li> <li>Select a control and place the tube. Repeat until there are no controls left on the screen.</li> <li>Press O.</li> </ol>
	Perform all interval QC mea- surements that are due.	Standby       10.12.2007 12:14         Overview       Workplace       Utilities         Overview>Add>0C Placement       P_NORM         P_NORM       P_PATH       Image: Comparison of the second sec	<ol> <li>Choose Overview &gt; Order &gt; ♥↓ or Workplace &gt; QC Status &gt; ♥.</li> <li>Press ♥.</li> <li>Press ♥.</li> <li>Select a control and place the tube. Repeat until there are no controls left on the screen.</li> <li>Press ♥.</li> </ol>
	Perform a single interval QC measurement.	Standby         23.06.2009 19:50         Image: Constraint of the state of the st	<ol> <li>Choose Workplace &gt; QC Status.</li> <li>Press ⊕.</li> <li>Select the test.</li> <li>Press ≫.</li> <li>Select a control and place the tube. Repeat until there are no controls left on the screen.</li> <li>Press Φ.</li> </ol>
	Validate the results.	Standby 23.06.2009 19:51 ( QC Status>Edit ! ASTL PNU Accept Ignore ()	<ol> <li>Choose Workplace &gt; QC Status.</li> <li>Press  P to look at result details. Close the details screen.</li> <li>Press  and press a button to validate the QC result.</li> <li>Remove the controls.</li> </ol>

The following table provides an overview of the steps that make up the QC process.

 Table B-7
 Steps for performing QC

# **Performing Default QC**

Performing **Default QC** follows a streamlined procedure whereby QC orders are automatically defined as soon as you identify a control. An order is defined for all tests for which this control is defined as the **Default QC**, provided the test is currently active on the system.

## ► To perform Default QC

- 1 Choose **Overview** > **Order**.
- 2 Press **S**.

A screen is displayed that contains a button for each of the controls that are required.

- **3** Prepare the controls.
- 4 Press a control button.

A screen is displayed, asking you to place the selected control.

**5** Place the control tube on any free position on the sample area.

The system registers the position and automatically defines an order for each test that has this control defined as its **Default QC**.

The screen with the control buttons is displayed again. The button for the control you just loaded is no longer active.

6 Press the next active control button.

A screen is displayed, asking you to place the control.

- 7 Place the control on a free position on the sample area.
- 8 Select and place the remaining controls as described in steps 6 through 7.

When all controls are placed, the screen for identifying samples is displayed.

**9** Press  $\bigotimes$  to close the screen.

**10** Press 0 to start processing the control orders.

# **Performing interval QC measurements**

## To define a QC order

- **1** To define a new order, do one of the following:
  - Choose **Overview** > **Order** > **O**. •
  - Choose Workplace > QC Status > (). ٠

A a test selection screen is displayed.

Standby 10.12.2007 12:13 ( ) QC Status>Add>Select QC							
ROUT*	STAT	DEV	DR	JG	5	6	
ALTL	ALTP	AS	TL	AS	TPL	CA	
СНО2А	GLU	2 <b>`</b> U	A2 💙				
							٦
							Ī
Ð		] [	R		I L	<b>&gt;&gt;</b>	ĩ
							/
Å							

A Press to select all tests with due QC. Press again to cancel the selection.

#### Figure B-24

Interpreting the display	The test is blocked for one of the following reasons:
	Calibration is required.
	The calibration failed.
	• For the reagent set, the number of available tests is 0, or a reagent bottle is
	missing (incomplete reagent set).
	A QC is due or its result has not yet been accepted.
i	The reagent set is on board and ready for use.
-	The reagent set is not on board.
-	
<b>2</b> T	o select the tests, do one of the following:

- - Press 🕑 to select all tests with due QC.

This selection also applies to controls of the type QC after Cal.

- Press a test button.
- 3 Press ».

A screen is displayed that contains a button for each of the controls that are required.

-Ò́-

If at this stage, the required controls are already loaded on the instrument, then the QC orders are automatically created.

4 Prepare the controls.

## **Roche Diagnostics**

**5** Press a control button.

A screen is displayed, asking you to place the selected control.

6 Place the control tube on any free position on the sample area.

The system registers the position and automatically defines an order.

If there are more controls to place, the screen with the control buttons is displayed again. The control you just placed is no longer active.

7 Select and place the remaining controls as described in steps 5 through 6.

When all controls have been selected and placed, the **Overview** tab is displayed.

**8** Press 0 to start processing the control orders.

# **Validating QC results**

Flagged QC results

If flags are generated, you must determine their cause and decide whether to accept the result or to ignore it.

The effect of flagged QC results depends on the configuration.

Configuration setting	Effect
QC Auto Accept: On	Results that do not contain a flag are auto- matically accepted. Results with flags that are marked in a predefined list of flags that should be ignored are also automatically accepted.
QC Auto Accept: Off	All results need to be manually accepted.
QC Acceptable Flags	Flags that are marked in this list will be ignored by the system.

 Table B-8
 QC flag configuration and its effect

See Result handling on page B-165.
 See Editing the acceptable flags list on page B-160.

## ► To validate the QC results

#### 1 Choose Workplace > QC Status.

A screen is displayed that lists the most recent QC results for each test that is installed on the system.

1	4 E	3	(	C	1	D		
S	andby	,			23.06.	2009	19:50 (	0
	Ove	rviev	v	W	'orkplace		Utilitie	s
Q	C Stat	us						
	3 Test	t	ID		Flag/Res	ult		
I	GLU	12	PPU		13 <b>.</b> 14mm	ol/L		
	ALT	PL	PNU	I .	52.0U/L			
l	AST	L	PNU		39.4U/L			
Г	AST	PL	PNU	l –	37.3U/L			1 🗸
	ALT	PL	PPU	I	R1(2.5s)			
	AST	L	PPU	I	R1(2.5s)			[⊗]
(	Д			F	»)⊕`		Ð	⊗


#### A Status

I: The result has not been accepted yet.

- I: The result was ignored.
- @: The result has not been transmitted yet.
- B Test name.
- C Control ID.

# Result, if no flag was generated. Flag with highest priority, if a flag was generated. Order status, if the control measurement has not been performed yet.

- Figure B-25
- **2** Select a result.
- **3** Press  $\swarrow$  to look at result details.

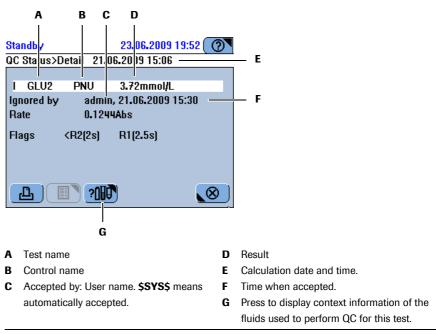



-Ò́-

**4** Press  $\bigotimes$  to close the screen.

If the QC result is outside the defined range, perform the QC again. If the results are still outside the range, check for other causes. If all fails, perform a calibration.

5 Press 🖉.

- **6** Do one of the following:
  - Press Accept to accept the QC results for the selected test.
  - Press **Ignore** to exclude the result from further QC result calculations such as QC history statistics.

When ignoring a result, the due status does not change.

In the QC results tables, ignored QC results are marked with "I".

Flags of ignored results are not inherited to dependent results.

• Press Retransmit to send the result again.

(This option is active if your instrument is connected to a host system.)

# Interpreting the QC history

The QC history provides—on individual screens—information on the QC results of the current and the previous calendar month as well as on results that were generated before the previous month.

A graphic representation of the results provides a convenient way for comparing results over a period of time.

## To interpret the QC history

1 Choose Workplace > QC History.

A screen is displayed that contains, one QC result entry for each lot of each test.


,	<b>A</b>	В		C	;	
Stand	ЬУ		2	3.06.2	)09 19:46	0
0	rview	,	Workpl	ace	Utiliti	es
QC H	story					
Test		D	Las	t meas	urement	
ALTI	PL İ	PNU	22.0	) <b>6.200</b>	9 14:13	
ALTI	PL I	PPU	22.0	) <b>6.200</b> :	9 14:13	
AST	LI	PNU	22.0	) <b>6.200</b> :	9 14:13	
AST	LI	PPU	22.0	) <b>6.200</b> :	9 14:13	
AST	PL I	PNU	22.0	) <b>6.200</b> :	9 14:13	
AST	PL I	PPU	22.0	) <b>6.2</b> 00	9 14:13	∎ <b>≥</b> ∣
ß	1		<u>P</u> (	Θ	7	$\otimes$
ΑΤ	est nam	ie.				С
B C	ontrol I	D.				

Date of most recent control. measurement. If the control does not belong to the current lot, **Previous Lot** is displayed instead of the date.

Figure B-27

**2** Press  $\square$ .

A graphic is displayed that shows either the QC results for the current or the previous month, or those generated before the previous month.

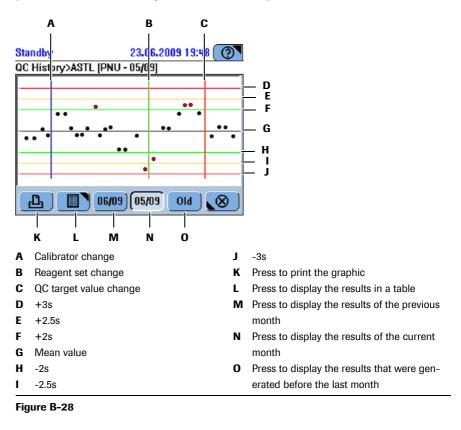




- The date format in the month selection buttons is always month/year.
- Ignored QC results are not displayed in the graph, but they are included in the QC History printout. Such results are marked with "I" at the end.
- Ignored QC results are not taken into account for QC history statistics.

**3** Press **1** to display the results in a table.

The results are sorted chronologically. An overview is followed by statistical data for all stored results and for the results of the current month. (If a QC target value change took place, only the results from this date onward are taken into account for both statistics.)

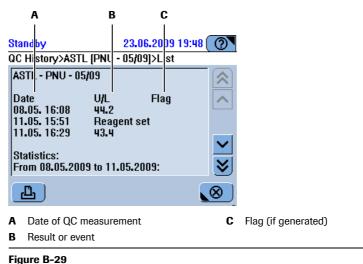

- **4** Press  $\bigotimes$  to close the screen.

# Performing reagent mixing

Reagents that contain Latex granules, for example D-Dimer, require periodic mixing. The mixing interval is part of the application definitions and cannot be changed by the user. The system checks every 30 minutes for reagent sets that require mixing.

- For mixing during the **Prepare** phase, see *Performing reagent mixing* on page B-29.
- To perform the mixing
  - 1 Choose Overview > III.
  - 2 Press the button of the reagent set you want to mix.

A screen is displayed that contains details on the set.

**3** Press **O**.

(This button is available if there is a mixing interval defined for this reagent set.)

Mixing starts. A screen is displayed informing you about the progress of the mixing action.

**4** When mixing is finished, press  $\bigotimes$  to close the screen.

Closing the dialog box while mixing is in progress stops the mixing action. Mixing would have to be performed anew.

The typical end of shift activities are organized in a single wizard. By performing the steps as suggested by this wizard, you put the instrument in a condition that allows you to hand over operation to another operator or to shut down the system and switch off the instrument.

# **Safety information**



Make sure that you have read and understood section *Safety information* on page B-5. The following safety messages in particular are relevant:

Warning messages:

- Injury through reagents and other working solutions on page B-5
- Infection by biohazardous materials on page B-5
- Injury through reagents and other working solutions on page B-5



#### Skin inflammation caused by reagents

Direct contact with reagents may cause skin irritation, inflammation, or burns.

When handling reagents, be sure to wear protective equipment and observe the cautions given in the package insert.

## **5 Daily operation**

Finishing the shift

# Short guide

The following table provides an overview of the steps that make up the end of shift process.

Ste	p		User action
1	Check for unfinished tasks.	Standby     10.12.2007 4:17       Result Review>Filter       All       Flagged       STAT       Not Accepted       Not Today       By ID	<ol> <li>Check for unfinished orders.</li> <li>Check for nonvalidated results.</li> <li>Check for results that were not transmitted.</li> </ol>
2	Start the <b>End Shift</b> wizard.	Standby     10.12.2007 13:46       Overview     Workplace     Utilities       1     2     3     4     5     6     7     8       0     Tests / 0     0     Tests / 0     1000     Tests     Tests     1000     Tests       0     Tests     admin     0     Tests	<ol> <li>On the Overview tab, press the End Shift but- ton.</li> </ol>
3	Perform data backup.	Standby 10.12.2007 5:35 End Shift>Backup  Press ⟨Backup> to copy the current database to the USB stick.	<ol> <li>Press .</li> <li>Insert the USB stick.</li> <li>Press Ø.</li> <li>Select the directory.</li> <li>Press Ø.</li> <li>Press Ø.</li> <li>Press D to proceed to the next stage in the End Shift wizard.</li> </ol>
4	Export support data.	Standby 10.12.2007 5:35 () End Shift>Export Press <export> to copy the full results to the USB stick (patient, QC, and calibration results).</export>	<ol> <li>Press ◆.</li> <li>Insert the USB stick, if you have removed it.</li> <li>Press ◆.</li> <li>Select the directory.</li> <li>Press ◆.</li> <li>Remove the USB stick.</li> <li>Press ▶ to proceed to the next stage in the End Shift wizard.</li> </ol>
5	Clean up the database.	Standby       10.12.2007 5:36       Image: Constraint of the standard s	<ol> <li>Press  .     <li>Press   to confirm the deletion.     <li>Press   to proceed to the next stage in the End Shift wizard.     </li> </li></li></ol>

 Table B-9
 Steps for finishing the shift

Ste	p		User action
6	Perform maintenance actions.	Standby       10.12.2007 4:23         End Shift>Check Maintenance         Clean Water and Waste       10.12.         Replace Photometer Lamp       10.12.         Clean Probe Manually       10.12.         Deproteinize Probe       11.12.         Clean Reagent Disk       07.01.         Replace Water Inlet Filter       25.02.         Clean Housing       02.06.	<ol> <li>Check which maintenance actions are due.</li> <li>Perform the maintenance actions. Perform at least all red maintenance actions.</li> <li>Press &gt;</li></ol>
7	Check the cuvette status.	Standby     10.12.2007 4:03       End Shift>Cuvettes       Select segment to replace       0     0       0     0       0     0       0     10       0     10	<ol> <li>Replace at least all red cuvette segments.</li> <li>Press the cuvette button.</li> <li>Open the main cover.</li> <li>Replace the cuvette segments.</li> <li>Close the main cover.</li> <li>Press &gt;</li></ol>
8	Remove the reagent disk.	Standby 10.12.2007 12:34 ( Overview>Disk 2 ALTL ASTL BARBT CA CHOL2 CRPLX D-DI GLUC2 My_MG THC UA2 CLEANT CLEAN NACL PYP I I III E M	<ol> <li>Choose Overview &gt; M.</li> <li>Press .</li> <li>Open the main cover.</li> <li>Lift the disk from the instrument.</li> <li>Place the disk in its container.</li> <li>Press .</li> <li>Close the main cover.</li> <li>Store the reagent disk in a cool place.</li> </ol>
9	Empty the waste container.	Standby       10.12.2007 12:28         End Shift>Check Bottles         Press fluid button when done.         Water       Waste         Full       Empty         Cleaner         100%         Standby       Standby	<ol> <li>Empty the waste container.</li> <li>Press the Waste button to confirm.</li> <li>Refill the water container.</li> <li>Press the Water button to confirm.</li> </ol>
10	Finish your shift.	Standby     10.12.2007 13:46 (2)       Overview     Workplace     Utilities       1     2     3     4     5     6     7     8       0     Tests / 0     Blocked     0     8       0     Tests / 0     To Validate       STAT     Order     U       Tests     admin     U       Prepare     End Shift     000,	<ol> <li>Log off the system.</li> <li>Shut down the system and switch off the main instrument. (If there is no other shift.)</li> </ol>

 Table B-9
 Steps for finishing the shift

# Checking for unfinished tasks

## Checking for unfinished orders

## To check for unfinished orders

- 1 Choose Workplace > Orders.
- **2** Choose  $\gamma$  > Not Finished.
- **3** Press  $\oslash$ .

A screen is displayed that contains all unfinished orders.

- 4 Do whatever is necessary to finish the orders.

## Checking for not accepted sample results



Results must be accepted before they can be printed or transmitted to the host.

#### To check for not accepted results

1 Choose Workplace > Result Review.

A screen is displayed that contains all sample results of the day.

- **2** Choose  $\gamma$  > Not Accepted.
- **3** Press 🔘.

A screen is displayed that contains all sample results that have not been accepted yet.

- 4 Select a result.
- **5** Press 🖉 to validate the result.

A screen is displayed for selecting your decision.

- See Accepting results on page B-57.
   See Repeating tests on page B-55.
   See Rerunning tests on page B-56.
- 6 Press one of the buttons.
- 7 Select and validate the remaining results as described in Steps 4 through 6.

If you did not accept all results, deal with the not accepted ones and check for not accepted results again.

#### Checking the transmission of results

This step is only relevant if you work with a host system.

Before deleting the results, you must make sure they were properly transmitted to the host.



Results must be accepted before they can be transmitted to the host.

See also the information on *Result handling* on page B-165, in particular on the **Auto Accept** settings.

#### ► To check for results that were not transmitted

1 Choose Workplace > Result Review.

A screen is displayed that contains all sample results of the day.

- **2** Choose  $\gamma$  > Not Sent to Host.
- **3** Press 🔘.

A screen is displayed that contains the sample results that have not been successfully transmitted to the host.

- 4 Select a result.
- 5 Press 🖉.

A screen for validating results is displayed.

6 Press Retransmit.

The list with the results that were not transmitted successfully is displayed again. The result you just retransmitted should no longer be present.

- 7 Select and retransmit the remaining results as described in Steps 4 through 6.
- 8 Choose  $\gamma$  > Not sent to Host.

The list should now be empty.

- **9** If the results were not successfully transmitted, contact the administrator of the host computer.

## Starting the End Shift wizard

The easiest way of performing the end of shift tasks is by following the **End Shift** wizard.

## To start the End Shift wizard

1 Choose Overview > End Shift.

A screen is displayed for performing daily backup.

# Performing the daily backup

The **Cedex Bio** instrument can store orders and result data for one working day. It is therefore necessary to export to an external medium all data that you need to keep.

During a database export, the full content of the database is copied to the USB stick.

The database data can be restored to the instrument if required (**Utilities** > **Import** > **Database**).

• For information on restoring the database, see *Importing a database* on page B-113.

## To perform the daily backup

1 Press 🖭.

A screen is displayed, asking you to insert the USB stick.

**2** Insert the USB stick.

A screen is displayed showing the directory structure of the USB stick.

- **3** Select the directory where you want the backup file to be copied.
- **4** Press 🔘 to confirm the selection.

The data are copied to the stick.

Database files names have the following name format: dba\_yyyymmddhhmmss.tgz.

**5** Press **>>** to proceed to the next stage in the **End Shift** wizard.

A screen is displayed for deleting sample results.





When performing backup without using the End Shift wizard: Choose Utilities > Export > Database.

# Exporting the result data

When exporting results, the full result data are exported. The following results are copied to the USB stick (each of the sets is contained in a separate file):

Data set	File name
Sample results	res_yyyymmddhhmmss.csv
QC results	qcs_yyyymmddhhmmss.csv
QC history results	qch_yyyymmddhhmmss.csv
Calibration results	cal_yyyymmddhhmmss.csv
Process event log	pev_yyyymmddhhmmss.csv

*File format* Result data are exported as comma-separated value (csv) files. The entries are separated by semicolons (;).

• For details on csv definitions, see *Exporting results* on page B-105.

These files can be processed with any spreadsheet program that can import csv files.

## ► To export the results

1 Press 📀.

A screen is displayed, asking you to insert the USB stick.

**2** Insert the USB stick.

A screen is displayed showing the directory structure of the USB stick.

- **3** Select the directory where you want the result files to be copied.
- **5** Press >>> to proceed to the next stage in the **End Shift** wizard.

A screen is displayed for deleting sample results.



When performing backup without using the End Shift wizard: Choose Utilities > Export > Full Results.

# Cleaning up the database

-χ-

Ò.

The **Cedex Bio** instrument can store orders and results data for one working day. It is therefore necessary to delete results and orders to ensure that there is enough space on the system for the next shift.

By deleting the sample results, you delete the corresponding orders as well.

## ► To clean up the database

1 Press  $\Theta$ .

The following results will be deleted:

- Accepted sample results.
- QC results that were accepted are deleted from the QC Status list (they remain in the QC history).
- QC results that were generated before the first day of the previous month are deleted from the QC history.
- Obsolete calibration results, provided there are more than five obsolete results for the same test.

(Results become obsolete if the empty set was removed or if the set was removed more than 30 days ago.)

**2** Press 🕑 to confirm the deletion.

The results are deleted.

**3** Press **>>** to proceed to the next stage in the **End Shift** wizard.

A screen for handling cuvettes is displayed.

# When deleting results outside the End Shift wizard:

You can delete accepted sample results, calibrations, orders, and accepted and ignored QC results by pressing  $\bigcirc$  in the respective result lists.

- To delete orders on page B-93
- To delete sample results on page B-94
- *To delete a calibration* on page B-95
- To delete QC results from the QC Status list on page B-97
- To delete QC results from the QC History on page B-97

#### Maintenance action Cleanup Database

The maintenance action **Cleanup Database** is intended for situations where the system does not work efficiently any more, it is not intended to clear up the data at the end of a shift.

### **Performing maintenance actions**

To ensure the smooth running of the system, you should perform all due maintenance actions. Performing these actions as part of the end of shift activities ensures that at the beginning of the next shift, when there might be many tests to be performed, the system is ready for processing quickly.

### ► To perform maintenance actions

- **1** Follow the instructions given in *To perform maintenance actions* on page B-19.
- **2** Press **>>** to proceed to the next stage in the **End Shift** wizard.

A screen is displayed that shows the status of the cuvettes.

`Q́`

-Ò́-

When performing maintenance actions without using the End Shift wizard: Choose Utilities > Maintenance.

### **Replacing cuvettes**

### To replace cuvettes

- 1 Follow the instructions given in *To prepare the cuvettes* on page B-28.
- **2** Press  $\gg$  to proceed to the next stage in the **End Shift** wizard.

A screen is displayed that shows the status of the external fluid bottles.

When replacing cuvettes without using the End Shift wizard: Choose Overview  $> \mathbb{T}$ . Finishing the shift

### Dealing with the external fluid bottles

#### Emptying the waste container



Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:

- Injury through reagents and other working solutions on page B-5.
- Infection by biohazardous materials on page B-5.
- Infection by waste solution on page B-5.

To prevent unpleasant smells and contamination of the environment, you need to empty the waste at the end of your shift.

### To empty the waste container

- **1** Remove the tubing adapter from the yellow waste container and insert it in the spare container.
- 2 Remove the bottle from the rack and place it on a firm and even surface.
- **3** Place the spare container on the rack.
- **4** Follow the instructions given in *To check the waste container* on page B-17.

Ù.

Roche recommend to refill the water container whenever you empty the waste container.

When emptying the waste without using the End Shift wizard:

Choose Overview > 🎯 > 🛍.

### Refilling the external water bottle



ò

#### Danger of poor measurement quality due to inadequate water quality

Inadequate water quality may lead to incorrect results. Always use purified water of the quality specified in section *Technical specifications*.

#### To refill the external water container

- 1 Remove the tubing adapter from the white water container and place it on a clean surface.
- **2** Follow the instructions given in *To check the water container* on page B-16.

Roche recommend to empty the waste container whenever you refill the water container.

#### When refilling the water without using the End Shift wizard:

Choose Overview > 🚳 > 🛍.

### Removing the reagent disk



During periods when you do not perform tests, the disk should be stored in a clean refrigerated place at temperatures in the range of +2 to  $+8^{\circ}$ C.

Make sure that you have read and understood section *Safety information* on page B-5. The following warning messages in particular are relevant:

- Injury through reagents and other working solutions on page B-5.
- Infection by biohazardous materials on page B-5.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page B-5.



#### Spillage through tipping reagent disk

The reagent disk container can slip off or tip over if it is not placed on an even horizontal surface.

When storing the reagent disk container, make sure to place it on a firm, even, horizontal surface that is easily accessible.

When handling the reagent disk, make sure not to tilt it.

### To remove the reagent disk

1 Choose Overview >  $\square$  >  $\square$ .

A screen is displayed, asking you to remove the reagent disk.

- **2** Open the main cover.
- **3** Remove the reagent disk.
- 4 Place the reagent disk in the reagent disk container.
- **5** Close the main cover.
- **6** Store the reagent disk container in a refrigerated place.

Logging off

## Logging off

You should log off the system before you hand it over to another operator. (Only one person can be logged on at any time.)

You can log off any time, even while the system is processing orders.

### To log off the system

1 Choose **Overview** > button with your user name.



-Ò́-

### Automatic logoff

You can configure the system to automatically log off the user after a configurable period of inactivity (Utilities > Configuration > System > Screen Saver Wait  $\neq$  0; Utilities > Configuration > System > Auto Log-off = On).

#### Viewing alarms while you are logged off

You can view alarms any time, even when you are logged off.

• For configuring automatic logoff, see *System* on page B-166.

If there is no other shift, you can now shut down the system and switch off the main instrument.

## Shutting down the system and switching off the instrument

When you switch off the system, reagent cooling stops. Therefore, you need to remove the reagent disk and store it in a cool place before switching off the system.

*Preconditions* The system must be in **Standby** status.

### ► To shut down the system and switch off the main instrument

- 1 Press 🞯
- 2 The Stop screen is displayed.
- **3** Press ① Shutdown to shut down the system.
  - Wait until the screen is black.
- 4 Press **O** on the power toggle switch.

## Using the barcode scanner

• For information on the barcodes, see *Barcode scanner* on page A-59.

Because the barcodes on reagent bottles and sample tubes are different, you use the barcode scanner in a different way. The various procedures are described in the following sections.



#### Loss of sight

The intense light of the LEDs may severely damage you eyes. Do not stare into the LEDs. Scanning equipment using LED technology is covered by the international standard IEC 60825-1 LED Safety: Class 1.

#### Injury through reagents and other working solutions

Direct contact with reagents, cleaning solutions, or other working solutions may cause personal injury. When handling reagents, exercise the precautions required for handling laboratory reagents, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics reagents and cleaning solutions.



### Malfunction due to spilled liquid

Any liquid spilled on the instrument may result in malfunction of the instrument. If liquid does spill on the instrument, wipe it up immediately and apply disinfectant.

Make sure you do not tilt the bottle or sample tube when scanning its barcode.

### **Reading reagent bottle barcodes**

### ► To use the barcode scanner for reading reagent bottle barcodes

- 1 When a screen is displayed, asking you to scan, use the barcode scanner and scan the barcode.
- **2** Make sure not to tilt the bottle while scanning its barcode.
- **3** Hold the scanner at a distance of approximately 20 cm (8 in) from the barcode, pull the trigger on the scanner and point the red light just outside the barcode.
- 4 Move the light slowly across the barcode.
- **5** Wait until you hear a beep and release the trigger.

If the reading was successful, a screen is displayed, asking you to place the bottle on the reagent disk.

### **Reading barcodes from sheets**

### ► To use the barcode scanner for reading barcodes from sheets

- 1 When a screen is displayed, asking you to scan, use the barcode scanner and scan the barcode.
- 2 Place the barcode transfer sheet on a flat surface and smooth it out.
- **3** Hold the scanner at a distance of approximately 20 cm (8 in) from the barcode, pull the trigger on the scanner and point the red light just outside the barcode.
- **4** Move the light slowly across the barcode.
- **5** Wait until you hear a beep and release the trigger.

If the reading was successful, a screen is displayed, asking you to place the item.

### **Reading sample tube barcodes**



For safety reasons, the barcode scanner is set to only read barcodes that contain a checksum.

### ► To use the barcode scanner for reading sample tube barcodes

- 1 When a screen is displayed, asking you to scan, use the barcode scanner and scan the barcode.
- 2 Make sure not to tilt the sample tube while scanning its barcode.
- **3** Hold the scanner at a distance of approximately 20 cm (8 in) from the barcode, pull the trigger on the scanner and point the red light at the barcode.
- 4 Wait until you hear a noise and release the trigger.

If the reading was successful, a screen for selecting tests is displayed.

# **Special operations**

The tasks you do not perform every day

In this chapter, you will find information on operator tasks that are not part of the daily routine of analyzing samples.

In this chapter	Chapter <b>6</b>
Deleting sample orders	B–93
Deleting sample results	B–94
Calibration	B–95
Deleting calibration results	B–95
Setting up your calibration schedule	B–96
Deleting QC results	B–96
Lot handling	B–98
Viewing the current lot definitions	B–98
Calibrator and QC lots	B–98
Reagent lots	B–99
Viewing the lot details of fluids that were used	B–100
Adding a new lot	
Exporting data	B–104
Exporting the database	B–104
Exporting results	B-105
Interpreting result data	B–106
Exporting log files	B–110
Importing data	B–111
Importing applications	B–111
Importing software	B–113
Importing a database	B-113
Importing certificates	B–115
Importing automatic mixing and extra wash cycle definitions	3B–115
Preparing a new disk	B–116
Assigning tests to test tabs	B–119
Deleting bottle sets from the Inventory list	B-120
Refilling printer paper	B–122
Removing condensation water from the reagent cooler	B-124

#### **Roche Diagnostics**

Replacing the probeB-1	125
Connecting and disconnecting the external fluid containers	
Adjusting the touchscreenB-1	131
Cleaning the touchscreenB-1	131

## **Deleting sample orders**

Sample orders are normally deleted as part of the daily end of shift activities. This section describes how to proceed if you need to delete sample orders outside the normal daily routine.

• For deleting orders at the end of a shift, see *Cleaning up the database* on page B-84.

*Preconditions for deleting* The orders should be fully processed before you delete them.

### To delete orders

1 Choose Workplace > Orders.

A screen is displayed that lists all currently defined sample orders.

#### **2** Do one of the following:

lf	Do this	
You want to delete an individual order:	1. Select the order entry in the list.	
	2. Press $\Theta$ .	
	3. Press Selected.	
You want to delete all or a specific group of	1. Press T.	
orders:	2. Press the appropriate filter button.	
	3. Press 🖉 to confirm the selection	
	4. Press $\Theta$ .	
	5. Press List [filter criterion].	

The orders are deleted. You cannot retrieve them any more from the system.



- · By deleting a sample order, you delete the associated results as well.
- You can delete a running order. After deleting the order, cuvettes and reagents that were not used are released again for future use

## **Deleting sample results**

Sample results are normally deleted as part of the daily end of shift activities. This section describes how to proceed if you need to delete sample results outside the normal daily routine.

• For deleting results at the end of a shift, see *Cleaning up the database* on page B-84.

*Preconditions for deleting* Sample results should be accepted and printed or sent to the host before deleting.

### ► To delete sample results

1 Choose Workplace > Result Review.

A screen is displayed that lists all results.

**2** Do one of the following:

lf	Do this		
You want to delete an individual sample	1. Select the result entry in the list.		
result:	2. Press $\Theta$ .		
	3. Press Selected.		
You want to delete all or a specific group	1. Press T.		
of sample results:	2. Press the appropriate filter button.		
-	3. Press 🖉 to confirm the selection.		
	4. Press $\Theta$ .		
	5. Press List [filter criterion].		

The results are deleted. You cannot retrieve them any more from the system.



By deleting all sample results of an order, you delete the corresponding order as well.

## Calibration

### **Deleting calibration results**

Obsolete calibration results are automatically deleted as part of the daily end of shift activities if there are more than five obsolete calibration results for the test. (Calibration results become obsolete when the empty set is removed or if it was removed more than 30 days ago.)

This section describes how to proceed if you need to delete results outside the normal daily routine.

For deleting calibration results at the end of a shift, see *Cleaning up the database* on page B-84.

### To delete a calibration

- 1 Choose Workplace > Calibrations.
- **2** Select the calibration entry.
- **3** Press  $\Theta$ .

A confirmation screen is displayed.

4 Press 🔘.

The order and the results are deleted. You cannot retrieve them any more from the system.

### Setting up your calibration schedule

The **Cedex Bio** instrument provides convenient software-supported procedures for the following major calibration scenarios:

- Performing all calibrations that are currently due
- Performing all calibrations that fall due within a configurable forecast period
- Performing individual calibrations

Performing calibration as part of the Prepare wizard includes performing all calibrations that are due and that will fall due during the forecast period.

The following table lists the configurable definitions and describes their effect.

Definition item	Configuration path	Effect
Forecast Hours	Utilities > Configuration > Calibration	Period of time, starting with the current date, within which calibration falls due. Defining a forecast period (value other than zero) allows to concentrate calibration per- formance to specific days.
Cal Auto Accept	Utilities > Configuration > Result Handling	If on, results that are not flagged are auto- matically accepted. Results with flags that are marked in a pre- defined list of acceptable flags would also be automatically accepted.
Cal Acceptable Flags	Utilities > Configuration > Result Handling	List of calibration relevant flags. Flags that are marked are automatically accepted if <b>Cal Auto Accept</b> is on.
Table B-10	Calibration definitions and their effect	

## **Deleting QC results**



- QC results are displayed in the QC Status list as soon as they are generated. A copy of the results is also placed in the QC History list. This list keeps the results of the previous and the current months.
- Deleting QC results, deletes the associated orders as well.

Accepted QC results are normally deleted from the **QC Status** list as part of the daily end of shift activities, at the same time, results that were generated before the first day of the previous month are deleted from the **QC History** list.

This section describes how to proceed if you need to delete QC results outside the normal daily routine.

● For deleting QC results at the end of a shift, see *Cleaning up the database* on page B-84.

Deleting QC results

### ► To delete QC results from the QC Status list

- 1 Choose Workplace > QC Status.
- **2** Press  $\Theta$ .

A screen for selecting results is displayed.

**3** Do one of the following:

lf	Do this		
You want to delete all QC results:	<ol> <li>Choose All.</li> <li>Press O to confirm the deletion.</li> </ol>		
You want to delete all accepted QC results:	<ol> <li>Choose All Accepted.</li> <li>Press O to confirm the deletion.</li> </ol>		
You want to delete the selected result:	1. Choose <b>Selected</b> .		

The results are deleted. You cannot retrieve them any more from the system.

### To delete QC results from the QC History

- 1 Choose Workplace > QC History.
- **2** Press  $\Theta$ .

A screen for selecting results is displayed.

**3** Do one of the following:

lf	Do this	
You want to delete all results in the list: (If a filter was applied, this list would contain only the results that fulfill the filter criterion, and the filter criterion would be part of the button name.)	<ol> <li>Choose List.</li> <li>Press O to confirm the deletion.</li> </ol>	
You want to delete all results that were generated before the first day of the previous month:	<ol> <li>Choose Older than Previous Month.</li> <li>Press O to confirm the deletion.</li> </ol>	
You want to delete the results of the selected control:	1. Choose <b>Selected</b> .	

The results are deleted. You cannot retrieve them any more from the system.

Lot handling

## Lot handling

Both the lot data and its handling depend on what item the lot refers to:

- Calibrators (calibrator lot)
- Controls (QC lot)
- Reagents, diluents (reagent lot)

### Viewing the current lot definitions

### **Calibrator and QC lots**

### To view calibrator or lot data

- 1 Choose Workplace > Lot Data.
- 2 Press either Controls or Calibrators.

A screen is displayed that lists all installed lots.



The following description shows an example of viewing a calibrator lot.

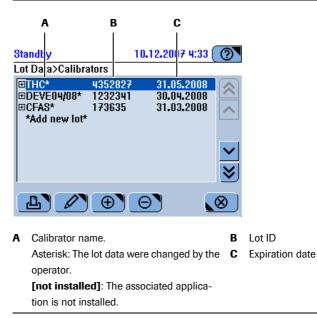


Figure B-30

**3** Select the calibrator and expand the entry.

A Standby 10.12.2007 4:36 0 Lot Data>Calibrators **⊡TH**C<sup>#</sup> 31.05.2008 30.04.2008 4352827  $\approx$ 1232341 173635 31.03.2008 ~ Material Code 401 В ⊟GLU2\* 10.5 1: mmol/L С 2: 0 mmol/L Last Is Water On **⊞ALTL**\* Ъ, Ð  $\Theta$  $\otimes$ Associated test А Calibrator name В Asterisk: The lot data were changed by the  $\ \ \boldsymbol{C}$ Lot values operator. [not installed]: The associated application is not installed.

**4** Select a test and expand the entry.

Figure B-31

Calibrators that were defined or changed manually are marked with an asterisk.

#### **Reagent lots**

You can view the lot number in the reagent set details.

#### To check the status of an individual reagent set

- 1 Choose **Overview** > **I**.
- 2 Press a reagent set icon.

A screen is displayed that contains information on the selected reagent set, including its lot number.

Lot handling

### ► To gain a general overview of the currently defined reagent sets

1 Choose Utilities > Inventory.

The **Bottle sets** list is displayed. It contains all bottle sets that are defined on any of the disks used on this **Cedex Bio** instrument, including sets that were removed from the disk, as long as they were not empty or their expiration date has not passed yet.

			A	В	5	
	andby ottle Se	ts	23.06.2009 20	:01	0	
A A A A A A A	et 1C-2 1C-2 LBT2 LP2S LTL LTL LTL	Lot 60165702 69409701 67505751 60915701 60915701 60630001	Expiration # 03.07.2009 100 05.06.2009 7 03.07.2009 100 01.12.2006 50 05.06.2009 79 08.06.2009 96 17.06.2009 75	8	× × ×	— c
A B		0	ft on which the set i	s	C	${\bf R}$ indicates that the set was removed from the disk.

#### Figure B-32

**2** Select a set and press  $\square$ .

A dialog box is displayed that contains detailed set information.

3 Press ?

A dialog box is displayed that contains set, calibration, and QC information.

- **4** Press 🛞 to close the dialog boxes until the **Utilities** tab is displayed.

### Viewing the lot details of fluids that were used

For sample, calibration, and QC results, you can check to which lot the used fluids belong (reagents, calibrators, controls).

### ► To check context information of a result

**1** Do one of the following:

lf	Do this		
You want to check the context information	<ol> <li>Choose Workplace &gt; Result Overview</li> <li>Select a result.</li> </ol>		
of a sample result:	2. Select a result. 3. Choose $\swarrow > \bowtie$ .		
You want to check the context information	1. Choose Workplace > QC Status.		
of a QC result:	2. Select a result.		
	3. Choose $\nearrow$ > MH.		
You want to check the context information	1. Choose Workplace > Calibrations.		
of a calibration result:	2. Select a result.		
	3. Choose $\nearrow$ > MHP.		

A screen is displayed that lists the lot IDs for the used reagent set, calibrators, and controls. (The following figure shows an example of sample-result lot-data.)

23.06.2009 19:53 Standby QC Status>Detail>Related Information Bottle set ALTL R1 60030501 4242  $\sim$ 1972 SR 18.06.2009 18:48 Calibrator CFAS 18144000 Control 1 PNU 18042500 டு  $\otimes$ 

- **2** Scroll to display the information, if required.
- **3** Press  $\bigotimes$  to close the screen.
- Adding a new lot

Lot data are usually entered manually from the package insert. With the exception of reagent and diluent lots: they can be defined by scanning their barcodes only.



**Defining lots manually requires utmost diligence** During lot definition you need to type data that are directly relevant to result generation.

Make sure that the controls and calibrators are assigned to the application.

Preparation tasks

• See Preparing applications on page B-140.

### To define a control lot

- 1 Choose Workplace > Lot Data.
- 2 Press Controls.

A screen is displayed that lists all installed control lots.

**3** Press 🕀.

A screen is displayed, asking you to scan the barcode transfer sheet or to type the data manually.

Lot handling

**4** Define the lot values.

Do one of the following:

lf	Do this
The lot data are	1. Scan the barcode.
available from a	The screen is displayed again that lists all installed control lots.
barcode	The new lot is selected.
You want to type	1. Press 📖.
the data manually	2. Type the name and press 🔊.
	3. Type the material code and press $\triangleright$ .
	You find this code in <b>Utilities</b> > <b>Applications</b> > select a test > <b>Control</b> .
	4. Type the Lot ID and press 🔊.
	5. Type the expiration date. Use the date format as indicated on
	the screen.
	6. Press 🞯 to confirm the definitions.
	If a lot with the same ID already exists on the system, a screen is displayed, asking you whether you want to replace the existing lot.
	Press 🔘 to confirm the definitions.
	7. Press *Add new test*.
	8. Press 🕀.
	A screen is displayed that contains all tests that use this control
	and that have not yet been assigned to this lot.
	9. Select a test and press <b>&gt;</b> .
	10. Type the value for the mean concentration and press $\gg$ .
	11. Type the value for the standard deviation.
	12. Press 🔘 to confirm the definitions.
	The screen for handling control lot data is displayed again.
	13. Perform steps 7 through 12 for all tests you want to use.

Lot handling

### ► To define a calibrator lot

1 Choose Workplace > Lot Data.

### 2 Press Calibrators.

A screen is displayed that lists all installed calibrator lots.

**3** Press ⊕.

A screen is displayed, asking you to scan the barcode or to type the data manually.

**4** Define the lot values. Do one of the following:

lf	Do th	his
The lot data are available from a bar- code	1	Scan the barcode. Fhe screen is displayed again that lists all installed calibrator lots Fhe new lot is selected.
You want to type the data manually	2. Т 3. Т	Press . Fype the name (up to 10 alphanumeric characters) and press Type the material code and press Refer to the package insert.
	4. Т	Fype the lot ID (up to nine alphanumeric characters) and press $\bigcirc$ Fype the expiration date and press $\bigcirc$ .
	τ	Use the date format as indicated on the screen.
	iı (	Fype the number of calibrators (cups) you need to place on the nstrument. Use the information given on the package insert. Exclude calibrators that use system water from this number. See step 14.)
	7. F	Press 🔘 to confirm the definitions.
		Press *Add new test*.
	9. F	$\operatorname{Press} \ \textcircled{\oplus} \ .$
		A screen is displayed that shows all tests that have this calibrator lefined for them and that are not yet assigned to this lot.
		Press one of the test buttons.
		Define the first calibration value (target value). Define the next calibration value.
		You can define up to six calibration values, and you must define hem in descending order.
	14. S	Select the value for Last is Water.
	b	Press <b>On</b> if you want to calibrate with system water as zero calibrator. (In this case, no cup needs to be placed on the system. (See step 7).
	b	Press <b>Off</b> if you want to use a special zero calibrator for the calibration. (The special zero calibrator needs to be placed on the sample area.)
	15. F	Press 🞯 to confirm the definitions.



Calibration values (target values) in method sheets may be defined in increasing order of concentration. Always define the values in *decreasing* order on the **Cedex Bio** instrument.

Exporting data

## **Exporting data**

→ Utilities > Export

### **Exporting the database**

The database is normally exported as part of the daily backup during the end of shift activities.

• See *Performing the daily backup* on page B-82.

During a database export, the full content of the database is copied to the USB stick.

The database data can be restored to the instrument if required.

• See *Importing a database* on page B-113.

### To export the database

1 Choose Utilities > Export > Database.

A screen is displayed, asking you to insert the USB stick.

- 2 Insert the USB stick.
- **3** Press 🔘.
- **4** Select the directory.
- **5** Press 🕖 to confirm the selection.

The data are copied to the USB stick.

Database files names have the following name format: dba\_yyyymmddhhmmss.tgz.

### **Exporting results**

When exporting results, the full result data are copied into an archive on the USB stick.

The archive has the name format csv\_yyyymmddhhmmss.tgz, and it contains the following files:

Data set	File name
Sample results	res_yyyymmddhhmmss.csv
QC results	qcs_yyyymmddhhmmss.csv
QC history results	qch_yyyymmddhhmmss.csv
Calibration results	cal_yyyymmddhhmmss.csv
Process event log	pev_yyyymmddhhmmss.csv

File format Result data are exported as line-oriented comma-separated value (csv) files.

The following definitions apply:

- Character set: ISO1LATIN1 ANSI ISO8859-1 ISO Latin 1, Western
- Separator: Semicolon (;)
- Element qualifier: Quote (") (ASCII-Code 34 (0x22)
- Empty element: Two quotes ("") (ASCII-Code 34 (0x22)
- Line terminator: Standard Win Style (CRLF) ASCII-Code 13+10 (0x0D + 0x0A)

These files can be processed with any spreadsheet program that can import csv files.

### To export the support data

1 Choose Utilities > Export > Full Results.

A screen is displayed, asking you to insert the USB stick.

- **2** Insert the USB stick.
- **3** Press 🔘.
- **4** Select the directory.

A screen is displayed for selecting a directory.

Press <\*.csv> to archive files only.

Press <\*.\*> to display all files and directories.

**5** Press 🔘 to confirm the selection.

The data are copied to the USB stick. The screen for exporting data is displayed again.

Exporting data

### Interpreting result data

The following tables list, for the sample, calibration, and QC result files, the column headings of the spreadsheet and provide information on the kind of information the columns contain.

res\_yyyymmddhhmmss.csv

Column Title	Description
Instr	Instrument ID (always 30 for Cedex Bio)
Msg	Message Type:
	7 Result for sample, control, or calibration
Index	166 ID of the data set used
App	Application code
Date	Date when the result was calculated
Time	Time when the result was calculated
SW-Version	Software version installed on the instrument
Serial	Serial number of the instrument
Test	Short name of the application
User	User name (\$SYS\$ if Auto Accept is on)
Sample	Sample ID
Order-Time	Time when the order was started
Result	Results of the measurement
Unit	Defined display unit
Flags	Flag name
Rates	Rate value calculated in Abs. (Last Calculation Point - First
	Calculation Point)
Raw1 - Raw40	Absorbance value of cycle 1 through 40 in Abs. (one cycle = 18 s)
isSTAT	Generated as part of a STAT order
Table 33	Explanations on test results

Exporting data

Instrument ID (always 30 for <b>Cedex Bio</b> ) Message Type: 7 Result for sample, control, or calibration						
7 Result for sample, control, or calibration						
151 Calibration data						
152 Std-1 value						
153 Std-2 value						
154 Std-3 value						
155 Std-4 value						
156 Std-5 value						
157 Std-6 value						
Application code						
Date when the result was generated						
Time when the result was calculated						
Software version installed on the instrument						
Serial number of the instrument						
Calibrator ID (C.f.a.s. = 401) and standard number User name ( <b>\$SYS\$</b> if <b>Auto Accept</b> is on) Lot number						
						Calibration type/Target value of corresponding calibrator
						L Lot calibration
S Set calibration						
Defined display unit						
Flag name						
Offset and factor of the calibration curve						
Kc parameter value or absorbance value of cycle 1 in Abs.						
A parameter value or absorbance value of cycle 2 in Abs.						
B parameter value or absorbance value of cycle 3 in Abs.						
C parameter value or absorbance value of cycle 4 in Abs.						
Absorbance value of cycle 5 through 40 in Abs. (one cycle = 18 s)						
Short name of the application						



If the calibration results were accepted using **Use Old**, "???" is displayed as the **Date** and **Time** values, and the corresponding values for **Raw1** - **Raw40** are missing.

### **6 Special operations**

Exporting data

qcs_yyyymmddhhmmss.csv	Column Title	Description						
	Instr	Instrument ID (always 30 for Cedex Bio)						
	Msg	Message Type:						
		7 Result for sample, control, or calibration						
	Index	161 Data values						
	Арр	Application code						
	Date	Date when the result was generated						
	Time	Time when the result was calculated						
	SW-Version	Software version installed on the instrument						
	Serial	Serial number of the instrument						
	Ctrl	Control ID						
	User	User name (\$SYS\$ if Auto Accept is on)						
	Lot	Lot ID						
	Mean	Mean value						
	Result	Result value						
	Unit	Defined display unit						
	Flags	Flag name						
	Rate	Rate value calculated in Abs. (First Calculation Point - Last Calculation Point)						
	Raw1 - Raw40	Absorbance value of cycle 1 through 40 in Abs. (one cycle = 18 s)						
	Test	Short name of the application						
	Ignored	"I" if the QC result was ignored						
	Table 35Explana	nations on QC results						
qch_yyyymmddhhmmss.csv	Column Title	Description						
	Instr	Instrument ID (always 30 for Cedex Bio)						
	Msg	Message Type:						
		8 QC history result						
	Index	151 Data values						
	Арр	Application code						
	Date	Date when the result was calculated						
	Time	Time when the result was calculated						
	SW-Version	Software version installed on the instrument						
	Serial	Serial number of the instrument						
	Ctrl	Control ID						
	User	User name (\$SYS\$ if Auto Accept is on)						
	Lot	Lot ID						
	Mean	Mean value						
	Result	Result value						
	Unit	Defined display unit						
	Flags	Flag name						
	<u>SD</u>	Standard deviation						
	Test	Short name of the application						
		Short name of the application "I" if the QC result was ignored						

Exporting data

pev_yyyymmddhhmmss.csv	
------------------------	--

Column Title	Description
Instr	Instrument ID (always 30 for Cedex Bio)
Msg	Message Type:
	9 Process event
Index	191 Reagent usage changed
	192 Calibrator assigned valued changed
	193 QC assigned valued changed
	199 Other process event
App	Application code
Date	Date when the event took place
Time	Time when the event took place
SW-Version	Software version installed on the instrument
Serial	Serial number of the instrument
Event	Description of the event
User	User name ( <b>\$SYS\$ "system user" = ???</b> if nobody is logged in)
Test	Short name of the application
Table 37	Explanations on process events

The following figure is an example of a test result file that was opened in a spreadsheet tool.

	A	B	C	D	E	F	G	H	1	J	K	L	M	N	0	P	Q	R	S	1
1	Instr	Msg	Index	Арр	Date	Time	SW-Version	Serial	Test	User	Sample	Order-Time	Result	Unit	Flags	Rate	Raw01	Raw02	Raw03	Ra
2	30	7	166	18	20071011	09:49	2.0.0.0708	50429	BILTS	\$SYS\$	1	09:38	1.	2 mg/dL		0.009	0.0029	0.0027	0.0006	
3	30	7	166	571	20071011	09:52	2.0.0.0708	50429	AMY-P	\$SYS\$	1	09:38	41.	1 U/L		0.00612	0.0072	0.0073	0.0006	0
4	30	7	166	685	20071011	09:57	2.0.0.0708	50429	ALTL	\$SYS\$	1	09:38	47.	3 U/L		0.00768	-0.0033	-0.0035	0.00272	0
5	30	7	166	18	20071011	09:49	2.0.0.0708	50429	BILTS	\$SYS\$	1	09:38	1.	2 mg/dL		0.00908	0.0026	0.0027	0.00067	0
6	30	7	166	571	20071011	09:54	2.0.0.0708	50429	AMY-P	\$SYS\$	1	09:38	41.	1 U/L		0.00611	0.0076	0.0075	0.00149	0
7	30	7	166	685	20071011	09:58	2.0.0.0708	50429	ALTL	\$SYS\$	1	09:38	4	6 U/L		0.00745	-0.0051	-0.005	0.00214	0
8	30	7	166	18	20071011	09:49	2.0.0.0708	50429	BILTS	\$SYS\$	1	09:38	1.	2 mg/dL		0.0091	0.0026	0.0029	0.00067	0
9	30	7	166	571	20071011	09:54	2.0.0.0708	50429	AMY-P	\$SYS\$	1	09:38	40.	8 U/L		0.00608	0.007	0.0067	0.00075	
10	30	7	166	685	20071011	09:59	2.0.0.0708	50429	ALTL	\$SYS\$	1	09:39	46.	8 U/L		0.0076	-0.004	-0.0039	0.0022	0
11	30	7	166	313	20071011	11:10	2.0.0.0708	50429	DCBLG	\$SYS\$	111	1 10:59	4.	2 mmol/1		0.121	-0.0027	-0.0028	0.81812	
12	30	7	166	313	20071011	11:11	2.0.0.0708	50429	DCBLG	\$SYS\$	222	2 11:00	6.8	2 mmol/1		0,1461	-0.0027	-0.0027	0.82014	0
12		res_2	007101	1111427	/					-		1								ЪЦ

	Т	U	V	W	X	Y	Z	AA	AB	AC	AD	AE	AF	AG	AH	AI	AJ	AK	AL	
1	Raw04	Raw05	Raw06	Raw07	Raw08	Raw09	Raw10	Raw11	Raw12	Raw13	Raw14	Raw15	Raw16	Raw17	Raw18	Raw19	Raw20	Raw21	Raw22	
2	0.0006	0.0006	0.0006	0.00128	0.00112	0.00112	0.0012	0.00096	0.0012	0.0012	0.00104	0.0012	0.0012	0.0074	0.0086	0.0091	0.0092	0.0096	0.0098	\$
3	0.00082	0.00082	0.00082	0.02161	0.02127	0.02127	0.02144	0.02135	0.02144	0.02135	0.0211	0.02152	0.02152	0.0407	0.0426	0.0442	0.0465	0.049	0.0514	Ε.
4	0.00256	0.00277	0.00261	0.03196	0.03043	0.02939	0.02859	0.0277	0.0269	0.02682	0.02658	0.02618	0.0257	0.9216	0.9199	0.9175	0.9151	0.913	0.9106	į –
5	0.00067	0.0006	0.0006	0.0012	0.00104	0.00112	0.00096	0.00112	0.00104	0.00104	0.00104	0.00112	0.00112	0.0073	0.009	0.009	0.0094	0.0098	0.0098	1
6	0.00157	0.00149	0.00164	0.02229	0.0222	0.02195	0.02229	0.02237	0.02203	0.02203	0.02178	0.02203	0.02212	0.0418	0.0429	0.0454	0.0475	0.0497	0.0518	4
7	0.00209	0.00214	0.0023	0.03132	0.02987	0.02883	0.02787	0.02674	0.02682	0.02642	0.02586	0.02546	0.02513	0.921	0.9191	0.9167	0.9146	0.9121	0.9099	ł.
8	0.00052	0.00075	0.00067	0.00112	0.00128	0.00112	0.0012	0.0012	0.00112	0.00104	0.00128	0.00104	0.0012	0.0073	0.0087	0.0092	0.0095	0.0098	0.0098	, –
9	0.0009	0.0009	0.00104	0.02161	0.02135	0.02144	0.02152	0.02135	0.02152	0.0211	0.02118	0.02127	0.02161	0.0405	0.0424	0.0444	0.0464	0.0487	0.0509	1
10	0.00225	0.0023	0.00214	0.03164	0.03043	0.02899	0.02795	0.02714	0.02674	0.02634	0.02578	0.02562	0.02522	0.9189	0.9168	0.9149	0.9126	0.9103	0.9078	1
11	0.8171	0.81643	0.81596	0.83628	0.83541	0.83476	0.8344	0.8336	0.83317	0.83309	0.83266	0.83252	0.83259	0.83273	0.83288	0.83338	0.83346	0.83295	0.83302	
12	0.81893	0.81825	0.81737	0.83454	0.83375	0.83302	0.8328	0.83252	0.8323	0.83194	0.83186	0.83179	0.83208	0.83259	0.83252	0.83208	0.83223	0.8323	0.83252	<u>.</u>
12	> H res	20071011	111427/									T							,	Ĩ
	AM	AN	AO	AP	AQ	AR	AS	AT	AU	AV	AW	AX	AY	AZ	BA	BB	BC	BD	BE	T
1	Raw23	Raw24	Raw25	Raw26	Raw27	Raw28	Raw29	Raw30	Raw31	Raw32	Raw33	Raw34	Raw35	Raw36	Raw37	Raw38	Raw39	Raw40	IsSTAT	
2	0.0099	0.01	0.0102	0.0102	0.0103	0.0102	0.0102	0.0104	0.0104	0.0105	0.0102	0.0103	0.0103	0.0106	0.0102	0.0103	0.0108	0.0102	0	1
3	0.0531	0.0554	0.0575	0.0591	0.0611	0.0629	0.065	0.0667	0.0687	0.0702	0.0723	0.0739	0.076	0.0778	0.0796	0.0812	0.0829	0.085	0	1
4	0.9083	0.9055	0.9036	0.9009	0.8986	0.8962	0.8945	0.8922	0.8896	0.8869	0.8849	0.8831	0.8806	0.879	0.8768	0.8741	0.8721	0.87	0	j .

5	0.01	0.01	0.0101	0.0102	0.0102	0.0104	0.0103	0.0106	0.0103	0.0105	0.0106	0.0103	0.0106	0.0101	0.0104	0.0104	0.0103	0.0105	0
6	0.0539	0.056	0.058	0.0599	0.0615	0.0633	0.0653	0.0672	0.069	0.0709	0.0726	0.0742	0.0761	0.078	0.08	0.0817	0.0836	0.0855	0
7	0.9067	0.9049	0.9024	0.8999	0.8979	0.8958	0.8937	0.8915	0.8886	0.8872	0.8855	0.8827	0.8807	0.8785	0.8759	0.8737	0.8722	0.8701	0
8	0.0101	0.0101	0.0103	0.0103	0.0102	0.0104	0.0104	0.0103	0.0104	0.0105	0.0103	0.0106	0.0103	0.0105	0.0104	0.0105	0.0105	0.0105	0
9	0.0531	0.055	0.0572	0.0588	0.0607	0.0625	0.0645	0.0663	0.0682	0.07	0.0714	0.0736	0.0753	0.0774	0.0789	0.081	0.0827	0.0846	0
10	0.9054	0.9031	0.9003	0.8977	0.896	0.8937	0.8915	0.8892	0.8869	0.8842	0.8821	0.8807	0.8789	0.8769	0.8743	0.8722	0.87	0.8681	0
11	0.83324	0.8377	0.8251	0.811	0.7959	0.7801	0.7641	0.7477	0.7317	0.7148	0.6982	0.6811	0.6642	0.6473	0.6304	0.6137	0.597	0.5805	0
12	0.83259	0.8362	0.8269	0.8182	0.8089	0.8	0.7905	0.7811	0.7712	0.7616	0.7521	0.7423	0.7329	0.723	0.7133	0.7032	0.6938	0.6837	0
12		00710111										14							- DIC
14 4	H\res_2	00/101111	1427/																,

Figure 3

Example of spreadsheet representation of a test result raw data file

#### **Roche Diagnostics**

### **Exporting log files**

During troubleshooting, you may be asked to export the log files and to send them to by the service representative for examination.

The log-files contain the alarm messages and system logs.

### ► To export the log files

1 Choose Utilities > Export > Log Files.

A screen is displayed, asking you to insert the USB stick.

- 2 Insert the USB stick.
- **3** Press 🔘 to confirm the insertion.

A screen is displayed for selecting a directory.

Press \*.tgz to display Cedex Bio system files only.

Press \*.\* to display all files and directories.

- **4** Select the directory.
- **5** Press 🕖 to confirm the definitions.

The data are copied to the USB stick. The screen for exporting data is displayed again.

Importing data

## **Importing data**

### → Utilities > Import

You can import the following kinds of data:

- Application data (application definitions)
- Software (software updates)
- Database (as exported using **Utilities** > **Export** > **Database**)
- Certificates (digital records that ensure the authenticity of reagent barcodes)
- Extra wash cycle and mixing rule definitions

### Importing applications

Importing an application consists of two steps:

- 1. Importing the data to the instrument by reading a barcode or a data file. This step stores the data on the instrument.
- 2. Installing the application. This step activates the application on the instrument and so makes it available for use.

### To import application data

- **1** Do one of the following:
  - Choose Utilities > Applications > Laboratory Parameters, then continue with step 2.
  - Choose Utilities > Import > Application, then continue with step 4.
- **2** Press ⊕.
- 3 Press Import Application.

A screen is displayed, asking you to scan the barcode or to import the data from the USB stick.

**4** Do one of the following:

lf	Do this
You intend to scan the barcode:	1. Scan the barcode from the barcode sheet.
You intend to import from the USB stick:	1. Press 📀.
	2. Insert the USB stick.
	3. Press 🖉.
	4. Select the directory that contains the application file.
	You recognize the application pack- ages by their file extension .tsb.
	5. Press 🔘 to confirm the selection.

The system checks whether there is already an application on the system with identical application code and short name.

#### Importing data

The following table lists the basic situations.

lf	This happens					
Both, the application code and the short name exist:	The laboratory parameters, for example the short name, are retained from the existing application.					
	The application definitions are replaced on the system by the new ones.					
The application code exists, and the short name does not exist for	The laboratory parameters of the new application are used, for example the short name.					
any application on the system:	The application definitions are replaced on the system by the new ones.					
The short name exists, but the application code does not:	A screen is displayed for changing the short name to a unique short name.					
	If you do not change the short name, the application is not imported.					



-`Ċ`-

If you want to replace both the laboratory parameters and the application definitions, you first need to uninstall and delete the existing application and then to import the new application.

- If **TRL Check** is *not* activated, the value for the lower limit of the measuring range is overwritten when re-importing the application.
- If **TRL Check** *is* activated, the value for the lower limit of the measuring range is *not* overwritten when re-importing the application. Check the relevant application documentation for the exact value, you may need to adjust it manually. See *Calculation definitions* on page B-142.
- There are applications where the lower limit of the measuring range cannot be changed, e.g. with ALBU2.
- **5** If you imported from a USB stick, wait until the copy action is complete before you remove the USB stick.

### 

### To install an application

- 1 Choose Utilities > Applications > Laboratory Parameters.
- 2 Select an application whose name is in square brackets, for example [GLU2].
- 3 Press ⊕.
- 4 Choose Install Application.

The system checks whether there is already an application installed on the system with identical application code or short name.

In the applications list, the application name is displayed without brackets and in the profile details the short name is displayed.

The profiles that use this application are included in the test selection screen, provided that all applications of the profiles are installed.

**5** Prepare the application.

• See *Preparing applications* on page B-140.

### Importing software

This function is usually used for installing software updates.

### To import system software

1 Choose Utilities > Import > Software.

A screen is displayed, asking you whether you want forced download of controller software, that is whether you want to overwrite the current controller firmware irrespective of its current status.

- **2** Do one of the following:
  - Press 🞯 to overwrite the controller firmware .
  - Press 🛞 to perform standard import.

A screen is displayed, asking you to insert the USB stick.

- **3** Insert the USB stick.
- 4 Press 🙆.
- **5** Select the software package.

You recognize the software packages by their file extension .tar.

**6** Press 🔘 to confirm the selection.

The software is installed on the system. A message will inform you when the installation is complete.

The system will automatically reboot.

When the import is complete, the **Overview** tab is displayed, and the system is in Standby status.

### Importing a database

Compatibility	the imported database wil	l be converted to the ver rted database is more rea	an the one on the target system, rsion on the system. cent than the one on the target				
Typical situations	Importing a database is main						
	<ul> <li>You want to restore your existing system to a previous status, for example from a backup version of the database that was generated on the same system.</li> </ul>						
	• You want to install a know	•	·				
	The following table illustrates	s the different effects in t	he two situations.				
	Data item	Identical source and	Different source and				
		target system	target systems				
	Installed applications	Overwrite	Write, overwrite if present				
	System configuration	Overwrite	Write, overwrite if present				

 Table B-11
 Effects on existing data when importing a database

Overwrite

Bottle set inventory

Write, overwrite if present

Data item	Identical source and target system	Different source and target systems
Bottle sets	Mark as removed	Mark as removed
Lot information for calibra- tors and controls	Overwrite	Write, overwrite if present
Sample orders and results	Overwrite	Delete if present
QC orders and results	Overwrite	Delete if present
QC history results	Overwrite	Delete if present
Calibration orders and results	Overwrite	Delete if present
Abs. Air/Water Calibration	Delete	Delete if present
Administrator password	Reset to default value	Reset to default value
Electrode definitions	Delete	Delete if present
ISE fluid bottle definitions	Delete	Delete if present

 Table B-11
 Effects on existing data when importing a database (Continued)

### To import the database

1 Remove all bottles of all bottle sets of all disks defined on the target system, using the appropriate software features.

• See *Preparing the reagents* on page B-22.

2 Choose Utilities > Import > Database.

A confirmation screen is displayed, informing you about the major effects of the import.

• For details see Table B-11 on page B-113.

**3** Press 🔘.

A screen is displayed, asking you to insert the USB stick.

- 4 Insert the USB stick.
- 5 Press 🙆.
- **6** Select the database file.

Typically, the file format looks like this: dba\_yyyymmddhhmmss.tgz.

Press \*.tgz to display Cedex Biosystem files only.

Press \*.\* to display all files and directories.

**7** Press 🞯 to confirm the selection.

When the data are imported, a message is displayed, asking you to restart the system.

**8** Choose  $\bigcirc$  > **Restart** to restart the system.

Importing data



### After the database installation

- Load all bottle sets that are defined, using the appropriate software features.
- Perform Abs. Air/Water Calibration maintenance action.
- If **TRL Check** is activated, the value for the lower limit of the measuring range is *not* overwritten when importing the database. Check the relevant application documentation for the exact value, you may need to adjust it manually. See *Calculation definitions* on page B-142. (Note that there are applications where the lower limit of the measuring range cannot be changed, e.g. ALBU2.)

### Importing certificates

Certificates are digital records that ensure the authenticity of reagent barcodes. On each **Cedex Bio** instrument there must be a certificate installed.

### To import certificates

1 Choose Utilities > Import > Certificate.

A screen is displayed, that lists the currently installed certificates.

**2** Press **♦**.

A screen is displayed, asking you to insert the USB stick.

- 3 Insert the USB stick.
- 4 Press 🔘.
- 5 Select the certificate.

You recognize a certificate by its file extension .prm.

**6** Press 🖉 to confirm the selection.

The certificates are installed on the system.

The screen is displayed again, that lists the currently installed certificates.

### Importing automatic mixing and extra wash cycle definitions

Mixing and extra wash cycle information is contained in a separate barcode on a barcode transfer sheet.

### ► To import mixing and extra wash cycle information

1 Choose Utilities > Import > EWC/Mixing.

A screen is displayed, asking you to scan the barcode.

**2** Scan the barcode.

The data are installed on the system. A message will inform you when the installation is complete. The screen for importing data is displayed again.

 For information on defining extra wash cycles see *Defining extra wash cycles* on page B-168.

Preparing a new disk

## Preparing a new disk



#### Injury through reagents and other working solutions

Direct contact with reagents, cleaning solutions, or other working solutions may cause personal injury.

When handling reagents, exercise the precautions required for handling laboratory reagents, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics reagents and cleaning solutions.

#### Infection by biohazardous materials

Contact with samples containing material of human origin may result in infection. All materials and mechanical components associated with samples of human origin are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

# Personal injury or damage to the analyzer due to contact with instrument mechanism

Contact with moving parts of the instrument may cause personal injuries, stop the instrument from processing, and cause damage to instrument parts.

- Do not touch moving parts during instrument operation.
- Keep all covers closed, operate them as instructed on the screen.


NOTICE

#### Damage to the reagent disk

The reagent disk is designed to handle reagents while it is loaded on the instrument. The cover is equipped with a locking mechanism.

Remove or insert bottles only when the reagent disk is loaded on the instrument and always use the software functions to perform these tasks.

You can use up to eight different reagent disks on one **Cedex Bio** instrument. Each reagent disk is equipped with numbered tabs. For automatic disk identification by the instrument, one—and only one—of these tabs is removed. The number of this removed tab is the disk ID. When you label the disk, make sure that the number on the label corresponds to that of the removed tab.

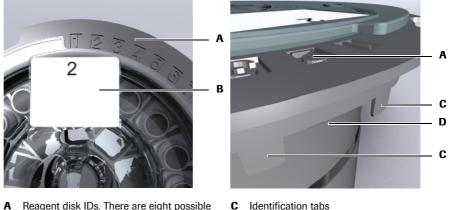

D

disk recognition

The tab has been removed for automatic

- A Reagent disk IDs. There are eight possible IDs
- **B** Disk label. The number must correspond to the reagent disk ID.

Figure B-38 Reagent disk ID

#### To prepare a new reagent disk

*Defining the disk ID* **1** Be sure to select an ID that is not used by any other reagent disk you intend to use on this **Cedex Bio** instrument.

You can choose a number between one and eight.

- 2 Print the number on a label and stick the label on the disk (A).
- **3** Using pliers, break off the ID tab (C, D) with the same number as the one printed on the label.
- *Loading the disk* **4** On the system, choose Overview > <u>()</u>.

A disk overview screen is displayed.

5 Press 🔃.

A screen is displayed, asking you to insert the disk.

- 6 Open the main cover.
- 7 Place the reagent disk in the reagent cooler.

Make sure the reagent port faces the front and align the cut-outs with their counterparts on the reagent cooler.

The system automatically detects that a disk was inserted.

A screen is displayed, asking you to close the main cover.

**8** Close the main cover.

At this stage, the system identifies the disk.

A screen is displayed that shows the status of the reagent sets.

Preparing a new disk

Loading reagent sets	9	Press	100
----------------------	---	-------	-----

A screen is displayed, asking you to scan the bottle barcode.

- **10** Scan the barcode on the bottle.
- **11** Remove the cap from the bottle and place it in the reagent bay.
- **12** Press 🔘 to confirm the insertion.

If you confirm without placing the bottle, the system assumes that the bottle is placed.

If you press 🛞 after placing the bottle, the position is deemed empty.

When the first reagent bottle of a set is loaded, the reagent set is defined. From this moment on, the reagents are handled as part of the set. You no longer handle them as individual reagents.

**13** Do one of the following:

lf	Do this		
There is another bottle	A screen is displayed, asking you to scan the bottle.		
belonging to the set:	1. Scan the bottle		
	2. Insert the bottle.		
	3. Press $\bigcirc$ to confirm the insertion.		
All bottles of the set are inserted:	1. Close the main cover.		
	On the screen with the reagent set buttons, the button for the new set is now present.		

14 Press 🛞.

**15** On the disk overview screen, press **10** to load the next reagent. Proceed as described in steps 10 through 15.

## Assigning tests to test tabs

The system provides two order modes: Easy and Full. When working in Easy mode, the test selection screen consists of one panel, if working in Full mode, the screen contains up to six panels, each is identified by a tab.

To make a test available for selection from the test selection screen, you need to assign it to a test tab. This is done in two steps:

- 1. Naming the tabs (if required).
- 2. Assigning the tests to the tabs.



You need Lab Manager or Administrator rights for assigning tests to tabs.

Assignment when importing applications

When you import applications, the tests are assigned to the tabs as follows:

- If you work in Easy mode, the tests are added to the Easy panel. (If the panel is full, they will not be displayed.)
- If you work in Full mode, the tests are added to the Easy panel (if there is space) and to the first Full mode tab. If there is not enough space on the first Full mode tab, the test is added to the next tab that has space available.

### To name a test tab

- 1 Choose Utilities > Configuration > Workflow.
- 2 Expand the Workflow entry, scroll down, and select Test Tab 1...6 Name.
- 3 Press 🖉.

A screen for typing text is displayed.

• For details on typing text see *Typing text* on page A-74.

4 Type up to four characters.

This is the name of the tab on the screen for assigning tests.

**5** Press 🔘.

The tab name is now available on the screen for assigning tests. See below.

### To assign a test to a tab

- 1 Choose Utilities > Applications > Laboratory Parameters.
- 2 Select the test you want to assign to a tab.
- 3 Press 🖉.
- **4** Select **General** and expand the entry.
- 5 Select Test Tabs.
- 6 Press 🖉.

A screen is displayed that provides a button for each of the possible tabs.

The **Easy** button is always available. It is the standard panel for working in **Easy** order mode. (You cannot rename it.)

- 7 Select the buttons for all the tabs where the test should be displayed.
- 8 Press 🔘.

The tests are now available on the corresponding tabs on the test selection screen.

### To remove a test from a tab

- 1 Choose Utilities > Applications > Laboratory Parameters.
- **2** Select the test you want to remove.
- 3 Press 🖉.
- 4 Select General and expand the entry.
- 5 Select Test Tabs.
- 6 Press 🖉.
- 7 Cancel the selection of the button of the tab to which the test is assigned to.
- **8** Press 🙆.

The test is no longer available from the corresponding tab on the test selection screen.

<u>`</u>\

To move a test from one tab to another, delete it from the original tab and assign it to the new tab.

## Deleting bottle sets from the Inventory list

The **Inventory** list serves to gain an overview on the status of all bottle sets that are defined on any of the reagent disks used on the **Cedex Bio** instrument.

You can delete bottle sets from the list; this would be appropriate if you removed a set from the disk and did not intend to use it again.



- Before you can delete a bottle set from the **Inventory** list you need to remove it from the disk.
- Bottle sets that have been removed from the disk and whose expiration date had passed more than 30 days ago are automatically deleted from the **Inventory** list, provided there is still a valid set for the same fluid on board. If this were not the case, the set that was loaded last would remain in the inventory.

#### ► To delete a bottle set from the Inventory list

Removing the bottle set

**1** Remove the bottle set from the disk.

• *Removing a reagent set* on page B-25.

If the bottle set that you want to delete is not on the currently installed disk, you need to change the disk first.

• *Removing the reagent disk* on page B-87, *Preparing the reagent disk* on page B-21.

#### *Deleting the bottle set* **2** Choose **Utilities** > **Inventory**.

The **Bottle sets** list is displayed. It contains all bottle sets that are defined on any of the disks used on this **Cedex Bio** instrument, including sets that were removed from the disk, as long as they were not empty or their expiration date has not passed yet.

Set	Lot	Expiration #	D	
A1C-2	60165702	03.07.2009 10	03	
A1C-2	60165702	05.06.2009 7	8	
ALBT2	69409701	03.07.2009 10	03	
ALP2S	67505751	01.12.2006 50	8	
ALTL	60915701	05.06.2009 79	R	V
ALTL	60915701	08.06.2009 96	R	
ALTL	60630001	17.06.2009 75	3	≫

#### Figure B-39

In the D column, removed bottle sets are marked with R (A).

- **3** Select the bottle set you want to delete.
- 4 Press ⊖

A confirmation dialog box is displayed. If the set cannot be deleted a message will inform you.

**5** Press 🔘 to confirm the deletion.



It is not possible to place a bottle set on a reagent disk again, once it was deleted from the **Bottle sets** list.

Refilling printer paper

# **Refilling printer paper**

<u>`</u>\.

You can refill the printer paper any time, provided printing is not in progress.

#### ► To check the printer status

1 On the **Overview** tab, press the System status button.

The printer button  $\square$  is red if the printer is out of paper.

- **2** Scroll to display the **Printer** entry and read the text.

#### • To replace the printer paper

- **1** Make sure the printer is not currently printing.
- **2** Open the printer panel.

Be sure to press down the release button (A) firmly *before* you pull the panel. The panel should open without resistance.




- **3** Lift the empty printer paper roll from its holder.
- **4** Remove the pin from the roll.
- **5** Insert the pin in the new roll.
- **6** Place the new roll on the holder.

Make sure the paper unrolls at the top and towards you.



7 Insert the paper in the slot in the printer panel and pull some through.

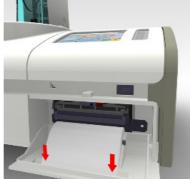

**8** Close the printer panel.

The system feeds some paper.

If the system ran out of paper during printing, it will resume printing.

**9** On the **Overview** tab, press the System status button.

The printer button 🕒 should now be green, and the status description should be **OK**.


# **Removing condensation water from the reagent cooler**

#### To remove water from the reagent cooler

- **1** Make sure the system is in **Standby** status.
- **2** Choose **Overview**  $> \square > \square$ .
- **3** Open the main cover.
- **4** Remove the reagent disk.

Be sure not to tilt it and place it on an even surface.

**5** Wipe the inside of the reagent cooler with a cloth or paper to remove the water.

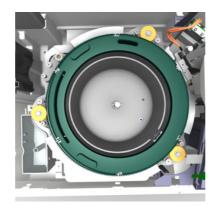

- **6** Insert the reagent disk.
- **7** Close the cover.

# **Replacing the probe**

If the probe is bent, broken, or corroded, you need to replace it.

Tools and materials required

- Probe and tubing set
- Tube with ISE Deproteinizer
- Tube with Activator
- Glass beaker



#### Injury through working solutions

Direct contact with cleaning solutions or other working solutions may cause personal injury. When handling such solutions, exercise the precautions required for handling them, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics cleaning solutions.

#### Infection by biohazardous materials

Contact with samples containing material of human origin may result in infection. All materials and mechanical components associated with samples of human origin are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- · If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

#### To remove the probe

- 1 Make sure the system is in **Standby** status.
- 2 Shut down the system and switch off the instrument.
- **3** Remove the transfer head cover.

Press the release buttons on both sides and lift.

4 Release the tube leading to the probe from all tubing clips.

Replacing the probe

**5** Remove the sensor tubing adapter (A).

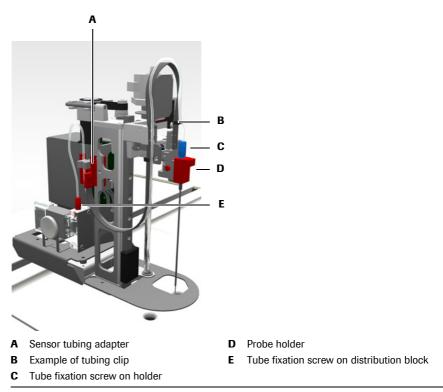


Figure B-40

**6** Remove the probe holder (D) from the carriage.

Press the release button on the side of the probe holder carriage and lift the probe holder with the probe.

- 7 Unscrew the probe from its holder and remove it from it.
- 8 Place the probe in the beaker and the probe holder on a clean surface.
- **9** Unscrew the tubing from the distribution block (E).
- 10 Lift the tubing and wait until all fluid has run into the beaker.
- 11 Dispose of the probe assembly. Treat it as biohazardous waste.

#### To install the new probe

- 1 Carefully insert the probe in the probe holder and fasten the screw that fixes the tubing to the holder (C).
- **2** Reinstall the probe holder with the probe.

Press the release button on the side of the probe holder carriage while you insert the probe holder. Release the button when the holder is inserted.

Push the holder firmly down until the release button latches on.

3 Reinstall the sensor tubing adapter.

Push down until the clips engage.

- **4** Screw the tubing to the distributor block.
- **5** Fix the tube to all tubing clips. Start near the probe holder.

Connecting and disconnecting the external fluid containers

**6** Install the transfer head cover.

Push down until the release buttons latch on.

**7** Switch on the system.

The system performs internal checks and routines.

The startup phase may take a few minutes.

When the instrument is ready for use, the **Overview** tab is displayed, and its status is **Standby**.

- 8 Log on the system.
- 9 Perform the **Deproteinize Probe** maintenance action.

• See *Deproteinize the probe* on page C-10.

# Connecting and disconnecting the external fluid containers

Before switching on the system, make sure that all external fluid containers are placed on the rack and properly connected.

#### To connect the water container

- 1 Place the white full water container on the tray. (Always use purified water of the quality specified in section *Technical specifications*.)
- **2** Connect the water tubing to the connector on the instrument by pushing the plug firmly in the socket.

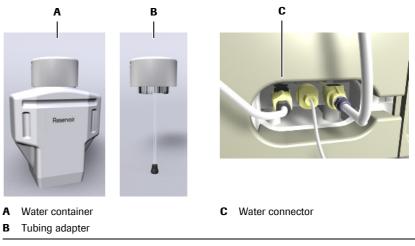



Figure B-41

**3** Insert the tubing adapter in the container and push it down firmly.

Connecting and disconnecting the external fluid containers

#### ► To connect the waste container

- **1** Place the empty yellow waste container on the tray.
- 2 Connect the waste tubing by pushing the plug firmly in the socket.

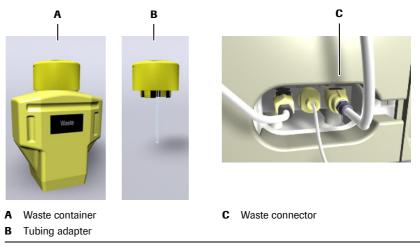




- **3** Insert the tubing adapter in the container and push it down firmly.

#### To connect the cleaner bottle


**1** Place the bottle on the tray.

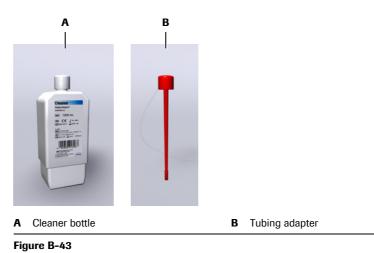

2 Thread the tube through the protective sleeve, make sure that the wider end of the sleeve points away from the instrument.

Connecting and disconnecting the external fluid containers


**3** Screw the cleaner tubing to the connector on the instrument (C). Do not fasten the screw too tight.

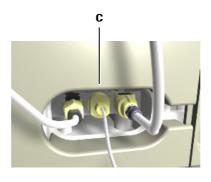

C Cleaner connector

#### Figure B-44

**4** Push the protective sleeve firmly on its base on the instrument (D).

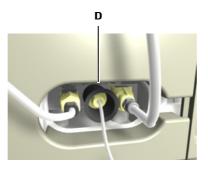

**D** Cleaner connector with protective sleeve


Figure B-45

- **5** Insert the tubing adapter in the container and push it down firmly.


#### ► To disconnect a tube from the instrument

- **1** Remove the right side panel.
- **2** Do one of the following:

lf	Do this		
You want to disconnect the water tubing:	Press down the release clamp (A) on the connecting socket and pull the connector away from the socket.		
You want to disconnect the cleaner tubing:	<ol> <li>Pull out the protective sleeve (B) and move it along the tube far enough that you can grip the connecting screw.</li> <li>Turn the connecting screw (D) counterclockwise until the connector is released.</li> </ol>		
You want to disconnect the waste tubing:	Press down the release clamp (C) on the connector and pull the connector away from the socket.		

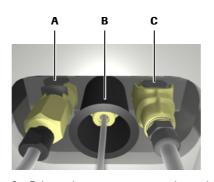

Cleaner connector with protective sleeve

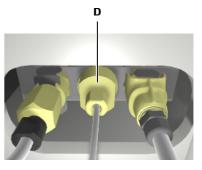

 $\textbf{A} \quad \text{Release clamp on water connecting socket} \quad \textbf{C} \quad \text{Waste connector release clamp}$ D Cleaner connector screw

Figure B-46

В

# Adjusting the touchscreen

With a touchscreen, it is important that the point where you press the screen corresponds exactly with its hardware equivalent. If this were not the case, pressing a screen item such as a button might not lead to the expected result.

#### To adjust the touchscreen

**1** Press the 🞯 global button.

A screen is displayed that offers several kinds of stopping. Ignore it.

**2** Press the 🞯 global button again.

A screen for adjusting the screen is displayed.

lack cross.

**3** Press exactly in the center of the black cross.

When you have touched it properly, it turns yellow and the next cross turns black.

- 4 Do the same for all crosses.
- 5 When all crosses are yellow, a message is displayed, informing you that you need to reboot the system before the new calibration values take effect.
- **6** Press 🔘 to close the calibration screen.
- **7** Restart the system.
  - Press 🕲 .
  - Press Restart.

The system shuts down and automatically restarts.

- Wait until the Overview tab is displayed and the system is in Standby status.

# **Cleaning the touchscreen**

Because the touchscreen gets easily soiled you should clean it regularly.

*Tools and materials required* 

- $\Box$  70% ethyl alcohol
- □ Tissues
- □ Protective gloves

#### **Roche Diagnostics**

Cleaning the touchscreen

#### ► To clean the touchscreen

- **1** Pour or spray some alcoholic solution on a tissue.
- **2** Wipe the screen.

Exert as little pressure as possible.

**3** Wipe the screen with a clean tissue.

# Configuration

# Integrating the system in your environment

In this chapter, you will find information on how to define the way you work with the **Cedex Bio** instrument.

In this chapter	Chapter 7
Introduction	B–135
Viewing values	B–135
Changing values	B–135
Applications	B–137
Installing applications	B–137
Activating and deactivating applications	B–139
Preparing applications	B-140
Uninstalling applications	B-143
Deleting applications	B-144
Defining the processing order	B–145
Defining ratio applications	B-146
Profiles	B-148
Defining profiles	B-148
Adding tests to a profile	B-148
Removing tests from a profile	B-149
Deleting profiles	B–149
Defining control definitions and lots	B–149
Defining calibrator definitions and lots	B–150
Short guide to application definitions	B–152
Version	B–152
Code	B–152
General application definitions	B–152
Calibration definitions	B–153
Control definitions	B–154
Calculation definitions	B–155
Calculation definitions (ratio applications)	B–155
Laboratory correlations (ratio applications)	B-156
Reference range definitions (ratio applications)	

Table of contents

Result conversion definitions	B–156
Configuration	B–157
Changing your password	B–157
Scheduling maintenance actions	B–158
User management	B–159
Defining a user	B–159
Deleting a user	B–159
Editing the acceptable flags list	B–160
Mapping the host codes	B–160
Short guide to configuration definitions	B–161
Workflow definitions	B–161
Host definitions	B–163
Result handling	B–165
Calibration	B–165
QC	B–165
System	B–166
Date and time	B–167
Maintenance	B–167
Abs adjustment	B–167
Defining extra wash cycles	B–168
Activating, deactivating and deleting extra wash cycles	B–169
Short guide to extra wash cycle definitions	B–170
Extra wash cycle definitions	B–170
Short guide to reagent mixing rules	B–171
Reagent mixing rule definitions	B–171
Printing mixing rules	B–171
Deleting mixing rules	B–172

Introduction

# Introduction

Ù.

It is the users' sole responsibility to determine whether they activate the **TRL Check** feature, and if they do so to define the lower measuring range in accordance with the country specific requirements. If the users activate the **TRL Check** feature, they are also responsible for validation of the ranges they define.

• See Calculation definitions on page B-155 and Installing applications on page B-137.


The process for viewing and changing configuration and application values is the same. The following procedures use the example of a calibration value.

## **Viewing values**

#### To view values

- 1 Choose Utilities > Applications > Laboratory Parameters.
- **2** Expand the list.
  - Select a list item marked with 🛨 .
  - Press 🔘.
  - Select a sublist item marked with 🛨.
  - Press 🔘.
  - Use the scrollbar, if required, to display the items you are interested in.

The definition items and their current values are displayed.

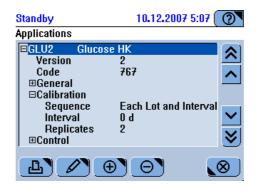

]

## **Changing values**

If the change of a value results in one or several other values needing to be changed, a wizard is started when you change the first value. You recognize a wizard by the presence of the  $\searrow$  and  $\bigotimes$  buttons.

#### ► To change a value

- 1 Choose Utilities > Applications > Laboratory Parameters.
- 2 Expand the list to display the items and their values.

#### **Roche Diagnostics**

Introduction

- **3** Select an item that shows a value.
- 4 Press 🖉.

A screen is displayed where you either can type a new value or select one by pressing its corresponding button.

Standby 10.12.2007 5:40 🤇	Standby	10.12.2007 5:40 🕐
Applications>Edit>Sequence	Applications>Edit>Interv	al
No Interval	24	h
L	789E	
Interval Only	4 5 6 -	
·	1 2 3 +	
	0 . C ←	
<u> </u>	) 🔍	) <b>» (@ (</b> & )

- **5** Type the new value or press a value button.
- **6** Do one of the following:

lf	Do this		
You are in a wizard and want to change a	1. Press ≫.		
further value:	A screen is displayed for changing the value.		
	2. Type the new value or press a value button.		
	3. Continue with step 6.		
	<ul> <li>4. When <i>≫</i> is no longer active, press</li> <li><i>⊘</i> to confirm the changes.</li> </ul>		
This was the only value you want to change:	1. Press 🖉 to confirm the changes.		

## **Installing applications**

To make available a new application on a **Cedex Bio** instrument, you need to perform the following steps:

1. Import the data to the instrument by reading its barcode or by reading them from a USB stick. This step saves the data on the instrument.

It is possible to import a new version of an application with identical application code, but you cannot install it.

- To import application data on page B-137.
- 2. Install the application. This step activates the application on the instrument and so makes it available for use.
  - *Installing applications* on page B-137.
- 3. Prepare the application. This step associates the application with lot data and auxiliary fluids such as calibrators and controls.
  - *To prepare an application* on page B-140.

#### To import application data

- **1** Do one of the following:
  - Choose Utilities > Applications > Laboratory Parameters, then continue with step 2.
  - Choose Utilities > Import > Application, then continue with step 4.
- **2** Press 🕀.
- 3 Press Import Application.

A screen is displayed, asking you to scan the barcode or to import the data from the USB stick.

**4** Do one of the following:

lf	Do this		
You intend to scan the barcode:	1. Scan the barcode from the barcode sheet.		
You intend to import from the USB stick:	<ol> <li>Press</li></ol>		
	<ul><li>You recognize the application packages by their file extension .tsb.</li><li>5. Press  to confirm the selection.</li></ul>		

The system checks whether there is already an application on the system with identical application code and short name.

The following table lists the basic situations.

lf	This happens
Both, the application code and the short name exist:	The laboratory parameters, for example the short name, are retained from the existing application.
	The application definitions are replaced on the system by the new ones.
The application code exists, and the short name does not exist for	The laboratory parameters of the new application are used, for example the short name.
any application on the system:	The application definitions are replaced on the system by the new ones.
The short name exists, but the application code does not:	A screen is displayed for changing the short name to a unique short name.
	If you do not change the short name, the application is not imported.



-`Ċ´-

If you want to replace both the laboratory parameters and the application definitions, you first need to uninstall and delete the existing application and then to import the new application.

- If **TRL Check** is *not* activated, the value for the lower limit of the measuring range is overwritten when re-importing the application.
- If **TRL Check** *is* activated, the value for the lower limit of the measuring range is *not* overwritten when re-importing the application. Check the relevant application documentation for the exact value, you may need to adjust it manually. See *Calculation definitions* on page B-142.
- There are applications where the lower limit of the measuring range cannot be changed, e.g. with ALBU2.
- **5** If you imported from a USB stick, wait until the copy action is complete before you remove the USB stick.

#### 

#### To install an application

- 1 Choose Utilities > Applications > Laboratory Parameters.
- 2 Select an application whose name is in square brackets, for example [GLU2].
- 3 Press ⊕.
- 4 Choose Install Application.

The system checks whether there is already an application installed on the system with identical application code or short name.

In the applications list, the application name is displayed without brackets and in the profile details the short name is displayed.

The profiles that use this application are included in the test selection screen, provided that all applications of the profiles are installed.

- **5** Prepare the applications.
  - See *Preparing applications* on page B-140.

## Activating and deactivating applications

When an application is installed on the system, it is automatically activated.

• *Installing applications* on page B-137.

You can deactivate an application to temporarily make unavailable the test on the test panel. All associated fluids and data remain unchanged. You can activate the application later again and continue using the test.

#### To activate or deactivate an application

- 1 Choose Utilities > Applications > Laboratory Parameters.
- **2** Expand the application entry.
- **3** Expand the **General** entry.
- 4 Select Active.
- 5 Press 🖉.
- 6 Choose **On** if you want to use the application.

Choose **Off** if you do not want to use the application.

- **7** Press  $\oslash$ .

## **Preparing applications**

Perform following steps for each new application. (Use the wizards to make the definitions.)

For the general procedure of making changes to definitions, see *To change a value* on page B-135.

#### ► To prepare an application

- 1 Choose Utilities > Applications > Laboratory Parameters.
- **2** Expand the application entry.
- *Code* The Roche application code identifies the application. It is used for host communication. If you work with a different code for a particular application, you need to map the Roche application code to your own code.

• See *Mapping the host codes* on page B-160.

*Version* The version uniquely identifies a particular set of application definitions.

The version number increases whenever there is a change to the application definitions, for example from version 1.0 to 1.1.

- *General definitions* **3** Expand the **General** entry, select **Short Name**, and press
  - 4 Change the short name (up to 5 alphanumeric characters) if required.

The short name affects the display of the test name in the user interface (for example test buttons, or test names in the orders and results lists, and in all printouts).

5 Press *≫* and change the long name (up to 30 alphanumeric characters) of the test, if required.

The long name is a telling description of the test. It is particularly useful if the short name is not generally familiar in the laboratory environment.

**6** Press  $\gg$  and then **On** to activate the application.

If an application is not active, its associated tests are not available for use. You cannot select them on the test selection screen. The application definitions remain on the system.

7 Press  $\gg$  and select the buttons of the tabs the test should be available from.

• For details, see *Assigning tests to test tabs* on page B-119.

**8** Press 🖉 to confirm the definitions.

Calibration definitions

For Roche calibrators, recommended definitions are automatically defined for each application. You can change the sequence, interval, and the replicate value, if required.



Roche recommend not to change the calibration definitions of Roche reagents.

9 Expand the Calibration entry, select Sequence, and press 🖉.

**10** Change the sequence, if required. (The sequence values define when calibration becomes due, and the system automatically informs you about due calibrations.)

*No Interval:* The system does not inform you about the due status. Use this value if the reagent is stable enough over the whole period until it is empty and you replace it. Calibration is due whenever a new reagent set is loaded on the instrument.

Interval only: Performed when the reagent interval has expired.

*Each Lot and Interval:* Performed whenever the first reagent of a new lot is loaded and then each time the interval has expired.

*Each set and Interval:* Performed whenever a reagent is loaded or when the reagent interval has expired.

- 11 Change the interval (number of days, hours for ISE applications), if required.
- **12** Change the number of replicates, if required.

Possible values: 1, 2, 3. (The default value is 2.)

**13** Press 🔘 to confirm the definitions.

Calibration definition changes require subsequent calibration of the tests.

Control definitions

-Ò́-

- *ns* **14** Expand the **Control** entry, select **Sequence**, and press **2**.
  - **15** Define the sequence.

Choose **No Interval** for cases when, for example, you intend to perform QC as part of standard sample testing and not as a separate task.

Choose **Interval Only** if QC should be performed whenever the reagent interval has expired. The system automatically informs you about due QCs.

**16** Define the interval in hours.

**17** For each control, perform the following steps:

• Define the material code.

Unique identifier for the controls. Check the package insert of the control.

- Define whether QC should be performed after calibration of the test (**QC After Cal**).
- Define whether the test should have QC performed as part of the **Default QC** function.

**Default QC** is an automated process for performing all QC measurements that are currently due. This is the ideal method if you want to perform QC periodically during routine operation

On: QC will be performed collectively for all tests that use a certain control.

**Off:** QC has to be ordered manually for each test.

**18** Press 🞯 to confirm the definitions.

Ratio Calculation definitions	<b>19</b> For applications that use a ratio, perform the following steps:
	• Expand the <b>Ratio Calculation</b> entry, select <b>Coefficient w</b> , and press 2.
	• Define one or two coefficients.
	• Define the formula.
	Incorrect results due to inappropriate formula
	The formula defines how the values of the applications and coefficients are mathemat-
WARNING	ically combined to generate a result.
	It is the responsibility of the user to ensure that the formula is appropriate for the appli- cation that is being defined.
	<b>20</b> Press 🖉 to confirm the definitions.
Calculation definitions	<b>21</b> Expand the <b>Calculation</b> entry, select <b>Factor</b> , and press
	<b>22</b> Change the factor and offset, if correlating two different methods.
	<b>23</b> Choose whether valuation (a reference range) should be employed or not.
	If you choose <b>Reference Range</b> , you need to define an upper and lower limit and decide which of them should be taken into account or whether both of them.
	<b>24</b> Define whether the lower limit is used, if required.
	<b>25</b> Define whether the higher limit is used, if required.
	<b>26</b> Change the lower limit value, if required.
	<b>27</b> Change the higher limit value, if required.
	<b>28</b> Define, for each control, whether sample results should be flagged that were generated using a test whose QC results were marked with a flag.
	<b>29</b> If <b>TRL Check</b> is on ( <b>Configuration</b> > <b>System</b> ), enter the TRL value (lower limit of measuring range) as defined in the relevant application documentation.
	See also System on page B-166.
	<b>30</b> Press 🖉 to confirm the definitions.
-ݣ-	If one of the initial tests for the ratio calculation did not produce a valid result (the result would be marked with a flag), the displayed ratio result would be <b>N/A</b> .
Result Conversion definitions	<b>31</b> Expand the <b>Result Conversion</b> entry, select <b>Laboratory Unit</b> , and press 🥒.
	<b>32</b> Change the laboratory unit if you intend to work with units that are different from the currently specified units.
	<b>33</b> Type the conversion factor, if required.
	This factor is required if lab units were defined.
	<b>34</b> Choose whether values should be displayed in standard or laboratory units.
	This definition affects displays on screens as well as printouts.
	<b>35</b> Change the decimal position, if required.
	This value corresponds to the number of digits after the decimal point that are displayed on the screens.
	<b>36</b> Press 💿 to confirm the definitions.

## **Uninstalling applications**

Uninstalling an application means making the application unavailable for use in testing.

You can reinstall an uninstalled application.

• See Installing applications on page B-137.

Before you can uninstall an application, you need to perform some preparation tasks:

#### ► To prepare uninstalling an application

- 1 Delete all associated sample results.
  - See *Deleting sample results* on page B-94.
- 2 Delete all associated QC results, both from the QC Status list and the QC History.
  - See *Deleting QC results* on page B-96.

#### To uninstall an application

- 1 Choose Utilities > Applications > Laboratory Parameters.
- **2** Select the application.

<u>`</u>Q́

Select an application that is not in square brackets. (Brackets indicate that the application is not currently installed.)

**3** Press  $\Theta$ .

A confirmation screen is displayed.

**4** Press 🕖 to confirm the action.

All associated calibration results and lot data are deleted.

In the applications list, the application name is enclosed in square brackets, for example [GLU2], and in the profile details, instead of the short name its application code is displayed.

The profiles that use this application are removed from the test selection screen. If you still want to use these profiles, you need to remove the application from them.

• See *Removing tests from a profile* on page B-149.

# **Deleting applications**

Deleting an application means removing the data from the system.

The process of removing an application from the system consists of the following steps:

- 1. Delete all associated sample results.
  - See Deleting sample results on page B-94.
- 2. Delete all associated QC results, both from the **QC Status** list and the **QC History**.

• See *Deleting QC results* on page B-96.

3. Uninstall the application. This process makes the application unavailable for use in testing.

• See Uninstalling applications on page B-143.

- 4. Remove all associated reagent sets from any of the reagent disks. (The disk may be off board, and you might need to load it in order to remove the reagents.)
  - See To load the reagent disk on page B-21 See Removing a reagent set on page B-25.
- 5. Delete the application. This process removes the application data from the system.

#### ► To delete an application

- 1 Choose Utilities > Applications > Laboratory Parameters.
- **2** Select the application.
- **3** Press  $\Theta$ .

The following table lists possible situations and how they affect the process of deleting.

Situation	Proceed as follows		
There is one entry for the application in the list, and the application is currently installed (not in brackets).	<ol> <li>Confirm to proceed with uninstalling.</li> <li>In the applications list, select the uninstalled application (it is now contained in brackets).</li> <li>Press .</li> <li>Confirm to proceed with deleting the application.</li> </ol>		
There is one entry for the application in the list, and the application is currently not installed (in brackets).	Confirm to proceed with deleting the appli- cation.		
There is a second entry for the application in the list. The selected application is installed (not in brackets)	Confirm to proceed with deleting the appli- cation. Note: You can only delete the installed application.		
The selected item is a profile.	Confirm to proceed with deleting the pro- file.		
There are still associated reagents on one of the disks.	Remove the associated reagent set from the disk. (The disk may be off board, and you might need to load it in order to remove the reagents.)		

A confirmation screen is displayed.

B-144

4 Press 🔘.


The application data are deleted.

## Defining the processing order

For a given sample, the tests are processed in the order defined by the time required to perform the tests (number of cycles), starting with the one that takes the longest. This order can be altered manually by defining a specific process sequence list.

#### ► To define the processing sequence

1 Choose Utilities > Applications > Process Sequence.

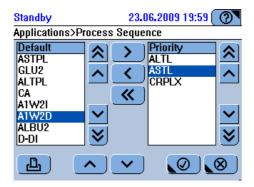

- 2 In the **Default** list, select a test you want to be processed with high priority.
- **3** Press > to move it to the **Priority** list.

The tests are processed in the order displayed in this list. When the tests contained in this list are processed, the rest of the tests are processed in the normal order.

- **4** Perform steps 2 and 3 to move the other tests you want to process with high priority.
- **5** To move a test from the **Priority** to the **Default** list, select the test in the **Priority** list and press ←. To move all tests to the **Default** lists, press ←.
- 6 In the **Priority** list, select a test and use and to move it up and down in the list.
- **7** Press 🔘 to save the processing sequence.

The processing sequence defined in the **Priority** list applies to all tests performed on the current **Cedex Bio** instrument.

ò

# **Defining ratio applications**

The user can manually define ratio applications. Such applications yield results that are mathematically arrived at, and are based on test results that were generated using up to four different applications.

In the results list, both the results of the test or tests as well as the calculated ratio result will be displayed.

#### ► To define a ratio application

- 1 Choose Utilities > Applications > Laboratory Parameters.
- 2 Press 🕀.
- 3 Press Add Ratio.
- **4** Type the code and press  $\gg$ .

The code is a number between 910 and 930 and identifies the application.

**5** Type the short name and press **>>**.

The short name must be unique on the **Cedex Bio** system and consists of up to 5 alphanumeric characters. It is used on the screens, for example on the test selection screen or the results list.

**6** Type the long name and press  $\gg$ .

The long name must be unique on the **Cedex Bio** system and consists of up to 30 alphanumeric characters. Use a name that is commonly known in your laboratory environment.

**7** Select the first application and press  $\gg$ .

You can define between one and four applications.

- **8** Define the other applications, if required.
- **9** Type the first coefficient and press **>>**.

You can define between one and four coefficients.

- **10** Define the other coefficients, if required.
- **11** Type the formula and press **>>**.

The formula relates the applications with the coefficients. You can use the following operands: +, -, \*, /, (, ), space, and application subindex 1 and 2.

Stand	iby			23.06.2009 20:02 🕐
Add Ratio>Calculation>Formula				
w*A-	ŀΖ			
A	В	C	D	1
w	x	У	z	2
I	*	-	+	
1	]		Clr	<b>—</b>
	<u> « » (0 (8</u>			



#### Incorrect results due to inappropriate formula

The formula defines how the values of the applications and coefficients are mathematically combined to generate a result.

It is the responsibility of the user to ensure that the formula is appropriate for the application that is being defined.

- 12 Type the factor to compensate for differences in measuring methods of different instruments, then press ≫.
- **13** Type the offset and press **>>**.
- 14 Select whether the offset/correlation factor is going to be used or not, then press ≫.
- **15** Select the valuation type, then press **>**.

No valuation means that no valuation will be performed.

**Reference range** means that the result will be compared with the defined reference ranges.

- **16** Select whether the low range will be used or not, then press  $\gg$ .
- 17 Select whether the high range will be used or not, then press  $\gg$ .
- **18** Type the low limit, then press  $\gg$ .
- **19** Type the high limit, then press **>>**.
- **20** Type the standard unit, then press  $\gg$ .
- **21** Type the laboratory unit, then press  $\gg$ .
- **22** Type the conversion factor, then press  $\gg$ .
- **23** Select whether the standard or the laboratory unit will be displayed for the final ratio result, then press **≫**.

(For the calculation of the individual results, the system uses standard units.)

- 24 Type how many decimal positions will be used in the displayed results.
- **25** Press 🔘.

#### 7 Configuration

Applications

# **Profiles**

#### **Defining profiles**

A profile is a set of tests that are frequently ordered in combination. When ordering a profile all tests of the profile are performed.

#### ► To define a profile

- 1 Choose Utilities > Applications > Laboratory Parameters.
- **2** Press 🕀.
- **3** Press Add Profile.
- **4** Type the profile code. (For customer defined profiles, use the numbers 910 through 930.)
- **5** Type the short name (up to 5 characters).

The short name will appear on the test selection screens.

- **6** Type the long name (up to 30 characters).
- 7 Select the tests that should be included in the profile.
- **8** Press 🕖 to confirm the selection.

#### Adding tests to a profile

#### ► To add tests to a profile

- 1 Choose Utilities > Applications > Laboratory Parameters.
- **2** Select the profile.
- **3** Expand the entry.
- 4 Expand the General entry.
- **5** Select Applications.
- 6 Press 🖉.

A test selection screen is displayed.

- **7** Select the tests.
- **8** Press 🕖 to confirm the selection.
- For information on assigning a profile to tabs on the test selection screen, see *Assigning tests to test tabs* on page B-119.

#### Removing tests from a profile

- To remove tests from a profile
  - 1 Choose Utilities > Applications > Laboratory Parameters.
  - **2** Select the profile.
  - **3** Expand the entry.
  - 4 Expand the General entry.
  - 5 Select Applications.
  - 6 Press 🖉.

A test selection screen is displayed.

- 7 Clear the selection of the tests you want to remove.
- **8** Press 🖉 to confirm the selection.

#### **Deleting profiles**

• See *Deleting applications* on page B-144.

## **Defining control definitions and lots**

→ Workplace > Lot Data

#### ► To define control definitions and lots

Perform the following steps for all controls.

- 1 Choose Workplace > Lot Data.
- 2 Press Controls.

A screen is displayed that lists all installed control lots.

3 Press ⊕.

A screen is displayed, asking you to scan the barcode transfer sheet or to type the data manually.

**4** Define the lot values.

Do one of the following:

lf	Do this
The lot data are	1. Scan the barcode.
available from a	The screen is displayed again that lists all installed control lots.
barcode	The new lot is selected.
You want to type	1. Press 📖.
the data manually	2. Type the name and press 🔊.
	3. Type the material code and press $\triangleright$ .
	You find this code in <b>Utilities</b> > <b>Applications</b> > select a test >
	4. Type the Lot ID and press $\gg$ .
	5. Type the expiration date. Use the date format as indicated on
	the screen.
	6. Press 🔘 to confirm the definitions.
	If a lot with the same ID already exists on the system, a screen i
	displayed, asking you whether you want to replace the existing lot.
	Press 🔘 to confirm the definitions.
	7. Press *Add new test*.
	8. Press 🕀.
	A screen is displayed that contains all tests that use this contro
	and that have not yet been assigned to this lot.
	9. Select a test and press 🔊.
	10. Type the value for the mean concentration and press $\triangleright$ .
	11. Type the value for the standard deviation.
	12. Press $\textcircled{O}$ to confirm the definitions.
	The screen for handling control lot data is displayed again.
	13. Perform steps 7 through 12 for all tests you want to use.

## Defining calibrator definitions and lots

You need to perform the definitions for all calibrators.

#### ► To define a calibrator lot

- 1 Choose Workplace > Lot Data.
- 2 Press Calibrators.

A screen is displayed that lists all installed calibrator lots.

**3** Press ⊕.

A screen is displayed, asking you to scan the barcode or to type the data manually.

If Do this		this
The lot data are	1.	Scan the barcode.
available from a bar- code		The screen is displayed again that lists all installed calibrator lot The new lot is selected.
You want to type the	1.	Press .
data manually	2. 3.	Type the name (up to 10 alphanumeric characters) and press $\searrow$ Type the material code and press $\bigcirc$ .
		Refer to the package insert.
	4. 5.	Type the lot ID (up to nine alphanumeric characters) and press $\bigcirc$ Type the expiration date and press $\bigcirc$ .
		Use the date format as indicated on the screen.
	6.	Type the number of calibrators (cups) you need to place on the instrument. Use the information given on the package insert. (Exclude calibrators that use system water from this number. So step 14.)
	7.	Press 🔘 to confirm the definitions.
		Press *Add new test*.
	9.	Press 🕒.
		A screen is displayed that shows all tests that have this calibrate defined for them and that are not yet assigned to this lot.
		Press one of the test buttons. Press 🔊.
		Define the first calibration value (target value). Define the next calibration value.
		You can define up to six calibration values, and you must defin them in descending order.
	14.	Select the value for Last is Water.
		Press <b>On</b> if you want to calibrate with system water as zero calibrator. (In this case, no cup needs to be placed on the system. (S step 7).
		Press <b>Off</b> if you want to use a special zero calibrator for the calibration. (The special zero calibrator needs to be placed on the sample area.)
	15.	Press 🔘 to confirm the definitions.

**4** Define the lot values. Do one of the following:



Calibration values (target values) in method sheets may be defined in increasing order of concentration. Always define the values in *decreasing* order on the **Cedex Bio** instrument.

# Short guide to application definitions

-`\$ <u>'</u> -		The following tables provide reference information on the meaning and use of application definitions the user might want to change.		
		The version uniquely identifies a particular set of application definitions.		
Code				
		The Roche application code identifies the application. If you work with a LIS you may need to map the Roche application code with your own code. All applications, profiles, and ratios must be identified with an application code. For profiles and ratios, use a number between 910 and 930 for these codes.		
		See <i>Mapping the host codes</i> on page B-160.		

#### General application definitions

ltem	Values	Comment
Short Name	Up to 5 alphanumeric characters	The short name is used on the screens, for exam- ple on the test selection screen or the results list.
Long Name	Up to 30 alphanu- meric characters	The long name is a telling description of the test. It is particularly useful if the short name is not generally familiar in the laboratory environ- ment.
Active	On, Off	Tests that not active are not available for use. You cannot see them on the screens. Their application definitions remain on the system.
Test Tabs	See Workflow definitions on	Choose one or several of the buttons. The test will be included in the corresponding test tabs.
	page B-161.	Choose Easy if you work with order mode Easy.
		Clear all selections to prevent the test from being selected in the test selection screens.
Table B-12	General application definitions	

#### **Calibration definitions**

-`Q́`-

ISE applications do not have calibration definitions.

ltem	Values	Comment		
Sequence	The sequence values define when calibration becomes due, and the system automatically informs you about due calibrations.			
	No Interval	The system does not inform you about the due status. Use this value if the reagent is stable enough over the whole period until it is empty and you replace it.		
		Calibration is due whenever a new reagent set is loaded on the instrument.		
	Interval Only	Due whenever the reagent interval has expired.		
	Each Set and Interval	Due whenever a new reagent is loaded and when the interval has expired.		
		The interval starts again whenever you place a new reagent set.		
		You can turn off the interval check by defining its duration as 0 (zero).		
	Each Lot and Interval	Due whenever the fist reagent of a new lot is loaded and then each time the interval has expired.		
		The interval (re)starts whenever you calibrate a reagent set.		
		You can turn off the interval check by defining its duration as 0 (zero).		
		If you install a reagent set of the same lot, no cal- ibration is required. (The system uses the Lot calibration result as the Set calibration result of the new reagent set.)		
Interval	Number of days.	Use 0 (zero) if the Sequence is set to No Inter-		
	(For ISE, 1 day fixed.)	val.		
Replicates	1, 2	Number of times each measurement should be repeated.		
Table B-13	Calibration definitions			

## **Control definitions**

statu	system does not inform you about the due
Use	
part arate	this value if you intend to perform QC as of standard sample testing and not as a sep- e task, and if QC should be handled by a host puter.
	e that an accepted QC result is required for a before it can be used.
The	whenever the reagent interval has expired. system automatically informs you about QCs.
Interval Number of days Use (hours for ISE applications)	0 (zero) if the Sequence is set to <b>No Interval</b>
For each of the controls:	
Mat. Code 0 999 Unio	que identification of the control material.
	nbers 801 through 999 are reserved for non ne controls.
You catio	need to define this initially for every appli- on.
	needs to be performed after each calibration e test.
	k the test to have QC performed as part of <b>Default QC</b> function.
	<b>nult QC</b> is an automated process for per- ning all QC measurements that are currently
due.	This is the ideal method if you want to per- a QC periodically during routine operation.

**Roche Diagnostics** 

#### **7** Configuration

Applications

#### **Calculation definitions**

ltem	Values	Comment
Factor		Used for sample and control results.
		Final result = (result * <b>Factor</b> ) + <b>Offset</b>
		The correlation factor is used to compensate for differences in measuring methods of different instruments. By applying a factor, the final results of the different instruments can be directly compared.
Offset		The offset is used to compensate for differences in measuring methods of different instruments. By adding the offset, the final results of the dif- ferent instruments can be directly compared.
Valuation	No Valuation	
	Reference Range	Use the limits defined as <b>Low Limit</b> and <b>High Limit</b> .
Low Used	On, Off	Refers to Low Limit.
High Used	On, Off	Refers to High Limit.
Low Limit		Results that are outside the limit are flagged.
High Limit		Results that are outside the limit are flagged.
13:Flag QC Fail	On, Off	Choose <b>On</b> to ensure that all sample results are flagged that were generated using a test whose QC results were marked with a flag.
		The initial number refers to the control that was used for performing the QC. (See also <i>Control definitions</i> .)
TRL		Threshold value below which the results are no longer displayed, instead it is indicated that the value is smaller than the TRL value and the <b><test b="" rng<=""> flag is added.</test></b>
		Note that you can only change this value if <b>TRL Check</b> is activated.

Table B-15

Calculation definitions

#### **Calculation definitions (ratio applications)**

ltem	Values	Comment
Application A through D	Up to 5 alphanumeric characters	Short names of the installed applications.
<b>Coefficient w</b> through z	000 999	Numeric variable for use in the formula.
Formula	Defined application and coefficient letters.	The formula relates the defined applications with the coefficients.
	Operands: +, -, *, /, (, ), and space	For example: $(A * w + B) * A / B$
Table B-16	Calculation definitions for ratio applications	

### Laboratory correlations (ratio applications)

ltem	Values	Comment
Factor		Used for sample and control results.
		Final result = (result * <b>Factor</b> ) + <b>Offset</b>
		The correlation factor is used to compensate for differences in measuring methods of different instruments. By applying a factor, the final results of the different instruments can be directly compared.
Offset		The offset is used to compensate for differences in measuring methods of different instruments. By adding the offset, the final results of the dif- ferent instruments can be directly compared.
Used	On, Off	Choose <b>On</b> to ensure that the defined factor and offset are applied the final ratio result.
Table B-17	Laboratory correlation definitions for ratio applications	

#### **Reference range definitions (ratio applications)**

ltem	Values	Comment
Valuation	No Valuation	No valuation is performed by the system.
	Reference Range	Use the limits defined as <b>Low Limit</b> and <b>High Limit</b> .
Low Used	On, Off	Refers to Low Limit.
High Used	On, Off	Refers to <b>High Limit</b> .
Low Limit	Values in standard units.	Results that are outside the limit are flagged.
High Limit	Values in standard units.	Results that are outside the limit are flagged.
Table B-18	Reference range definitions for ratio applications	

#### **Result conversion definitions**

ltem	Values	Comment
Standard Unit		
Laboratory Unit	Alphanumeric characters	Type the unit.
Factor		Conversion factor for converting results from one unit to another, for example from mg/dL to mmol/L.
Displayed Unit	Standard Unit	
	Laboratory Unit	
Dec. Position	0 4	Number of digits after decimal point that are displayed on the screens.
Table B-19	Result conversion definitions	

# **Changing your password**

NOTICE

# Damage to data and programs due to unauthorized access to the Cedex Bio instrument

User access to the **Cedex Bio** instrument is controlled by a login name and password. Users are administered by a system administrator.

Roche recommend to periodically change the password of the system administrator and to store the current password in a safe place with adequate access control.

<u>ک</u>

User name and password are case sensitive. This means that for example Admin and admin are two different names.

# To change you password

**1** Do one of the following:

lf	Do this
You are logged off:	1. Press Log On.
You are logged on:	1. Press the button showing your user
	name.
	2. Press Log On.

**2** Type your user name.

As soon as you type the first characters of your user name, the system looks for a name that starts with these letters and, if it finds one, displays the complete name.

- **3** Press 🔘.
- 4 Press Change Password.
- **5** Type your old password.
- 6 Press 🔘.
- 7 Type the new password
- 8 Press 🔘.
- **9** Type the new password a second time.
- **10** Press 🔘.

# Scheduling maintenance actions

The **Cedex Bio** instrument facilitates performing the maintenance actions in bundles at times that suit your laboratory work processes. To that purpose, you can define one day of the week as your maintenance day.

*Concept of the main* The main maintenance day is the day in the week when you want to perform the maintenance day majority of the maintenance actions.

Let us suppose that you chose Monday as your main maintenance day. On the Maintenance screen, actions that are due on Monday or an earlier day are marked in red, actions that will fall due between Tuesday and Sunday will be marked in yellow.

The idea is to perform on the main maintenance day all red and yellow maintenance actions. This way, between Tuesday and Sunday, you only need to perform maintenance that need to be performed daily.

### To define the main maintenance day

- 1 Choose Utilities > Configuration > Maintenance.
- 2 Expand the main entry and select Major Maint. Day.
- **3** Press 🥢.

A screen is displayed for selecting a day of the week.

- 4 Press one of the weekday buttons.
- **5** Press 🔘.



The definition of the minor maintenance day has currently no practical influence on the maintenance scheduling, because there are no maintenance actions with intervals of between two and six days.

# **User management**

You need Administrator user rights to define or delete a user.

### **Defining a user**

# To define a user

1 Choose Utilities > Users.

A screen is displayed that lists all currently defined users.

- **2** Press ⊕.
- **3** Type the user name.

Make sure it is unique within your laboratory.

- 4 Press ≫.
- **5** Type the password.
- 6 Press ≫.
- 7 Press the appropriate user level button.
  - Press Operator for users who should perform routine tests.
  - Press **Supervisor** for users who should perform routine tests as well as calibration and QC.
  - Press Lab Manager for users who should perform all test actions and who should be able to update or add applications and perform diagnostics tasks.
  - Press **Administrator** to define the administrator of the **Cedex Bio** instrument. The administrator can, in addition to all **Lab Manager** actions, perform user administration.
- **8** Press 🔘.

### **Deleting a user**

### ► To delete a user

1 Choose Utilities > Users.

A screen is displayed that lists all currently defined users.

- **2** Select the user you want to delete.
- **3** Press  $\Theta$ .

A confirmation screen is displayed.

- **4** Press  $\swarrow$  to confirm the deletion.

# Editing the acceptable flags list

When working with automatic accepting of results, results that are not flagged are automatically accepted. Results with flags that are marked in an acceptable flags list are accepted as well. There is such a list for sample, QC, and calibration results.

# ► To add or remove a flag from an acceptable flags list

- 1 Choose Utilities > Configuration > Result Handling.
- 2 Expand the list.
- **3** Select one of the following entries:
  - Sample Accept. Flags
  - QC Acceptable Flags
  - Cal Acceptable Flags
- 4 Press 🖉.

A list of relevant flags is displayed.

**5** Press the square to the left of a flag name to select it or to cancel its selection.

A selected flag shows a check mark ( $\checkmark$ ).

- 6 Press 🔘.

# Mapping the host codes

Applications are delivered with a unique ID called code. Laboratories often use their own codes for applications. By mapping the codes used by the laboratory (host codes) to those defined by the application manufacturer, the laboratory can work with their own accustomed codes. The **Cedex Bio** instrument provides a mapping table that can easily be edited.

Editing a mapping table involves the following steps:

- 1. Exporting the mapping table to an USB stick. (The table is written as a text file.)
- 2. Editing the file on a computer and save it as a text file to the stick.
- 3. Importing the file to the **Cedex Bio** instrument.

# To export the mapping table

- 1 Choose Utilities > Applications > Host Codes.
- **2** Press **♦**.

A screen is displayed, asking you to insert the USB stick.

- **3** Insert the USB stick.
- 4 Press 🖉.
- **5** Select the directory where you want to save the mapping file.
- **6** Press 🖉 to confirm the selection.

The file is exported as a text file of the name format hct\_yyyymmddhhmmss.txt.

### ► To re-import the mapping table

- 1 Choose Utilities > Host Codes.
- **2** Press **♦**.

A screen is displayed, asking you to insert the USB stick.

- 3 Insert the USB stick.
- 4 Press 🔘.
- **5** Select the mapping file.

Typically, mapping files have the name format hct\_yyyymmddhhmmss.txt.

Press \*.txt to display text files only.

Press \*.\* to display all files and directories.

**6** Press 🕖 to confirm the selection.

The file is installed on the system. The screen with the mapping table is displayed again.

### To use the manufacturer's application codes

- 1 Choose Utilities > Host Codes.
- 2 Press 11.

The host codes are set to values that are identical to the application codes.

By pressing 11 again, you can revert to the original values.

# Short guide to configuration definitions

# Workflow definitions

ltem	Values	Comment
Working Mode		Defines the way in which you select tests during order definition.
	Manual	Use this definition if the <b>Cedex Bio</b> instrument is used as a stand-alone system.
		During order definition, the test selection screen is displayed, allowing you to select the tests or to make changes to the selection.
	Host	Use this definition if the <b>Cedex Bio</b> instrument is connected to a host computer.
		The orders are automatically defined, the test selec- tion screen is not displayed. After identifying the sample, you are asked to place the sample.
Table B-20	Workflow defin	itions

ltem	Values	Comment
Sample Barcode	On, Off	You can work with or without barcodes on sample tubes. Use <b>Off</b> only if you exclusively work without barcodes.
		If you use <b>On</b> , during order definition the option to type the sample ID manually is skipped
Order Mode	Easy	You work with one test panel on the test selection screen, containing up to 25 tests.
		Typical situation for this mode: You work with one or two reagent disks, manual order identification, identical sample and order IDs, possibly but not nec- essarily with host connection.
	Full	You can work with up to six test tabs on the test selec- tion screen, each containing up to 20 tests.
		Typical situation for this mode: You distribute the reagents across several (up to eight) reagent disks and you predominantly work with specific groups of tests. You also work with a host connection with automatic order identification and result upload.
Sample ID Handling	Order ID = Sample ID	The order ID is automatically defined when you iden- tify the sample.
	Independent IDs	You define the sample ID and order ID indepen- dently of each other.
	Grouped Sample	Order and sample ID are identical.
	ID	For a given sample, the same sample ID is used all the time. A three digit number is appended and automatically increased with each new order.
		Examples: 789-001 for the first order, 789-002 for the second order, and so on.
Sample ID Separator	One alphanumeric character	This definition is used in combination with <b>Grouped</b> <b>Sample ID</b> . It defines the character that separates the sample ID from the automatically appended number.
		Examples: 789-001, 789.001, 789_001.
Auto Order-ID	On, Off	If you use <b>On</b> , the system automatically increments the order ID number by one whenever you define a new order.
		If you use <b>On</b> , choose <b>Off</b> for <b>Order ID</b> = <b>Sample ID</b> .
Auto Print Results	On, Off	If you use <b>On</b> , sample results are printed as soon as all results of the order are available and accepted.
Test Tab 1 6	Up to four	Display name for the test tab 1 6
Name	alphanumeric characters	Test tabs are used to group tests that are usually used together.
		You can assign a test to more than one tab.
Test Tab 1 6	On, Off	Use <b>On</b> to make available the tab on the test selection
Visible		screen.

# Host definitions

ltem	Values	Comment
Baud Rate		Data transmission rate for host interface communica-
		tion.
		Select one of the predefined values.
Handshake	Off	An exchange of signals between two devices when
	Hardware	communication begins in order to ensure synchroni-
	XON/XOFF	zation.
Line Mode		Number of bits and stop bits used during data trans-
		mission.
		Select one of the predefined values.
Parity Check	On, Off	An error detection technique that tests the integrity
		of digital data in memory or on disk.
Communication	Off	Download to the Cedex Bio instrument is possible,
Mode		but upload to host is not.
	Batch	Only manual upload to the host is possible.
	Realtime	Results and queries are sent during processing.
Checksum	On, Off	A value used to ensure that data are stored or trans-
		mitted without error. It is stored with the data. When
		the data are retrieved from memory or received at the
		other end of a network, a new checksum is computed
		and matched against the existing checksum. A non-
		match indicates an error.
Order Query		Query mode allows real-time communication during
Mode		order definition of regular, rerun, and repeated tests.
	Off	No queries are sent.
		Test orders have to be defined either manually or by
		downloading them from the host.
	Once	A query is sent only when the sample is identified the
		first time.
	Always	A query is sent every time the sample is identified.
Timeout	10 300 s	Period of time the system will wait for a response
		from the host when working in Order Query Mode.
		This definition applies to routine, STAT and rerun
		sample tests.
Send Result	Off	The results are sent to the host manually.
Mode	Complete	The results are automatically sent to the host when all
	_	results of the order are available and accepted.
	Immediate	Each result is automatically sent to the host as soon as
		it has been accepted.
Send Raw Data	On, Off	Raw data include all measurement results made dur-
		ing the measuring cycles, not just the calculated test
		results.
Table B-21	Host definitions	

ltem	Values	Comment
Trace	On, Off	The Trace function records the content of the com- munication with the host in a trace file.
		This is a troubleshooting function. Do not work with it during normal daily operation.
		The trace file can be exported as a log file ( <b>Utilities</b> > <b>Export</b> > <b>Log Files</b> ; file format: htryyyymmd-dhhmmss.log.)
System Device ID	Up to 10 alphanumeric characters.	The ID is used on reports and for communication between the <b>Cedex Bio</b> instrument and the host.
Host Device ID	Up to 10 alphanumeric characters.	The ID is used on reports and for communication between the <b>Cedex Bio</b> instrument and the host.
Table B-21	Host definitions	

Configuration

# **Result handling**

ltem	Values	Comment
Sample Auto Accept	On, Off	Results that do not contain a flag are automatically accepted. Results with flags that are marked in a list of acceptable flags are accepted as well.
Sample Accept. Flags		List of possible flags. Results with marked flags are automatically accepted if <b>Sample Auto Accept</b> is on.
QC Auto Accept	On, Off	Results that do not contain a flag are automatically accepted. Results with flags that are marked in a list of acceptable flags are accepted as well.
QC Acceptable Flags		List of possible flags. Results with marked flags are automatically accepted if <b>QC Auto Accept</b> is on.
Cal Auto Accept	On, Off	Results that do not contain a flag are automatically accepted. Results with flags that are marked in a list of acceptable flags are accepted as well.
Cal Acceptable Flags		List of possible flags. Results with marked flags are automatically accepted if <b>Cal Auto Accept</b> is on.
Table B-22	Result handling	g definitions

# Calibration

ltem	Values	Comment
Forecast Hours	0 999 h	Calibrations that fall due within this period will be performed collectively.
		0 means working without this feature.
		The default value is 8.
Table B-23	Calibration defin	itions

QC

Values	Comment
	Rule: A flag is generated if one result is outside a limit.
	s : Standard deviation. The value for the standard deviation is defined in the QC lot.
Off	
2.5s	
3s	
	Rule: A flag is generated if two consecutive results are outside the same limit.
	s : Standard deviation. The value for the standard deviation is defined in the QC lot.
Off	
2s	
QC definitions	
	2.5s 3s Off 2s

Configuration

# System

ltem	Values	Comment
Language	Language ID according to ISO 639-1.	The language English is always included.
Date Format	dd.mm.yyyy dd-mm-yyyy yyyy.mm.dd yyyy-mm-dd mm/dd/yyyy	Examples for June 14, 2006: • dd.mm.yyyy: 14.06.2006 • mm/dd/yyyy: 06/14/2006
Time Format	24 Hours 12 Hours	
Screen Saver Wait	0 120 min	Time after which the screen saver feature becomes active.
		0 (zero) means that the screen saver function is off. The screen saver time is linked to the <b>Auto Log-off</b> feature. For example, if you define a Screen Saver Wait time of 5 minutes, then after 5 minutes of inac- tivity, the screen saver feature becomes active and the user is automatically logged off, provided <b>Auto Log-</b> <b>off</b> is on.
Auto Log-off	On, Off	If you work with <b>On</b> the user is automatically logged off after a certain number minutes of inactivity on the instrument. This number is defined in the <b>Screen</b> <b>Saver Wait</b> definition.
Sound Effects	On, Off	Acoustic signals when handling instrument items, e.g. clicking sound when typing. (Does not affect the acoustic signals for alarms and warnings.)
Volume	0 100	Volume of <i>any</i> acoustic signal generated by the instrument, including alarm and warning signals. 0 (zero) means silent.
Host Server	On	The host interface of the instrument is enabled, allowing it to communicate with a host.
	Off	The host interface of the instrument is disabled. Communication with a host computer is not possible. Host definitions have no effect.
TRL Check	Off	The actual results for the lower limit of the measuring range are displayed.
		You cannot change the lower limit of the measuring range ( <b>Low Limit</b> ). See <i>Calculation definitions</i> on page B-142.
	On	If you choose <b>On</b> , you must, for each test, check the TRL value (lower limit of the measuring range) in the relevant application documentation. You may have to manually adjust it. See <i>Calculation definitions</i> on page B-142.
		If a measured value is below this TRL threshold, the actual results are not displayed, instead a value such as <0.5 is displayed, where 0.5 is an example of the TRL value defined for this test.
Table B-25	System definitions	

**Roche Diagnostics** 

# Date and time

ltem	Values	Comment
Date	Format: dd.mm.yyyy	Date that is used for time records, for example with results. Example: 17.02.2006 for February 17, 2006. This format is for entry purposes only. It does not influence the format of displayed dates.
Time	Format: hh:mm	Time that is used for time records, for example with results. Example: 08:32. This format is for entry pur- poses only. It does not influence the format of dis- played times.
Table B-26	Date and time defi	nitions

# Maintenance

ltem	Values	Comment
Major Maint. Day		Day in the week when you usually perform mainte- nance actions.
Minor Maint. Day		This definition has currently no effect.
Table B-27	Maintenance of	definitions

# Abs adjustment

The absorbance adjustment values cannot be changed.

ltem	Values	Comment
Adjustment Valid	On, Off	
H2O Signal		For each wave length, the adjustment value for the cuvette filled with water is given.
H2O Cuv Signal		For each wave length, the adjustment value for the empty cuvette is given.

 Table B-28
 Absorbance adjustment definitions

Roche Diagnostics

# Defining extra wash cycles

Extra wash cycles are used to minimize carry-over.

You can define an extra wash cycle either by reading a barcode or by defining the values manually.

# ► To define an extra wash cycle

1 Choose Utilities > Applications > Extra Wash Cycles.

A screen is displayed that lists the currently defined extra wash cycles.

- 2 Select Add new rule.
- **3** Press ⊕.
- **4** Do one of the following.

lf	Do this	
The definitions are available	1.	Read the barcode.
on a barcode:		The definitions will be installed on the system.
You want to define the	1.	Press 📖.
values manually:	2.	Choose when—in relation to the pipetting action from
		the bottle defined in Trigger Bottle Code and Pipet-
		ting Type—the extra wash cycle should be performed
		(action), then press 🔊 .
	3.	Type the trigger-bottle code and press $\triangleright$ .
		The trigger-bottle code is the reagent set code. You
		find it on the barcode sheet.
	4.	Choose the pipetting type. This defines the individual
		bottle (bottle of the set or the sample tube) whose
		pipetting triggers performance of the extra wash cycle.
	5.	Define whether the wash cycle should be active or not.
	6.	Type the cleaner bottle code to define which cleaner
		should be used.
	7.	Type the volume of cleaner in $\mu L$ that will be pipetted.
	8.	Press 🔘 to confirm the definitions.

The new wash cycle is listed in the **Extra Wash Cycles** list. Its name is made up of the action, the short name of the trigger bottle and the pipetting type.

# Activating, deactivating and deleting extra wash cycles

You can deactivate an extra wash cycle if you do not want to use it. (The definitions remain on the system. You can activate the extra wash cycle later again.)

# ► To activate or deactivate an extra wash cycle

1 Choose Utilities > Applications > Extra Wash Cycles.

A screen is displayed that lists the currently defined extra wash cycles.

- **2** Select and expand the extra wash cycle entry.
- 3 Press 🖉.
- 4 Choose **On** to activate the extra wash cycle.

Choose Off to deactivate the extra wash cycle.

5 Press 🔘.

A screen is displayed for entering your password.

- **6** Type your password and confirm by pressing **O**.

### To delete an extra wash cycle

1 Choose Utilities > Applications > Extra Wash Cycles.

A screen is displayed that lists the currently defined extra wash cycles.

- **2** Select the extra wash cycle entry.
- **3** Press  $\Theta$ .

A screen is displayed for entering your password.

- 4 Type your password and confirm by pressing 
   O

   The extra wash cycle definitions are deleted.

# Short guide to extra wash cycle definitions

→ Utilities > Applications > Extra Wash Cycles > 🖉

# Extra wash cycle definitions

ltem	Values	Comment
Active	On, Off	Defines whether the extra wash cycle can be applied or not.
Action	None Before After	Defines when, in relation to the pipetting action from the bottle defined in <b>Trigger Bottle Code</b> and <b>Pipet-</b> <b>ting Type</b> , the extra wash cycle should be performed.
Trigger-Bottle Code	0 999	Bottle set code of the fluid that triggers performance of the extra wash cycle.
		If you use the pipetting type <b>Sample</b> , it defines the bottle whose fluid will be prevented from being carried over to the sample by using the extra wash cycle.
Pipetting Type	Reagent 1 Reagent 2 Start Reagent Sample	Defines the individual bottle (bottle of the set or a sample tube) whose pipetting triggers performance of the extra wash cycle.
Cleaner-Bottle Code	0 999	Material code of the cleaner.
Volume	μL	Amount of cleaner that will be pipetted during the extra wash cycle.
Table B-29	Extra wash cycle c	definitions

# Short guide to reagent mixing rules

→ Utilities > Applications > Reagent Mixing > 🖉

You can import and delete mixing rules, but cannot change them.

◆ For importing reagent mixing rules, see *Importing automatic mixing and extra wash cycle definitions* on page B-115.

### **Reagent mixing rule definitions**

You cannot change these definitions.

ltem	Values	Comment
Active		Defines whether this rule can be applied to an application or not.
Initial Mixing Time		Defines the duration of the mixing for the first mix- ing.
Mixing Interval		Defines the time after which mixing must be per- formed again.
Periodic Mixing Time		Defines the duration of the mixing for the second or subsequent mixing.
Relaxing Time		Defines the time after mixing before the reagent can be used.
Table B-30	Reagent mixing rule	e definitions

# **Printing mixing rules**

You can print mixing rules.

# ► To print mixing rules

- 1 Choose Utilities > Applications > Reagent Mixing.
- **2** Select the mixing rule you want to print.
- **3** Press 🔘.

A printing screen is displayed.

4 Select whether you want to print the selected rule set or all defined rule sets. The rules are printed on the built-in printer.

<sup>`̈́</sup>Q́

# **Deleting mixing rules**

*Prerequisites* You need Lab Manager or Administrator user rights to perform this task.

# ► To delete a mixing rule

1 Choose Utilities > Applications > Reagent Mixing.

A screen is displayed that lists the currently defined mixing rules.

- **2** Select the mixing rule.
- **3** Press  $\Theta$ .

A screen is displayed for entering your password.

- 4 Type your password and confirm by pressing ②.The mixing rule definitions are deleted.

# Maintenance **C**

8 General maintenance C-3	8	General maintenance	C-3
---------------------------	---	---------------------	-----

# **General maintenance**

# Keeping the instrument going

In this chapter, you will find step-by-step instructions of the maintenance actions that you must perform to keep the instrument running smoothly and efficiently.

In this chapter	Chapter	8
Overview		C-5
Performing maintenance actions		C-6
Maintenance actions and their intervals		C-7
Maintenance actions		C-8
Safety information		
Deproteinize the probe		. C-10
Prime the fluid system		
Clean the probe manually		
Clean reagent disk and sample area		
Clean the water and waste containers		. C–16
Replace water inlet filter		. C-19
Clean and disinfect housing		
Clean up the database		
Replace photometer lamp		
Contact service representative		
Perform an air/water calibration		
Initialize degasser fluid sensor		
Clean rotor and heating channel		

### 8 General maintenance

Table of contents

Overview

# **Overview**

	Completing maintenance actions correctly and on time helps to ensure smooth and uninterrupted operation of your instrument.
Maintenance scheduling	The <b>Cedex Bio</b> instrument facilitates performing the maintenance actions in bundles at the times that suit your laboratory work processes. To that purpose, you can define in the configuration settings one day of the week as your maintenance day.
	For information on scheduling maintenance actions, see <i>Scheduling maintenance actions</i> on page B-158.
	All maintenance actions can be performed any time.
Interval	For most maintenance actions a fixed maintenance interval is defined. (You cannot change this interval.) This is the basis upon which the system calculates the date when the actions need to be performed.
	The interval timers and counters are reset whenever you confirm that the maintenance action has been performed.
	Maintenance actions without predefined intervals are performed whenever necessary, or they are triggered by another maintenance action.
Due date	The due date is the last possible maintenance day. This is the date you see when you check the status of maintenance actions.
Ensuring smooth operation	Performing all due maintenance actions during the daily <b>Prepare</b> or <b>End Shift</b> phase ensures that routine operation does not have to be interrupted for performing maintenance actions.

Overview

# Performing maintenance actions

►	То	perform maintenance actions	
Checking when maintenance actions are due	1	Choose Utilities > Maintenance	
		A B Standby 10.12.2007 4: 22 ⑦ Maintenance Clean Water and Waste 10.12  Replace Photometer Lamp 10.12  Clean Probe Manually 10.12  Deproteinize Probe 11.12  Clean Reagent Disk 07.01  Replace Water Inlet Filter 25.02  Clean Housing 02.06  V	— C — D
		A     Maintenance action name.     D       B     Due dates, in ascending order.     D       C     This maintenance action is selected.	Use the scroll bar to display the maintenance actions that are currently hidden.
Interpreting the colors		The defined maintenance interval h action now. This maintenance action should be	as expired. Perform this maintenance
		maintenance day.	
		No action is currently required.	
		This maintenance action is selected	
Preparing the maintenance actions		The maintenance actions are sorted accord performed. Use these dates for planning the ordering the required materials.	• • •
Performing the maintenance	2	Select the maintenance action you want to	perform.
action		The selected line turns blue.	
	3	Press 💽.	
		The first screen of the wizard for performin displayed. Usually, this is the maintenance	•

Press 🝙 if you want to print the instructions before you start.

4 Follow the instructions on the screen.

A message will inform you when the maintenance action is complete. In addition, with maintenance actions that take some time, an acoustic signal is sounded. This allows you to do work away from the instrument while the maintenance action is in progress.

- For information on how to perform individual maintenance actions, see *Maintenance actions* on page C-8.



### Incorrect results or processing stop due to incomplete maintenance actions

You can cancel a maintenance action any time by pressing  $\bigotimes$ .

If you interrupt the performing of a maintenance action that was due, its status will remain due, and you need to fully re-perform the action later.

If at all possible, complete a maintenance action without interrupting it.

# Maintenance actions and their intervals

The following table lists the maintenance actions and shows how frequently they need to be performed.



Every effort has been made to ensure that all the information contained in this table is correct at the time of publication. However, Roche Diagnostics GmbH reserves the right to make any changes necessary without notice as part of ongoing product development.

Maintenance actions with no defined interval need to be performed in particular situations, for example during troubleshooting.

Interval	Maintenance action
None	Prime Fluid System
None	Abs. Air/Water Calibration
None	Clean Rotor And Heating Channel
None	Cleanup Database
Daily	Deproteinize Probe
Weekly	Clean Probe Manually
Monthly	Clean Reagent Disk
Three-monthly	Clean Water and Waste
Three-monthly	Replace Water Inlet Filter
Six-monthly	Clean Housing
Six-monthly	Replace Photometer Lamp
Yearly	Call Service Rep.
Yearly	Initialize Degasser Fluid Sensor
Table C-1	Periodicity of maintenance actions

Table C-1

Periodicity of maintenance actions

# **Maintenance actions**

The following sections describe the maintenance actions of the main instrument.

● For specific ISE maintenance actions, see ISE maintenance actions on page E-42.

# **Safety information**



### Injury through working solutions

Direct contact with cleaning solutions or other working solutions may cause personal injury. When handling such solutions, exercise the precautions required for handling them, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics cleaning solutions.

# Personal injury or damage to the analyzer due to contact with instrument mechanism

Do not touch moving parts during instrument operation.

### Infection by biohazardous materials

Contact with samples containing material of human origin may result in infection. All materials and mechanical components associated with samples of human origin are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

### Incorrect results due to build-up of contaminants

During use, contaminants may adhere to the probe. As a result, traces of analytes or reagents may be carried over to the next. Make sure to perform the probe maintenance actions as soon as they are due in order to prevent potentially false results.

### Incorrect results or damage to the analyzer due to dust and soiling

The user can leave the main cover open while the system is in **Standby** status or while the instrument is shut down. This can cause dust and dirt being collected in the heating channel, which in turn might decrease the quality of the cuvettes.

Keep all covers closed. Open them only to perform operation actions.

### Incorrect results or processing stop due to skipping maintenance actions

Not performing maintenance actions that are due may lead to situations where the system cannot continue processing orders, or it may lead to incorrect results. If at all possible, perform the maintenance actions when they are due.



#### Incorrect results or damage to the analyzer due to wrong operation

Operators are required to have a profound knowledge of relevant guidelines and norms as well as the information and procedures contained in the Operator's Manual.

- Do not carry out operation and maintenance unless you have been trained by Roche Diagnostics.
- Start all maintenance actions on the screen. Do not perform maintenance actions without the assistance of the user interface.
- Carefully follow the procedures specified in the Operator's Manual for the operation and maintenance of the system.
- Leave maintenance that is not described in the Operator's Manual to a trained service representative.
- · Follow standard laboratory practices, especially when working with biohazard material.

#### Incorrect results or processing stop due to incomplete maintenance actions

You can cancel a maintenance action any time by pressing 🛞 at the end of a step. Using 🛞 while the system is performing a maintenance action may lead to unexpected errors. If you interrupt the performing of a maintenance action that was due, its status will remain due, and you need to fully re-perform the action later.

- If at all possible, complete a maintenance action without interrupting it.
- If you need to interrupt the maintenance action, wait until you see confirmation on screen that the current step has been completed.

#### Danger of poor measurement quality due to inadequate water quality

Inadequate water quality may lead to incorrect results. Always use purified water of the quality specified in section *Technical specifications*.

#### Infection by waste solution

Contact with waste solution may result in infection. All materials and mechanical components associated with the waste systems are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.



Whenever the term "purified water" is used in this document, water of at least the quality specified in chapter *Technical specifications* must be used.



### Malfunction due to inappropriate interrupt

Using (③) while the system is performing a maintenance action may lead to unexpected errors.

• Do not use the 🛞 button while the system performs maintenance actions. If you need to interrupt the maintenance process, wait until you see confirmation on screen that the current step has been completed.

Tools a

# Deproteinize the probe

Utilities > Maintenance > Deproteinize Probe	
To maintain the efficiency of the instrument, you must clean and deproteinize the probe regularly to prevent the build-up of contaminants.	
Approximately 5 minutes.	
The system must be in <b>Standby</b> status.	
<ul><li>□ ISE Deproteinizer</li><li>□ Activator</li></ul>	
	To maintain the efficiency of the instrument, you must clean and deproteinize the probe regularly to prevent the build-up of contaminants. Approximately 5 minutes. The system must be in <b>Standby</b> status.



Make sure that you have read and understood section *Safety information* on page C-8. The following warning messages in particular are relevant:

- Injury through working solutions on page C-8.
- Incorrect results due to build-up of contaminants on page C-8.

### ► To deproteinize the probe

- 1 Prepare one tube with ISE Deproteinizer and one with Activator.
- 2 Select the maintenance action Deproteinize Probe.
- 3 Press 💽.

The maintenance definition screen is displayed.

4 Press 🔘.

A placement list is displayed.

- **5** Place the ISE Deproteinizer and Activator on the positions indicated on the screen.
- **6** Press 🕖 to confirm the placing.

The system first deproteinizes and then activates the probe.

A message will inform you when the maintenance action is complete.

- **7** Remove the tubes.
- **8** Press 🛞.

# Prime the fluid system

### → Utilities > Maintenance > Prime Fluid System

Priming flushes the fluid system with water to ensure the paths in the fluid system (including pipette, tubing, and probe) are filled with fluid and are free of air bubbles. If air bubbles are present in the fluid system, your test results will be inaccurate.

Operator time Approximately 2 minutes.

*Prerequisites* The system must be in **Standby** status.

*Tools and materials required*  $\Box$  Purified water.

# ► To prime the fluid system

- 1 Check the external water and waste containers. If required, refill the water and empty the waste container.
- 2 Select the maintenance action **Prime Fluid System**.
- **3** Press **()**.

The maintenance definition screen is displayed.

4 Press 🔘.

The system performs the priming actions automatically.

A message will inform you when the maintenance action is complete.

- 5 Press 🛞.

Tools

# Clean the probe manually

$\rightarrow$	<b>Utilities &gt; Maintenance &gt; Clean Probe Manually</b>
	You must clean the probe regularly by hand to prevent any build up of deposits that may affect results.
Operator time	Approximately 1 minute.
Prerequisites	The system must be in <b>Standby</b> status.
s and materials required	□ 70% ethyl alcohol
	□ Tissues
	Make aura that you have read and understand partian Cafety information on page C. 0



Make sure that you have read and understood section *Safety information* on page C-8. The following warning messages in particular are relevant:

- Injury through working solutions on page C-8.
- Incorrect results due to build-up of contaminants on page C-8.
- Infection by biohazardous materials on page C-8.

# To clean the probe

- 1 Select the maintenance action Clean Probe Manually.
- **2** Press **()**.

The maintenance definition screen is displayed.


**3** Press *⊘*.

The system deactivates the transfer motors.

**4** Remove the transfer head cover.

Press the release buttons on both sides and lift.

- **5** Remove the probe with its holder.
  - Remove the tube from at least the first two tubing clips.
  - Press the release button on the side of the probe holder carriage and lift the probe holder with the probe. Be careful not to pull the tubing.

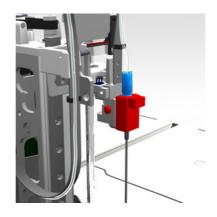

NOTICE

6 Hold the assembly by the probe holder and clean the probe using a tissue moistened with 70% ethyl alcohol.


Wipe from top to bottom several times.

# Probe damage and instrument malfunction due to inappropriate probe handling

Applying unilateral side pressure on the probe can lead to its deformation and consequently to instrument malfunction.

Make sure to apply equal pressure on both sides and to move exactly in the direction of the probe when wiping it.

7 Wipe the side of the probe holder that faces the level detection contact. Use a tissue moistened with 70% ethyl alcohol.

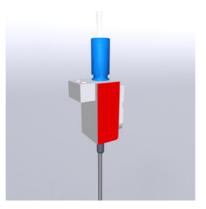

**8** Reinstall the probe holder with the probe.

Press the release button on the side of the probe holder carriage before you insert the probe holder. Release the button when the holder is inserted.

Push down firmly until the release button latches on.

- **9** Fix the tube to the tubing clips.
- **10** Reinstall the transfer head cover.

Press down until the release buttons latch on.

11 Press 🔘.

The system automatically initializes the transfer head and performs the **Prime Fluid System** maintenance action.

A message will inform you when the maintenance action is complete.

**12** Press 🛞.

# Clean reagent disk and sample area

→ Utilities > Maintenance > Clean Reagent Disk

To remove possible spillages you must clean the reagent disk and sample area regularly.

This maintenance action involves the following steps:

- Cleaning the reagent disk
- Cleaning the sample area

Operator time Approximately 10 minutes.

*Prerequisites* The system must be in **Standby** status.

Tools and materials required

□ Alcoholic cleansing tissues

- □ Tissues
- $\hfill\square$  Commercial nonabrasive detergent

Make sure that you have read and understood section *Safety information* on page C-8. The following warning messages in particular are relevant:

- Injury through working solutions on page C-8.
- Infection by biohazardous materials on page C-8.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page C-8.


### To clean the reagent disk

- 1 Select the maintenance action Clean Reagent Disk.
- **2** Press **()**.

The maintenance definition screen is displayed.

- **3** Press  $\oslash$ .
- 4 Open the main cover.
- 5 Remove the reagent disk.
- 6 Remove spillages and stains on the disk with alcoholic cleansing tissues.
- 7 Wipe the inside of the reagent cooler with a cloth dampened with soapy water.
- 8 Reinsert the reagent disk.
  - For information on inserting the reagent disk, see *Preparing the reagent disk* on page B-21.
- **9** Open the right service flap.

**10** Lift the sample tray from its fixture.

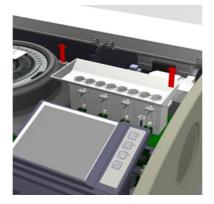


11 Wash the sample tray using detergent. (You can wash the tray in a dish washer.)



- Do not disassemble the sample tray.
- **12** Leave the sample tray to dry or dry it using tissues.
- **13** Reinsert the sample tray.

Make sure to align the guides with the pins.

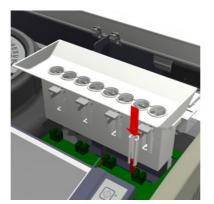

Press down firmly.

- 14 Close the right service flap and the main cover.
- **15** Press to confirm that you have performed the maintenance action.The system initializes automatically.
  - A message will inform you when the maintenance action is complete.
- **16** Press 🛞.

# Clean the water and waste containers

→ Utilities > Maintenance > Clean Water and Waste

You must clean the external water and fluid waste containers to prevent the build-up of contaminants, which can affect the water quality and therefore the quality of the results produced.

This maintenance action consists of the following steps:

- Cleaning the water container
- Cleaning the waste container

The water and waste containers are placed on the rack next to the **Cedex Bio** instrument.

*Operator time* Approximately 10 minutes.

*Prerequisites* The system must be in **Standby** status.

Tools and materials required

□ Spare waste container

- □ Purified water
- □ Hypochlorite solution (0.6%)
- $\hfill\square$  Alcoholic cleansing issues



Make sure that you have read and understood section *Safety information* on page C-8. The following warning messages in particular are relevant:

• Injury through working solutions on page C-8.

- Infection by waste solution on page C-9.
- Infection by biohazardous materials on page C-8.

### To clean the water container

- 1 Select the maintenance action Clean Water and Waste.
- **2** Press **()**.

The maintenance definition screen is displayed.

- **3** Press 🖉.
- **4** Lift the tubing adapter from the white water container.



**5** Place the tubing adapter on a clean surface.


- 6 Dispose of any remaining water.
- 7 Pour 500 mL hypochlorite solution in the water container.
- 8 Clean the water container with a brush.
- **9** Dispose of the solution.
- **10** Clean the interior of the water container manually using tissues. Remove any residue.
- 11 Rinse the water container with tap water at least four times.
- **12** Rinse the water container with purified water.
- **13** Wipe the suction pipe with a tissue.
- 14 Fill the water container with purified water.
- 15 Reinsert the tubing adapter into the water container. Press down firmly.

### To clean the waste container

1 Have the spare waste container ready.

The system periodically performs wash actions. Therefore, an external waste container must be connected at all times.

**2** Remove the tubing adapter from the yellow waste container and insert it in the spare container.

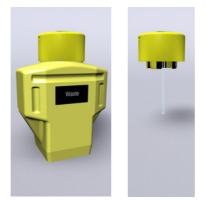

- **3** Remove the bottle from the rack and place it on a firm and even surface.
- 4 Place the spare container on the rack.
- **5** Check that the tubing adapter is properly inserted.
- 6 Empty the removed waste container. Treat the fluid as biohazardous waste.
- 7 Pour 500 mL hypochlorite solution in the waste container.
- 8 Clean the container with a brush.
- **9** Dispose of the fluid. Treat it as biohazardous waste.
- **10** Clean the interior of the waste container manually using tissues. Remove any residue.
- **11** Rinse the waste container with tap water.
- **12** Empty the container completely.

### **Roche Diagnostics**

**13** Wipe the waste inlet with a tissue.

**14** Reinsert the tubing adapter into the waste container. Press down firmly.

15 Press 🔘.

The maintenance action **Prime Fluid System** is automatically performed.

A message will inform you when the maintenance action is complete.

**16** Press 🛞.

# **Replace water inlet filter**

# → Utilities > Maintenance > Replace Water Inlet Filter

The water inlet filter in the external water container removes particulate matter from the water supply. You must replace this filter according to the maintenance action schedule.

The water container is placed on the rack next to the Cedex Bio instrument.

Operator time Approximately 10 minutes.

Prerequisites None

Tools and materials required

- □ Purified water
- $\Box$  Hypochlorite solution (0.6%)
- □ Tissues
- $\Box$  Water inlet filter



Make sure that you have read and understood section *Safety information* on page C-8. The following warning messages in particular are relevant:

• Injury through working solutions on page C-8.

# To replace the water inlet filter

- 1 Select the maintenance action Replace Water Inlet Filter.
- **2** Press **()**.

The maintenance definition screen is displayed.

- **3** Press  $\oslash$ .
- **4** Remove the tubing adapter from the water container.
- **5** Place the tubing adapter on a clean surface.

**6** Replace the water inlet filter.

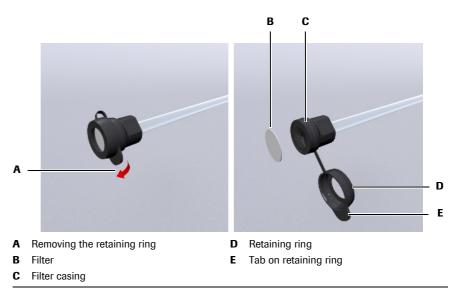


Figure C-2

- Remove the retaining ring by pulling it by its tab away from the filter casing (A, D, E).
- Remove the filter (B) and replace it with a new one.
- Place the retaining ring.

Press it firmly on the filter casing by placing your thumb across it an applying equal pressure all around.

- 7 Reinsert the tubing adapter into the water container. Press down firmly.
- 8 Press 🔘.

The maintenance action **Prime Fluid System** is automatically performed.

A message will inform you when the maintenance action is complete.

- **9** Press 🛞.

#### **Clean and disinfect housing**

#### → Utilities > Maintenance > Clean Housing

You must clean the housing regularly to ensure that contaminants do not build up and affect the efficiency of the instrument as a measuring system.

Operator time	Approximately 10 minutes.
---------------	---------------------------

Prerequisites Instrument is switched off.

Tools and materials required

Ethyl alcoholTissues



Make sure that you have read and understood section *Safety information* on page C-8. The following warning messages in particular are relevant:

- Injury through working solutions on page C-8.
- Infection by biohazardous materials on page C-8.
- To find out what steps are required, you can choose Utilities > Maintenance > Clean Housing and then press (). (The maintenance definition screen is displayed.)
- Do not press is at this stage. You do this only when you have completed the maintenance action. Instead, press is to close the screen.

#### To clean the housing

- 1 Make sure the system is in **Standby** status.
- 2 Shut down the system and switch off the instrument.
- **3** Wipe all the outside of the instrument with alcoholic cleansing tissues.
- **4** Open the main cover and the right service flap.
- 5 Wipe all the inside of the cover and flap with alcoholic cleansing tissues.
- **6** Close the right service flap and then the main cover.
- **7** Push the transfer head to the rightmost position.
- 8 Open the main cover and the left service flap.
- **9** Wipe all the inside of the flap with alcoholic cleansing tissues.
- **10** Close the left service flap and the main cover.
- **11** Push the transfer head to its home position.
- **12** Remove the transfer head cover.

(Press the release buttons on both sides and lift.)

- 13 Wipe all the inside of the cover with alcoholic cleansing tissues.
- **14** Reinstall the transfer head cover.

(Push down until the release buttons latch on.)

- **15** Remove the left and right side panels.
- 16 Wipe the panels with alcoholic cleansing tissues.
- **17** Reinstall the left and right side panels.

#### **Roche Diagnostics**

**18** Switch on the system.

The system performs internal checks and routines.

The startup phase may take a few minutes.

When the instrument is ready for use, the **Overview** tab is displayed, and its status is **Standby**.

- **19** Log on the system.
- **20** Choose Utilities > Maintenance > Clean Housing.
- **21** Press 💽.

The maintenance definition screen is displayed.

**22** Press 🔘.

A detailed maintenance definition screen is displayed.

**23** Press 🕖 to confirm the completion of the maintenance action.

#### Clean up the database

#### → Utilities > Maintenance > Cleanup Database

This maintenance action restores the database to the condition where the instrument can work at its optimum. You typically use this action when you think that the instrument suddenly works slow, for example when displaying information on the screen takes long or when it takes long until an action starts after you have initiated it.

Operator time Approximately 10 minutes.

*Prerequisites* The system must be in **Standby** status.

Tools and materials required None

#### ► To clean up the database

- 1 Select the maintenance action **Cleanup Database**.
- **2** Press **()**.

The maintenance definition screen is displayed.

**3** Press 🔘.

An instructions screen is displayed while the database cleanup is performed.

- **4** When the maintenance action is complete, press 🕖 to confirm the completion.

#### **Replace photometer lamp**

#### → Utilities > Maintenance > Replace Photometer Lamp

It is essential that the photometer lamp maintains a constant intensity in successive absorbance measurements. After a certain time of use, the intensity of the light emitted by the lamp deteriorates, and the accuracy of the measurements may no longer meet the required standards. Therefore, the lamp must be periodically replaced. The system alerts you when the lamp should be replaced by changing the maintenance action status to due (entry turns red).

Operator time Approximately 20 minutes.

*Prerequisites* The system must be in **Standby** status.

- D Photometer lamp assembly (Lamp, socket and cable are pre-assembled.)
- □ Cuvette segments
- $\Box$  Clean cloth



Make sure that you have read and understood section *Safety information* on page C-8. The following warning messages in particular are relevant:

- Injury through working solutions on page C-8.
- Infection by biohazardous materials on page C-8.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page C-8.

#### To replace the photometer lamp


- 1 Select the maintenance action **Replace Photometer Lamp**.
- **2** Press **()**.

The maintenance definition screen is displayed.

**3** Press 🔘.

The system switches off the lamp, and the transfer head moves to the home position.

- *Preparing the instrument* **4** Open the main cover and the right service flap.
  - **5** Remove the sample tray by lifting it.

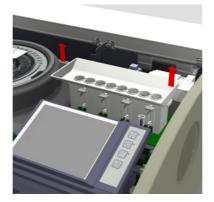




#### Injury through hot parts

Direct contact with hot parts may result in injury.

Do not touch the lamp assembly with bare hands. Let the lamp cool down for about ten minutes before touching any part of the lamp housing.

*Removing the lamp* 6

Unplug the power connector (A).

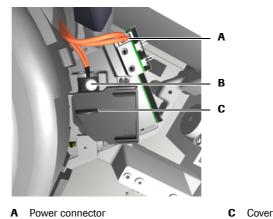

B Assembly screw

#### Figure C-3

- 7 Lift the photometer lamp housing cover (C).
- 8 Unscrew (B) the lamp assembly and remove it.


WARNING

Installing the new lamp

#### **9** Insert the new lamp assembly.

#### Incorrect results through soiled lamp


Touching the lamp with bare fingers reduces the life of the bulb and may affect the consistency of measurements made with the absorbance photometer. Hold the lamp assembly by its screw.

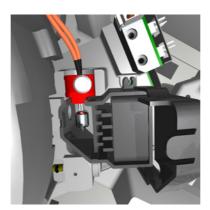

Hold the assembly by the screw and fit the lamp socket in the guides.

- **10** Fasten the screw of the lamp assembly.
- **11** Close the photometer lamp housing cover.
- **12** Plug the power connector into its socket.

#### *Preparing the instrument* **13** Reinsert the sample tray.

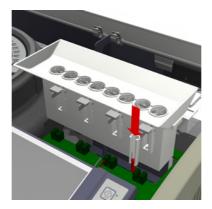

- 14 Close the right service flap and the main cover.
- 15 Press 🔘.

The system automatically initializes the instrument.

A message will inform you when the maintenance action is complete.

- **16** Press 🛞.
- **17** Perform an air/water calibration.

You cannot process orders unless you perform this action.

• See *Perform an air/water calibration* on page C-28.

#### **Contact service representative**

→ Utilities > Maintenance > Call Service Rep.

To ensure smooth operation of the instrument, a service representative periodically needs to perform some preventive maintenance tasks.

The system displays a message when this maintenance action is due.

Operator time Approximately 2 minutes.

(A service representative takes about 4 hours to perform preventive maintenance.)

#### Prerequisites None

#### ► To initiate preventive maintenance

- 1 A message screen reminds you that you should contact your service representative.
- 2 Select the maintenance action Call Service Rep.
- **3** Press **()**.

The maintenance definition screen is displayed.

4 Press 🔘.

An instructions screen is displayed.

- **5** Contact your service representative to arrange for preventive maintenance.
- 6 Press 🖉.

#### Perform an air/water calibration

#### → Utilities > Maintenance > Abs. Air/Water Calibration

The absorbance photometer calibration determines the air/water correction values that are needed for result calculation. In addition, the obtained absorbance values for an empty cuvette at each wavelength are used as a reference for the cuvette blank check.

Operator time Approximately 5 minutes.

*Prerequisites* The system must be in **Standby** status.

*Tools and materials required* 



Make sure that you have read and understood section Safety information on page C-8.

The following warning messages in particular are relevant:

• Injury through working solutions on page C-8.

 $\Box$  2 cuvette segments with empty cuvettes

- Infection by biohazardous materials on page C-8.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page C-8.

#### To perform an air/water calibration

- 1 Choose Utilities > Maintenance > Abs. Air/Water Calibration.
- **2** Press **()**.

Your are asked to insert a cuvette segment with empty cuvettes.

- **3** Open the main cover.
- **4** If there is a segment in the cuvette ring remove it and check whether there are used cuvettes.

If all cuvettes of this segment are empty, you can reinsert it.

- 5 Place a cuvette segment with empty cuvettes on the cuvette ring.
- 6 Press 🔘.

Your are asked to insert a second cuvette segment with empty cuvettes.

7 If there is a segment on in the cuvette ring remove it and check whether there are used cuvettes.

If all cuvettes of this segment are empty, you can reinsert it.

- 8 Place the cuvette segment with empty cuvettes on the cuvette ring.
- **9** Close the main cover.
- **10** Press 🔘.

The system performs the calibration measurements. (It first measures all cuvettes empty, pipettes water in the cuvettes, and measures them again.)

A message will inform you when the measurements are complete.

- **11** Scroll to display the results, if required.
- **12** Check the **Outlier statistics**.

**13** Press **(2)** to accept the results.

`Q́

If at this stage, you press (③), the results are not stored and the maintenance action is deemed not performed.

14 Remove the used cuvettes or replace them with empty ones.

#### Initialize degasser fluid sensor

$\rightarrow$	Utilities > Maintenance > Initialize Degasser Fluid Sensor
	The degasser fluid sensor checks whether there is air in the water supply to the syringe assembly. Air in the water supply may lead to inaccurate pipetting and ultimately to incorrect results. Therefore, the sensor needs to be initialized to ensure its proper functioning.
Operator time	Approximately 5 minutes.
Prerequisites	The system must be in <b>Standby</b> status.
	All covers must be closed.
Tools and materials required	□ Full external water container
	□ Clean cloth

#### ► To initialize the degasser fluid sensor

- **1** Make sure that all covers on the instrument are closed.
- 2 Choose Utilities > Maintenance > Initialize Degasser Fluid Sensor.

An instructions screen is displayed.

**3** Press 🔘.

The maintenance definition screen is displayed.

**4** Check whether the tube connecting the external water container with the instrument is full of water or contains air bubbles.




**5** Do one of the following:

lf	Do	o this	
You cannot see any air bub- bles	1.	Continue with step 6.	
You can see air bubbles or the tube is empty		Press 🛞 to abort this maintenance action. Perform the maintenance action <b>Prime Fluid Sys</b> -	
	3.	tem. Restart the maintenance action Initialize Degasser Fluid Sensor.	

**Roche Diagnostics** 

**6** Lift the tubing adapter from the white water container.

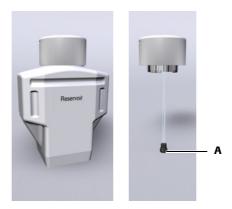

**7** Place the tubing adapter on a clean surface.

Make sure the filter (A) is not obstructed and air can get to it.

**8** Press 🖉 to initialize the fluid sensor.

The system aspirates air through the tubing adapter and initializes the fluid sensor.

When initialization is complete, you will be asked to insert the tubing adapter.

- **9** Insert the tubing adapter in the water container.
- **10** Press 🔘.

The fluid system is primed.

A message will inform you when the maintenance action is complete.

- **11** Press 🛞.

#### **Clean rotor and heating channel**

Perform this maintenance action whenever a message asks you to do so.

Operator	time	

Tools and materials required

□ 95% ethyl alcohol□ Lint free tissues

Approximately 15 minutes.



Make sure that you have read and understood section *Safety information* on page C-8. The following warning messages in particular are relevant:

- Injury through working solutions on page C-8.
- Infection by biohazardous materials on page C-8.

#### ► To clean the rotor and heating channel

- **1** Remove the reagent disk.
  - Choose **Overview** > 🔟 > 🔃.

A screen is displayed, asking you to remove the reagent disk.

- Open the main cover.
- Remove the reagent disk.
- Close the main cover.
- **2** Remove all cuvettes.
  - Choose **Overview** > **T**.
  - Open the main cover.
  - Press a segment button.
  - Remove the segment.
  - Press **T** to confirm the removal.


Remove all segments as described above.

- Close the main cover.
- **3** Shut down the system and switch off the instrument.
- **4** Move the transfer head to the sample area.
- **5** Open and remove the main cover and left service flap.




6 Move the transfer head to its leftmost position.

- **7** Open and remove the right service flap.
- **8** Remove the sample tray.

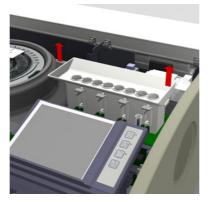

Removing the rotor

**9** Loosen the display fixation screw (A).

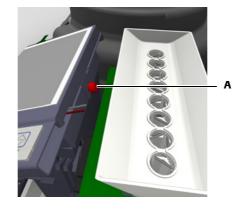

- **10** Tilt forward the display.
- **11** Move transfer head to its rightmost position.
- **12** Remove the upper rotor shell.

**13** For each of the two rotor bearings to the right, loosen slightly one of the fixation screws and completely unscrew the other. (Use a screw driver to loosen the screws, if required.) Then pull away from the rotor the side of the bearing that you completely unscrewed.

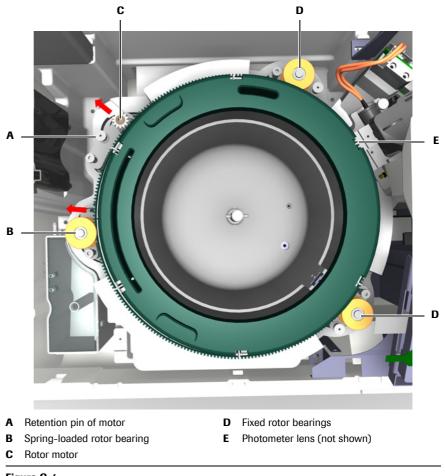



Figure C-4

**14** From the rotor motor, slightly lift the retention pin, move the motor away from the rotor and engage the pin to fix the motor in the removed position.

Make sure to use the hole nearer to the front of the instrument for arresting the bearing with the pin.

**15** With one hand, release the spring-loaded bearing from the rotor, and with the other hand lift the rotor.

*Cleaning the rotor* **16** Rinse the cuvette ring (A) under running water.

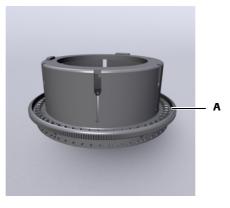


NOTICE

Do not use any cloth to clean or dry the rotor. This could leave dust and lint on the Abs. inspection holes and so prevent proper measurement.

- **17** Leave the rotor to dry.
- *Cleaning the heating channel* **18** Clean the heating channel (B) with a lint free tissue or a cloth moistened with ethyl alcohol.

Make sure not to touch the photometer lens (A).

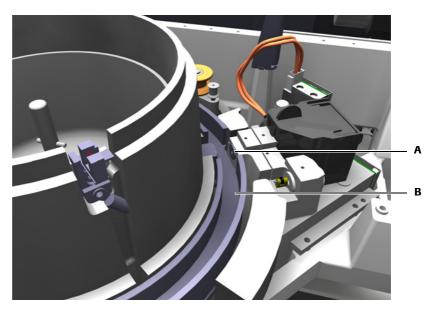


Reinstallation

- ion 19 With one hand, pull the spring-loaded bearing away from the rotor position, and with the other hand insert the rotor.
  - 20 Engage the right rotor bearings and fasten the fixation screws.
  - **21** Move the rotor to test its smooth running.
  - **22** Hold the spring-loaded rotor motor in position, lift the retention pin and release the motor gently.
  - 23 Move the transfer head to its leftmost position.
  - 24 Fit the upper rotor shell.
  - **25** Fold down the display and fasten the fixation screw.

#### **Roche Diagnostics**

**26** Install the sample tray.

Make sure to align the guides with the pins.

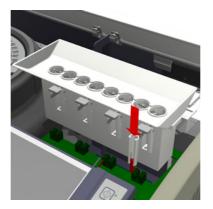

Press down firmly.

- **27** Install and close the right service flap.
- **28** Move the transfer head to its rightmost position.
- **29** Install and close the left service flap and main cover.

*Finishing the maintenance action* 

- **30** Switch on the instrument.
- **31** Wait for initialization to finish.
- 32 Log on to the system.
- **33** Insert the reagent disk.
  - See Preparing the reagent disk on page B-21

# D

## Troubleshooting

9	Messages and alarms	D-3
10	Result flags	D-23
11	Troubleshooting	D-39

## How to make the most of the available information

In this chapter, you will find information on messages generated by the **Cedex Bio** instrument, and on how to use them and react to them.

In this chapter	Chapter	9
About messages		D–5
Message screen		D–5
Acoustic signals		D–6
Alarm monitor		D–6
Reacting to alarm messages		D–7
Printing alarm messages		D–8
Confirming alarm messages		D–9
Exporting the alarm message log		D-9
List of alarm messages		.D–10

Table of contents

About messages

## **About messages**

The system performs numerous checks. When a certain event has occurred or when an irregularity is detected, a message is generated. Messages are displayed in two ways:

- Immediate feedback on user actions is displayed in a pop-up message screen.
- Information concerning a problem that occurred during operation is reported as *alarm messages* in the *alarm monitor*.

Alarm messages are stored in a log file, which you can export.

• See *Exporting the alarm message log* on page D-9.

### Message screen

Message screens are displayed automatically as soon as the message is generated.

Standby	10.12.2007 13:16 🕐
Invalid Data	
	insert all values.
	<b>S</b>



Read the message and press  $\textcircled{\otimes}$  to close the screen.

## **Acoustic signals**

The operator is alerted to certain events by an acoustic signal.

The following events trigger an acoustic signal.

- An alarm was generated.
- The Main cover is open while the system is in **Operating** status.
- A run is finished. (The system status has changed from **Operating** to **Standby**.)
- A maintenance action is complete (provided it is a longer maintenance action where no user intervention is required).

You can adjust the volume of all generated acoustic signals (**Utilities** > **Configuration** > **System** > **Volume**).

You can also turn on and off acoustic signals:

- Choose Utilities > Configuration > System > Sound Effects to turn on and off the acoustic signals other than alarm and warning signals.
- Choose Utilities > Configuration > System > Volume and set the value to zero to turn off all acoustic signals.
- System on page B-166.

## **Alarm monitor**

Messages concerning an irregularity that occurred during operation can be viewed in the alarm monitor. The alarm LED alerts you when such messages are generated.

#### Alarm button and LED

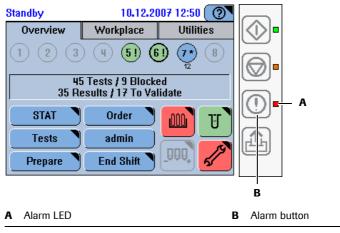


Figure D-2 Alarm LED

The alarm button ① is always active, even if you are not logged on the system.

#### **Reacting to alarm messages**


#### To react to an alarm message

#### **1** Observe the alarm LED.

□ No color Off	There are no unconfirmed alarm messages.
Yellow	There is at least one unconfirmed alarm message. You need to deal with it as soon as possible.
Red	There is at least one unconfirmed alarm message. You need to deal with it immediately, processing may not be able to continue unless you do so.
An acoustic signal is sounded when an alarm is generated. You can adjust the volume (Utilities > Configuration > System > Volume).	

#### **2** Press the **Alarm** button **①**.

The most recently generated alarm message in this list is displayed. Which messages are included in this list depends on the filter criteria that are applied to it.

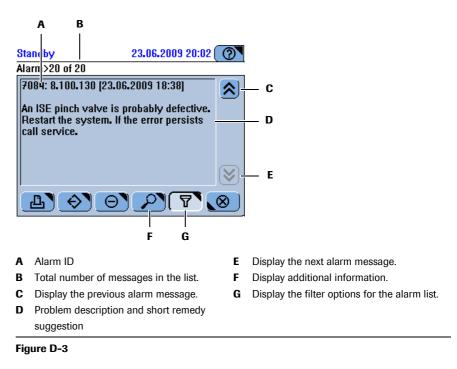



The alarm list can contain up to 250 messages. When this number is reached, the oldest message is automatically deleted when a new one is generated.

**3** To limit the number of messages that are contained in the list, press **Y**.

A list of filter options is displayed.

- Select **Not confirmed** to include in the list all alarm messages that have not yet been confirmed by the user.
- Select **Today's alarms** to include in the list all alarm messages that so far have been generated on the current day.
- Select **Complete list** to include in the list all alarm messages that are stored in the alarm database.

- 4 Select the message you are interested in. Use ▲ and ¥ to display the previous and next message respectively.
- **5** Do one of the following:
  - Press *P* to display more detailed information.

(If this button is not available there is no additional information.)

- Press 🕒 to print alarm messages.
- Press 📀 to export the log file.
  - See *Exporting the alarm message log* on page D-9.
- 6 When you have dealt with the issue, press ⊖ to confirm that you have dealt with the issue. The message is deleted from the **Not confirmed** list but it remains in the other lists and the alarm messages log, which you can export.
  - See *Confirming alarm messages* on page D-9.

#### **Printing alarm messages**

When you print alarm messages, their details are included in the printout.

#### To print alarm messages

- 1 Press ①.
- 2 Press 🔔.

A screen is displayed for defining which alarm messages should be printed.

- **3** Do one of the following:
  - Press All to print all messages that have not been dealt with.
  - Press **Selected** to print the current message.

Alarm monitor

#### **Confirming alarm messages**

Confirming a message means that you have dealt with the issue mentioned in it.

- ► To confirm a message
  - 1 Press ①.
  - 2 Select the message. (Use ▲ and ♥ to display the previous and next message respectively.)
  - **3** Press  $\Theta$ .

A confirmation dialog box is displayed.

- Press All to remove all messages from the list.
- Press Selected to remove the currently displayed message.

The message or messages are removed from the **Not confirmed** list, but they remain in the other lists and the alarm messages log, which you can export.

When all messages are confirmed the Alarm LED is turned off (not lit).

#### Exporting the alarm message log

#### ► To export the message log file

- 1 Press ①.
- **2** Press 📀.

A screen is displayed asking you to insert a USB stick.

- **3** Insert the USB stick.
- **4** Press 🔘 to confirm the insertion.

A screen is displayed for selecting a directory.

Press \*.tgz to display log files only.

Press \*.\* to display all files and directories.

- **5** Select a directory for the error log.
- **6** Press 🖉 to confirm the selection.

All log files are added to a file with the name format log\_yyyymmddhhmmss.tgz and copied to the stick. The error log file has the name format err\_yyyymmddhhmmss.log.

List of alarm messages

## List of alarm messages



In the following table, the variables %s and %d are used, they represent text and numbers.

Unknown message, id=%d! A fatal system exception software error (%d) occurred.	
A fatal system exception software error (%d) occurred	
A latar system exception software error (%u) occurred.	Restart the system. If the error recurs export the log files while the system stays in the error condition. If the error persists call service.
An internal software error (%d) occurred while updating the configuration.	Restart the system. If the error recurs export the log files while the system stays in the error condition. If the error persists call service.
A software error occurred. Order (%d, %d) could not be processed.	Delete the orders that are not processed yet. If the error recurs export the log files while the system stays in the error condition. If the error persists call service.
A software error occurred. The unit is not ready, order (%d) rejected.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
A mathematical exception arose.	Repeat the order. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
Last calculation cycle (%d) too short.	Repeat the order. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
Failed to calculate linear regression.	Repeat the order. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
The cycle range (%d%d) is not valid.	Repeat the order. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
The reaction direction (%d) is invalid.	Repeat the order. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
An internal communication error occurred. The received IPC message (%d, %d) could not be interpreted.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
A software error occurred. The status of an order is invalid (%d.%d).	Delete and redefine the order. If the error recurs export the log files with the system remaining in this error condi- tion. If the error persists call service.
The installed language version is out of date (%s, %s).	Install the language version that belongs to the currently installed system software version. If the error persists call service.
The current system date and time are invalid.	Correct the date and time settings in Configuration. If the error persists call service.
Printer controller revision %d is not supported.	Call service.
	the configuration. A software error occurred. Order (%d, %d) could not be processed. A software error occurred. The unit is not ready, order (%d) rejected. A mathematical exception arose. Last calculation cycle (%d) too short. Failed to calculate linear regression. The cycle range (%d%d) is not valid. The reaction direction (%d) is invalid. An internal communication error occurred. The received IPC message (%d, %d) could not be interpreted. A software error occurred. The status of an order is invalid (%d.%d). The installed language version is out of date (%s, %s). The current system date and time are invalid.

ID	Message	Comment
0020	This reagent set has more bottles than there is space available on the disk.	Load another disk or remove unused reagents.
0021	Disk space problem.	
0022	You are about to delete results that have been printed or sent to host!	
0023	Enter the A1c part.	
0024	Now enter the A1c part.	
5001	Air detected in the fluid system. Water container or cleaner bottle is empty.	
5002	[%d] Air detected in the fluid system.	Refill the water container. Check the filling level of the cleaner bottle. Replace the cleaner bottle if required. Perform <prime fluid="" system="">. If the error persists call service.</prime>
5003	[%s] An internal software error occurred. A parameter is not valid.	Restart the system. If the error persists call service.
5004	[%s] Error syringe motor. The syringe motor could not be started.	Restart the system. If the error persists call service.
5010	[%s] Initialization of syringe failed.	Switch off instrument. Remove right side panel. Check for syringe mechanism jamming or blockage. Ensure correct syringe installation. Restart the system. If the error per- sists call service.
5013	[%s] Error mixer motor. The mixer motor could not be started.	Restart the system. If the error persists call service.
5019	[%s] Initialization of the mixer motor failed.	Switch off the instrument. Remove the transfer head cover. Check if the mixing mechanism is jamming or blocked. Restart the system. If the error persists call ser- vice.
5022	[%s] Error temperature sensor analyzer heating system. The sensor does not react anymore.	Restart the system. If the error persists call service.
5024	[%s] Error temperature sensor analyzer cooling system. The sensor does not react anymore.	Restart the system. If the error persists call service.
5026	[%s] Error temperature sensor cooling system. The heat sink temperature sensor does not react anymore.	Restart the system. If the error persists call service.
5028	[%s] Error temperature sensor instrument. The ambient temperature sensor does not react anymore.	Restart the system. If the error persists call service.
5030	[%s] Error wash pump P1. The wash pump could not be started.	Restart the system. If the error persists call service.
5032	[%s] Error waste pump P2. The waste pump could not be started.	Restart the system. If the error persists call service.
5034	[%s] Error wash valve V1. The valve does not react anymore.	Restart the system. If the error persists call service.
5036	[%s] Error cleaner valve V2. The valve does not react anymore.	Restart the system. If the error persists call service.
5038	[%s] Error analyzer heating system. The heating system does not react anymore.	Restart the system. If the error persists call service.
5040	[%s] Error cooling system. One of the Peltier elements does not react anymore.	Restart the system. If the error persists call service.
Table D-1	List of alarm messages (Continued)	

ID	Message	Comment
5044	[%s] Error main board fan. The fan could not be started or the fan speed is not correct.	Restart the system. If the error persists call service.
5047	[%s] Error cooling fan. The fan could not be started.	Restart the system. If the error persists call service.
5049	[%s] Error sample area fan. Switch off the instrument.	Remove the right side panel and check if the fan is jam- ming or blocked. If possible eliminate the problem. Restart the system. If the error persists call service.
5052	[%s] Error cooling system. The cooling heat sink overheated.	Ensure unrestricted cooling air circulation. Check the cooling fan and make sure the external vent is not blocked. Restart the system. If the error persists call ser- vice.
5054	[%s] The analyzer heating system overheated.	Switch off the instrument and wait 15 minutes to let cool down the analyzer. Restart the system. If the error persists call service.
5055	[%s] The analyzer temperature is outside the range.	Restart the system. If the error persists call service.
5056	[%s] The calibration of the transfer head fluid sensor failed.	Ensure that the sample probe tube is correctly installed. Ensure that the syringe is correctly installed. Restart the system. If the error persists call service.
5058	[%s] Error fluid sensor transfer head. The sensor does not react anymore.	Restart the system. If the problem persists, call service.
5060	[%s] A not specified hardware error (%d) was reported by the Multi-Slave control.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
5061	[%d] Air detected in the fluid system.	Refill the water container. Check the filling level of the cleaner bottle. Replace the cleaner bottle if required. Per- form <prime fluid="" system="">. If the error persists call ser- vice.</prime>
5062	The degasser fluid sensor is not initialized.	Perform <initialize degasser="" fluid="" sensor="">. Follow the instructions provided on the display. If the error persists call service.</initialize>
5063	[%s] Air bleed valve error. The air bleed valve does not react anymore.	Restart the system. If the error persists call service.
5064	[%s] Error fluid sensor degasser. The sensor does not react anymore.	Restart the system. If the error persists call service.
6001	[%s] Photometer error. Signal is invalid or unstable.	Switch off the instrument. Remove and clean the rotor. In case of a recurrence perform <replace photometer<br="">Lamp&gt; and <abs air="" calibration="" water="">. If the error persists call service.</abs></replace>
6002	[%s] Photometer error. The photometer has a wrong status because initialization has not been successfully performed yet.	Restart the system. If the error persists call service.
6018	[%s] A not specified photometer hardware error (%d) was reported by the photometer.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
6019	[%s] Photometer error. The photometer signals were outside the range while performing a photometer initialization or a photometer measurement.	Restart the system. If the error persists call service.
Table D-1	List of alarm messages (Continued)	

ID	Message	Comment
6020	[%s] Photometer lamp error.	Perform <clean and="" channel="" heating="" rotor="">. If the error persists, replace the photometer lamp. Perform first the maintenance action <replace lamp="" photometer=""> and then <abs air="" calibration="" water="">.</abs></replace></clean>
6021	[%s] Photometer error. The lamp voltage is outside the range.	Restart the system. If the error persists call service.
6022	[%s] Photometer error. The internal photometer 15 V power supply failed.	Restart the system. If the error persists call service.
6023	[%s] Photometer error. The lamp current is outside the range.	Perform first the maintenance action <replace photome-<br="">ter Lamp&gt; and then <abs air="" calibration="" water="">. Restart the system. If the error persists call service.</abs></replace>
6024	[%s] Photometer error. Either the lamp fan is not running or its speed is not correct.	Switch off the instrument. Open the main cover and check if the lamp fan is jamming or blocked. Restart the system. If the error persists call service.
6025	[%s] Photometer error. Either the dark shutter motor does not work or the shutter mechanism is jamming or blocked.	Restart the system. If the error persists call service.
6026	[%s] Photometer error. A problem caused by the internal circuitry has been detected.	Restart the system. If the error persists call service.
6027	[%s] Photometer error. The photometer signals are outside the range while performing a dark measurement.	Restart the system. If the error persists call service.
6029	[%s] Photometer error. A problem caused by the internal circuitry has been detected.	Restart the system. If the error persists call service.
6030	[%s] Initialization of the rotor failed.	Switch off the instrument. Remove the rotor and clean all rotor measurement holes. Reinstall the rotor. Restart the system. If the error persists call service.
6031	[%s] Photometer error. The photometer dark signal is not stable.	Restart the system. If the error persists call service.
7001	[%d] An Instrument Control software exception error occurred.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7002	[%d] A software error occurred.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7004	[%d] An internal communication software error occurred.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7006	[%s] A not specified hardware error (%d) was reported.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7007	[%d] An internal communication error (%d) occurred. A control unit could not be connected.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7008	[%d] A software error occurred that was caused by an unexpected system situation.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
Table D-1	List of alarm messages (Continued)	

List of alarm messages

ID	Message	Comment
7009	The LAN connection to the IC software failed.	Check the external Ethernet cable. Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7010	Instrument initialization failed.	Check for previous errors. Switch off instr. Check transfer mechanism. Restart system. If error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7011	Failed to aspirate sample from tube on position %d. This is probably a consequential error.	Ensure that there is enough sample in the tube.
7012	Failed to identify the reagent disk. This is probably a consequential error.	Check for previous errors and eliminate the cause. Remove the reagent disk and check the ID tabs. Reinsert the reagent disk correctly.
7013	[%s] Position error X-transfer. Switch off the instrument.	Check if the X-transfer is blocked or jamming. Eliminate the cause. Restart the system. If the error persists call ser- vice.
7014	[%s] %s electrode is not stable during sample measurement.	Perform restart. Perform <electrode service="">. Check ISE Ref solution tubing for obstructions. Replace electrode.</electrode>
7015	[%s] %s electrode is not stable during calibrator measurement.	Perform restart. Perform <electrode service="">. Check ISE Ref solution tubing for obstructions. Replace electrode.</electrode>
7016	[%s] The ISE sample sensor could not detect ISE Deproteinizer.	Check the ISE Deproteinizer level, proper electrode mounting, tube leakages and obstructions.
7017	[%s] The ISE sample sensor could not detect ISE Etcher.	Check the ISE tower and the ISE tubing for obstructions and leaks. Ensure that the tubing is correctly inserted into the sample sensor. If the error persists call service.
7018	[%d] The probe is not straight or not properly mounted.	Check that the probe is properly mounted. Perform <replace probe="">.</replace>
7019	[%d] Transfer initialization failed due to a level detection problem.	Clean the transfer init plate. Ensure that the probe is not bent and installed correctly. Restart the system. If the error persists call service.
7020	[%d] Transfer initialization failed due to an open main cover.	Ensure that all covers are closed. Restart the system.
7021	[%d] Instrument failed to initialize due to dirty init plate or probe.	Perform <clean manually="" probe=""> and restart the system. If the problem persists, call service.</clean>
7022	[%s] Motor %d movement impeded.	Perform restart. If the problem persists call service.
7033	[%s] Position error Z-transfer.	Switch off the instrument. Check if the Z-transfer is blocked or jamming. Eliminate the problem. Restart the system. If the error persists call service.
7034	[%s] Either a rotor or an X-transfer or a Z-transfer motion control error occurred.	Restart the system. If the error persists call service.
7036	[%s] The level detection frequency is outside the range.	Clean the probe and probe holder. Replace the water inlet filter. Ensure that the system water meets the specifica- tions. Restart the system. If the error persists call service.
7037	[%s] The level detection frequency is outside the range.	Clean the probe and probe holder. Check the level detec- tion cable for damage. Ensure that the cable is connected correctly. Restart the system. If the error persists call ser- vice.
Table D-1	List of alarm messages (Continued)	

Table D-1

List of alarm messages (Continued)

ID	Message	Comment
7038	[%s] Motor %d error (overheated).	Switch off the instrument. Check for smooth transfer movements. Restart the system. If the problem persists call service.
7039	[%s] Power supply error. One or more internally used voltages are outside the range.	Switch off the instrument. Check the fuses F1 through F4 and replace blown fuses. Restart the system. If the error persists call service.
7042	[%s] Fuse F1 blown.	Switch off the instrument. Check fuse F1. If necessary replace fuse F1. Restart the system. If the error persists call service.
7043	[%s] Fuse F2 blown.	Switch off the instrument. Check fuse F2. If necessary replace fuse F2. Restart the system. If the error persists call service.
7044	[%s] Fuse F3 blown.	Switch off the instrument. Check fuse F3. If necessary replace fuse F3. Restart the system. If the error persists call service.
7045	[%s] Fuse F4 blown.	Switch off the instrument. Check fuse F4. If necessary replace fuse F4. Restart the system. If the error persists call service.
7046	[%s] The initialization of the rotor failed.	Switch off the instrument. Check if the rotor is jamming or blocked. Clean the rotor initialization light barrier. Restart the system. If the error persists call service.
7051	[%s] Error rotor motor. The motor could not be started.	Restart the system. If the error persists call service.
7052	[%s] Error motor X-transfer. The motor could not be started.	Restart the system. If the error persists call service.
7053	[%s] Error motor Z-transfer. The motor could not be started.	Restart the system. If the error persists call service.
7057	[%s] The ISE sample sensor is not calibrated or no calibration data is available.	Perform <initialize ise="" unit="">. If the error persists call service.</initialize>
7058	[%s] ISE sample sensor could not detect fluid.	Check the ISE tower and the ISE tubing for obstructions and leaks. Ensure that the electrodes are installed cor- rectly. If the error persists call service.
7059	[%s] The ISE sample sensor could not detect the end of the transported fluid segment.	Check the ISE tower and the ISE tubing for obstructions and leaks. If the error persists call service.
7060	[%s] The calibration of the ISE sample sensor failed.	Check the filling level of the ISE Cal bottle. Check the ISE tubing for obstructions and leaks. Perform <initialize ise="" unit="">. If the error persists call service.</initialize>
7061	[%s] The calibration of the ISE sample sensor failed.	Check the filling level of the ISE Cal bottle. Check the ISE tubing for obstructions and leaks. Perform <initialize ise="" unit="">. If the error persists call service.</initialize>
7062	[%s] The ISE sample sensor could not detect calibrator.	Check the filling level of the ISE Cal bottle. Check the ISE tubing for obstructions and leaks. Perform <initialize ise="" unit="">. If the error persists call service.</initialize>
7063	[%s] The ISE sample sensor could not detect the end of the calibrator segment.	Check the ISE tubing for obstructions and leaks. Perform <initialize ise="" unit="">. If the error persists call service.</initialize>
7064	[%s] The ISE sample sensor could not detect Activator during initialization.	Check the ISE tower and ISE tubing for obstructions and leaks. Perform <initialize ise="" unit="">. If the error persists call service.</initialize>
7065	[%s] Calibration of the ISE reference sensor failed, no valid calibration data are available.	Check the filling level of ISE Ref. Perform <initialize ise<br="">Reference Sensor&gt;. If the error persists call service.</initialize>
Table D-1	List of alarm messages (Continued)	
Roche Diag		

List of alarm messages

ID	Message	Comment
7066	An internal control unit (%d) reported a wrong firmware version.	Reinstall the system software.
7067	An internal control unit (%d) reported a checksum or an attach error.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7068	[%d] Wash aborted due to open cover.	Close the cover and restart the system.
7069	[%d] Transfer initialization failed due to an open main cover.	Ensure that all covers are closed. Restart the system.
7070	The available reagent bottle volume in bottle set %s is not sufficient. The bottle set has been marked as empty.	Replace the bottle set with a new one.
7071	Photometer error. Signal is invalid or unstable.	Switch off the instrument. Remove and clean the rotor. If the error recurs perform <replace lamp="" photometer=""> and <abs air="" calibration="" water="">. If the error persists call service.</abs></replace>
7072	[%d] The reagent mixing failed due to a level detection problem.	Check if the probe is obstructed. Check the filling level of the associated reagent bottle. If the error persists call ser- vice.
7073	[%s] Peristaltic pump speed is outside the range.	Check ISE tubing for obstructions and leaks. Ensure that the electrode tension lever and the peristaltic pump cover are closed. Perform <initialize ise="" unit="">. If the error per- sists call service.</initialize>
7074	[%s] The detected sample or calibrator segment is too short.	Check the ISE tubing for obstructions and leaks. Ensure that the electrodes are installed correctly. If the error per- sists call service.
7075	[%s] A timeout error occurred while performing an ISE measurement.	Restart the system and repeat the measurement. If the error persists call service.
7076	[%s] Initialization of the ISE unit failed due to a fluid transport problem.	Check the ISE tower and the ISE tubing for obstructions and leaks. Check the electrodes. Perform <initialize ise<br="">Unit&gt;. If the error persists call service.</initialize>
7083	[%s] The ISE bypass tubing is blocked.	Perform <replace ise="" tubing="" unit="">. Follow the instruc- tions in the Operator's Manual. If the error persists call service.</replace>
7084	[%s] An ISE pinch valve is probably defective.	Restart the system. If the error persists call service.
7086	[%s] The ISE air pressure is outside the range.	Check the air tubing underneath the ISE tower for obstructions. If the error persists call service.
7087	[%s] ISE air pressure sensor detected that the pressure is too low.	Restart the system. If the problem persists, call service.
7088	[%s] A failure of the internal ISE power supply has been detected.	Restart the system. If the error persists call service.
7099	[%s] Weight sensors of the ISE Ref and ISE Cal bottle are not calibrated correctly.	In <diagnostics> perform <adjust bottle="" ise="" ref="" sensor=""> and <adjust bottle="" cal="" ise="" sensor="">. If the error persists call service.</adjust></adjust></diagnostics>
7100	A power fail has been detected and an automatic system restart was performed.	
7101	[%s] Rotor position error. Switch off the instrument.	Check if the rotor is blocked or jamming. Eliminate the cause. Restart the system. If the error persists call service.
7102	[%s] Transfer bumper init failed. Perform restart.	
Table D-1	List of alarm messages (Continued)	

#### **Roche Diagnostics**

List of alarm messages

ID	Message	Comment
7103	[%s] Adjustment of the ISE Ref bottle or ISE Cal bottle weight sensor failed.	Repeat the adjustment: In <diagnostics> perform <adjust ISE Ref Bottle Sensor&gt; and <adjust bottle="" cal="" ise="" sen-<br="">sor&gt;. If the error persists call service.</adjust></adjust </diagnostics>
7107	[%s] A not specified hardware error (%d) was reported by the ISE Unit.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7108	[%s] A not specified hardware error (%d) was reported by the DC-Slave control.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7109	[%d] The sample probe is bent or not properly aligned.	Check the probe and replace it if necessary. Restart the system. If the error persists call service.
7110	[%d] A timing error occurred while fluid is expelled into a cuvette. Probably the error was caused by a previous hardware problem.	Check for previous errors. Restart the system. If the error persists call service.
7111	[%d] Transfer movement failed due to open cover.	Keep covers closed while the system is in Operating status.
7112	[%s] The ISE reference sensor detected air when ISE Ref solution was transported.	Check the ISE Ref bottle filling level. Perform <prime ise<br="">Reference and Calib.&gt; and check the fluid flow. If the error persists call service.</prime>
7113	[%d] A prime cycle failed due to an open cover.	Ensure that all covers are closed while the system is being initialized, in the Operating or Maintenance status.
7114	[%d] An initialization error occurred. The X-transfer zero positions are invalid.	Check the X-transfer for obstructions. Ensure that all cov- ers are closed. Restart the system. If the error persists call service.
7115	[%d] Dispensing into the ISE tower stopped because the ISE tower was not empty.	Check the ISE tower waste tubing. Check the ISE tower outlet for obstructions. Check the pinch valves. If the error persists call service.
7116	[%d] The needle crashed while the system was trying to dispense fluid.	Switch off the instrument. Check if the transfer or probe movement is obstructed. Restart the system. If the error persists call service.
7117	[%d] Consequential error caused by a previous hardware or software problem.	Check previous errors. Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7118	[%s] Rotor initialization error. The light barrier did not react.	Switch off the instrument. Remove the rotor and clean the light barrier. Reinstall the rotor. Restart the system. If the error persists call service.
7119	[%d] Consequential error caused by a previous hardware or software problem.	Check previous errors. Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7120	[%s] Either a rotor, an X-transfer or a Z-transfer position error occurred. Maybe one of these devices has been moved manually by accident.	Restart the system. If the error persists call service.
7121	[%d] Consequential error caused by an internal communication problem.	Check previous errors. Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7122	[%d] The Prepare action has been stopped due to an open cover.	Ensure that all covers are closed while the system is being initialized, in the Operating or Maintenance status.
7123	Extra wash cycle after reagent pipetting failed.	Perform <deproteinize probe=""> before next run.</deproteinize>
7124	Analyzer temperature outside the range.	
Table D-1	List of alarm messages (Continued)	

#### **Roche Diagnostics**

ID	Message	Comment
7125	[%s] Motion control error in rotor motor.	Restart the system to correct the problem. If the error per- sists call service.
7126	[%s] Motion control error X-transfer motor.	Restart the system to correct the problem. If the error per- sists call service.
7127	[%s] Motion control error Z-transfer motor.	Restart the system. If the error persists call service.
7128	[%s] Initialization error X-transfer.	Switch off the instrument. Ensure that all covers are closed. Check if the X-transfer is blocked or jamming. Restart the system. If the error persists call service.
7129	[%s] Initialization error Z-transfer.	Switch off the instrument. Check if the Z-transfer is blocked or jamming. Check the seating of the transfer head cover. Restart the system. If the error persists call service.
7130	[%s] Rotor position error.	Switch off the instrument and check if the rotor is blocked or jamming. If the error recurs call service.
7131	[%s] Position error X-transfer.	Switch off the instrument and check if the X-transfer is blocked or jamming. If the error recurs call service.
7132	[%s] Initialization error Z-transfer.	Switch off the instrument. Check if the Z-transfer is blocked or jamming. Check the seating of the transfer head cover. Restart the system. If the error persists call service.
7133	[%s] Rotor motor overheated.	Switch off the instrument. Check if the rotor is blocked or jamming. Restart the system. If the error persists call service.
7134	[%s] Motor X-transfer overheated.	Switch off the instrument. Check if the X-transfer is blocked or jamming. Ensure that the X-transfer can be moved smoothly. Restart the system. If the error persists call service.
7135	[%s] Motor Z-transfer overheated.	Switch off the instrument. Check if the Z-transfer is blocked or jamming. Ensure that the Z-transfer can be moved smoothly. Restart the system. If the error persists call service.
7136	[%s] Tube bottom detector is blocked or jamming.	Switch off the instrument. Remove the transfer head cover. Check the tube bottom detector mechanism. Install the transfer head cover. Restart the system. If the error persists call service.
7137	[%s] Na sample measurement unstable.	Check the expiry date of the Ref and Na electrodes. Replace expired electrodes. Check the ISE tubing for obstructions and leaks. Perform <electrode service="">. If the error persists call service.</electrode>
7138	[%s] Cl sample measurement unstable.	Check the expiry date of the Ref and Cl electrodes. Replace expired electrodes. Check the ISE tubing for obstructions and leaks. Perform <electrode service="">. If the error persists call service.</electrode>
7139	[%s] K sample measurement unstable.	Check the expiry date of the Ref and K electrodes. Replace expired electrodes. Check the ISE tubing for obstructions and leaks. Perform <activate electrodes="">. If the error per- sists call service.</activate>
7140	[%s] Na calibrator measurement unstable.	Perform <electrode service=""> and <prime ise="" reference<br="">and Calib.&gt;. Check ISE Cal and Ref flow. Check the ISE tubing for obstructions. If the error persists call service.</prime></electrode>
Table D-1	List of alarm messages (Continued)	
Roche Diag		
	1051105	

ID	Message	Comment
7141	[%s] Cl calibrator measurement unstable.	Perform <electrode service=""> and <prime ise="" reference<br="">and Calib.&gt;. Check ISE Cal and Ref flow. Check the ISE tubing for obstructions. If the error persists call service.</prime></electrode>
7142	[%s] K calibrator measurement unstable.	Perform <electrode service=""> and <prime ise="" reference<br="">and Calib.&gt;. Check ISE Cal and Ref flow. Check the ISE tubing for obstructions. If the error persists call service.</prime></electrode>
7143	[%d] An internal software problem (%d) occurred. The access to the internal EEPROM failed.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
7144	[%d] The instrument battery is low or not connected. The date and time may no longer be valid. A replacement of the battery is probably required.	Call service.
7145	[%d] The action <clean ise=""> has been stopped due to an open cover.</clean>	Ensure that all covers are closed while the system is being initialized, or in the Operating or Maintenance status.
7146	[%d] The initialization of the system has been stopped due to an open cover.	Ensure that all covers are closed while the system is being initialized, or in the Operating or Maintenance status.
7147	[%d] The initialization of the ISE Unit has been stopped due to an open cover.	Ensure that all covers are closed while the system is being initialized, or in the Operating or Maintenance status.
7148	[%d] Initialization of the ISE Unit has been stopped due to a blocked transfer.	Switch off the instrument. Check if the transfer is blocked or jamming. Ensure that all covers are closed. Restart the system. If the error persists call service.
7149	[%d] Initialization of the rotor failed. The calculated offset is outside the range.	Switch off the instrument. Check if the rotor is blocked or jamming. Restart the system. If the error persists call ser- vice.
7150	[%d] Initialization of the transfer failed due a level detection problem.	Switch off the system. Clean the probe and probe holder and ensure that they are installed correctly. Restart the system. If the error persists call service.
7151	[%d] Dispensing fluid into the ISE tower failed due to an incorrect X-transfer or probe position.	Switch off the instrument. Check if the X- transfer and Z- transfer are blocked or jamming. Restart the system. If the error persists call service.
7152	[%d] A pipetting or dispensing cycle failed due to an incorrect X-transfer or probe position.	Switch off the instrument. Check if the X- transfer and Z- transfer are blocked or jamming. Restart the system. If the error persists call service.
7153	[%d] A reagent mixing cycle failed because the probe crashed.	Switch off the instrument. Check if the probe and probe holder are correctly installed. Replace a bent probe. Restart the system. If the error persists call service.
7154	[%d] A probe wash cycle failed because the probe crashed.	Switch off the instrument. Check if the probe and probe holder are correctly installed. Replace a bent probe. Restart the system. If the error persists call service.
7155	[%d] Initialization of the transfer failed because the probe crashed.	Switch off the instrument. Check if the probe and probe holder are correctly installed. Replace a bent probe. Restart the system. If the error persists call service.
7156	[%d] A prime cycle failed because the probe crashed.	Switch off the instrument. Check if the probe and probe holder are correctly installed. Replace a bent probe. Restart the system. If the error persists call service.
7157	[%d] Initialization of the Z-transfer failed due to a level detection problem.	Switch off the instrument. Check if the probe and probe holder are correctly installed. Clean the init plate. Restart the system. If the error persists call service.

ID	Message	Comment
7158	[%d] A tube bottom sensor problem was detected.	Switch off the instrument. Check if the tube bottom sen- sor mechanism can be moved freely. Restart the system. If the error persists call service.
7159	[%d] Movement stopped because the probe crashed.	Switch off the instrument. Check if the probe and probe holder are correctly installed. Replace a bent probe. Check the magnetic slider. Restart the system. If the error per- sists call service.
7160	Sample pipetting on position %d was aborted due to an unstable LD signal%s.	
7161	Reagent pipetting of bottle set %s was aborted due to an unstable level detection signal%s. The test is automatically repeated.	
7162	The probe crashed due to a jamming main cover slider.	Check the slider and ensure that it is operating correctly.
7163	The run aborted due to a jamming main cover slider.	Check the slider and ensure that it is operating correctly.
7164	An internal communication error occurred. Controller %s could not be connected.	Restart the system. If the error persists call service.
12001	The host computer did not answer the order query for sample %s within the time limit.	Check the communication and existing orders on the host.
12002	The communication port is not accessible. Communication to the host server is not possible.	Check the Host Settings and compare them with the host server settings.
12003	Sending a message to the host computer failed.	Enable the host trace file in the Host Settings in order to record HIF communication activities.
12004	Receiving or processing a message from the host computer failed.	Enable the host trace file in the Host Settings in order to record HIF communication activities.
12005	Host server could not send message to the host.	Select result and perform Retransmit. If the problem recurs enable the host trace file in the Host Settings. Retransmit again and export the log files. If the error per- sists call service.
12006	Changes of the host communication settings could not be applied.	Restart the system. If the error persists call service.
13001	A software error (%d) occurred. The database could not be opened.	Restart the system. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
13002	A software error occurred. The calculated result for the order (%d) could not be stored in the database.	Export the log files. If the error persists call service.
13003	A software error occurred. Data of the order (%d) could not be retrieved from the database.	Export the log files. If the error persists call service.
13004	A software error occurred. The required parameters for application (%d) could not be found.	Export the log files. If the error persists call service.
13005	A software error occurred. The calibration result for order (%d) could not be stored in the database.	Export the log files. If the error persists call service.
13006	A software error occurred. A required parameter set for a certain application (%d) could not be found.	Export the log files. If the error persists call service.
13007	Reagent disc %d was replaced by disc %d while the system was switched off.	Ensure that the information provided in <disk status=""> corresponds to the actually loaded reagent bottle sets.</disk>
13008	The import of the database failed. The database version does not correspond to the currently installed software version. An old database version was provided (detected V %d ; expected V %d ).	
Table D-1	List of alarm messages (Continued)	
Roche Diag	nostics	

#### 9 Messages and alarms

List of alarm messages

ID	Message	Comment
13009	A software error occurred. An invalid database entry (%d) for a control was detected.	Export the log files. If the error persists call service.
13010	A software error occurred. An invalid database entry (%d) for a control was detected.	Export the log files. If the error persists call service.
13011	Validation of Development Channels failed. The number of installed DC applications does not match the number of available channels. (%d DC applications found, %d channels installed).	
13012	Auxiliary reagent (%d) not found.	Ensure that the reagent bottle has been loaded. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
13013	Reagent bottle %d not found.	Ensure that the reagent bottle has been loaded. If the error recurs export the log files with the system remaining in this error condition. If the error persists call service.
13014	Calibrator lot definition for application %d inconsistent. Check number of standards used with number of calibrator cups and Last Is Water definition.	
20001	<replace filter="" inlet="" water=""> completed.</replace>	
20002	<clean and="" waste="" water=""> completed.</clean>	
20003	<replace lamp="" photometer=""> completed.</replace>	
20004	<clean manually="" probe=""> completed.</clean>	
20005	<prime fluid="" system=""> completed.</prime>	
20006	<clean housing=""> completed.</clean>	
20007	<deproteinize probe=""> completed.</deproteinize>	
20008	<clean disk="" reagent=""> completed.</clean>	
20009	<database backup=""> completed.</database>	
20010	<call rep.="" service=""> completed.</call>	
20011	<clean ise="" manually="" tower=""> completed.</clean>	
20012	<replace ise="" tubing="" unit=""> completed.</replace>	
20013	<replace ise="" pump="" tubing=""> completed.</replace>	
20014	<replace %s="" electrode="" ise=""> completed.</replace>	
20015	<initialize ise="" reference="" sensor=""> completed.</initialize>	
20016	<clean automatically="" ise="" tower=""> completed.</clean>	
20017	<initialize ise="" unit=""> completed.</initialize>	
20019	<electrode service=""> completed.</electrode>	
20020	<activate electrodes=""> completed.</activate>	
20021	<prime and="" calibrator="" ise="" reference=""> completed.</prime>	
20022	<condition ise="" tubing=""> completed.</condition>	
20023	An unknown action %d was completed successfully.	
20024	Installation completed. The intervals of the maintenance actions were reset successfully (except daily maintenance actions).	
20025	<abs. air="" calibration="" water=""> completed.</abs.>	
20026	The external water reservoir was refilled.	
20027	The external waste container was emptied.	
Table D-1	List of alarm messages (Continued)	

#### **Roche Diagnostics**

#### 9 Messages and alarms

List of alarm messages

ID	Message	Comment
20028	<daily actions="" prepare=""> completed.</daily>	
20029	Database cleanup completed.	
20030	<initialize degaser="" fluid="" sensor=""> completed.</initialize>	
Table D-1	List of alarm messages (Continued)	

# **Result flags**

In this chapter, you will find a list of the flags generated by the **Cedex Bio** instrument, their associated error messages, and the possible user actions.

In this chapter	Chapter	10
About flags		D-25
Safety	••••••	D-27
List of flags		D-28
>		D-28
V		D-28
Ag Excess		D-28
? Cal		D-29
Cal Error		D-29
Calc Error		D-29
Curv Dir	••••••	D-30
Ep Unstab	••••••	D-30
High Abs	••••••	D-31
High Act		D-31
< Kin Read	••••••	D-32
Low Act	••••••	D-32
Non Linear		D-33
Non Mono		D-33
Out of Rng		D-34
? QC		D-34
R 1(2.5s)	••••••	D-34
R 1(3s)		D-35
>R 2(2s)		D-35
<r 2(2s)<="" td=""><td></td><td> D-35</td></r>		D-35
>Reag Rng		D-36
<reag rng<="" td=""><td></td><td> D-36</td></reag>		D-36
> Repl Dev		D–36
> RR		D-37
< RR		D-37

**Roche Diagnostics** 

Table of contents

> Std Dev	D-37
> Test Rng	D-38
< Test Rng	D-38

## **About flags**

Flags are automatically generated with results if during processing certain technical checks were not passed or if the result exceeds or does not reach predefined limits.

Measurements that did not generate flags can be considered technically correct.

Flags are displayed and printed with the results.

#### To display flag information


1 Display the result.

Do one of the following:

lf	Do this
You want to look at sample results:	Choose Workplace > Result Review
You want to look at calibration results:	Choose Workplace > Calibrations
You want to look at QC results:	Choose Workplace > QC Status

A screen is displayed that lists the results. If flags were generated, the flag with the highest priority is displayed. For ratio results, the flag with the highest priority of all constituent test results is displayed.

- **2** Select the flagged result.
- **3** Press  $\nearrow$  to display all flags that were generated for this test (A). They are sorted according to their priority.

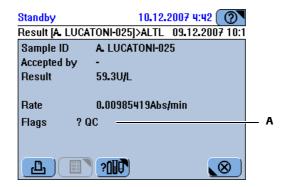

**4** Look up the flag in section *List of flags* on page D-28 for detailed information on the flag.

#### ► To display flag information with the sample still on board

1 Choose **Overview** > tube button.

A screen is displayed that shows the result, together with the flag of the highest priority.

To view all flags that were possibly generated see *To display flag information* on page D-25.

About flags

Flags and error messages	Some flags trigger error messages if they appear in consecutive measurements. Each flag has its own counter. If one measurement does not generate the flag, the counter is reset to zero.
Flag priority	If the conditions in the <b>Cedex Bio</b> instrument are such that multiple flags were generated for a single measurement, only the flag with the highest priority is displayed in the results lists.
Flags and user actions	With each flag description contained in this chapter, the recommended user actions are given.
	• See <i>List of flags</i> on page D-28.

## Safety

Before you start troubleshooting, it is essential that you both read and understand the safety information listed below.

Read carefully all safety notices given in instructions and make sure you understand them.



#### Injury through working solutions

Direct contact with cleaning solutions or other working solutions may cause personal injury. When handling such solutions, exercise the precautions required for handling them, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics cleaning solutions.

#### Infection by biohazardous materials

Contact with samples containing material of human origin may result in infection. All materials and mechanical components associated with samples of human origin are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

## Personal injury or damage to the analyzer due to contact with instrument mechanism

Do not touch moving parts during instrument operation.

#### Incorrect results or damage to the analyzer due to wrong operation

Operators are required to have a profound knowledge of relevant guidelines and norms as well as the information and procedures contained in the Operator's Manual.

- Do not carry out operation and maintenance unless you have been trained by Roche Diagnostics.
- Start all maintenance actions on the screen. Do not perform maintenance actions without the assistance of the user interface.
- Carefully follow the procedures specified in the Operator's Manual for the operation and maintenance of the system.
- Leave maintenance that is not described in the Operator's Manual to trained service representatives.
- Follow standard laboratory practices, especially when working with biohazard material.



Where a flag description gives a list of recommended actions, work through each step in turn until you have solved the problem. Proceed to the next step only after you have completed the previous one.

	This section lists the general flags and those that concern absorbance measurements only, and it provides information on each of the flags.
	• For information on ISE-specific flags, see <i>List of ISE flags</i> on page E-71.
Meaning	Result generated by re-performing the test with the same running parameters (repeated).
Message ID	120
Priority	32
Possible cause	The user initiated the repeat.
Recommended actions	No action required.

V

>

Meaning	A rerun result with diluted sample.
Message ID	121
Priority	33
Possible cause	The user initiated the rerun.
Recommended actions	No action required.

## Ag Excess

Meaning	Antigen excess.	
	The sample contains an excess antigen and a valid result cannot be calculated.	
Message ID	6	
Priority	19	
Recommended actions	Rerun the test with postdilution.	

## ? Cal

Meaning	A result where its calibration is flagged.
Message ID	110
Priority	29
Possible cause	The calibration was in question and has been flagged, however the calibration could be used to calculate the result (compare with Cal Error where the calibration did not provide a usable result).
Recommended actions	Check and redo calibration if needed.

### **Cal Error**

Meaning	No valid calibration data available.
Message ID	43
Priority	24
Possible cause	Caused by an alarm that occurred during calibration because the calibration could not provide a usable result.
Recommended actions	Check the flags of the calibration and proceed to deal with the calibration first.

## **Calc Error**

Meaning	Calculation error.	
	General calculation error.	
	• Slope or nonlinear standard curve cannot be calculated due to a calibration error.	
Message ID	16	
Priority	13	
Possible cause	Calibrator outdated or deteriorated.	
	Misplacement of the calibrator tubes.	
Recommended actions	Repeat calibration with fresh calibrators.	
	• If the calibrators were incorrectly positioned, replace the calibrators and repeat the calibration.	

#### 10 Result flags

List of flags

## **Curv Dir**

Meaning	Curve direction.
	The direction of the calibration curve is incorrect or not as expected.
Message ID	14
Priority	10
Possible cause	Misplacement of calibrator tubes
Recommended actions	<ol> <li>Check the details of the calibration.</li> <li>If the calibrators were incorrectly positioned, replace the calibrators correctly and rerun the calibration.</li> </ol>

## Ep Unstab

Meaning	End point unstable.
	The absorbance values near the endpoint show unacceptable fluctuation.
Message ID	9
Priority	8
Recommended actions	CAL
	1. Repeat with fresh calibrators if only calibration is flagged.
	2. If the flag is on control, sample, or on calibration repeat results, replace the reagent set.
	CNTL
	1. If only one control is flagged, repeat with fresh control.
	2. If more than one control and sample result is flagged, replace the reagent set.
	TEST
	1. Rerun diluted.
	2. If only one sample result is flagged, repeat with fresh sample.
	3. If more than one sample result is flagged or a repeat result, replace the reagent set.
	4. If a test result is flagged, repeat with dilution.

## **High Abs**

Meaning	Excessive absorbance.
	The absorbance value to be used for calculation and checks is $> 2.0$ Abs.
Message ID	5
Priority	1
Recommended actions	CAL
	Repeat with fresh calibrator.
	CNTL
	• If a single control is flagged, rerun with a fresh control.
	• If more than one control is flagged, perform a calibration.
	TEST
	• If a single sample result is flagged, rerun with dilution.
	• If the problem persists, replace the reagent set.
High Act	
Meaning	High activity.
	• The absorbance change during measurement is above the accepted limit, i.e. the reaction has taken off too fast.
Message ID	56

Priority 9

Recommended actions CAL

- 1. Check if the correct calibrators are placed on the instrument.
- 2. Repeat with fresh calibrators.

CNTL

1. Check control. Repeat with fresh control.

#### TEST

1. Rerun with dilution.

## < Kin Read

Meaning	Insufficient kinetic reading points.
Message ID	11
Priority	11
Possible cause	Not enough kinetic readings.
	• Insufficient absorbance readings were found in the linear part of the reaction to be able to calculate a rate.
	• Not enough readings in kinetic range for calc mode Kinetic.
Recommended actions	CAL
	1. Check reaction plot.
	2. Repeat with fresh calibrator.
	CNTL
	1. Check reaction plot.
	2. Repeat with fresh control.
	TEST
	1. Check reaction plot.
	2. Rerun with dilution or concentration if necessary.

#### Low Act

Meaning	Low activity.
	This check is used to identify zero pipetting for samples and reagents.
	• The absorbance change during measurement is below the accepted limit, i.e. the reaction has taken off too slowly. Possibility of substrate exhaustion.
	• The absorbance change during the measurement is lower than the defined limit.
Message ID	56
Priority	9
Recommended actions	CAL
	1. Check if the correct calibrators are placed on the calibrator rack.
	2. Repeat with fresh calibrators.
	CNTL
	1. Check the control. Repeat with fresh control.
	TEST
	1. Rerun with dilution.

## **Non Linear**

Meaning	Abnormal linearity.
Message ID	11
Priority	11
Possible cause	• The slope changes between the first and last parts of the reaction curve.
Recommended actions	CAL
	1. Check reaction plot.
	2. Repeat with fresh calibrator.
	CNTL
	1. Check reaction plot.
	2. Repeat with fresh control.
	TEST
	1. Check reaction plot.
	2. Rerun with dilution or concentration if necessary.

#### Non Mono

Meaning	Curve not monotonic.
	No calibration curve could be calculated because the rates of the calibrators were non-monotonic.
Message ID	80
Priority	14
Possible cause	Misplacement of a series of calibrators, automatic dilution series failed.
Recommended actions	1. If the calibrators were incorrectly positioned, re-place the calibrators correctly and rerun the calibration.
	2. Check the fluid system
	3. Repeat the calibration

## Out of Rng

Meaning	Curve out of range.
	The calibration curve is outside the acceptable/ programmed range.
Message ID	84
Priority	16
Possible cause	Deteriorated or outdated calibrator or reagent.
Recommended actions	<ol> <li>Repeat with fresh calibrators.</li> <li>If the calibrators were incorrectly positioned, re-place the calibrators correctly and repeat the calibration.</li> <li>If flag reoccurs, repeat with fresh reagent.</li> </ol>

## ? QC

Meaning	A result where the quality control measurement is flagged.
Message ID	111
Priority	30
Possible cause	The last QC measurement before this measurement was flagged. The dependency is derived from the time sequence.
Recommended actions	Check QC measurement.

## R 1(2.5s)

Meaning	One control value is above 2.5 standard deviation or below -2.5 standard deviation.
Message ID	36
Priority	28
Possible cause	<ul><li>The reagent has deteriorated (linearity of working curve degraded).</li><li>One control is concentrated or has deteriorated.</li></ul>
Recommended actions	<ol> <li>Repeat with fresh control.</li> <li>Check that the correct control material has been used.</li> <li>If the flag reappears, check the calibration and reagent status. Repeat with fresh reagent.</li> </ol>

## R 1(3s)

Meaning	One control value is above 3 standard deviation or below -3 standard deviation.
Message ID	35
Priority	27
Possible cause	<ul><li>Improper control is set.</li><li>Proper control values (mean value, standard deviation) are not specified.</li></ul>
Recommended actions	<ol> <li>Repeat with fresh control.</li> <li>Check that the correct control material has been used.</li> <li>If the flag reappears, check the calibration and reagent status. Repeat with fresh reagent.</li> </ol>

## >R 2(2s)

Meaning	Two sequential control measurements are above the 2 standard deviation.
Message ID	29
Priority	26
Possible cause	<ul><li>The controls are not properly prepared.</li><li>The controls are not properly positioned on the instrument.</li></ul>
Recommended actions	<ol> <li>Repeat with fresh control.</li> <li>Check that the correct control material has been used.</li> <li>If the flag reappears, check the calibration and reagent status. Repeat with fresh reagent.</li> </ol>

## <R 2(2s)

Meaning	Two sequential control measurements are below the -2 standard deviation.		
Message ID	29		
Priority	26		
Possible cause	<ul><li>The controls are not properly prepared.</li><li>The controls are not properly positioned on the instrument.</li></ul>		
Recommended actions	<ol> <li>Repeat with fresh control.</li> <li>Check that the correct control material has been used.</li> <li>If the flag reappears, check the calibration and reagent status. Repeat with fresh reagent.</li> </ol>		

## >Reag Rng

Meaning	Absorbance above defined reagent range.	
	• During calibration, the absorbance value of the lowest calibrator is above the upper defined range.	
Message ID	12	
Priority	7	
Possible cause	Deteriorated or outdated calibrator or reagent.	
Recommended actions	<ol> <li>Check the calibration and reagent expiry dates.</li> <li>Repeat with fresh lowest calibrator.</li> <li>If the flag reappears, replace reagent set (a reagent may be contaminated).</li> </ol>	

## <Reag Rng

Meaning	Absorbance below defined reagent range.
	• During calibration, the absorbance value of the lowest calibrator is below the lower defined range.
Message ID	12
Priority	7
Possible cause	Deteriorated or outdated calibrator or reagent.
Recommended actions	<ol> <li>Check the calibration and reagent expiry dates.</li> <li>Repeat with fresh lowest calibrator.</li> <li>If the flag reappears, replace reagent set (a reagent may be contaminated).</li> </ol>

## > Repl Dev

Meaning	Greater than replicate deviation.		
	Deviation between replicates exceed programmed limit. One or more replicates are erroneous.		
Message ID	13		
Priority	12		
Possible cause	Problems in air/water system, air bubbles, clogged probe.		
Recommended actions	<ol> <li>Check the calibrators for air bubbles and repeat the calibration.</li> <li>If the flag reappears, repeat with fresh calibrators.</li> <li>Check the probe.</li> <li>Check the fluid system.</li> </ol>		

## > RR

Meaning	Above reference range.
	The calculated concentration is greater than the upper limit of the expected value range.
	nnnnn is the actual value that was checked against.
Message ID	40
Priority	22
Recommended actions	No action required.

#### < RR

Meaning	Below reference range.
	The calculated concentration is less than the lower limit of the expected value range.
	nnnnn is the actual value that was checked against.
Message ID	41
Priority	23
Recommended actions	No action required.

#### > Std Dev

Meaning	Greater than standard deviation.		
	A point in the calibration fell outside the defined limits.		
Message ID	17		
Priority	17		
Recommended actions	<ol> <li>Check calibration curve.</li> <li>Repeat with fresh calibrators.</li> <li>If the calibrators were incorrectly positioned, replace the calibrators correctly.</li> </ol>		

#### 10 Result flags

List of flags

## > Test Rng

Meaning	PANIC value over (upper) Technical Limit.	
	The result is higher than the upper limit for the test.	
Message ID	26	
Priority	20	

## < Test Rng

Meaning	PANIC value below (lower) Technical Limit.		
	The result is lower than the lowest limit for the test.		
Message ID	27		
Priority	21		

# Troubleshooting

# Dealing with exceptional situations

In this chapter, you will find information on how to deal with selected exceptional situations.

In this chapter	Chapter	11
Introduction		D–41
Dealing with exceptional situations		D–42
Reacting to messages		D–44
Detailed procedures		D–46
Checking the printer status		D–46
Clearing a paper jam		D–46
Changing the mains fuses		
Changing the low voltage fuses		D–49
Realigning the reagent disk		D–50
Exporting diagnostics screens to USB		D–51

#### **11 Troubleshooting**

Table of contents

Introduction

	Effective troubleshooting requires a good understanding of all basic operating procedures for the <b>Cedex Bio</b> instrument and the system software.
Preventive maintenance	You can avoid most unforeseen situations if you observe the recommended procedures at all times and if the operating environment is effectively controlled by performing all maintenance actions when they become due in the manner described in the Operator's Manual.
	• See <i>Maintenance actions</i> on page C-8.
Information about the status of the system	The system performs numerous checks. When a certain event has occurred or when an irregularity is detected, a message is generated. Messages are displayed in two ways:
	• Immediate feedback on user actions is displayed in a <i>pop-up message screen</i> .
	• See Message screen on page D-5.
	• Information concerning a problem that occurred during operation is reported as <i>alarm messages</i> in the <i>alarm monitor</i> .
	• See <i>Alarm monitor</i> on page D-6.

Dealing with exceptional situations

## **Dealing with exceptional situations**

The following table list possible exceptional situations, provides information on the possible causes, and suggests ways of remedying the situation. (Subsequent sections contain detailed procedures for selected remedies.)

Situation	Affected items	Possible causes	Comments
The fluid connectors need to be disconnected from the instrument.	External fluid connectors.	<ul><li>Moving the instrument to another location.</li><li>Removing blockages.</li></ul>	<ul> <li>To get easier access to the release clamps on the connectors, remove the side panel before removing the connectors.</li> <li>See <i>To disconnect a tube from the instrument</i> on page B-130.</li> </ul>
A service representative asks you to perform diag- nostics tasks.	Diagnostics features.	These features are usually used at the request of a service repre- sentative.	<ul> <li>Follow the instructions given by the service representative.</li> <li>See <i>Exporting diagnostics screens to USB</i> on page D-51.</li> </ul>
The system does not work or does not switch on and the power switch light is off.	Instrument.	No electricity.	<ol> <li>Check that the instrument is connected to the mains power supply.</li> <li>If the instrument is properly connected replace the mains fuses.</li> <li>See <i>Changing the mains fuses</i> on page D-47.</li> </ol>
No printing.	Printer.	Printer is out of paper.	<ul><li>Load printer paper.</li><li><i>Refilling printer paper</i> on page B-122.</li></ul>
		Printer jam.	To clear the paper jam on page D-46.
Reagent disk cannot be inserted in the reagent cooler.	Reagent disk.	The bottle storage assembly in the reagent disk was moved while the latter was outside the instrument.	The reagent disk is designed to handle reagents while it is loaded on the instrument. The cover is equipped with a locking mechanism.
			Always remove and load reagents while the reagent disk is on the instrument and by using the software sup- ported procedures.
			1. Align the bottle storage assembly in the reagent disk.
			$\odot$ See Realigning the reagent disk on page D-50.
			2. Insert the reagent disk in the cooler.

Figure D-4 Troubleshooting overview

Dealing with exceptional situations

Situation	Affected	Possible causes	Comments
	items		
The system does not go	System status	There are not enough free cuvettes on the instrument.	Load empty cuvette segments.
to <b>Operating</b> status.			• See <i>Preparing cuvettes</i> on page B-28.
		Air/water calibration is required.	Perform Abs. Air/Water Calibration maintenance
			action.
			• See Perform an air/water calibration on page C-28.
		In the corresponding calibrator lot definitions of at least one of the tests you want to perform, the value for <b>Last Is Water</b> is <b>Off</b> .	The <b>Last Is Water</b> definition defines what is used as the zero calibrator. It is either system water ( <b>On</b> ) or a standard, usually water ( <b>Off</b> ), which would be placed in a tube like any other calibrator.
			With some tests, calibration does not work with special zero calibrators, they need the <b>Last Is Water</b> definition on.
			Possible remedies:
			<ol> <li>Delete the current calibration order and redefine it.</li> <li>If the Last Is Water definition is Off, make sure to place a special zero calibrator on the instrument when calibrating. (This is usually water.)</li> </ol>
			If this does not help change the <b>Last Is Water</b> value to <b>On</b> . (Calibration for this test does not work with a special zero calibrator.)
			<ol> <li>Choose Workplace &gt; Lot Data &gt; Calibrators.</li> <li>Select the entry for the test and expand the entry.</li> <li>Select the Last Is Water entry and press Ø.</li> <li>Press On.</li> <li>Press Ø.</li> <li>Delete the current calibration order and redefine it.</li> </ol>
		A hardware error had occurred.	Some hardware conditions prevent the system from
			performing tests.
			1. Press 🛈 to check the alarm messages.
			2. Follow the advise given in the messages.
			3. Choose $\bigcirc$ > <b>Restart</b> to restart the system.

Figure D-4 Troubleshooting overview (Continued)

Reacting to messages

## **Reacting to messages**

A **Cedex Bio** error message starts with a message ID, which consists of a number followed by the date and time, for example:

7009.19.19.32.1402 [25.08.2006 8:53]

Providing information to service<br/>representativesWhenever you are asked to provide message information to a service representative,<br/>provide the error codes contained in the explanatory text, not the message ID of the<br/>first line. If there are no error codes in the text, supply the complete message text.

Also keep in mind that an event may trigger several messages, therefore it is important to provide the information for all messages that were generated within about one minute.

Basically there are three kinds of error message texts:

- Messages that contain text only.
- Messages that contain an error code of the format a.bbb.xxx, for example [3.000.121].
- Messages that contain a nine digit error code, for example [104000551].

The following table lists selected messages and provides more detailed information on how to react to them.

<u>`</u>ل

In the "Message or error code" column of the following table, the most relevant information is the number, the texts may be different on the actual instrument.

Message or error code	Affected items	Possible causes	Comments
1.xxx.81 1.xxx.82	Instrument.	Fuse F1, F2, or F4 is defective.	Replace the low voltage fuse that is mentioned in the alarm message.
1.xxx.84			See Changing the low voltage fuses on page D-49.
[109000572] code raised by IC software caused by an unexpected handling or sys- tem situation.	Instrument.	The probe is bent or the trans- fer head is poorly adjusted.	<ol> <li>Check that the probe is mounted properly and replace it if necessary.</li> <li>See <i>Replacing the probe</i> on page B-125.</li> <li>If the problem persists, contact your service representative.</li> </ol>
No connection to Instrument Control.	Instrument.	Fuse <b>F3</b> is defective.	<ul> <li>Check the LED of fuse F3. Replace the fuse if necessary. (If the fuse has blown, the LED is off.)</li> <li>See <i>Changing the low voltage fuses</i> on page D-49.</li> </ul>
-		The LAN cable is not or not properly connected.	<ol> <li>Check that the LAN cable is properly connected:</li> <li>Switch off the instrument.</li> <li>If the cable was not connected, connect it. If the cable was connected, remove it at both ends and reconnect it.</li> <li>Start the instrument.</li> </ol>
		Electronic problem	If the above measures are not successful, contact your service representative.



Troubleshooting with the help of messages

#### **11 Troubleshooting**

Reacting to messages

Message or error code	Affected items	Possible causes	Comments
[1.xxx.73] Level detection error (conductivity).	Level detection.	<ul> <li>Poor water quality.</li> <li>Contamination with cleaner (valve is leaking)</li> </ul>	<ol> <li>Make sure you use the correct water quality.</li> <li>See <i>Technical specifications</i> on page A-65.</li> <li>Perform the <b>Prime Fluid System</b> maintenance action.</li> <li>See <i>Prime the fluid system</i> on page C-11.</li> <li>If the problem persists, contact your service representative.</li> </ol>
[4.xxx.41] through [4.xxx.44] Various texts related to photometer Gc-0166807.	Photometer.	<ul><li>Lamp defective.</li><li>Photometer lens soiled.</li></ul>	<ul><li>Do not attempt to clean the photometer lenses.</li><li>1. Replace the photometer lamp.</li><li>2. If the problem persists, contact your service representative.</li></ul>

Figure D-5 Troubleshooting with the help of messages (Continued)

Detailed procedures

## **Detailed procedures**

#### **Checking the printer status**

#### To check the printer status

- 1 On the **Overview** tab, press the **System status** button.
  - The printer button  $\square$  is red.
- **2** Scroll to display the Printer entry.


#### **Clearing a paper jam**

#### To clear the paper jam

**1** Choose 🔘 > 🕒 **Stop Printing**, if required.

Press down the release button (A) firmly before you pull the panel. The panel should

**2** Open the printer panel.

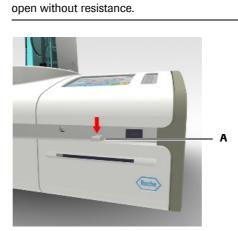

- **3** Lift the printer paper roll from its holder.
- 4 Remove all loose paper. (Cut or tear it off).
- **5** Place the roll back on the holder.

Make sure the paper unrolls at the top and towards you.



- **6** Insert the paper in the slot in the printer panel and pull some through.

**7** Close the printer panel.

The system feeds some paper and then automatically resumes printing.

- 8 Choose **Overview** > 🛃.
- **9** The printer button should now be green, and the status description should be **OK**.

#### Changing the mains fuses

#### To change the mains fuses

- 1 Press 🞯
- 2 The Stop screen is displayed.
- **3** Press ① Shutdown to shut down the system.

Wait until the screen is black.

- 4 Press O on the main power toggle switch to switch off the instrument.
- **5** Disconnect the mains cable from the instrument.




Detailed procedures


6 Using a screwdriver disengage the clips on both sides of the fuse box cover.



**7** Remove the holder.

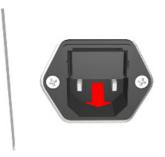

**8** Remove the two fuses from the holder.

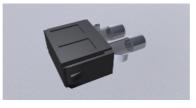

**9** Replace both fuses.



**Damage to the analyzer due to use of wrong fuses** Always replace fuses with new ones of the same type and specifications. See *Technical specifications* on page A-65.

**10** Insert the fuse holder in the fuse box.

Press it in firmly until the clips engage.

- **11** Connect the mains cable to the instrument.
- **12** The instrument switch should be illuminated.

#### Changing the low voltage fuses

#### To change a low voltage fuse

- 1 Shut down the system and switch off the instrument.
- **2** Using a screwdriver size 2 or 3, remove the fuse indicated in the alarm message.

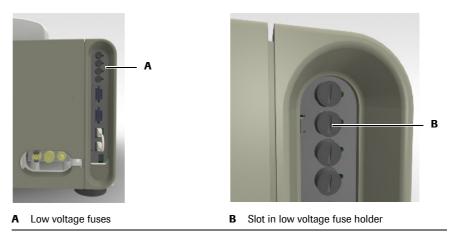


Figure D-6

- i
- Insert the screwdriver in the slot and turn anticlockwise until the holder disengages.
- Remove the holder with the fuse.



- **3** Remove the fuse from its holder.
- 4 Insert the new fuse in the holder.

NOTICE

#### Damage to the analyzer due to use of wrong fuses

Always replace fuses with new ones of the same type and specifications. See *Technical specifications* on page A-65.

- **5** Insert the holder in the socket.
- **6** Using the screwdriver, lock the holder in position by pressing it in and turning the screwdriver clockwise until the slot is vertical.
- **7** Switch on the instrument.

Detailed procedures

#### Realigning the reagent disk




#### To realign the reagent disk


1 Insert your fingers in the holes in the reagent disk cover an lift the reagent disk.

**Spillage through tipping reagent disk** When handling the reagent disk, make sure not to tilt it.

- **2** Insert a pencil in the opening in the center of the underside of the reagent disk.
- **3** Turn the reagent disk until the arrows of the reagent disk and the cover face each other.

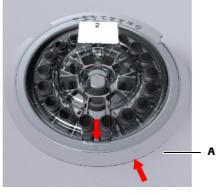

A Incorrect position: The two arrows do not face each other.

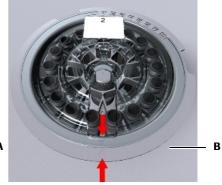

**B** Correct position: The two arrows face each other.

Figure D-7

#### **Exporting diagnostics screens to USB**

When performing diagnostics actions you can export the content of the screen to the USB stick.

*Prerequisites* You need Lab Administrator or Administrator user rights to perform diagnostics actions, and the system must be in Standby or Stopped status.

#### ► To export the content of the screen

- **1** Press **()** to start the diagnostics action.
- **2** Press **(**
- 3 Press Send to File.

As screen is displayed, asking you to insert the USB stick.

- 4 Insert the USB stick.
- **5** Press 🔘.
- 6 Select the directory.
- **7** Press 🔘 to confirm the selection.

The data are copied to the USB stick. (File name format: prt\_yyyymmddhhmmss.txt.)

#### **11 Troubleshooting**

Detailed procedures

# ise E

12	ISE description	<i>E-3</i>
13	ISE operation	<i>E-15</i>
14	ISE maintenance	E-39
15	ISE troubleshooting	E-67

# **ISE description**

Overview of the ISE unit

In this chapter, you will find a general overview of the **Cedex Bio** ISE unit. You also will find a description of the main components and their principle operation.

In this chapter	Chapter	12
Overview		E–5
Abbreviations	••••••	E-5
Measuring modes		E–6
Principles of operation	••••••	E–7
ISE solutions	••••••	E–8
Hardware	••••••	E–9
Panels		E–9
ISE LED	••••••	E–10
Main components	••••••	E–10
Fluid containers	••••••	E–11
Basic operation		E–12
Samples	••••••	E–12
Calibration	•••••••••••••••••••••••	E-12
ISE Standby		E-12
Maintenance actions		E-13
Technical specifications		E–13

#### 12 ISE description

Table of contents

Overview

## **Overview**

The Ion-Selective Electrode (ISE) module is a measurement system for electrolytes that runs independently of the main instrument. The ISE unit uses the same samples and sample tubes that are used for photometric measurements, and the samples are transferred using the same probe. The ISE unit uses specific solutions, some of which are kept in bottles on the ISE unit itself, the others are transferred to sample tubes and placed on the sample area.

## **Abbreviations**

Abbreviation	Definition
F	Solution 1 factor
K-I	Potassium indirect
mV	Measured voltage in millivolt
NA-I	Sodium indirect
S	Slope in mV/dec
Std 1/1	First measurement of ISE Solution 1
Std 1/2	Second measurement of ISE Solution 1
Std 2	Measurement of ISE Solution 2

The following ISE specific abbreviations are used:

Overview

## **Measuring modes**

	The ISE unit makes, in aqueous solutions, quantitative determinations of the following electrolytes:
	• Sodium (Na <sup>+</sup> )
	• Potassium (K <sup>+</sup> )
	Measurements are done using <i>indirect</i> mode. The samples, controls, and standard solutions are diluted with system water 1:6 (1+5). The dilution and mixing are performed automatically in the ISE tower.
Pipetting volumes	Sample: 15 µL
	Diluent (H2O): 75 µL
ISE measurements	ISE measurements and photometric measurements can be carried out at the same time and are independent of each other. (ISE measurements do not use cuvettes, the reagent rotor, or the photometer.) All requested ISE measurements on a specific sample, control, or standard solution are performed in parallel, that is, measurements are made at each electrode at the same time.

Overview

## **Principles of operation**

The ISE unit uses flow-through ion-selective electrodes and a reference electrode with an open liquid junction. Each electrode has a membrane that is sensitive to a particular type of ion.

*Measuring process* 1. The ISE maintenance and standard solutions (ISE Deproteinizer, ISE Etcher, Activator, and ISE Solution 1 and 2) are pipetted from sample tubes on the sample area to the ISE tower as required.

- 2. The sample is pipetted from the sample tube (located on the sample area) into the ISE tower. The sample is diluted with system water. Mixing is performed with four air jets arranged in a circle. These jets blow air into the tower to produce a homogenous mixture.
- 3. The sample is divided into segments with the aid of a special arrangement of valves. The first (shorter) segments are used for cleaning, these are followed by a longer segment, on which the measurements are made.
- 4. The sample is passed to the ion-selective electrodes by the action of the peristaltic pump.

The exact positioning of segments is ensured by the ISE sample sensor.

- 5. In the meantime, the ISE tower is washed with distilled water and dried.
- 6. ISE Reference Solution is passed through the ISE Reference Electrode and into the measuring channel downstream of the electrodes. The ISE Reference Solution completes the electrical circuits for each electrode so that measurements can be made. While the measurements are made, the sample and ISE Reference Solution are stationary.
- 7. A one-point calibration is performed after each sample measurement using the ISE Calibrator Indirect, which is located on ISE unit.
- 8. The electrolyte concentration of the sample is calculated.

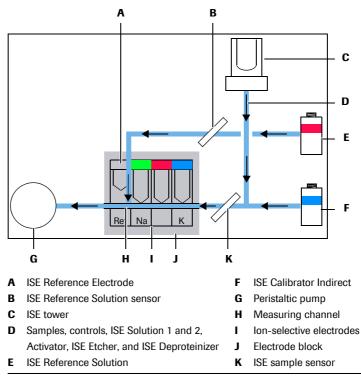


Figure E-1 Main parts of the ISE measuring system

## **ISE** solutions

The ISE unit requires several specific solutions for performing the analyses and maintaining the system. They are listed in the following table.

Solution summary	Solution	Location	Use
	ISE Solution 1	Sample area	Used in the two-point calibration of sodium and potas- sium.
	ISE Solution 2	Sample area	Used in the two-point calibration of sodium and potas- sium.
	ISE Calibrator Indirect	ISE unit	Used in the one-point calibration after each measure- ment. It is also used during ISE initialization and standby, and for maintenance purposes.
	ISE Reference Solution	ISE unit	Used in all ISE measurements of sodium, potassium, and chloride.
			Also used during ISE standby.
	ISE Etcher	Sample area	Used as a cleaning solution for cleaning the ISE Sodium Electrode during ISE maintenance.
	ISE Deproteinizer	Sample area	Used as a cleaning solution for cleaning the probe, ion- selective electrodes, ISE tower, and tubing during ISE maintenance.
	Activator	Sample area	Activates the electrodes, tubing, and the ISE tower dur- ing ISE maintenance. It is also used for the initialization of the ISE unit and for activating the probe.
			Roche recommend using their Activator on the <b>Cedex Bio</b> instrument. (For details see the package insert of the Activator.)
	Water	External water container	Used for diluting the samples, controls, and standard solutions, and for cleaning purposes. It is also used during ISE initialization.
	Sample	Sample area	Sample tubes are placed on the sample area of the main instrument and pipetted by the main instrument probe.
	Waste	Internal waste tank	The waste is first pumped into the internal waste tank of the main instrument and then to the external waste container.

Fluid stability

question.



Incorrect results due to changes in fluids

The chemical composition of ISE fluids changes over time. The assigned on-board stability is the interval within which the quality of the fluid remains within the prescribed tolerances. Using fluids whose interval has expired may lead to incorrect results.

For information on the stability of fluids see the package insert of the fluid in


Always exchange the fluids when an interval has expired. Follow the instructions given on the screen.

Hardware

## Hardware

**Panels** 

The ISE cover has two joints. The front cover provides access to the parts you may need to handle during daily operation. Opening the back cover as well provides access to the peristaltic pump and its tubing.

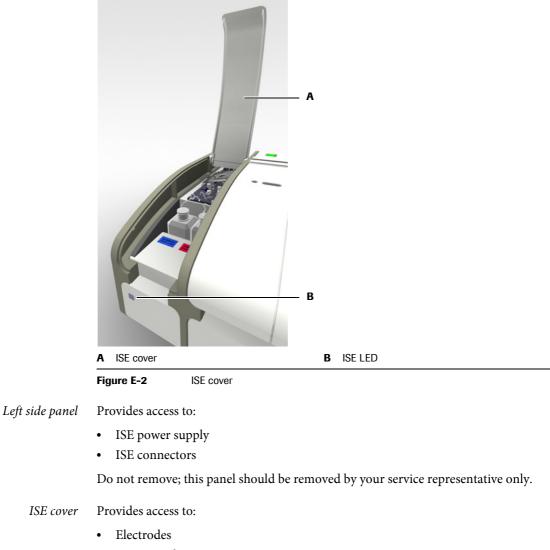

- ISE unit tubing
- ISE fluid bottles

Do not open this cover when the system is in **Operating** status.

*Left service flap* Provides access to:

• ISE tower


Remove this flap as instructed during maintenance actions, or shut down the system and switch off the instrument before you remove it.

## **ISE LED**

### The ISE LED is color coded.

	Power off (e.g. power supply disconnected.)
	The ISE unit is OK and ready for use.
Blinking	The fluid level in the ISE Reference or ISE Calibrator Indirect bottle is low.
	Error, processing stopped.

## **Main components**

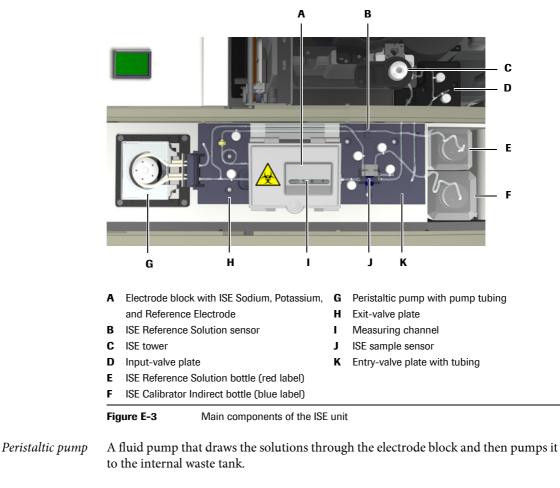

*Input-valve plate* A set of valves to create cleaning segments and to control the flow of waste fluids.

- *Entry-valve plate* A set of valves to control and monitor the flow of fluids.
  - *Exit-valve plate* A set of valves to control the aspiration action of the peristaltic pump.
  - *Electrode block* A block that contains the electrodes.

ISE tower	A hollow, transparent cylinder with multiple inlets and outlets for air and solutions.
	In the ISE tower, samples and diluents are mixed using an air stream.
Tubing	Tubing is used for transporting the fluids. It connects the ISE unit with the water and waste system, and also with the ISE Reference Solution and ISE Calibrator indirect/urine bottles, which are placed on the ISE unit.
Sensors	The ISE sample sensor detects the presence of liquid (sample, control, standard solution, or ISE Calibrator indirect/urine) or air and enables the correct placing of segments for measurement and cleaning.
	The ISE Reference Solution sensor detects the presence of ISE Reference Solution.
ISE fluid bottles	Two bottles are placed on the ISE unit:
	• ISE Calibrator indirect/urine bottle (blue label)
	• ISE Reference Solution bottle (red label)

Level monitoring is performed on the basis of weight, the bottles are placed on scales.

## **Fluid containers**

The following table shows which container is used for which fluid:

Fluid	Container(s)	Position
Samples	Tube	Sample area
Controls	Tube	Sample area
ISE Solution 1 and 2	Tube	Sample area
Cleaner	External cleaner bottle	External fluid rack
ISE Deproteinizer	Tube	Sample area
Activator	Tube	Sample area
ISE Etcher	Tube	Sample area
ISE Reference Solution	Bottle	ISE unit
ISE Calibrator indirect/urine	Bottle	ISE unit
Water	External water container	External fluid rack
Waste	External waste container	External fluid rack

The term tube includes all kinds of tubes, as long as their dimensions lay within prescribed limits. It also includes secondary tubes (cups). See *Tubes* on page A-53.

<u>ک</u>

**Basic operation** 

**Cedex Bio System** 

•	
	Operation of the ISE unit is integrated in the operation of the main instrument.
	The following sections contain some ISE specific information on operation-related issues.
	For general information on operating the main instrument, see the corresponding sections in Chapter 2 Introduction to the instrument and Chapter 5 Daily operation. For specific ISE tasks, see Chapter 13 ISE operation.
Samples	
	The sample tubes are placed on the sample area of the main instrument.
Calibration	
	The ISE unit requires frequent calibration to ensure the accuracy of the test results obtained.
Main calibration	The electrodes are calibrated with a two-point calibration using ISE Solution 1 and 2.
	The main calibration is typically performed during the Prepare phase.
One-point calibration	The electrodes are calibrated after each ISE measurement using the on-board ISE Calibrator indirect/urine. This calibration is an integral part of each ISE measurement and is performed automatically.
ISE Standby	
	If there is no measurement for more than three minutes, the ISE unit switches to <b>Standby</b> status. In this status, ISE Calibrator indirect/urine and ISE Reference Solution are pumped into the measuring channel and moved a short distance at regular intervals. Also, every two hours the ISE tubing is primed with ISE Calibrator indirect/urine and ISE Reference Solution. This is done to prevent the following problems:
	• Flow of ISE Reference Solution backwards into the measuring channel, which can damage the electrodes because of the high ion concentration of the ISE Reference Solution.
	• Crystallization of salts in the tubing, causing blockages.
	The ISE unit has a separate power supply. Therefore, the <b>Standby</b> status is maintained even if the main instrument is switched off.
NOTICE	Damage to electrodes and possible tubing blockage
	<ul> <li>Do not unplug or switch off the ISE power supply. Periodic flow of solutions must be performed at all times.</li> </ul>
	If you intend not to use the ISE unit for more than one week, you should deactivate it. This saves ISE fluids, reduces wear and tear of the tubing, and lowers the likelihood of malfunctions and errors occurring. (See <i>Deactivating the ISE unit</i> on page E-64.)

## **Maintenance actions**

Maintenance of the ISE unit is integrated in the maintenance of the main instrument. The system software guides you through the maintenance procedures.

Maintenance actions are performed after a certain event or after a defined interval has expired.

● For specific ISE maintenance actions, see Chapter 14 *ISE maintenance*.

## **Technical specifications**

• See *Technical specifications* on page A-65.

## 12 ISE description

Technical specifications

**ISE** operation

In this chapter, you will find information on performing the routine tasks that are required for processing tests using the ISE unit.

In this chapter	Chapter	13
Daily operation		E-17
Safety information		E–19
Preparing the system		E-20
Short guide		E-20
Starting the Prepare wizard		E-21
Checking the status of the external fluid containers		E-21
Checking the ISE status		E-21
Performing maintenance actions		E-22
Completing the preparation tasks		E-23
Performing calibrations		E-23
Defining orders		E-23
Monitoring the progress		E-24
Validating sample results		E-25
Main calibration		E-26
Performing QC		E-26
Finishing the shift		E-27
Short guide		E-28
Replacing ISE fluid bottles		E-30
Replacing electrodes		E-32
Cleaning the ISE tower off the instrument		E-36

## 13 ISE operation

Table of contents

Daily operation of the ISE unit is integrated in the daily operation of the main instrument.

See Chapter 5 Daily operation.

The following table gives an overview of the tasks you might need to perform during daily operation.

	Task	Steps		Navigation
			With wizard	As individual steps
1	Starting the system	1. Switch on the system.		
2	Logging on the system			Overview > Logon
3	Preparing the system	Start the <b>Prepare</b> wizard.	Overview > ]	Prepare
		1. Check the external fluid containers.		Overview > 🚱 > 🛍
		2. Check the ISE status.		Overview >
		3. Perform Daily Prepare Actions.		Utilities > Maintenance
4	Defining orders	Start the Orders wizard.	Overview >	Order (or Overview > STAT)
		1. Identify the sample.		n/a
		2. Select the tests.		n/a
		3. Place the sample.		n/a
		4. Start the run.		$\Diamond$
5	Monitoring the progress		n/a	Overview
				$\odot$
6	Validating results	1. View results.	n/a	Workplace > Result Review
		2. Handle flagged results.	n/a	Workplace > Result Review > 🖉
				> Repeat
				> <b>Rerun</b>
		3. Accept results.	n/a	Workplace > Result Review > 🖉 >
				Accept
7	Performing calibrations			
	Performing individual	1. Start the wizard.	Workplace >	> Calibrations > 🕀
	calibrations	2. Select the test.		n/a
		3. Prepare and place the calibrators.		n/a
		4. Start the calibration.		$\bigcirc$
		5. Validate the calibration results.		Workplace > Calibrations >
		6. Remove the calibrators.		
	Performing	1. Start the wizard.	Workplace >	> Calibrations > 🛞
	all due calibrations	2. Select all tests with due calibrations.		n/a
		or	*	
		Select all tests with calibration due within the forecast period.	****	
		3. Prepare and place the calibrators.	_	n/a
		<ol> <li>Frepare and place the calibrators.</li> <li>Start the calibration.</li> </ol>		
		<ol> <li>Start the calibration.</li> <li>Validate the calibration results.</li> </ol>		Workplace > Calibrations > 🖉
		<ol> <li>Valuate the calibration results.</li> <li>Remove the calibrators.</li> </ol>		·····

### **Roche Diagnostics**

## 13 ISE operation

Daily operation

Task	Ste	ps		Navigation
			With wizard	As individual steps
Performing controls				
Performing			Overview > 0	Order > 📲
Default QC	2.	Select a control and place the tube. Repeat until there are no controls left on the screen.		n/a
	3.	Start the QC measurement.		$\bigcirc$
		Validate the QC results.		Workplace > QC Status >
	5.	Remove the controls.		n/a
Performing an individual	1.	Start the wizard.	Workplace >	• QC Status > 🕀
QC measurement	2.	Select a test.		n/a
	3.	Select a control and place the tube. Repeat until there are no controls left on the screen.	-	n/a
	4.	Start the QC measurement.		$\odot$
	5.	Validate the QC results.		Workplace > QC Status >
	6.	Remove the control.		
Performing all	1.	Start the wizard.	Overview > 0	Order > 🕑
due QC measurements	2.	Select a control and place the tube. Repeat until there are no controls left on the screen.	_	n/a
	3.	Start the QC measurement.		$\bigcirc$
	4.	Validate the QC results.		Workplace > QC Status > 🖉
	5.	Remove the controls.		
Finishing the shift	1.	Check for unfinished orders.		Workplace > Orders Choose 🝸 > Not Finished
	2.	Check for non-validated results.		Workplace > Result Review Choose Y > Not Accepted
	3.	Check for non-transmitted results.		Workplace > Result Review
		(If working with a host system only.)		Choose Y > Not Sent to Host
	4.	Start the <b>End Shift</b> wizard.	Overview > 1	
	5.	Perform the daily backup.	-	Utilities > Export > Database
	6.	Export the full results.	-	Utilities > Export > Results
		Clean up the database.	-	Workplace > Orders > Θ
		-		Workplace > Result Review> Θ
				Workplace > QC Status > $\Theta$
				Workplace > QC History > Θ
				Workplace > Calibrations > $\Theta$
	8.	Perform the maintenance actions that are due.	-	Utilities > Maintenance
	9.	Check the external fluid containers.	_	Overview > 🔄 > 🛍
	10.	Log off the system.		<b>Overview</b> > button with your user nam
	11.	Shut down the system and switch off the instrument (if last shift).	n/a	n/a

Table E-3

Overview of the daily operation tasks

The following sections describe the specific ISE operation tasks.

## Safety information



#### Injury through reagents and other working solutions

Direct contact with reagents, cleaning solutions, or other working solutions may cause personal injury.

When handling reagents, exercise the precautions required for handling laboratory reagents, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics reagents and cleaning solutions.

#### Infection by biohazardous materials

Contact with samples containing material of human origin may result in infection. All materials and mechanical components associated with samples of human origin are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

## Personal injury or damage to the analyzer due to contact with instrument mechanism

Do not touch any parts of the instrument other than those specified. During operation and maintenance of the instrument, proceed according to the instructions.

Keep all covers closed, operate them as instructed on the screen.

#### Incorrect results or damage to the analyzer due to wrong operation

Operators are required to have a profound knowledge of relevant guidelines and norms as well as the information and procedures contained in the Operator's Manual.

- Do not carry out operation and maintenance unless you have been trained by Roche Diagnostics.
- Start all maintenance actions on the screen. Do not perform maintenance actions without the assistance of the user interface.
- Carefully follow the procedures specified in the Operator's Manual for the operation and maintenance of the system.
- Leave maintenance that is not described in the Operator's Manual to trained Roche support representative personnel.
- · Follow standard laboratory practices, especially when working with biohazard material.

## Preparing the system

## Short guide

The following table provides an overview on the steps that make up the preparation process.

Ste	p		Us	er action
1	Start the Prepare wizard.	Standby     10.12.2007 13:46 (?)       Overview     Workplace     Utilities       1     2     3     4     5     6     7     8       0     Tests / 0     Blocked     0     7     8       0     Tests / 0     To Validate       STAT     Order     0     0       Tests     admin     0     0       Prepare     End Shift     000     0	1.	On the <b>Overview</b> tab, press <b>Prepare</b> .
2	Check the status of the external containers.	Standby 10.12.2007 12:27 () Check Bottles Press fluid button when done. Water Waste Cleaner 100% ()	2. 3.	Refill the water container, if required, and con- firm by pressing the water button. (If you refill the water, empty the waste as well.) Empty the waste container, if required, and confirm by pressing the waste button. Replace the cleaner bottle, if required, and con firm by pressing the cleaner button. Press >>>>>>>>>>>>>>>>>>>>>>>>>>>>>>>>>>>
3	Check the onboard ISE fluids.	Standby 10.12.2007 3:23 ( Prep3BE Status ISE Unit: Ready Ref Na K Ref 100% Cai 100% Cai	1. 2. 3.	Replace the ISE fluid bottles, if required. Replace the electrodes, if required. Press D to proceed to the next stage in the <b>Prepare</b> wizard.
4	Perform the maintenance actions that are due.	Standby       10.12.2007 3:23         Prep2Check Maintenance         Replace ISE Unit Tubing       10.12.         Replace ISE Pump Tubing       10.12.         Older Probe Manually       10.12.         Daily Prepare Actions       11.12.         Electrode Service       11.12.         Deproteinize Probe       11.12.         Clean ISE Tower Automatically       17.12.		Perform the <b>Daily Prepare Actions</b> mainte- nance action. Perform the other ISE maintenance actions that are due. Perform at least all red maintenance actions.

 Table E-4
 Steps for preparing the system

• For non-ISE-specific tasks, see *Preparing the system* on page B-12.

Starting the Prepare wizard

- ► To start the preparation process
  - 1 Choose Overview > Prepare.

#### Checking the status of the external fluid containers

#### To check the status of the external fluid containers

- **1** Follow the instructions given in *Checking the status of the external fluid containers* on page B-15.

A screen is displayed that shows the status of the electrodes and the on-board ISE fluids.



When checking external fluid containers without using the Prepare wizard: Choose Overview > ( ) > ( )

### **Checking the ISE status**



#### Injury through working solutions

Direct contact with cleaning solutions or other working solutions may cause personal injury.

When handling such solutions, exercise the precautions required for handling them, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics cleaning solutions.

#### Replacing electrodes

#### To replace an electrode

1 Follow the instructions given in *Replacing electrodes* on page E-32.



- After replacing an electrode, the Electrode Service maintenance action needs to be performed. This action will be performed as part of the maintenance procedure of the Prepare wizard.
- After replacing an electrode, a main calibration needs to be performed. This action will be performed as part of the **Daily Prepare Actions** maintenance action.

*Replacing ISE fluid bottles* A bottle with a blue label for ISE Calibrator indirect/urine and a bottle with a red label for ISE Reference Solution are located on the ISE unit. The procedure for exchanging these bottles is the same.

## ► To replace an on-board ISE fluid bottle

- 1 Follow the instructions given in *Replacing ISE fluid bottles* on page E-30.
  - Allow at least 15 seconds for priming the ISE bottle before continuing with any other activity, e.g. replacing the other ISE bottle.
  - After replacing an ISE fluid bottle, the **Prime ISE Reference and Calib.** maintenance action needs to be performed. This action will be performed as part of the maintenance procedure of the **Prepare** wizard.
  - After replacing an ISE fluid bottle, a main calibration needs to be performed. This action will be performed as part of the **Daily Prepare Actions**.
- 2 When you have finished replacing electrodes and ISE fluid bottles, press ≫ to proceed to the next stage in the **Prepare** wizard.

A screen is displayed that lists the maintenance actions.

### **Performing maintenance actions**

-`()´-

The routine maintenance actions that are relevant for working with the ISE unit are integrated in one single maintenance action: **Daily Prepare Actions**.

For general information on performing maintenance actions, see *Performing maintenance actions* on page B-19.
 For details on how to perform individual ISE maintenance actions, see *ISE maintenance actions* on page E-42.

### To perform maintenance actions

- 1 Select Daily Prepare Actions.
- **2** Press **()**.

The maintenance definition screen is displayed.

- **3** Follow the instructions given on the screen.
  - See *Daily prepare actions* on page E-44.
- 4 Perform the non-routine ISE maintenance actions.
  - Follow the instructions given on the screen.





When performing maintenance actions without using the Prepare wizard: Choose Utilities > Maintenance.

#### Completing the preparation tasks

1 When you have finished performing maintenance actions, press *▶* to proceed to the next stage in the **Prepare** wizard.

A screen is displayed that shows the status of the reagents on the reagent disk. This is not relevant for ISE operation.

**2** Press **>>**.

A screen is displayed that shows the status of the cuvettes that are on board. This is not relevant for ISE operation.

**3** Press ≫.

A screen is displayed that shows the tests that need calibrating. Because the calibrations have already been performed as part of the **Daily Prepare Actions** maintenance action, this step is not relevant for ISE operation.

4 Press ≫.

The Overview tab is displayed. The preparation tasks are completed.

#### **Performing calibrations**

The ISE calibrations are integrated in the Daily Prepare Actions maintenance action.

• See *Daily prepare actions* on page E-44.

## **Defining orders**

ISE sample analysis is integrated in the operation of the main instrument. The process of defining orders is the same.

### ► To define orders

1 Follow the instructions given in *Defining orders* on page B-38.

ISE tests are defined and requested as profiles. (There is hardly ever the need for electrolytic measurements of just one electrode.)

Ù.


## Monitoring the progress

### ► To monitor the progress

- 1 Follow the instructions given in *Monitoring the analysis progress* on page B-44.

## To check the instrument status

1 Press the **Overview** tab.

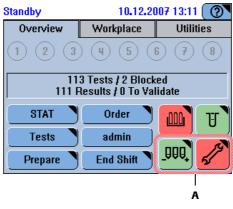

A Relevant instrument status buttons

#### Figure E-4

Interpreting the instrument status buttons

lcon		Meaning		ssible action
The		The ISE unit is in standby.	No	action is required.
		The expiration date of an electrode has passed or an on-board ISE fluid level is low.	1. 2.	Check the details on the dedi- cated <b>System Status</b> screen. Replace the electrode or ISE fluid bottle as soon as possible.
		The ISE unit is blocked.	1. 2.	Check the details on the dedi- cated ISE status screen. Act accordingly.

lcon	Meaning	Possible action	
System Status	buttons of the underlying <b>System Statu</b> tized by color, first priority being red, fo	System Status button displays both the icon and the color of one of the tons of the underlying System Status screen. (The icons are first priori- d by color, first priority being red, followed by yellow and green, and the ording to the sequence in which they are listed below.)	
	Analyzer (main cover)	Press the button and on the <b>Sys-</b> <b>tem Status</b> screen check the text about the status of the main cover	
1	Reagent cooler and cuvette ring temperature	Press the button and on the <b>Sys-</b> <b>tem Status</b> screen check the text about the temperature.	
<b>4</b>	Sample area ventilation	Press the button and on the Systen Status screen check the text about the ventilation status.	
<u>ÔÔ</u> )	External fluid containers	Press the button and on the <b>Sys-</b> tem Status press it again to display the screen for handling the exter- nal fluid containers.	
J.	Maintenance	Press the button and on the <b>Sys-</b> tem Status press it again to display the maintenance actions list.	
Ŀ	Printer	Press the button and on the <b>Sys-</b> <b>tem Status</b> screen check the text about the printer status.	

Acoustic signal

An acoustic signal informs you of the fact that all tests are finished and the system status has changed to **Standby**.

## Validating sample results

ISE result handling is integrated in the operation of the main instrument. The process of reviewing results is the same.

### To validate sample results

- 1 Follow the instructions given in *Validating sample results* on page B-51.
  - For information on ISE flags, see *List of ISE flags* on page E-71.

## Main calibration

The main calibration of the electrodes must be performed periodically. It also needs to be performed after certain maintenance actions.

For the indirect method, the sodium (NA-I) and potassium (K-I) tests are calibrated with a two-point calibration using ISE Solution 1 and 2.

The main calibration is typically performed during the **Prepare** phase and, if there follows a second shift, at the end of a shift. It is integrated in the **Daily Prepare Actions** maintenance action.

For details on performing the Daily Prepare Actions maintenance action, see Daily prepare actions on page E-44.
 For details on performing individual calibrations, see Process of performing calibrations on page B-61.

## **Performing QC**

ISE QC handling is integrated in the operation of the main instrument.

## ► To perform QC

• Follow the instructions given in *Performing QC* on page B-68.

## Finishing the shift

ISE Standby status	If there is no measurement for more than three minutes, the ISE unit switches to Standby status. In this status, ISE Calibrator indirect/urine and ISE Reference Solution are pumped into the measuring channel and moved a short distance at regular intervals. Also, every two hours, the ISE tubing is primed with ISE Calibrator indirect/urine and ISE Reference Solution. This is done to prevent the following:				
	• Flow of ISE Reference Solution backwards into the measuring channel, which can damage the electrodes because of the high ion concentration of the ISE Reference Solution.				
	Crystallization of salts in the tubing, which can cause blockages.				
	The ISE unit has a separate power supply. Therefore, the Standby status is maintained even if the main instrument is switched off.				
NOTICE	Damage to electrodes and possible tubing blockage				
	Do not unplug or switch off the ISE power supply. Periodic flow of solutions must be per- formed at all times.				
<b>End Shift</b> wizard	The end of shift activities are organized in a single wizard. By performing the steps as suggested by this wizard, you put the system in a condition that allows you to hand over operation to another operator or to shut down the system and switch off the instrument.				
	The following table provides an overview on the steps that make up the end of shift process.				
	• For details on performing the individual tasks, see the instructions in <i>Finishing the shift</i> on page B-77.				

## 13 ISE operation

Daily operation

## Short guide

Step			User action	
1	Check for unfinished tasks.	Standby     10.12.2007 4:17       Result Review>Filter       All       Flagged       STAT       Not Accepted       Not Today       Not Sent to Host       By ID	1. 2. 3.	Check for nonvalidated results.
2	Start the <b>End Shift</b> wizard.	Standby     10.12.2007 13:46 (?)       Overview     Workplace     Utilities       1     2     3     4     5     6     7     8       0     Tests / 0     Blocked     0     Results / 0     To Validate       STAT     Order     Image: Constant of the second se	1.	On the <b>Overview</b> tab, press the <b>End Shift</b> button.
3	Perform data backup.	Standby 10.12.2007 5:35 () End Shift>Backup Press <backup> to copy the current database to the USB stick.</backup>	2. 3. 4. 5.	Press       □         Insert the USB stick.         Press       □         Select the directory.         Press       □         Press       □         Press       □         Press       □         Press       ■         Press       ■         to proceed to the next stage in the         End Shift wizard.
4	Export support data.	Standby       10.12.2007 5:35 ()         End Shift>Export         End Shift>Export         Image: Standback of the	2. 3. 4. 5. 6.	Press €.         Insert the USB stick, if you have removed it.         Press ∅.         Select the directory.         Press ∅.         Remove the USB stick.         Press Ď to proceed to the next stage in the         End Shift wizard.
5	Clean up the database.	Standby       10.12.2007 5:36 ()         End Shift>Clean Up         Fress <delete> to clean up results:         - Accepted patient results (sent to host and/or printed)         - Accepted QC results         - Obsolete calibration results         - Obsolete calibration results         - Oc history older than previous month</delete>	2.	Press

**Table E-5**Steps for finishing the shift

### 13 ISE operation

Daily operation

Step		User action		
5	Perform maintenance actions.	Standby       10.12.2007 3:23         Prep>Check Maintenance         Replace ISE Unit Tubing       10.12.         Replace ISE Pump Tubing       10.12.         Clean Probe Manually       10.12.         Daily Prepare Actions       11.12.         Electrode Service       11.12.         Clean ISE Tower Automatically       17.12.         Clean ISE Tower Automatically       17.12.		Check which maintenance actions are due. Perform the maintenance actions. Perform at least all red maintenance actions. Press 🔊 to proceed to the next stage in the End Shift wizard.
7	Check the cuvette status. (Not relevant for ISE operation.)	Standby       10.12.2007 4:03         End Shift>Cuvettes         Select segment to replace         0       0	1.	Press <b>&gt;&gt;</b> to proceed to the next stage in the <b>End Shift</b> wizard.
;	Remove the reagent disk. (Not relevant for ISE operation.)	Standby     10.12.2007 12:34       Overview>Disk 2       ALTL     ASTL       BARBH     CA       CHOL2       CRPLX     D-DI       GLUC2     My_MG       THC       UA2     CLEAN       CLEAN     ALCL       PYP       Image: Standard St		
1	Empty the waste container.	Standby       10.12.2007 12:28         End Shift>Check Bottles         Press fluid button when done.         Water       Waste         Full       Waste         (20)         (20)       (20)	1. 2. 3. 4.	Empty the waste container. Press the <b>Waste</b> button to confirm. Refill the water container. Press the <b>Water</b> button to confirm.
0	Finish your shift.	Standby     10.12.2007 13:46       Overview     Workplace     Utilities       1     2     3     4     5     6     7     8       0     Tests / 0     Blocked     0     0     8       0     Tests / 0     To Validate       Tests     admin       Prepare     End Shitt     PRO	1. 2.	Log off the system. Shut down the system and switch off the main instrument. (If there is no other shift.)

<u>ک</u>

If you intend not to use the ISE unit for more than one week, you should deactivate it. This saves ISE fluids, reduces wear and tear of the tubing, and lowers the likelihood of malfunctions and errors occurring. (See *Deactivating the ISE unit* on page E-64.)

**Roche Diagnostics** 

Replacing ISE fluid bottles

## **Replacing ISE fluid bottles**



ISE fluid bottles are supplied with a barcode that contains information on their expiration date, which is monitored by the system.

Roche recommend replacing ISE fluid bottles as soon as their expiration date has passed. (Their icon on the screen turns yellow.)

For information on the stability of fluids, see their package inserts.

ISE Calibrator indirect/urine or ISE Reference Solution bottle

Tools and materials required

•

Tissues




## Injury through working solutions

Direct contact with cleaning solutions or other working solutions may cause personal injury.

When handling such solutions, exercise the precautions required for handling them, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics cleaning solutions.

### ► To replace an on-board ISE fluid bottle

- 1 Check that the system is in **Standby** status.
- 2 Choose Overview > []].
- **3** Check the status of the ISE bottles (A).

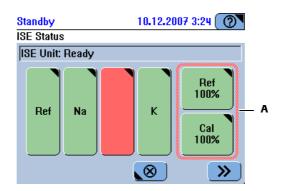

Interpreting the bottle status screen	No fluid registered by an ISE fluid sensor. (Operation has stopped.)You need to replace the bottle now.
	The fluid level in the bottle is low. (Operation will proceed until one of the sensors detects that there is no fluid.)
	The fluid's expiration date has passed.
	You should replace the bottle as soon as possible.
	No action is required.

### 4 Press Cal or Ref.

A screen is displayed that shows details about the bottle status.

#### **Roche Diagnostics**

Replacing ISE fluid bottles

5 Press Ito.

A screen is displayed, asking you to scan the bottle barcode or to type it manually.

- 6 Scan the bottle barcode.
- **7** Press  $\oslash$ .
- 8 Observe the messages on screen and react accordingly.

If the barcode cannot be read, type it manually.

When the scanning process is finished, a message is displayed, asking you to replace the bottle.

- **9** Open the ISE cover.
- 10 Remove the tubing adapter by lifting it and placing it on a clean lint-free tissue.
- **11** Remove the bottle.
- **12** Remove the cap of the new bottle.
- **13** Place the new bottle.



Make sure to place a full bottle.

- 14 Insert the tubing adapter.
- **15** Close the ISE cover.
- **16** On the screen, press 🞯 to confirm the placement.

The system performs the Prime ISE Reference and Calib. maintenance action.

**17** Press  $\textcircled{\otimes}$  to close the screen for reading barcodes.

Allow at least 15 seconds for priming the ISE bottle before continuing with any other activity.



You need to perform a main calibration before you can process orders.

**18** On the ISE status screen, press **>>**.

A placement list is displayed, telling you where to place ISE Solution 1 and 2.

- **19** Place ISE Solution 1 and ISE Solution 2 on the positions indicated on the screen.
- 20 Press 🔘.

The **Overview** tab is displayed.

- **21** Press 0 to perform the main calibration.
- **22** When the system is in **Standby** status, remove ISE Solution 1, and ISE Solution 2 from the sample area.
- **23** Press 🛞.

Replacing electrodes

## **Replacing electrodes**

For information on the stability of the electrodes, see their package inserts.		
There are up to three ion-selective electrodes and one ISE Reference Electrode on the ISE unit. The replacement procedure is the same for all of them.		
As part of replacing electrodes, the <b>Electrode Service</b> maintenance action and a main calibration need to be performed.		
Replacing an electrode takes about 15 minutes.		
• ISE Calibrator indirect/urine (on board)		
ISE Reference Solution (on board)		
• ISE Deproteinizer		
• ISE Etcher		
• Activator		
• ISE Solution 1		
• ISE Solution 2		
• ISE Sodium Electrode (if you intend to replace this type of electrode)		
• ISE Potassium Electrode (if you intend to replace this type of electrode)		
• ISE Reference Electrode (if you intend to replace this type of electrode)		
• ISE Dummy Electrode (if you intend to replace this type of electrode or if you no longer intend to use one of the ion-selective electrodes)		

- Injury through reagents and other working solutions on page E-19.
- Injury through reagents and other working solutions on page E-19.

Replacing electrodes

#### ► To replace an electrode

- 1 Check that the system is in **Standby** status.
- 2 Choose Overview > III.

The ISE status is displayed.

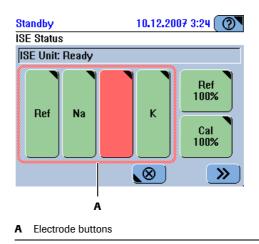


Figure E-5

3 Press the button of the electrode you want to replace.


Details of the electrode status are displayed.

4 Press 10 to start the exchange wizard.

The system rinses the electrodes with ISE Calibrator indirect/urine to remove any sample residue and then drains all electrodes.

You are asked to remove the electrode. (Make sure to remove the electrode that is marked on the screen.)

- **5** Open the ISE cover.
- **6** Open the electrode block lid.
- **7** Release the tension lever (A).

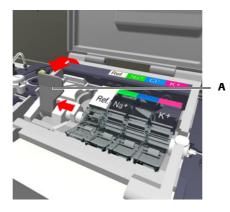


A Tension lever

#### Figure E-6

Turning the tension lever releases the electrodes. You may have to push the anchoring piece towards the tension lever to create enough space for removing the electrodes.

**8** Remove the electrode.

Lift it at the front to disengage, then remove it.

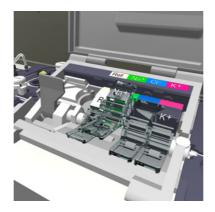

**9** On the screen, press the electrode button to confirm that you have removed the electrode.

You are asked to scan the barcode of the new electrode or to type its ID manually.

**10** Scan the electrode barcode on the package insert or type the ID manually.

(If you want to install a dummy electrode scan its barcode.)

You are asked to insert the new electrode.



### **Fixed electrode positions**

Always replace an electrode with the same type or with an ISE Dummy Electrode. (The ISE Reference Electrode must always be replaced with another ISE Reference Electrode. ISE Dummy Electrode is not allowed on this position.)

#### Keeping the barcode

Keep the electrode barcode in a safe place. You will need it again in the case of a database import.

**11** Insert the new electrode.

Push it back towards the contacts and press it down until it clicks into place.

- **12** Fasten the tension lever.
- **13** Close the electrode block lid.
- **14** On the screen, press the electrode button to confirm that you inserted the electrode.

(If you inserted a dummy electrode, press the button of the electrode you replaced with a dummy electrode.)

A screen is displayed that shows detailed information on the electrode.

15 Press 🛞.



At this stage, if you want to replace another electrode, press its button and continue with step 7.

- **16** Close the ISE cover.
- 17 On the ISE status screen, press  $\gg$  to continue the process.
  - A placement list is displayed.

- **18** Place the Activator, ISE Deproteinizer, ISE Etcher, ISE Solution 1, and ISE Solution 2 on the sample area positions indicated on the screen.
- **19** Press O to confirm the placement and to start the action.

The system performs the **Electrode Service** maintenance action.

**20** Press 0 to perform the main calibration.

A message will inform you when the action is complete.

- **21** When the system is in **Standby** status, remove the Activator, ISE Deproteinizer, ISE Etcher, ISE Solution 1, and ISE Solution 2 from the sample area.
- **22** Press 🛞.

## **Cleaning the ISE tower off the instrument**

If performing the **Clean ISE Tower Manually** maintenance action did not lead to the desired results, you need to remove the ISE tower to clean it.

Cleaning the ISE tower off the instrument consists of the following steps:

- 1. Removing the ISE tower
- 2. Soaking the ISE tower in ISE Deproteinizer
- 3. Cleaning and drying the ISE tower
- 4. Installing the ISE tower
- 5. Performing the maintenance action Clean ISE Tower Automatically.

Tools and materials required

- Cotton swabs
- ISE Deproteinizer
- Small glass beaker

Make sure that you have read and understood section *Safety information* on page E-19. The following warning messages in particular are relevant:

- Infection by biohazardous materials on page E-19.
- Injury through reagents and other working solutions on page E-19.

## To clean the ISE tower

Removing the ISE tower

1 Make sure the system is in **Standby** status.

- 2 Shut down the system and switch off the instrument.
- **3** Move the transfer head to its rightmost position.
- 4 Open the main cover and the left service flap.
- **5** Remove the ISE overflow collector by pulling it upwards and turning it from side to side as you do so.


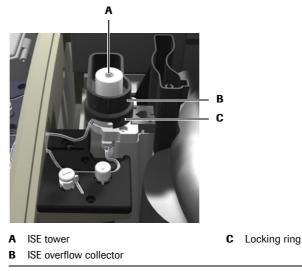


Figure E-7

Place the collector behind the wash station.

- **6** Remove the locking ring from the base of the ISE tower by turning it counterclockwise one quarter of a turn and then lifting it over the ISE tower.
- 7 Lift the ISE tower and carefully disconnect the tubing.

Check whether any O-ring is stuck to the ISE tower. If so, remove it and place it back in its position on the ISE tower base.

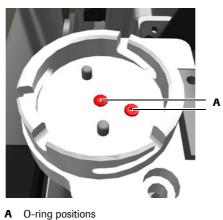

• o mg poo

Figure E-8

- *Cleaning the ISE tower* **8** Clean the inside of the ISE tower with a cotton swab to remove any visible clots.
  - **9** Pour ISE Deproteinizer in the beaker.


10 Soak the ISE tower in the ISE Deproteinizer for about 5 minutes.(If there are clots in the tower, you can extend the soaking time to 30 minutes.)

11 Rinse the ISE tower with deionized water and leave it to dry.

*Reinstalling the ISE tower* 

- **12** Reconnect the tubing to the ISE tower.
- **13** Make sure that both O-rings are present and seated on the tower base.
- **14** Place the ISE tower on its base.

Align the two pins on the base with the two holes in the ISE tower.

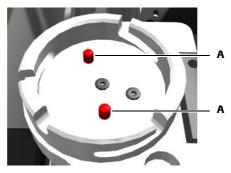

A Alignment pins



**15** Reinstall the locking ring.

Make sure it is properly engaged and turn it firmly clockwise.

- **16** Reinstall the overflow collector.
- **17** Close the left service flap and the main cover.

### **Roche Diagnostics**

Cleaning the ISE tower off the instrument

<b>18</b> Switch on the instrument.
<b>19</b> Wait for initialization to finish.
<b>20</b> Log on to the system.
<b>21</b> Perform the maintenance action <b>Clean ISE Tower Automatically</b> .

• See *Clean ISE tower automatically* on page E-50.

Table of contents

# **ISE** maintenance

In this chapter, you will find step-by-step instructions of the specific ISE maintenance actions that you must perform to keep the instrument running smoothly and efficiently.

In this chapter	Chapter <b>14</b>
ntroduction	
ISE maintenance actions and their intervals	E-41
ISE maintenance actions	E-42
Safety information	E-42
Daily prepare actions	E-44
Activate electrodes	E-45
Condition ISE tubing	E-46
Initialize ISE Reference Solution sensor	E-47
Prime ISE reference and calibrator	E-48
Electrode service	E-49
Clean ISE tower automatically	E-50
Clean ISE tower manually	E-51
Initialize ISE unit	E-53
Replace ISE pump tubing	E-54
Replace ISE unit tubing	E–57
Deactivating the ISE unit	E-64

## 14 ISE maintenance

Table of contents

# Introduction

The maintenance actions of the ISE unit are integrated in the maintenance actions of the main instrument.

● For general information on performing maintenance actions, see Overview on page C-5.

This chapter describes the specific ISE maintenance actions.

## ISE maintenance actions and their intervals

-Ò

The following table lists the ISE maintenance actions and shows how frequently they need to be performed.

Every effort has been made to ensure that all the information contained in this table is correct at the time of publication. However, Roche Diagnostics GmbH reserves the right to make any changes necessary without notice as part of ongoing product development.

Maintenance actions with no defined interval need to be performed in particular situations, for example during troubleshooting.

Interval	Maintenance action
None (Initiated by other maintenance action)	Activate Electrodes
None (Initiated by other maintenance action)	Condition ISE Tubing
None (Initiated by other maintenance action)	Initialize ISE Reference Sensor
None (Initiated by other maintenance action)	Prime ISE Reference and Calib.
Daily	Daily Prepare Actions
Daily	Electrode Service
Weekly	Clean ISE Tower Automatically
Monthly	Clean ISE Tower Manually
Monthly	Initialize ISE Unit
Yearly	Replace ISE Unit Tubing
Yearly	Replace ISE Pump Tubing
Table E-6         Periodici	ty of maintenance actions



For the expiration dates of electrodes, see their package inserts.

# **ISE** maintenance actions

## **Safety information**



#### Injury through working solutions

Direct contact with cleaning solutions or other working solutions may cause personal injury. When handling such solutions, exercise the precautions required for handling them, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics cleaning solutions.

#### Personal injury or damage to the analyzer due to contact with instrument mechanism

Do not touch moving parts during instrument operation.

#### Infection by biohazardous materials

Contact with samples containing material of human origin may result in infection. All materials and mechanical components associated with samples of human origin are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- · If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

#### Incorrect results due to build-up of contaminants

During use, contaminants may adhere to the probe. As a result, traces of analytes or reagents may be carried over to the next. Make sure to perform the probe maintenance actions as soon as they are due in order to prevent potentially false results.

#### Incorrect results or damage to the analyzer due to dust and soiling

The user can leave the main cover open while the system is in **Standby** status or while the instrument is shut down. This can cause dust and dirt being collected in the heating channel, which in turn might decrease the quality of the cuvettes.

Keep all covers closed. Open them only to perform operation actions.

#### Incorrect results or processing stop due to skipping maintenance actions

Not performing maintenance actions that are due may lead to situations where the system cannot continue processing orders, or it may lead to incorrect results. If at all possible, perform the maintenance actions when they are due.

#### Incorrect results or processing stop due to incomplete maintenance actions

You can cancel a maintenance action any time by pressing 🛞 at the end of a step. Using 🛞 while the system is performing a maintenance action may lead to unexpected errors.

If you interrupt the performing of a maintenance action that was due, its status will remain due, and you need to fully re-perform the action later.

- If at all possible, complete a maintenance action without interrupting it.
- If you need to interrupt the maintenance action, wait until you see confirmation on screen that the current step has been completed.



#### Incorrect results or damage to the analyzer due to wrong operation

Operators are required to have a profound knowledge of relevant guidelines and norms as well as the information and procedures contained in the Operator's Manual.

- Do not carry out operation and maintenance unless you have been trained by Roche Diagnostics.
- Start all maintenance actions on the screen. Do not perform maintenance actions without the assistance of the user interface.
- Carefully follow the procedures specified in the Operator's Manual for the operation and maintenance of the system.
- Leave maintenance that is not described in the Operator's Manual to a trained service representative.
- · Follow standard laboratory practices, especially when working with biohazard material.

#### Danger of poor measurement quality due to inadequate water quality

Inadequate water quality may lead to incorrect results. Always use purified water of the quality specified in section *Technical specifications*.

#### Infection by waste solution

Contact with waste solution may result in infection. All materials and mechanical components associated with the waste systems are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.



Whenever the term "purified water" is used in this document, water of at least the quality specified in chapter *Technical specifications* must be used.

**Daily prepare actions** 

# **Utilities > Maintenance > Daily Prepare Actions**

 $\rightarrow$ 

To make it easier and more efficient for the operator, some daily maintenance actions that may take quite some time are grouped into one maintenance action. This way the operator does not have to remain near the instrument while the actions are

The following maintenance actions are included if due:

- **Initialize ISE Unit**
- **Electrode Service**
- **Clean ISE Tower Automatically**
- **Deproteinize** Probe

In addition, a main calibration is performed.

**Operator** time Approximately 25 minutes.

performed.

**Prerequisites** The system must be in Standby status.

Tools and materials required

#### □ ISE Deproteinizer

- □ Activator
- □ ISE Etcher
- □ ISE Solution 1
- □ ISE Solution 2



Make sure that you have read and understood section Safety information on page E-42.

The following warning messages in particular are relevant:

- Injury through working solutions on page E-42.
- Incorrect results due to build-up of contaminants on page E-42.

#### To perform the daily prepare actions

- 1 Prepare a tube with ISE Deproteinizer, ISE Etcher, Activator, ISE Solution 1, and ISE Solution 2 each.
- 2 Select the maintenance action Daily Prepare Actions.
- 3 Press 💽.

The maintenance definition screen is displayed.

4 Press ⊘.

A placement list is displayed.

- **5** Place the fluids on the positions indicated on the screen.
- **6** Press 🔘 to confirm the placing.

The system performs the maintenance actions.

- 7 When the system is in **Standby** status, remove the tubes.
- 8 Press 🛞.

## **Activate electrodes**

#### → Utilities > Maintenance > Activate Electrodes

The electrode surfaces must be activated with fresh Activator so that the correct potentials are measured. Electrodes may be damaged if they are deprived of regular contact with Activator.

Operator time Approximately 4 minutes.

*Prerequisites* The system must be in **Standby** status.

#### *Tools and materials required* $\Box$ Activator



Make sure that you have read and understood section *Safety information* on page E-42. The following warning messages in particular are relevant:

- Injury through working solutions on page E-42.
- Infection by biohazardous materials on page E-42.

#### To activate the electrodes

- 1 Select the maintenance action Activate Electrodes.
- **2** Press **()**.

The maintenance definition screen is displayed.

**3** Press 🖉 to start the action.

A placement list is displayed.

- 4 Place the Activator on the sample area position indicated on the screen.
- **5** Press 🔘 to confirm the placement and to start the action.

- **6** Remove the Activator from the sample area.
- **7** Press  $\bigotimes$ .

Tools

# **Condition ISE tubing**

<b>→</b>	Utilities > Maintenance > Condition ISE Tubing
	Conditioning the tubing with Activator ensures the proper flow of fluids.
Operator time	Approximately 3 minutes.
Prerequisites	The system must be in <b>Standby</b> status.
s and materials required	□ Activator
	Make sure that you have read and understood section <i>Safety information</i> on page E-42. The following warning messages in particular are relevant:
	Injury through working solutions on page E-42.
	Infection by biohazardous materials on page E-42.
To condition the ISE tubing	
	<b>1</b> Select the maintenance action <b>Condition ISE Tubing</b> .
	<b>2</b> Press <b>()</b> .
	The maintenance definition screen is displayed.
	<b>3</b> Press 🔘 to start the action.
	A placement list is displayed.
	<b>4</b> Place the Activator on the sample area position indicated on the screen.
	<b>5</b> Press 🔘 to confirm the placement and to start the action.

- **6** Remove the Activator from the sample area.
- 7 Press  $\otimes$ .

## **Initialize ISE Reference Solution sensor**

→ Utilities > Maintenance > Initialize ISE Reference Sensor

Initializing the ISE Reference Solution sensor ensures proper detection of possible air in the ISE tubing.

This maintenance action forms part of the **Replace ISE Unit Tubing** maintenance action.

Operator time Approximately 1 minute.

□ Tissues

*Prerequisites* The system must be in **Standby** status.

Tools and materials required



Make sure that you have read and understood section *Safety information* on page E-42. The following warning messages in particular are relevant:

- Injury through working solutions on page E-42.
- Infection by biohazardous materials on page E-42.

#### ► To initialize the ISE Reference Solution sensor

- 1 Select the maintenance action Initialize ISE Reference Sensor.
- **2** Press **()**.

The maintenance definition screen is displayed.

**3** Press 🔘 to start the action.

The instrument pumps ISE Reference Solution through the ISE tubing.

You are asked to remove the tubing adapter from the ISE Reference Solution bottle.

- 4 Open the ISE cover.
- **5** Remove the tubing adapter from of the ISE Reference Solution bottle.
  - Have a clean lint-free tissue ready.
  - Lift the tubing adapter and place it on the tissue.
- **6** Press 🕖 to confirm completion of the action.

The system now starts to initialize the ISE Reference Solution sensor.

You are asked to insert the tubing adapter in the ISE Reference Solution bottle and to close the ISE cover.

- 7 Insert the tubing adapter in the ISE Reference Solution bottle.
- 8 Close the ISE cover.
- **9** Press 🔘.

The instrument pumps ISE Reference Solution through the ISE tubing.

- 10 Press 🛞.

## Prime ISE reference and calibrator

$\rightarrow$	Utilities > Maintenance > Prime ISE Reference and Calib.
	This maintenance action ensures that no traces of a previous ISE Calibrator indirect/urine and ISE Reference Solution are left in the tubing and so prevents drift in the results.
Operator time	Approximately 1 minute.
Prerequisites	The system must be in <b>Standby</b> status.
Tools and materials required	None
Hazards and precautions	None
	To prime the ISE calibrator and ISE Reference Solution

### ► To prime the ISE calibrator and ISE Reference Solution

- 1 Select the maintenance action **Prime ISE Reference and Calib.**
- **2** Press **()**.

The maintenance definition screen is displayed.

**3** Press 🔘 to start the action.

The instrument pumps first ISE Reference Solution and then ISE Calibrator indirect/urine through the ISE tubing and past the electrodes.

- 4 Press 🛞.

## **Electrode service**

→ Utilities > Maintenance > Electrode Service	Maintenance > Electrode Service	e
---	---------------------------------	---

The electrodes must be cleaned regularly to prevent the buildup of deposits and so to maintain the efficiency of the ISE unit.

During this action, the ISE Sodium Electrode is etched and all electrodes are deproteinized and activated, and the tubing is conditioned.

Operator time Approximately 9 minutes.

*Prerequisites* The system must be in **Standby** status.

Tools and materials required

- □ ISE Etcher
- □ Activator
- □ ISE Deproteinizer

Make sure that you have read and understood section *Safety information* on page E-42. The following warning messages in particular are relevant:

- Injury through working solutions on page E-42.
- Infection by biohazardous materials on page E-42.

#### To perform electrode service

- 1 Select the maintenance action Electrode Service.
- 2 Press 💽.

The maintenance definition screen is displayed.

**3** Press 🔘 to start the action.

A placement list is displayed.

- **4** Place the Activator, ISE Deproteinizer, and ISE Etcher on the sample area positions indicated on the screen.
- **5** Press 🔘 to confirm the placement and to start the action.

A message will inform you when the maintenance action is complete.

- 6 Remove the Activator, ISE Deproteinizer, and ISE Etcher from the sample area.
- 7 Press ⊗.



You need to perform a main calibration before you can process orders.

## **Clean ISE tower automatically**

#### → Utilities > Maintenance > Clean ISE Tower Automatically

The ISE tower must be cleaned regularly to ensure the proper functioning of the ISE unit. Cleaning includes deproteinizing and activating the ISE tower.

Operator time Approximately 3 minutes.

*Prerequisites* The system must be in **Standby** status.

Tools and materials required

□ ISE Deproteinizer

□ Activator



Make sure that you have read and understood section *Safety information* on page E-42. The following warning messages in particular are relevant:

- Injury through working solutions on page E-42.
- Infection by biohazardous materials on page E-42.

#### ► To clean the ISE tower automatically

- 1 Select the maintenance action Clean ISE Tower Automatically.
- 2 Press 💽.

The maintenance definition screen is displayed.

**3** Press 🔘 to start the action.

A placement list is displayed.

- **4** Place the Activator and the ISE Deproteinizer on the sample area positions indicated on the screen.
- **5** Press 🔘 to confirm the placement and to start the action.

- 6 Remove the Activator and ISE Deproteinizer from the sample area.
- 7 Press 🛞.

Tools

## **Clean ISE tower manually**

$\rightarrow$	<b>Utilities &gt; Maintenance &gt; Clean ISE Tower Manually</b>
-\\$-	When the ISE tower is blocked, perform first the <b>Clean ISE Tower Automatically</b> mainte- nance action (See <i>Clean ISE tower automatically</i> on page E-50). If this does not help, per- form the <b>Clean ISE Tower Manually</b> maintenance action. If the ISE tower is still blocked, clean the ISE tower off the instrument. (See <i>Cleaning the ISE tower off the instrument</i> on page E-36.)
Operator time	Approximately 6 minutes.
Prerequisites	The system must be in <b>Standby</b> status.
s and materials required	□ ISE Deproteinizer
	□ Activator
	□ Cotton swabs
	Make sure that you have read and understood section Safety information on page E-42.
<u> </u>	The following warning messages in particular are relevant:

- Injury through working solutions on page E-42.
- Infection by biohazardous materials on page E-42.

#### To clean the ISE tower

- 1 Select the maintenance action Clean ISE Tower Manually.
- **2** Press **()**.

The maintenance definition screen is displayed.

**3** Press 🔘 to start the action.

A placement list is displayed.

- **4** Place the Activator and ISE Deproteinizer on the sample area positions indicated on the screen.
- **5** Press 🔘 to confirm the placement and to start the action.


The ISE tower is first washed.

The transfer head moves to the right side of the instrument.

You are asked to clean the ISE tower.

**6** Open the main cover and then the left service flap.

7 Clean the inside of the ISE tower with a cotton swab to remove any residual of ISE Deproteinizer.

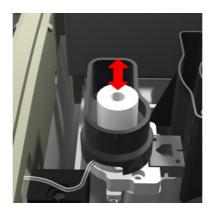

- **8** Close the left service flap and then the main cover.
- 9 Press Ø to confirm the cleaning of the ISE tower.The ISE tower is first washed.

The transfer head moves to the right side of the instrument.

You are asked to clean the ISE tower.

- **10** Open the main cover and then the left service flap.
- **11** Clean the inside of the ISE tower with a cotton swab to remove any residual water.
- **12** Close the left service flap and then the main cover.
- **13** Press 🔘 to confirm the cleaning of the ISE tower.

The ISE tower is first washed and dried and then conditioned with Activator.

- A message will inform you when the maintenance action is complete.
- 14 Remove the Activator and ISE Deproteinizer from the sample area.
- 15 Press 🛞.

# Initialize ISE unit

#### → Utilities > Maintenance > Initialize ISE Unit

Initializing the ISE unit uses Activator to condition the electrodes and to adjust the sensors and the pump action. This ensures proper transportation of fluids through the ISE tubing.

Operator time Approximately 3 minutes.

*Prerequisites* The system must be in **Standby** status.

*Tools and materials required*  $\Box$  Activator



Make sure that you have read and understood section *Safety information* on page E-42. The following warning messages in particular are relevant:

• Infection by biohazardous materials on page E-42.

### To initialize the ISE unit

- 1 Select the maintenance action Initialize ISE Unit.
- **2** Press **()**.

The maintenance definition screen is displayed.

**3** Press 🖉 to start the action.

A placement list is displayed.

- 4 Place the Activator on the sample area positions indicated on the screen.
- **5** Press Ø to confirm the placement and to start the action.

- **6** Remove the Activator from the sample area.
- 7 Press  $\bigotimes$ .

## **Replace ISE pump tubing**

Replacing the pump tubing ensures proper ISE pump action.

During the **Replace ISE Pump Tubing** maintenance action, the following maintenance actions are automatically performed:

- Prime ISE Reference and Calib.
- Initialize ISE Unit

Operator time Approximately 10 minutes.

*Prerequisites* The system must be in **Standby** status.

□ Activator

Tools and materials required



□ Roche approved ISE pump tubing set

Make sure that you have read and understood section *Safety information* on page E-42. The following warning messages in particular are relevant:

- Injury through working solutions on page E-42.
- Infection by biohazardous materials on page E-42.
- Infection by waste solution on page E-43.

#### To replace the ISE pump tubing

- 1 Select the maintenance action **Replace ISE Pump Tubing**.
- **2** Press **()**.

The maintenance definition screen is displayed.

**3** Press 🔘 to start the action.

The placement screen is displayed.

- 4 Place the Activator on the sample area position indicated on the screen.
- **5** Press 🔘 to confirm the placement.

The tubing is emptied.

Your are asked to remove the tubing.

**6** Open the ISE cover.


*Removing the tubing* **7** Open the cover of the peristaltic pump.



A Cover of peristaltic pump

#### Figure E-10


- 8 Push the pump towards the back of the instrument.
- **9** Disengage the tube connector plate.

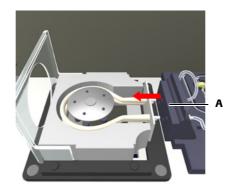

A Connector plate

Figure E-11

**10** Remove the ISE pump tubing set.

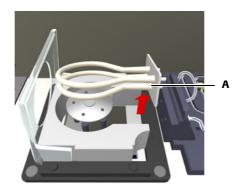

A ISE pump tubing set


Figure E-12

Installing the tubing

*tubing* Install the tubing in the reverse order of removing it.

**11** Pull the new tubing set over the pump head, with the connector plate facing the electrode block.

**12** Align the tubing connector plate with the tube connectors on the ISE unit.

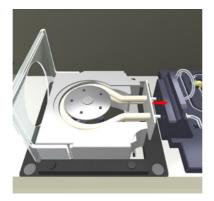

**13** Carefully close the peristaltic pump lid.

The tube connector plate is automatically pushed into the tube connectors of the ISE unit.

- **14** Close the ISE cover.
- **15** Press 🕖 to confirm completion of the action.

The system first performs the **Prime ISE Reference and Calib.** maintenance action, then **Initialize ISE Unit**.

A message will inform you when the maintenance actions are complete.

**16** Remove the Activator from the sample area.

**17** Press 🛞.

# **Replace ISE unit tubing**

## → Utilities > Maintenance > Replace ISE Unit Tubing

Due to the abrasion by the ISE valve action, the ISE tubing needs to be replaced periodically. Replacing the ISE tubing ensures the proper functioning of the ISE unit.

During the **Replace ISE Unit Tubing** maintenance action, the following maintenance actions are automatically performed:

- Prime ISE Reference and Calib.
- Initialize ISE Reference Sensor
- Condition ISE Tubing
- Initialize ISE Unit



Always replace the ISE tubing with the help of the **Replace ISE Unit Tubing** maintenance action.

Operator time Approximately 15 minutes.

*Prerequisites* The system must be in **Standby** status.

Tools and materials required

## □ Activator

- $\hfill\square$  Roche approved ISE unit tubing set
- □ Tissues

Make sure that you have read and understood section *Safety information* on page E-42. The following warning messages in particular are relevant:

- Injury through working solutions on page E-42.
- Infection by biohazardous materials on page E-42.
- Infection by waste solution on page E-43.
- Personal injury or damage to the analyzer due to contact with instrument mechanism on page E-42.

#### To remove the ISE unit tubing

- 1 Select the maintenance action **Replace ISE Unit Tubing**.
- 2 Press 💽.

The maintenance definition screen is displayed.

**3** Press 🔘 to start the action.

The placement screen is displayed.

- 4 Place the Activator on the sample area position indicated on the screen.
- **5** Press 🕖 to confirm the placement.

The tubing is emptied.

Your are asked to open the covers.

**6** Open the main cover, then the left service flap and the ISE cover.

Your are asked to remove the tubing adapters from the ISE solution bottles.

#### 14 ISE maintenance


ISE maintenance actions

- *Disconnecting the ISE solution* **7** Remove the tubing adapters from ISE Calibrator indirect/urine and ISE Reference Solution bottles.
  - **8** Press 🞯 to confirm the removal.

The tubing is drained.

**9** Remove the tubing from the adapters.

The two tubing adapters are slightly different, but the process of removing the tube is the same.

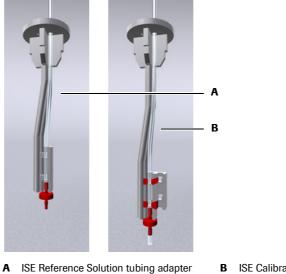


**B** ISE Calibrator indirect/urine tubing adapter

#### Figure E-13

- Have a tissue ready.
- Remove the tubing adapter by lifting it and place it on the clean lint-free tissue.
- Remove the tube from its clips.
- Remove the nozzle from the tube by pulling it off.
- Pull through the tube from the top of the adapter.


Place the adapter together with its nozzle on the tissue.

**10** Remove the pinch valve caps and clamps of the input (B), entry (C), and exit (A) valve-plates.

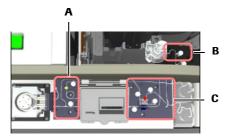


Removing the pinch valve caps and clamps

To remove a pinch valve cap and clamp:

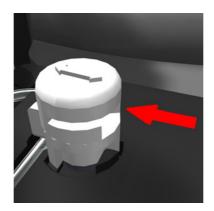

- Push the cap horizontally in the direction of the arrow, away from the cut-out.
- Lift and put it aside.
- Lift the valve clamp and put it aside.

Disconnecting the tubing from ISE tower **11** Disconnect the tubing from the ISE tower

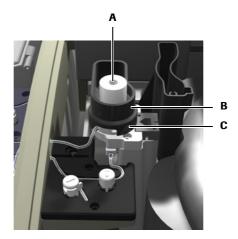

• Remove the ISE overflow collector (B) by pulling it upwards and turning it from side to side as you do so.


Place the collector behind the wash station.

- Remove the locking ring (C) from the base of the ISE tower by turning it counter-clockwise one quarter of a turn and then lifting it over the ISE tower.
- Lift the ISE tower (A). Check whether any O-ring is stuck to the ISE tower. If so, remove it and place it back in its position on the ISE tower base.
- Disconnect the tubing.



Do not remove the piece of tubing at the underside of the mixing tower. It will be changed when a service representative performs periodic maintenance.


**12** Slide off the lid of the fluid distributor.

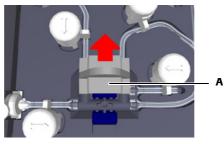

A Lid of fluid distributor

#### Figure E-14

**13** Disconnect the tubing from the electrode block.

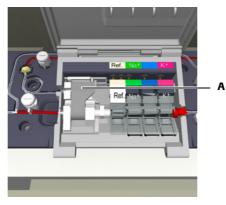

#### A Tension lever

#### Figure E-15

- Open the electrode block lid.
- Release the tension lever.
- Lift the connector on the side of the K electrode.
- Disconnect the tube on the side of the tension lever.
- Lift the connector to the ISE Reference Electrode and disconnect the tube leading to the ISE Reference Solution bottle.

**14** Disconnect the tubing connecting to the peristaltic pump.

Removing the tubing	<b>15</b> Lift the tubing from its routing, follow carefully the individual tubes.
	<b>16</b> Dispose of the tubing. Treat it as biohazardous waste.
stalling the new tuhing	<b>17</b> Install and connect the tubing in the order recommended on the illustrat

*Installing the new tubing* **17** Install and connect the tubing in the order recommended on the illustration on the ISE tubing set cover.

When installing the single items, perform the individual steps in reverse order of removing the items.

#### ► To install the new ISE unit tubing

1 Remove the new tubing set from its packaging.

<u>`</u>\

Make sure to keep the piece of tubing designed to fit at the underside of the mixing tower. It will be required when a service representative performs periodic maintenance.

**2** Place the tubing roughly in position on the ISE unit. Refer to the illustration on the ISE tubing set cover.

Take care not to cut the tubing, avoid any sharp objects.


**3** Press the tubes into their routing.

Take care not to compress the tubes or to kink them.

Make sure to press the tubing firmly in its routing around the ISE Reference Solution sensor.

- 4 Connect the tubing to the ISE tower and then install the tower.
- **5** After connecting the tubing to the electrode block, make sure the electrodes are properly mounted when closing the tension lever. Close the electrode block cover.
- **6** When connecting the tubing to the peristaltic pump connector block, take care not to introduce sharp angles in the tubing. (Sharp angles may lead to flow problems.)


Always insert the tubing connectors in the lower set of holes in the pump connector block, and make sure to fully insert the tubing in the connectors.

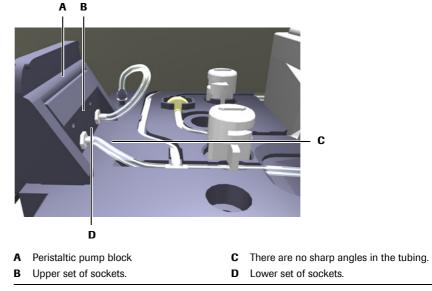

- Figure E-16
- 7 Install all clamps and pinch valve caps. Be sure to place the clamps in the correct position, covering the tubing.
- **8** Install the lid of the fluid distributor.

**9** Install the tubing on the tubing adapters of the ISE Calibrator indirect/urine and ISE Reference Solution bottles.

The two tubing adapters are slightly different, but the process of installing the tube is the same.

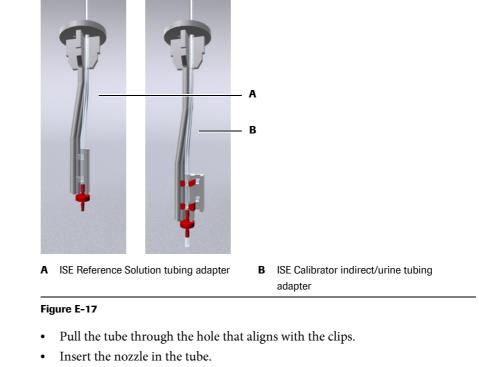

- At the top of the adapter, pull through the tube until the nozzle is flush with the end of the adapter. There should be no slack.
- Fix the tube to the clips.

With the adapter that offers two clip positions, use the inner position.

- Insert the tubing adapter in the appropriate bottle.
- **10** Close the ISE cover, then the left service flap and the main cover.
- **11** Press 🕖 to confirm completion of the action.

*Conditioning the tubing* The system now performs the maintenance action **Prime ISE Reference and Calib.** 

A message will inform you when the maintenance action is complete.

The system now performs the maintenance action **Initialize ISE Reference Sensor**.

You are asked to remove the tubing adapter from of the ISE Reference Solution bottle.

**12** Open the ISE cover.

**13** Remove the tubing adapter from of the ISE Reference Solution bottle.

- Have a clean lint-free tissue ready.
- Lift the tubing adapter and place it on the tissue.

**14** Press 🕖 to confirm completion of the action.

The system now starts to initialize the ISE Reference Solution sensor.

You are asked to insert the tubing adapter in the ISE Reference Solution bottle and to close the ISE cover.

- **15** Insert the tubing adapter in the ISE Reference Solution bottle.
- **16** Close the ISE cover.
- **17** Press 🔘.

The system first pumps ISE Reference Solution through the ISE tubing. Then it performs the **Condition ISE Tubing** maintenance action, and finally **Initialize ISE Unit**.

- **18** Remove the Activator from the sample area.
- **19** Press 🛞.

## **Deactivating the ISE unit**

If you intend not to use the ISE unit for more than one week, you should deactivate it. This saves ISE fluids and reduces wear and tear of the tubing, and lowers the likelihood of malfunctions and errors occurring.

This situation may occur, for example, if you need close the laboratory temporarily, or if you want to relocate the **Cedex Bio** instrument with the ISE unit.

Tools and materials required

- □ Glass beaker
- $\hfill\square$  Deionized water
- □ Tissues



Make sure that you have read and understood section *Safety information* on page E-42. The following warning message in particular is relevant:

- Injury through working solutions on page E-42.
- Infection by biohazardous materials on page E-42

#### To prepare the ISE data

- 1 Delete all ISE sample, calibration, and control results.
  - See Deleting sample results on page B-94, Deleting calibration results on page B-95, Deleting QC results on page B-96.
- **2** Disable all ISE applications.
  - See Activating and deactivating applications on page B-139.
- **3** Delete the ISE lot data.
  - See *Lot handling* on page B-98.

#### To deactivate the ISE unit

- **1** Open the ISE cover.
- 2 Remove the tubing adapters from the ISE solution bottles.
- **3** Place the adapters in a beaker filled with deionized water.
- 4 Remove bottles and screw on their lids.
- **5** Place the beaker in the position where your removed the ISE solution bottles from.
- 6 Choose Utilities > Maintenance > Prime ISE Reference and Calib.
- 7 Press 💽.

The maintenance definition screen is displayed.

**8** Press 🔘 to start the action.


The ISE unit tubing is primed with water.

A message will inform you when the maintenance action is complete.

- 9 Repeat steps 6 through 8.
- **10** Remove the adapters from the beaker and place them on a lint-free tissue.
- **11** Repeat steps 6 through 8.

The water is drained from the ISE unit tubing.


**12** Open the cover of the peristaltic pump.

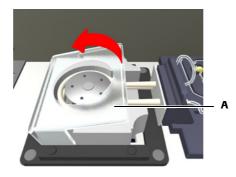

A Cover of peristaltic pump

#### Figure E-18

- **13** Push the pump towards the back of the instrument.
- **14** Disengage the tube connector plate.

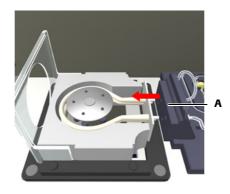

A Connector plate

#### Figure E-19

- **15** Remove all pinch valve caps and clamps, store them in secure place.
  - See *Removing the pinch valve caps and clamps* on page E-58.
- **16** Close the ISE cover.
- 17 Shut down the system and switch off instrument.
- **18** Disconnect the ISE power supply from the wall socket.
- **19** Switch on instrument.
  - On the **Overview** tab, the ISE icon  $\bigcirc$  is deactivated (grey).

### 14 ISE maintenance

ISE maintenance actions

# **ISE troubleshooting**

# How to deal with exceptional ISE situations

In this chapter, you will find information on the result flags that may be generated with ISE measurements, and also guidance for general troubleshooting procedures and on how to deal with error messages.

## Chapter In this chapter Introduction ......E-69 Messages ......E-69 List of ISE flags ......E-71 Air Fluid ......E-71 Calc Error ......E-73 Ise Unstab ...... E-73 Out of Rng ......E-75 Seg Fluid ......E-77 Seg Isecal ..... E-78 Reacting to error messages ..... E-80

**Roche Diagnostics** 

## 15 ISE troubleshooting

Table of contents

# Introduction

Effective troubleshooting on the ISE unit requires a good understanding of all basic operating procedures. However, you can avoid most of the common problems if you observe the recommended procedures at all times and if the operating environment is effectively controlled.

• For information on ISE maintenance, see ISE maintenance actions on page E-42.

## Information about the ISE unit status

The user interface keeps you informed about the status of the ISE unit as a whole, and about particular hardware, software, and chemistry events as they arise. It does this in the following ways:

• Color coded LEDs on the instrument inform you when to open covers or place sample tubes.

• See *Color concept* on page A-72.

• The colors of buttons tell you whether you need to intervene.

You can check the meaning of a button and its color using the Online Help 2.

- Buttons on the **Overview** tab lead to detailed information on the status of selected processes and hardware items.
- Messages on the screen provide information on individual tasks and events.
- The text in the Status line provides information on the status and activities of the analyzer unit.
- Flags with results from samples, calibrations, and controls are automatically generated if during processing certain technical checks were not passed or if the result exceeds or does not reach predefined limits.

## Messages

The system performs numerous checks. When an irregularity is detected, an alarm message is generated. Alarm messages are displayed in two ways:

- Immediate feedback on user actions is displayed in a *pop-up message screen*.
- Information concerning a problem that occurred during operation is reported in the *Alarm Monitor*.
- For details on alarm messages, see *Alarm monitor* on page D-6.
   For details on instrument messages, see *Message screen* on page D-5.

# Safety

Before you start troubleshooting, it is essential that you both read and understand the safety information listed below.

Read carefully all safety notices given in instructions and make sure you understand them.



#### Injury through working solutions

Direct contact with cleaning solutions or other working solutions may cause personal injury. When handling such solutions, exercise the precautions required for handling them, observe the cautions given in the package insert, and observe the information given in the Safety Data Sheets available for Roche Diagnostics cleaning solutions.

#### Infection by biohazardous materials

Contact with samples containing material of human origin may result in infection. All materials and mechanical components associated with samples of human origin are potentially biohazardous.

- Be sure to wear protective equipment. Take extra care when working with protective gloves; these can easily be pierced or cut, which can lead to infection.
- If any biohazardous material is spilled, wipe it up immediately and apply disinfectant.
- If waste solution contacts your skin, wash it off immediately with water and apply a disinfectant. Consult a physician.

#### Injury or damage to the analyzer due to contact with instrument mechanism

Do not touch any parts of the instrument other than those specified. During operation and maintenance of the instrument, proceed according to the instructions.

#### Incorrect results or damage to the analyzer due to wrong operation

Operators are required to have a profound knowledge of relevant guidelines and norms as well as the information and procedures contained in the Operator's Manual.

- Do not carry out operation and maintenance unless you have been trained by Roche Diagnostics.
- Start all maintenance actions on the screen. Do not perform maintenance actions without the assistance of the user interface.
- Carefully follow the procedures specified in the Operator's Manual for the operation and maintenance of the system.
- Leave maintenance that is not described in the Operator's Manual to trained Roche Technical Support personnel.
- · Follow standard laboratory practices, especially when working with biohazard material.

# List of ISE flags

<u>`</u>0́-

This section lists the general and ISE-specific flags in alphabetic order and provides information on each of the flags.

• For information on flags that concern absorbance measurements only, see *List of flags* on page D-28.

Where a flag description gives a list of recommended actions, work through each step in turn until you have solved the problem. Proceed to the next step only after you have completed the previous one.

Meaning	Result generated by re-performing the test with the same running parameters (repeated).
Message ID	120
Priority	32
Possible cause	The user initiated the repeat.
Recommended actions	No action required.

## **Air Fluid**

Meaning	<ul><li>The length of the sample segment was too short.</li><li>The end of the sample segment was not detected. It takes too long for a sample segment to reach the electrodes.</li></ul>
Message ID	19
Priority	4
Possible cause	<ul> <li>Air bubbles are present.</li> <li>The ISE tower is clogged.</li> <li>The dosing from the sample probes is inaccurate.</li> <li>The ISE tubing is leaking, blocked, or not conditioned.</li> <li>Mixing is too vigorous.</li> </ul>
Recommended actions	<ul><li>Check the quality of the sample.</li><li>Check the sample tubes.</li><li>Check the ISE tower and ISE tubing for leaks and blockages.</li><li>Repeat the measurement.</li></ul>

## 15 ISE troubleshooting

List of ISE flags

# **Air Isecal**

Meaning	<ul><li>The length of the calibrator segment was too short.</li><li>The end of the calibrator segment was not detected. It takes too long for a calibrator segment to reach the electrodes.</li></ul>
Message ID	91
Priority	5
Possible cause	<ul><li>Air bubbles are present.</li><li>The ISE tubing is leaking or blocked.</li></ul>
Recommended actions	• Ensure that the ISE Calibrator Indirect on the ISE unit is available and check the level of fluid; replace if necessary.
	<ul><li>Repeat the measurement.</li><li>Check the ISE tubing for leaks and blockages.</li></ul>

## ? Cal

Meaning	A result where its calibration is flagged.
Message ID	110
Priority	29
Possible cause	The calibration was in question and has been flagged, however the calibration could be used to calculate the result (compare with Cal Error where the calibration did not provide a usable result).
Recommended actions	Check and redo calibration if needed.

## **Cal Error**

Meaning	No valid calibration data available.
Message ID	43
Priority	24
Possible cause	Caused by an alarm that occurred during calibration because the calibration could not provide a usable result.
Recommended actions	Check the flags of the calibration and proceed to deal with the calibration first.

# **Calc Error**

Meaning	Calculation error.
	General calculation error.
	• Slope or nonlinear standard curve cannot be calculated due to a calibration error.
Message ID	16
Priority	13
Possible cause	Calibrator outdated or deteriorated.
	Misplacement of the calibrator tubes.
Recommended actions	Repeat calibration with fresh calibrators.
	• If the calibrators were incorrectly positioned, replace the calibrators and repeat the calibration.

## Ise Unstab

Meaning	The signal from the electrode(s) was not stable during the measurement.
Message ID	18
Priority	6
Possible cause	<ul> <li>There are air bubbles in the measuring segment.</li> <li>The concentration of sample is too low (for example, dialyzed samples).</li> <li>The ISE Reference Solution tubing is blocked or leaking.</li> <li>The pH of the sample is too low (for example, pH &lt; 5.5).</li> <li>One or more O-rings are missing or damaged.</li> <li>There are salt deposits or there is an obstruction at the adapter of the ISE Reference Solution bottle.</li> <li>The electrodes have not been serviced.</li> <li>The expiration date of the electrode has expired or the electrode is damaged.</li> <li>The electrodes and/or the electrode block is not dry.</li> <li>The preamplifier is defective.</li> </ul>
Recommended actions	<ol> <li>Perform the maintenance action Electrode Service and repeat the measurement.</li> <li>Repeat the measurement with a suitable sample (where pH &gt; 5.5 and sample concentration within the test range) to establish that the ISE module is OK.</li> <li>Check the position and condition of the O-rings and replace if necessary.</li> <li>Check the adapter of the ISE Reference Solution bottle for salt deposits and obstruction.</li> </ol>

# No Fluid

Meaning	The ISE sample or ISE Reference Solution sensor could not detect any fluid. It takes too long for a sample segment to reach the electrodes.
Message ID	68
Priority	2
Possible cause	<ul> <li>The fluid was not pipetted.</li> <li>The ISE tower is clogged.</li> <li>The ISE tubing is blocked or leaking.</li> <li>The ISE sample sensor is defective.</li> <li>The ISE Electrodes are blocked or not properly mounted.</li> </ul>
Recommended actions	<ol> <li>Repeat the measurement.</li> <li>Check the sample tubing for clots.</li> <li>Check ISE Electrodes and O-Rings for proper mounting.</li> <li>Check the ISE tower and ISE tubing for leaks and blockages.</li> <li>Ensure that the ISE bottles on the ISE unit are available and check the level of fluid; replace if necessary. (Allow at least 15 seconds for priming the ISE bottle before continuing with any other activity, e.g. replacing the other ISE bottle.)</li> </ol>

# No Isecal

Meaning	The ISE sample sensor could not detect any calibrator. It takes too long for a calibrator segment to reach the electrodes.
Message ID	90
Priority	3
Possible cause	<ul><li>ISE Calibrator Indirect bottle is not available.</li><li>The ISE tubing is leaking or blocked.</li><li>The ISE sample sensor is defect.</li></ul>
Recommended actions	<ol> <li>Repeat the measurement.</li> <li>Ensure that the ISE Calibrator Indirect on the ISE unit is available and check the level of fluid; replace if necessary</li> </ol>

# Out of Rng

Meaning	Slope out of range.
	The slope for the specified electrode is outside the acceptable/programmed range.
Message ID	20
Priority	15
Possible cause	<ul> <li>This flag appears with calibrations because:</li> <li>There are problems with the ISE solutions on the instrument.</li> <li>There are problems with the electrodes or the ISE unit was not properly serviced.</li> <li>The ISE Reference Solution tubing is blocked or leaking.</li> </ul>
	<ul><li>The dosing from the sample probe is not accurate.</li><li>The preamplifier is not working properly.</li></ul>
Recommended actions	<ol> <li>If only one calibration is flagged:         <ol> <li>Repeat the calibration.</li> <li>Repeat the calibration with fresh ISE Solutions 1 and 2, and check visually the correct flow of all ISE solutions and the ISE Reference Solution.</li> <li>If successive calibration with fresh ISE Solutions 1 and 2, and check visually the correct flow of all ISE solutions and the ISE Reference Solution.</li> <li>Repeat the calibration with fresh ISE Solutions 1 and 2, and check visually the correct flow of all ISE solutions and the ISE Reference Solution.</li> <li>Check whether the electrode expiration date has expired. Replace the electrodes, if necessary.</li> <li>Perform the maintenance action Electrode Service.</li> <li>Perform the maintenance action Deproteinize Probes.</li> <li>Check ISE Reference Solution tubing for leaks and blockages.</li> <li>If the slopes of all electrodes are zero, replace the ISE Reference Electrode.</li> <li>Contact your service representative.</li> </ol> </li> </ol>
? QC	
Meaning	A result where the quality control measurement is flagged.
Message ID	111
Priority	30
Possible cause	The last QC measurement before this measurement was flagged. The dependency is derived from the time sequence.
Recommended actions	Check OC measurement.

# R 1(2.5s)

Meaning	One control value is above 2.5 standard deviation or below -2.5 standard deviation.
Message ID	36
Priority	28
Possible cause	<ul><li>The reagent has deteriorated (linearity of working curve degraded).</li><li>One control is concentrated or has deteriorated.</li></ul>
Recommended actions	<ol> <li>Repeat with fresh control.</li> <li>Check that the correct control material has been used.</li> <li>If the flag reappears, check the calibration and reagent status. Repeat with fresh reagent.</li> </ol>

# R 1(3s)

Meaning	One control value is above 3 standard deviation or below -3 standard deviation.
Message ID	35
Priority	27
Possible cause	<ul><li>Improper control is set.</li><li>Proper control values (mean value, standard deviation) are not specified.</li></ul>
Recommended actions	<ol> <li>Repeat with fresh control.</li> <li>Check that the correct control material has been used.</li> <li>If the flag reappears, check the calibration and reagent status. Repeat with fresh reagent.</li> </ol>

# >R 2(2s)

Meaning	Two sequential control measurements are above the 2 standard deviation.
Message ID	29
Priority	26
Possible cause	<ul><li>The controls are not properly prepared.</li><li>The controls are not properly positioned on the instrument.</li></ul>
Recommended actions	<ol> <li>Repeat with fresh control.</li> <li>Check that the correct control material has been used.</li> <li>If the flag reappears, check the calibration and reagent status. Repeat with fresh reagent.</li> </ol>

List of ISE flags

# <R 2(2s)

Meaning	Two sequential control measurements are below the -2 standard deviation.
Message ID	29
Priority	26
Possible cause	<ul><li>The controls are not properly prepared.</li><li>The controls are not properly positioned on the instrument.</li></ul>
Recommended actions	<ol> <li>Repeat with fresh control.</li> <li>Check that the correct control material has been used.</li> <li>If the flag reappears, check the calibration and reagent status. Repeat with fresh reagent.</li> </ol>

# > RR

Meaning	Above reference range.		
	The calculated concentration is greater than the upper limit of the expected variance.		
	nnnnn is the actual value that was checked against.		
Message ID	40		
Priority	22		
Recommended actions	No action required.		

# < RR

Meaning	Below reference range.		
	The calculated concentration is less than the lower limit of the expected value range.		
	nnnnn is the actual value that was checked against.		
Message ID	41		
Priority	23		
Recommended actions	No action required.		

# Seg Fluid

• See Air Fluid.

### 15 ISE troubleshooting

List of ISE flags

# Seg Isecal

• See Air Isecal.

# Sol1 F Dev

Meaning	The ISE Solution 1 factor was not within the expected range.		
Message ID	22		
Priority	18		
Possible cause	<ul> <li>This flag occurs with calibrations because:</li> <li>The dosing from the sample probe is not accurate.</li> <li>The ISE tower is blocked or the mixing was not sufficient.</li> <li>There may be problems with the ISE unit tubing, the electrodes, or the ISE unit was not serviced properly.</li> <li>The ISE Calibrator Indirect has expired.</li> <li>ISE Solution 1 has expired.</li> </ul>		
Recommended actions	<ul> <li>If only one calibration is flagged:</li> <li>Repeat the calibration.</li> <li>If successive calibrations are flagged:</li> <li>Check whether the ISE Calibrator Indirect is in the correct position.</li> <li>Check whether the ISE Calibrator Indirect has expired.</li> <li>Check whether the ISE Solution 1 has expired.</li> <li>Check whether the ISE Solution 1 has expired.</li> <li>Perform the maintenance action Clean ISE Tower Automatically.</li> <li>Perform the maintenance action Deproteinize Probe.</li> <li>Perform the maintenance action Prime ISE Calibrator &amp; Reference.</li> <li>Check whether the electrode expiration date has passed. Replace the expired electrode, if necessary, and perform the maintenance action Electrode Service.</li> <li>Check the ISE unit tubing for leaks and blockages.</li> <li>Contact your service representative.</li> </ul>		

List of ISE flags

# > Test Rng

Meaning	PANIC value over (upper) Technical Limit.		
	The result is higher than the upper limit for the test.		
Message ID	26		
Priority	20		

# < Test Rng


Meaning	PANIC value below (lower) Technical Limit.			
	The result is lower than the lowest limit for the test.			
Message ID	27			
Priority	21			

# **Reacting to error messages**

For general information on error messages and troubleshooting, see *Reacting to messages* on page D-44.

The following table lists selected messages and provides more detailed information on how to react to them.

Message or error code Affected items Possible causes		Comments		
[8.xxx.56] The ISE reference sensor has detected too much	ISE unit.	ISE Reference Solution is empty.	<ul><li>Replace the ISE Reference Solution bottle.</li><li>See <i>Replacing ISE fluid bottles</i> on page E-30.</li></ul>	
air during ISE Reference Solution transport.		Leakage in the tubing or at elec- trodes.	<ul> <li>Check that the ISE unit tubing is properly connected.</li> <li>Check the ISE unit tubing for leakages and blockages.</li> <li>See <i>Replace ISE unit tubing</i> on page E-57.</li> <li>Check that the electrodes are mounted properly.</li> <li>See <i>Replacing electrodes</i> on page E-32.</li> </ul>	
		The cover of the peristaltic pump is open or not properly closed.	Close the cover.	
	erence Solution	The tubing around the ISE Ref- erence Solution sensor is not properly installed.	Make sure that the tubing is properly pressed into the groves and that the tubing itself is not compressed.	
[109000567] code raised by IC software caused by an unexpected handling or sys- tem situation.	ISE unit.	This is a flow problem within the ISE unit, the ISE tower is not empty. The system checks whether the tower is empty before dispensing.	<ul> <li>Perform the following checks:</li> <li>Check that the ISE unit tubing is properl connected.</li> <li>Check the ISE unit tubing for leakages ar blockages.</li> <li>Check that the electrodes are mounted p erly.</li> <li>Check that the pinch valve caps and clam are properly installed.</li> <li>Check that the peristaltic pump cover is properly closed.</li> </ul>	
			<ul> <li>See Replace ISE unit tubing on page E-57.</li> </ul>	

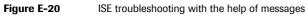

# Glossary and Index

Glossary	F-3
Index	F-11

# Glossary

This glossary is a compendium in which to look up the meaning of technical terms used in conjunction with the **Cedex Bio** instrument.

### Numbers

**2-dimensional barcode** A type of barcode on reagent bottles, calibrators, and control barcode sheets. These matrix barcodes, which use PDF417 symbology, contain more information than traditional linear barcodes.

# Α

**absorbance photometer** A device used to make absorbance photometry measurements. It consists of a halogen light source, optical groupings, a monochromator (diffraction grating), a preamplifier, filters, and a diode array. See also photometer.

**accuracy** The absolute deviation of a result from a predefined target value in percent or absolute units.

**activate** The process of conditioning electrodes with Activator so that the correct potentials are measured.

**Activator** Used to coat ISE electrodes, tubing and the sample probe after cleaning procedures. (The **Cedex Bio** Activator is based on human sample material. For further details refer to the Activator package insert.)

**alarm** A visual or audible operator notification of any system irregularity.

**Alarm button** A button used to display the alarm messages.

**alarm level** A level that identifies the source and severity of a problem. There are two levels: warning, and alarm.

**alarm message** A message that reports a system irregularity. It can be viewed in the alarm monitor screen. See Alarm button.

**aliquot** Portion of sample material pipetted into any secondary cup.

**alphanumeric sorting** The listing of information, in a printout or on a screen, in a pre-defined order by letters or numbers.

**analyte** The constituent in the sample that is to be determined.

**analytical instrument** A device or a combination of devices used to carry out an analytical process.

**analyzer unit** Central analyzing unit including cooling assembly, heating channel, rotor and insulation. Does not include the photometer unit.

**application** Chemistry and method to determine a parameter. See also test.

**assay** 1. A specific test.2. The process of measuring a substance.

# В

**backup** The saving of data onto supplementary storage media such as disks or tape. If such data is required again but is no longer available from the main storage (instrument hard disk), it can be restored from a backup copy.

**barcode** A numeric or alphanumeric code used on sample tubes and reagent bottles to identify the samples and reagents. Different barcode standards are available. See also barcode type.

**barcode scan** The process of reading barcode information into the memory of an instrument.

**barcode scanner** The device that reads the code from sample or reagent barcode labels.

**barcode transfer sheet (BTS)** A sheet bearing a barcode that encodes test-related (or calibrator or control) information for an analytical instrument.

**barcode type** Typical sample barcode types commonly used are Code39, NW7 (Codabar), ITF, and Code 128.

**biohazardous** A classification used to identify material that poses a health threat, for example something contaminated with biological material.

**button** A button is found on a screen or pop-up window. It can be touched either to initiate an action or to move to a different screen.

### Glossary

calculated result - diluent (DIL)

# С

**calculated result** A test result calculated from different individual analytical methods with a given formula such as ratio A/B.

**calibration** The set of operations that establish, under specified conditions, the relationship between values indicated by the analytical instrument and the corresponding known values of an analyte.

**calibration interval** A specified interval at which an assay should be calibrated. Typically found on reagent package inserts.

**calibration type** Describes the event that triggers the calibration, for example lot change, and interval expiration.

calibrator 1. A material of known composition or properties that can be presented to the analytical instrument for calibration purposes.2. The test portion or test solution used for calibration of

an analytical procedure.

**carryover** A process by which materials are carried into a reaction mixture where they do not belong.

**check digit** A verification number used in barcodes and software.

**check sum** The result of a mathematical procedure to validate the integrity of a set of data.

**cleaner** A solution used to wash the probe.

**clot detection** 1. A device built into the pipetting system to detect clots and to avoid false pipetting.2. The procedure of detecting a clot.

**cobas** A modular range of in vitro diagnostics systems from Roche Diagnostics/Hitachi High-Technologies.

**communication** The exchange of data between different computers.

**conditioning** The process of letting protein containing liquid flow through the flow path before electrolyte measurement.

**consumables** A generic term for items that are used during test processing and must be replaced on a regular basis by the operator. Examples of consumables include printer paper and cuvette segments.

**control** See control material.

**Roche Diagnostics** 

**control ID** The abbreviated name for a control material. Control IDs are used on screens where limited space prevents the use of longer names.

**control material** A material used to assess the performance of an analytical procedure or part of an analytical procedure. Also called the control sample.

**control name** The name of a control material.

**CPU** The Central Processing Unit of the system or computer.

**cup** See sample cup.

**cup-on-tube** The placement of a smaller secondary sample container (for example a Hitachi Cup) on top of a primary sample tube.

**cuvette** Disposable plastic container into which sample, reagents, diluent, and water is pipetted, and where the sample pre-dilution and reaction takes place.

**cuvette port** Section of the analyzer unit where cuvette segments are inserted and removed.

**cuvette segment** Cuvette holder containing 10 cuvettes.

**cycle** The instrument time interval during which pipetting or measurement can be carried out.

# D

**database** A defined section of the computer memory where all instrument- and assay-relevant data are processed and stored.

**DB** Abbreviation for database.

**DBMS** Abbreviation for database management system.

**dead volume** The amount of residual sample material in a sample tube or cup.

**default QC** An automated process for performing multiple QC measurements at the time when you define the QC orders.

**default value** A set value registered in advance (initial setting).

**diluent (DIL)** A liquid agent used to reduce the concentration of a sample.

**dilution factor** A preset dilution ratio that is used by the analyzer to perform a requested dilution.

**disposable** Typically a plastic tip, vessel, or cuvette that is discarded after a single use.

**dynamic range** The reportable range of an assay. This range extends from the lower detection limit to the limit of linearity.

# E

**E-stopped** A status indicating that the system has performed an emergency stop. This could be due to hardware failure or because any of the safety devices have requested an emergency stop.

**expected range** The pre-defined range of test result values expected for a defined group of materials. Also known as *normal* range or *reference* range.

**expected value** A value for a test result that can be considered as a *normal* result.

**expiration date** Also called the expiry date. The end of a period until which Roche Diagnostics guarantees product claims for its reagents, calibrators, and controls.

**extended dynamic range** The measuring range for an assay at its highest dilution.

# F

**filter** A process that sorts data for viewing, documenting, or printing according to pre-defined criteria

**flag** An identifier used to call an operator's attention to a result.

# G

global action button See global button.

**global button** A button that allows access to the global software screens and that can be used at any time. For example, <Start>, <Stop>, or <Alarm>.

H

**home position** The position to which a certain part of the instrument returns on reset. The start position of a mechanism.

host computer 1. A computer used for overall management and control of the computer network.2. A laboratory computer that stores and processes requests and results. A host is able to communicate with analytical instruments.

**host communication** Data exchange with a laboratory information system (LIS).

**host interface protocol** A technical description that defines data transfer between a host computer and an analytical system.

# 

**Initialization** Also known as Initializing. The operational mode of an analyzer that occurs immediately after power on and during which the instrument prepares itself for operation.

**installing applications** Process of making the application available for testing, as opposed to importing the application data, which just saves the data on the system.

**interval QC** A streamlined process of performing (all) due QC measurements (for QC that needs to be performed periodically.)

**ISE** Abbreviation for ion-selective electrode.

**ISE Reference Electrode** The electrode through which the ISE Reference Solution flows to set the electronic baseline to zero for ISE measurement.

**ISE Reference Solution** A KCl solution pulled through the ISE Reference Electrode to set the electronic baseline to zero for ISE measurement. Also known as the REF or reference electrode solution.

### Glossary

level detection - order ID

### **Cedex Bio System**

# L

**level detection** A check for the availability of sufficient liquid in a container.

**linear barcode** A conventional one-dimensional barcode with limited data capacity.

**liquid level detection (LLD)** The ability of an analytical instrument to sense liquid by using the probe.

**log file** A set of data, typically stored in the control unit, that traces instrument-related or operator-related activities such as maintenance.

**log off** The procedure of terminating access to a system. Also known as log out or logoff. The reverse procedure is known as log on or log in.

**Log off button** A button used to terminate access to a system. See also log off.

**log on** The procedure of gaining access to a system by entering a user name and, if required, a password. Also known as log in or logon. The reverse procedure is known as log off or log out.

**Log on button** A button used to gain access to a system. See also log on.

**lot calibration** A mandatory calibration when a new lot of reagents is introduced to an analytical instrument.

# Μ

**main calibration** Two-point calibration of the ISE tests with ISE Solution 1 and 2. (As opposed to the calibration with ISE Calibrator Indirect, which is part of each ISE measurement.)

**maintenance action** A maintenance procedure performed by the system or the operator; and that must be performed on a regular basis (for example daily, weekly, or monthly) to secure reliable operation of the analyzer.

maintenance procedure See maintenance action.

**manual dilution** An off-system, pre-analytical step performed by laboratory staff to reduce the analyte concentration in a sample.

**mean** The sum of values divided by the number of values.

**measure point** Time at which absorbance reading is taken and used to calculate results.

measuring range See reportable range.

**message** In computing, a defined set of alphanumeric data that transfers information from computer to computer or from an analytical instrument to the operator.

**Microcup** A secondary sample cup made by Hitachi with a small dead (residual) volume.

**minimum sample volume** The amount of residual sample material plus the volume required for assaying all requested tests to ensure faultless sample aspiration.

**monochromatic** Absorbance measurement at one (primary) wavelength.

**multi-wavelength spectrophotometer** A spectrophotometer in which detectors are placed at multi-wavelength positions to enable simultaneous light reception.

# 0

**on-board** 1. A technical device or function that is part of the analytical instrument and can be used by the instrument at any time.

2. The availability of reagents and consumables on an analytical instrument for use at any time.

**online help** On-screen documentation that a user can request in a context-sensitive manner.

**operating system** A software program that controls all basic functions of a computer (for example, Windows, Linux, and Palm OS).

**operator** The person who uses and controls the analytical instrument or a computer system. See also user.

**operator ID** An alphanumeric ID that a system uses to identify a particular operator.

**order** Collection of all tests that are defined to be performed for a particular sample.

**order ID** A set of alphanumeric data that unmistakably identifies a particular sample order in the lab and is often manually written on the sample label.

# Р

**parameters** A set of criteria or definitions used to establish how an assay is performed. Examples of parameters include sample and reagent volumes and incubation times and temperatures. Such information is typically contained in the application definitions and cannot be changed by the operator.

**password** A form of authentication that uses secret data to control access to a resource.

**photometer** A device that measures the intensity of light.

**photometer unit** Complete photometer assembly, including optical components, power supply, preamplifyer and absorbance controller.

**pipette (to** ...**)** The process of aspirating and dispensing sample and reagents, performed by an appropriate probe.

**pop-up screen** A screen that contains additional information or additional options required for making entries or decisions.

**potentiometric assay** An assay in which analytes (for example Na or K) are measured by ion-selective electrodes.

**precision** The closeness of agreement between independent test results obtained under prescribed conditions.

**pre-dilution** A dilution step performed before samples are analytically processed on the analyzer.

**preventive maintenance** A series of actions, suggested by the system, that the operator should perform to ensure smooth and uninterrupted operation of the instrument (for example, emptying the waste container or replenishing reagents).

**primary tube** The original tube containing the sample that has been drawn. See sample tube.

**prime** The process of flushing a system, for example the tubing, with ISE Calibrator indirect/urine and water to clear it of possible obstacles, traces of fluid, and of air bubbles.

**profile** A user-defined set of tests.

**protocol** 1. A convention or standard that controls or enables the connection, communication, and data transfer between two computing end points. Protocols can be implemented by hardware, software, or a combination of the two.

2. A set of rules that guides how an activity should be performed.

7		1
L	•	

**QC** Abbreviation for quality control. The operational techniques and activities that are used to fulfill requirements for quality.

**QC measurement** The process of performing a test with control material.

**query download** A communication process between instrument PC and LIS by which a pre-defined data set is transmitted upon request of the analytical instrument.

# R

**RAM** Abbreviation for random access memory. Semiconductor memory devices used in computers. RAM content is lost when a computer is switched off.

**ratio** A test result calculated from different individual analytical methods with a given formula such as A/B.

**raw data** The unprocessed values obtained during the analytical process on an instrument (for example mVolt or absorbance).

**reagent** A composition of chemicals used to determine the concentration of substances.

**reagent bottle** Standard plastic bottle for photometric reagents.

**reagent disk** The removable device into which the reagent bottles are placed.

**reagent disk position** One of the multiple positions on the reagent disk.

**reagent port** Section of the analyzer unit where reagent bottles are inserted and removed.

reference electrode See ISE Reference Electrode.

reference range See expected range.

reference solution See ISE Reference Solution.

### **Cedex Bio System**

### Glossary

repeat - SW

**repeat** The performance of the same test on a sample again with identical dilution.

**reportable range** The range of results that can be reported for the assay. It stretches from the lower detection limit to the high end of the calibration curve.

### request See order.

**rerun** The performance of the same test on a sample again with a different dilution.

**result** The value reported by an analytical device during or after the assay of a sample or control.

**ROM** Abbreviation for read-only memory. Semiconductor memory devices used in computers. ROM content remains when a computer is switched off.

**rotor** Assembly consisting of cuvette ring and plastic frame.

# S

**sample area** Area on the instrument where sample tubes are placed.

**sample cup** A small container that is used for samples and also for calibrator and control material. A sample cup can be placed either directly on the sample area, or on sample tubes. Compared to a sample tube, a sample cup allows the use of smaller liquid volumes and so reduces the dead volume.

**sample ID** A set of alphanumeric data that unmistakably identifies a particular sample within a whole organization. It is used on barcodes and is the unique identifier used to communicate with a laboratory information system.

**sample splitting** The act of making one or more aliquots from a primary or secondary specimen.

**sample tube** A glass or plastic container for liquid samples to be used with the system. It may or may not have a bar code label, which may be used for positive sample identification. A sample tube contains sample of one specific specimen (sample) type.

**sample type** A type of sample that can be analyzed.

**sampling stop** An alarm level that indicates a problem with the sampling system. See also S.Stop.

**Roche Diagnostics** 

scan See barcode scan.

**scroll** The action of moving through text or graphics (up, down, left, or right) to see parts of the file or list that cannot fit on the screen.

**scroll arrow** An arrow on either end of a scroll bar that you use to scroll through the contents of the screen or list box.

**scroll bar** A bar that appears at the bottom or right edge of a screen whose contents are not entirely visible.

**SD** Abbreviation for standard deviation.

**secondary tube** A sample container of variable size into which aliquots are transferred.

**shutdown** The process of powering off an instrument.

**software** A computer-operated program that processes data in a defined manner. Software is usually intellectual property of the software supplier or its licensee.

**specimen** Generic term for a calibrator, control and sample.

**standard** Traceable reference material used to create the master calibration curve.

**standard deviation** A statistic used as a measure of the dispersion or variation in a distribution of data.

**Start button** A button used to start system operation (Operating status) and begin the pipetting of samples, measurement, and the result calculation process.

**STAT** Abbreviation for Short Turn Around Time.

**status** A general term used to refer to the current status of the system. Explicit terminology may be used to address sub-parts of the system status (such as analyzer status and printer status).

**Stop button** A button used to display options for different kinds of processing stop.

**SW** Abbreviation for software.

# Т

**target range** The allowed range of recovery for an analyte in a control material.

**target value** The mean of all participant responses after removal of outlying values.

**TAS** Abbreviation for Test Application Software.

**test** Running an analysis of a specific parameter for a sample. See also application.

**test ID** The abbreviated name for a test. This code is displayed on test buttons shown on screens.

**test sheet** A document that lists all the information necessary to perform a specific assay or test on an instrument.

**touchscreen** An input device that allows the user to interact with the computer by touching the display screen.

**turn-around-time** 1. Time between receiving a sample and sending out the result.

# U

**unit** A chosen reference quantity of an analyte used to compare quantities of the same dimension (for example, mol/L, g/L, or U/L).

**user** The person who operates and controls the analytical instrument or a computer system. See also operator.

**user interface** The part of a system exposed to a user. In a computer system, the user typically interacts with an operating system or with application software. With these the user interacts by using menus, icons, keystrokes, mouse clicks, and similar means.

# V

**validation** The process of checking results or data against defined rules. Validation can be against technical criteria.

# W

**waste** Anything discarded by the analyzer; waste can be liquid or solid.

**worklist** A report generated by an analytical instrument. A worklist aids a user by listing calibrators, controls, and samples currently loaded on the sample area.

### Glossary

worklist - worklist

# Index

# A

abbreviations, 7 abbreviations, ISE, E-5 aborting the process, A-116 Abs. air/water calibration, maintenance action, C-28 absorbance adjustment, definitions, B-167 absorbance measurements, process, A-19 absorbance photometer, A-61 acceptable flags list, editing, B-160 accepting - calibration results, B-66 - QC results, B-74 - sample results, B-57 acoustic signals, A-81, D-6 Activate electrodes, maintenance action, E-45 adding lots, B-101 adjusting, touchscreen, B-131 air/water calibration, C-28 alarm button, A-80 alarm LED, A-80 alarm messages - checking, A-81 - confirming, D-9 - exporting log, D-9 - overview, A-81 - printing, D-8 - reacting to, D-7, D-44 alarm monitor - overview, A-80, D-6 alarms - checking, A-81 -LED, A-80 - LED colors, A-81 - monitor, A-80 analyzer - See, instrument. analyzing - checking the progress, B-44 - samples, B-34 application codes, mapping, B-160 applications, B-137 - activating, B-139 - calculation definitions, B-155 - calibration definitions, B-153 - control definitions, B-154 - deactivating, B-139 - definitions, B-152 - deleting, B-144 - general definitions, B-152 - importing, B-111, B-137 - installing, B-112, B-138

### **Roche Diagnostics**

preparing, B-140
result conversion definitions, B-156
screen, A-109
uninstalling, B-143
applications, screen, A-108
approved parts, A-8
assemblies, syringe, A-46
auto accept

- QC results, B-64, B-72

- sample results, B-51

# в

backup, B-82 barcode scanner, A-59, B-89 barcode sheets, reading, B-90 barcodes - reading from reagent bottles, B-89 - reading from sample tubes, B-90 - reading from sheets, B-90 - reagent bottles, A-60 - samples, A-60 - scanner, A-59 biohazardous materials, safety, A-6 bottle sets - See, reagent sets bottles - cleaner, A-52 - external cleaner, B-18 - external fluid containers, B-15 - external waste container, B-17 - external water container, B-16, B-86 - ISE fluids, E-11 - reagents, A-53 - reagents, specifications, A-66 - replacing ISE, E-22, E-30 bottom detection, A-49 buttons - alarms, A-80 - global action, A-71 - overview, A-119

# С

calculation, B-155 - ratios, B-155

calibration - checking if due, B-31 - definitions, B-153, B-165 - deleting orders, B-63 - deleting results, B-95 - during Prepare phase, B-30 - forecast period, A-29 - introduction, A-26 - ISE, E-12, E-26 - main ISE calibration, E-12 - overview, A-26 - performing, B-59 - reserve, A-29 - schedule, B-96 - selecting tests, B-62 - short guide, B-60 - validating results, B-64, B-65 calibration list, screen, A-101 calibration results - auto accept, B-64 - obsolete, A-28 calibrator lot, defining, B-103 calibrators - placing, B-32, B-62, B-63 - specifications, A-66 certificates, importing, B-115 changing - configuration values, B-135 -low voltage fuses, D-49 - mains fuses, D-47 - orders, B-42 - password, B-157 - user definitions, B-159 characters, special, A-76 checking – alarm messages, B–48 - bottle sets, B-23, B-100 - cleanerbottle, B-18 - external fluid containers, B-15 - instrument status, B-46 - ISE status, E-21 - printer status, D-46 - sample status, B-44 -wastecontainer, B-17 - watercontainer, B-16, B-86 classification of safety notices, A-5 Clean and disinfect housing, maintenance action, C-21 Clean ISE tower automatically, maintenance action, E-50 Clean ISE tower manually, maintenance action, E-51 Clean probe manually, maintenance action, C-12 Clean reagent disk and sample area, maintenance action, C-14 Clean rotor and heating channel, maintenance action, C-10

Clean the water and waste containers, maintenance action, C-16 cleaner - checking external bottle, B-18 - replacing bottle, B-18 cleaner bottle, A-52 cleaning - ISE tower off the instrument, E-36 - ISE tower on the instrument, E-51 - waste container, C-17 - water container, C-16 cleaning up - database, C-23 - results, B-84 Cleanup database, maintenance action, C-23 clearing paper jam, D-46 collapsing lists, A-74 colors - alarm LED, A-81, B-48 - cleaner button, B-18 - concept, A-69, A-72 - cuvette buttons, B-28 - instrument buttons, B-47 – ISE bottle buttons, E–30 - ISE LED, E-10 - LEDs, A-41, A-117 - maintenance actions, B-19 - reagent buttons, B-25 - sample buttons, B-44 - test buttons, B-31 - waste button, B-17 - water button, B-16 components - disposal, A-11 - instrument, A-43 – ISE unit, E–10 concepts - calibration, A-26 - colors in screens, A-69, A-72 - wizards, A-69 condensation, removing from the reagent cooler, B-124 Condition ISE tubing, maintenance action, E-46 configuration, B-133, B-157 - Abs. adjustment, B-167 - calibration definitions, B-165 - changing values, B-135 - date and time, B-167 - extra wash definitions, B-168, B-169, B-170, B-171 - host definitions, B-163 - maintenance days, B-167 - QC definitions, B-165 - result handling, B-165 - screen, A-106 - system definitions, B-166 - viewing values, B-135 - workflow definitions, B-161

confirming, alarm messages, D-9 connecting - external cleaner bottle, B-128 - external waste container, B-128 - external water container, B-127 connectivity, host, A-32 connectors - fluids, A-47 - overview, A-63 contact addresses, 2 Contact service representative, maintenance action, C-27 containers - fluids, A-50 - reagent disks, A-58 -waste, A-50 -water, A-51 control lot, defining, B-101 controls - definitions, B-154 - placing, B-70 conventions used in manual, 6 cooling, reagents, A-58 copyrights, 2 correlations, B-156 covers - ISE, E-9 - overview, A-39 cross-contamination, A-9 csv files, B-83 cups, A-53 cuvette ring, A-59 cuvette status, screen, A-89 cuvettes, A-54 - checking the status, B-28 - handling, A-55 - replacing, B-29, B-85 - ring, A-59 - segments, A-55 - specifications, A-66

# D

daily backup, B-82 daily operation, B-3 - introduction, A-22 - overview table, A-22 Daily prepare actions, maintenance action, E-44 data security, A-10 database - cleaning up, C-23 - exporting, B-104 - importing, B-113 - restoring, B-113 date and time, definitions, B-167

### **Roche Diagnostics**

deactivating ISE unit, E-64 Default OC - overview, B-68 - performing, B-70 defining - calibrator lots, B-103, B-150 - control lot, B-101 - control lots, B-149 -lots, B-101 - orders, B-38 - ratios, B-146 - users, B-159 definitions, applications, B-152 degasser, A-44 deleting - applications, B-144 - bottle sets, B-120 - calibration orders, B-63 - calibration results, B-95 - mixing rules, B-172 - orders, B-93 - profiles, B-149 - QC results, B-96 - results at end of shift, B-84 - sample results, B-94 - users, B-159 Deproteinize probe, maintenance action, C-10 detection - fluid levels, A-49 - tube bottom, A-49 diagnostics data, exporting, D-51 diagnostics, screen, A-112 dilutions, A-31 dimensions, specifications, A-65 disconnecting tubes from instrument, B-130 disinfecting housing, C-21 display items, A-71 disposing - instrument, A-11 - waste, B-86 document information, 2

# E

edition notice, 2 electrical safety, A–5 electrode block, E–10 Electrode service, maintenance action, E–49 electrodes – activating, E–45 – maintenance, E–49 – position in block, E–34 – replacing, E–32 electromagnetic devices, safety, A–8 emptying, waste container, B–17, B–86

End of Shift short guide, B-78 environment, specifications, A-65 error messages, reacting to, D-44, E-80 exceptional situations - ISE unit, E-80 - main instrument, D-42 exchanging reagents, B-25 expanding lists, A-74 export, screen, A-115 exporting - alarm message log, D-9 - database, B-104 - diagnostics data, D-51 -log files, B-110 - results, B-83, B-105 external bottle status, screen, A-93 external waste container, A-51 extra wash cycles - activating, B-169 - deactivating, B-169 - defining, B-168 - definitions, B-170 - deleting, B-169 - importing, B-115 - screen, A-110

### F

features of instrument, A-17 filtering, A-76 finishing the shift, B-77 flags - Abs. results, D-28 - ISE results, E-71 - overview, D-25 - priorities, D-26 - viewing details, D-25 flaps - instrument, A-39 – ISE unit, E–9 fluid system - overview, A-45 - priming, C-11 fluids -bottles, A-50 - bottles for ISE, E-11 - checking external containers, B-15 - connecting external cleaner bottle, B-128 - connecting waste container, B-128 - connecting water container, B-127 - external connectors, A-47 -ISE, E-8- replacing ISE bottles, E-22, E-30 forecast period, A-29

fuses, A–64 – changing for mains, D–47 – changing low voltage, D–49

### G

global action buttons, A-71

### Η

handling - cuvettes, A-55 -lots, B-98 - orders, A-31 - reagents, A-24, A-56 - results, A-33 - samples, A-30 - software data, A-66 hardware - overview, A-37 - specifications, A-65 heads, transfer, A-48 heating channel, cleaning, C-10 host - configuration, B-163 - connectivity, A-32 - definitions, B-163 host codes, B-160 - screen, A-110 housing, cleaning, C-21

### I

ignoring QC results, B-74 import, screen, A-114 importing - applications, B-111, B-137 - certificates, B-115 - database, B-113 - extra wash cycles, B-115 - mixing information, B-115 - software, B-113 Initialize ISE reference sensor, maintenance action, E-47 Initialize ISE unit, maintenance action, E-53 installation, safety, A-7 installing - applications, B-112, B-138 - ISE pump tubing, E-54 – ISE unit tubing, E–61

instrument - disposal, A-11 - features, A-17 - introduction, A-15 - main components, A-43 - overview, A-17 - switching on, B-10 - where to use, A-7 intended use - instrument, 5 - of operator's manual, 2 interface - See, user interface interfaces, specifications, A-66 internal waste tank, A-50 interpreting, QC History, B-74 Interval QC - overview, B-68 - performing, B-71 intervals - ISE maintenance actions, E-41 - maintenance actions, C-7 introduction - See, overview. inventory - checking bottle sets, B-23, B-100 - deleting bottle sets, B-120 - screen, A-113 ISE - calibration, E-12 - calibrator, priming, E-48 - checking status, E-21 - cleaning tower automatically, E-50 - cleaning tower manually, E-51 - cleaning tower off the instrument, E-36 - conditioning tubing, E-46 -flags, E-71 - fluid containers, E-11 - installing pump tubing, E-54 - installing tubing, E-61 -list of flags, E-71 - main calibration, E-26 - maintenance, E-39 - measurement process, E-7 - measuring modes, E-6 - monitoring status, E-24 - operation, E-17 - overview, E-5 - preparing system, E-20 - priming calibrator, E-48 - principles of measurement, E-7 - pump tubing, E-54 - replacing pump tubing, E-54 - replacing tubing, E-57 - sensors, E-11

# solutions, E-8 standby, E-12 status, E-69 tower, E-36, E-50, E-51 troubleshooting, E-67 tubing, E-11, E-61 ISE reference sensor, initializing, E-47 ISE status, screen, A-91 ISE tower, E-11 cleaning automatically, E-50 cleaning manually, E-51 cleaning off the instrument, E-36 ISE unit, E-9 deactivating, E-64 main components, E-10

- specifications, A-66

### Κ

keyboards, A-74

# L

laboratory correlations, B-156 layout of screens, A-70 LEDs - alarms, A-81 - color table, A-41, A-117 - ISE, E-10 - overview, A-41 level detection, A-49 lids - ISE, E-9 - overview, A-39 lists - Abs. flags, D-28 - alarm messages, D-10 - expanding and collapsing, A-74 - ISE flags, E-71 -lots, A-103 - orders, A-95 loading - reagent disk, B-21, D-50 - reagents, B-26 loadlist, screen, A-104 log files, exporting, B-110 logging off, B-88 logging on, B-11 logs, alarm messages, D-9 lot data, screen, A-102 lot list, A-103

lots - adding, B-101 - defining, B-101 - defining for calibrators, B-150 - defining for controls, B-149 - handling, B-98 - list, A-103 - viewing definitions, B-98, B-99 - viewing details of used fluids, B-100 lower-case letters, A-75

### Μ

main calibration, E-12, E-26 maintenance - instrument, C-5 - ISE, E-39 - overview, C-5 - screen, A-107 maintenance actions - Abs. air/water calibration, C-28 - Activate electrodes, E-45 - Clean housing, C-21 – Clean ISE tower automatically, E–50 - Clean ISE tower manually, E-51 - Clean probe manually, C-12 - Clean reagent disk, C-14 - Clean rotor and heating channel, C-10 - Clean rotor heating channel, C-32 - Clean water and waste, C-16 - Cleanup database, C-23 - colors of entries, B-19 - Condition ISE tubing, E-46 - Contact service representative, C-27 - Daily prepare actions, E-44 - Deproteinize probe, C-10 - Electrode service, E-49 - Initialize ISE reference sensor, E-47 - Initialize ISE unit, E-53 - intervals, C-7 - intervals for ISE, E-41 - performing, C-6 - performing during End Shift phase, B-85 - performing during Prepare phase, B-19 - Prime fluid system, C-11 – Prime ISE calibrator and reference, E–48 - Replace ISE pump tubing, E-54 - Replace ISE unit tubing, E-57 - Replace photometer lamp, C-24 - Replace water inlet filter, C-19 maintenance days, B-167 manual, how to use, 5 mapping, application codes, B-160 measurements, specifications, A-65 measuring modes, ISE, E-6

messages - alarms, A-81, D-10 -list, D-10 - overview, A-80, D-5 - screen, A-80, D-5 mixing motor, A-49 mixing reagents, B-76 - during Prepare phase, B-29 mixing rules, B-171 - definitions, B-171 - deleting, B-172 - importing, B-115 - printing, B-171 monitor - See, alarm monitor. monitoring - analysis process, B-44 - ISE status, E-24

# Ν

navigation, through screens, A-72

# 0

obsolete calibration, A-28 online Help, using, 5, A-79 operation - overview, B-7 - principles, A-19 - routine, B-3 - routine ISE, E-17 - safety, A-6 operators - qualifications, A-9 - regular breaks, A-9 optical safety, A-6 order of processing, B-44, B-145 orders - changing, B-42 - changing with sample on board, B-42 - checking for unfinished, B-80 - defining, B-38, B-39 - defining profiles, B-148 - deleting, B-93 - handling, A-31 - IDs, A-31 - introduction to handling, A-31 -list, A-95 - order of processing, B-44 - overview, A-31 - STAT, B-41 - workflow setup, B-37

overview -buttons, A-119 - calibration, A-26 - daily operation, A-22 -flags, D-25 - instrument, A-15, A-17 – ISE, E–5 - ISE operation, E-17 - maintenance, C-5 - messages, A-80, D-5 - order handling, A-31 - reagent handling, A-24 - result handling, A-33 - routine operation, B-7 - sample handling, A-30 - user interface, A-20, A-69 -wizards, A-21 Overview tab, A-82

# Ρ

panels - ISE, E-9 - overview, A-39 - See also, test panels. paper jam, clearing, D-46 paper, refilling, B-122 password - changing, B-157 - typing, B-11 performing - calibration, B-59 - calibration during Prepare phase, B-30 - Default QC, B-70 – Interval QC, B–71 - maintenance actions, C-6 – QC, B–68 - reagent mixing, B-76 - reagent mixing during Prepare phase, B-29 - tests, B-34 peristaltic pump, E-10 photometer - overview, A-61 - specifications, A-66 photometer lamp, replacing, C-24 power - interruption, A-8 - requirements, A-65 precalibration, A-29 Prepare phase, short guide, B-13 Prepare wizard, B-15

preparing - applications, B-140 - ISE unit, E-20 - new reagent disks, B-116 - reagents, B-22, B-25 - system, B-12 Prime ISE reference and calibrator, maintenance action, E-48 Prime the fluid system, maintenance action, C-11 principles - absorbance measurements, A-19 - calibration, A-26 - ISE measurements, E-7 - specifications, A-65 printer - clearing paper jam, D-46 - overview, A-61 - refilling paper, B-122 - specifications, A-66 printer status, checking, D-46 printing alarm messages, D–8 - mixing rules, B-171 - overview, A-78 – sample results, B–58 - stopping, A-78, B-58 priority, flags, D-26 probe - cleaning manually, C-12 - deproteinizing, C-10 - description, A-49 – hardware, A–46 - replacing, B-125 process sequence - screen, A-111 processing - defining sequence, B-145 processing, stop, B-49 profiles, B-148 - adding tests, B-148, B-149 - deleting, B-149 - removing tests, B-149 pumps, ISE, E-10

# Q

OC – Default QC, B–68 - definitions, B-165 - interpreting history, B-74 - Interval QC, B-68 -list, A-98 - overview, A-29 - performing, B-68 - performing Default QC, B-70 - performing Interval QC, B-71 - short guide, B-69 – status list, A–98 - when due, A-29 QC history – graph, B–75 - interpreting, B-74 - screen, A-99 QC history graph, screen, A-100 QC results - auto accept, B-72 - deleting, B-96 -validating, B-72 QC status list, screen, A-98 qualifications, of operators, A-9 quality control - See, QC.

# R

ratios - calculation definitions, B-155 - defining, B-146 reagent bottles, A-53 - reading barcodes, B-89 - specifications, A-66 reagent cooler, removing condensation, B-124 reagent disks - cleaning, C-14 - ID, A-58 -loading, B-21 - overview, A-57 - preparing, B-116 - realigning, D-50 - removing, B-87 - storage container, A-58 reagent handling, overview, A-24 reagent lots - viewing, B-99 reagent mixing -screen, A-111 reagent sets, A-25 - checking status, B-23 - deleting from inventory, B-120

reagent status, screen, A-90 reagents - barcodes, A-60 - bottles, A-53 - cooling, A-58 - exchanging, B-25 - handling, A-56 - introduction to handling, A-24 -loading, B-26 - preparing, B-22, B-25 - removing, B-25 - replacing, B-25 - safety, A-7 - sets, A-25 reference ranges, ratios, B-156 refilling – printer paper, B–122 - water container, B-16, B-86 removing, B-50 - reagent disk, B-87 - reagents, B-25 - sample tubes, B-50 repeating, tests, B-55 Replace ISE pump tubing, maintenance action, E-54 Replace ISE unit tubing, maintenance action, E-57 Replace photometer lamp, maintenance action, C-24 Replace water inlet filter, maintenance action, C-19 replacing - cleaner bottle, B-18 - cuvettes, B-29, B-85 - electrodes, E-32 - ISE bottles, E-22, E-30 - printer paper, B-122 - probe, B-125 - reagents, B-25 - water container, B-16, B-86 rerunning, tests, B-56 restarting the system, A-116 restoring database, B-113 result conversion, B-156 result handling - definitions, B-165 - overview, A-33 result list, screen, A-96 results - calibration, B-65 - checking for not accepted sample results, B-80 checking for transmission, B–81 - cleaning up at end of shift, B-84 - deleting of calibration, B-95 - deleting of QC, B-96 - deleting of samples, B-94 - exporting, B-83, B-105 - flags, D-23 - handling, A-33 - introduction to handling, A-33

results - list, A-96 - printing of samples, B-58 - samples, B-51 - samples, if sample on board, B-52 - samples, if sample removed, B-53 - validating for QC, B-72 - validating of samples, B-51 revision history, 2 rotor - cleaning, C-10 - overview, A-56

### S

safety, A-3 - approved parts, A-8 - biohazardous materials, A-6 - calibration, B-59 - classification, A-5 - cross-contamination, A-9 - data security, A-10 - disposal of instrument, A-11 - during End of Shift phase, B-77 – during Prepare phase, B–12 - during routine operation, B-5 - electrical, A-5 - electromagnetic devices, A-8 - insoluble contaminants, A-9 - installation, A-7 - instrument covers, A-6 - ISE operation, E-19 - operation and maintenance, A-6 - operator breaks, A-9 - operator qualifications, A-9 - optical, A-6 - performing ISE maintenance, E-42 - performing QC, B-68 - power supply, A-8 - reagents and working solutions, A-7 - spillages, A-9 - third-party software, A-8 - transport, A-5 -waste, A-7 - when analyzing samples, B-34 - when exchanging external bottles, B-15 - when handling cuvettes, B-28 - when handling external bottles, B-15 - when handling reagents, B-22 - when removing reagent disk, B-87 - when troubleshooting ISE, E-70 - when using the barcode scanner, B-89 - with flags, D-27 - with maintenance actions, C-8

sample area - cleaning, C-14 - overview, A-44 sample handling, overview, A-30 sample ID, A-31 sample overview, screen, A-84 sample results - accepting, B-57 - auto accept, B-51 - checking for not accepted, B-80 - deleting, B-94 - printing, B-58 - validating, B-51 - validating if sample on board, B-52 - validating if sample removed, B-53 samples - analyzing, B-34 - barcodes, A-60 - checking the status, B-44 - handling, A-30 - introduction to handling, A-30 - overview, B-50 - reading barcodes, B-90 – short guide, B–35 - specifications, A-65 sampling stop, A-116 scanners, A-59 scheduling, calibration, B-96 screens - alarms, A-80 - applications, A-108 - applications list, A-109 - calibration list, A-101 - configuration, A-106 - cuvette status, A-89 - diagnostics, A-112 - display items, overview, A-71 - expanding and collapsing lists, A-74 - export, A-115 - external bottle status, A-93 - extra wash cycles, A-110 - host codes, A-110 - import, A-114 - inventory, A-113 - ISE status, A-91 - keyboards, A-74 -layout, A-70 - Loadlist, A-104 -lot data, A-102 -lot list, A-103 - maintenance, A-107 - messages, A-80, D-5 - navigation, A-72 - orders list, A-95

screens - QC history, A-99 – QC history graph, A–100 - QC list, A-98 - OC status, A-98 - reagent mixing, A-111 - reagent status, A-90 - result list, A-96 - sample overview, A-84 - scrolling, A-73 - stopping processing, A-116 - system status, A-92 - test selection, A-86 - test status, A-87 - users, A-115 -worklist, A-104 scrolling, A-73 segments of cuvettes, A-55 selecting groups of list items - See, filtering. selecting, tests to be calibrated, B-62 sensors, ISE, E-11 sequence of processing, B-145 service representative, contacting, C-27 setting up, order definition, B-37 sheets, barcodes, B-90 shift - finishing, B-77 - starting, B-10 short guides - analyzing samples, B-35 - calibration, B-60 - End of Shift, B-78 – ISE Prepare phase, E–20 - Prepare phase, B-13 – QC, B–69 shutting down the system - end of shift, B-88 - stop option, A-116 software - importing, B-113 - updating, B-113 software data handling, specifications, A-66 solutions, ISE, E-8 special characters, A-76 specifications, technical, A-65 spillages, safety, A-9 standby, ISE status, E-12 starting – a run, B–44 - processing, B-44 - shift, B-10 STAT orders, B-41

status - cuvettes, B-28 - instrument, B-46 - ISE unit, E-69 - reagent sets, B-23 - sample tubes, B-44 - standby for ISE, E-12 - system, A-34 stop, screen, A-116 stopping - printing, A-78, B-58 - processing, B-49 – run, B–49 storage facilities, specifications, A-66 switching - between upper and lower-case, A-75 – off, B–88 - on, B-10 symbols, -6 syringe - assembly, A-46 - hardware, A-46 system - definitions, B-166 - logging on, B-11 - preparing, B-12 - preparing ISE, E-20 - See also, instrument. system status - overview, A-34 - screen, A-92 systems - fluid, A-45 - See also, units, assemblies.

# Т

tabs, B–119 – assigning tests, B–119 – Overview, A–82 – removing tests, B–120 – Utilities, A–105 – Workplace, A–94 technical specifications, A–65 test panels – assigning tests, B–119 test selection, screen, A–86 test status, screen, A–87

tests - adding to profile, B-148 – assigning to tabs, B–119 - defining profiles, B-148 - performing, B-34 - removing from profile, B-149 - removing from tabs, B-120 - repeating, B-55 - rerunning, B-56 text, typing, A-74 third-party software, A-8 throughput, specifications, A-65 touchscreen, adjusting, B-131 trademarks, 2 transfer head, A-48 transfer unit, A-48 transmission, checking of results, B-81 transport, safety, A-5 TRL check, B-52, B-53 - after database installation, B-115 - installing application, B-137 troubleshooting - ISE, E-67 - main instrument, D-39 tube bottom detection, A-49 tubes, A-53 - disconnecting from instrument, B-130 tubing - conditioning, E-46 - connecting to instrument, B-127 - disconnecting from instrument, B-130 - installing for ISE, E-61 - ISE, E-11 - replacing of ISE, E-57 - replacing of ISE pump, E-54

typing – password, B–11 – text, A–74

– user name, B–11

# U

unfinished orders, B–80 uninstalling applications, B–143 units – ISE, E–9 – laboratory, B–156 – transfer, A–48 updating software, B–113 upper-case letters, A–75 user interface, overview, A–20, A–69 user name, typing, B–11 users – defining, B–159 – deleting, B–159 – screen, A–115 using – manual, 5 – online Help, 5, A–79 Utilities tab, A–105

# V

validating
calibration results, B-64
QC results, B-72
sample results, B-51
viewing
cal and QC lot definitions, B-98
configuration values, B-135
flag details, D-25
lot details of used fluids, B-100
reagent lot definitions, B-99

# W

wash cycles - See, extra wash cycles wash station, A-47 waste - checking external container, B-17 - cleaning container, C-17 - emptying, B-17, B-86 - external container, A-51 - internal tank, A-50 - safety, A-7 water - checking externalcontainer, B-16, B-86 - cleaning container, C-16 - container, A-51 - refilling, B-16, B-86 - specifications, A-66 water filter, changing, C-19 wavelengths, A-62 wizards - concept, A-69 - End Shift, B-81 - introduction, A-21 - orders, B-35 - Prepare, B-15 workflow, definitions, B-161 working solutions, safety, A-7 worklist, A-104 Workplace tab, A-94



Table of contents

# Revisions

This chapter contains amendments to the Cedex Bio System Operator's Manual.

In this chapter	Chapter	18
Changes in Cedex Bio System Operator's Manual 3.1		G-5
Structure of a change entry		G-6
Change 1: Warning statement added to inform about sharps,	rough edges, a	nd/or
moving parts		G-7
Change 2: Safety statement about data security updated		G-8

# Changes in *Cedex Bio System* Operator's Manual 3.1

Manual	Manual Version	Software Version	<b>Revision Date</b>	Changes
<i>Cedex Bio System</i> Operator's Manual	3.1.1	2.1	October 2021	Updated data security, updated safety information regarding sharp edges

In order to ensure that the readers of the *Cedex Bio System* Operator's Manual use the up-to-date content, Roche recommends to manually mark the affected passage in the original manual.

### ► To update the *Cedex Bio System* Operator's Manual


- 1 In the current chapter *Revisions*, go to the first change.
- 2 Familiarize yourself with the items that changed.
- **3** Identify the page number affected in the original manual.
- **4** Go to the appropriate page in the original and manually mark the changes. For example cross out text that is no longer applicable or add a reference where to find new and changed content.



If you print these pages, Roche recommends printing them single-sided. This way you can easily insert them in the existing manual.

Changes in Cedex Bio System Operator's Manual 3.1

### Structure of a change entry

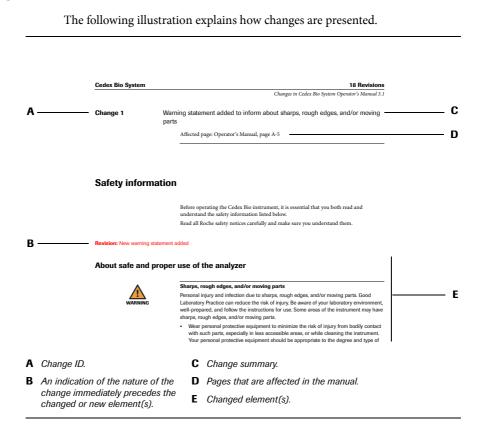

 Change 1
 Warning statement added to inform about sharps, rough edges, and/or moving parts

Affected page: Operator's Manual, page A-5

# **Safety information**

Before operating the Cedex Bio instrument, it is essential that you both read and understand the safety information listed below.

Read all Roche safety notices carefully and make sure you understand them.

Revision: New warning statement added

# About safe and proper use of the analyzer



### Sharps, rough edges, and/or moving parts

Personal injury and infection due to sharps, rough edges, and/or moving parts. Good Laboratory Practice can reduce the risk of injury. Be aware of your laboratory environment, well-prepared, and follow the instructions for use. Some areas of the instrument may have sharps, rough edges, and/or moving parts.

 Wear personal protective equipment to minimize the risk of injury from bodily contact with such parts, especially in less accessible areas, or while cleaning the instrument.
 Your personal protective equipment should be appropriate to the degree and type of potential hazard, e.g. suitable lab gloves, eye protection, lab coat, and footwear. Changes in Cedex Bio System Operator's Manual 3.1

### Change 2

### Safety statement about data security updated

Affected page: Operator's Manual, page A-10

Revision: Section updated

# **Data security**

Data security	<ul> <li>The physical access to all hardware and servers (including analyzer, computer systems, cables, network equipment, etc.) should be protected.</li> <li>Use the software's operator management functionality as intended.</li> <li>All users must exit the software or lock their computer before leaving their workplace.</li> <li>Do not use shared user accounts for laboratory staff.</li> <li>It is recommended to have individual accounts with appropriate operator level, to access the software and the laboratory's credentials.</li> <li>It is recommended to delete any obsolete user accounts from the Roche systems.</li> </ul>
Network security	<ul> <li>The laboratory network and any laboratory infrastructure must be secured against unauthorized access, malicious software, and cyber-attacks.</li> <li>All devices and services used in the customer laboratory infrastructure must be safe and secured against malicious software and unauthorized access.</li> <li>A secure network environment which is resilient to traffic redirection and eavesdropping must be installed.</li> </ul>
Data transfer and data entry on analyzers	<ul> <li>Data transfer using any host protocol (e.g. HL7, ASTM) is not encrypted; data is transferred as plain text and readable with IT tools (e.g. sniffer).</li> <li>Do not download from any host system (e.g. LIS, middleware) any privacy-sensitive data to the analyzer.</li> </ul>
Secure data storage	<ul> <li>Any data backup and data archive that has been exported from the analyzer must be securely stored and protected from any unauthorized access.</li> <li>Data access to any data backup or data archive is permitted to authorized persons only.</li> <li>Make sure that system backups and archive files are protected from any unauthorized access and disaster. Do not take any identification data outside of the laboratory.</li> <li>Implement controls for handling and disposal of media and assets (e.g., hard disks). Media and assets containing data from Roche products must be securely destroyed.</li> </ul>
Cybersecurity and Privacy awareness	<ul> <li>Perform regular security and privacy awareness training for staff handling personal data and according the data privacy principles that are mandatory according to regulations.</li> <li>Check your analyzer for suspicious activity and report any suspected compromise immediately to your Roche representative.</li> </ul>

- Always leave software updates to a Roche Service representative, or perform such updates with their assistance. New software updates should be implemented as soon as possible.
- Read carefully and follow the recommendations and advice written in the safety chapter.
- Tor further information, contact your Roche Service representative

Changes in Cedex Bio System Operator's Manual 3.1